



Medical & Drug • IVD & POCT • Food Safety • Smart Wear

China Headquarter- Shenzhen

Address : R777, Hangcheng Wisdom Science Park,
Hangcheng street, Bao'an District Shenzhen 518128, China
Phone : +86 0755-23006699
Email: leo@reagen.us
Web : www.reagen.cn

USA Headquarter- California

Address : 7098 Miratech Dr Ste 110
San Diego CA 92121 USA
Phone : +1 619-3898888
Email: max@reagen.us
Web : www.reagen.us

Europe Headquarter-Norway

Address : Vollveien 7, 1433 Ås,
Norway
Phone : +47 46393984
Email: michael@reagen.us
Web : www.reagen.eu

REAGEN®

REAGEN Aflatoxin B1 ELISA Test Kit

RNM98007

Declaration

This kit can detect Aflatoxin B1 in fish and meat, Please use this extraction step to process the sample:

- 1) Weigh out 1mL of the ground sample and place into a suitable container.
- 2) Add 5mL of **70% Methanol** and vortex vigorously for 3 minutes manually or 10 minutes using a multi-vortexer.
- 3) Centrifuge the samples for 5 minutes at 4,000 x g at room temperature (20-25°C / 68–77°F). Dilute 200 µL of the obtained supernatant with 800 µL of **1X Sample Dilution Buffer**. Vortex for 1 minute.
- 4) Use 50 µL of the diluted sample per well in the test.

Note: Dilution Factor = 25

Hereby declared.



REAGEN