Novobiocin Supplement

Instructions For Use
ENGLISH

Selective supplement for detection of motile *Salmonella* according to ISO 6579.

DESCRIPTION

A lyophilised supplement for use with MSRV Medium Base (Ref. 610018 / 620018) for the selective enrichment of motile *Salmonellae* from food and animal feed, environmental samples and samples from the primary production stage such as animal faeces, dust and swabs.

COMPOSITION

Ingredient	Content per vial
Novobiocin Sodium Salt	10 mg

Each vial is sufficient for 1 litre of medium.

METHOD PRINCIPLE

Novobiocin is an antibiotic active mostly against Gram-positive bacteria.

DIRECTIONS

1. Suspend 31.6 g of MSRV Medium Base in 1 litre of purified water.
2. Aseptically reconstitute the content of 1 vial of Novobiocin Supplement with 5 ml sterile distilled water.
3. Mix to complete dissolution and add to 1 litre of MSRV Medium Base, boiled and cooled to 45-50°C.
4. Mix well and distribute into sterile Petri dishes.

TECHNIQUE AND INTERPRETATION OF THE RESULTS

Refer to the technical sheet of the medium being prepared.

STORAGE

2-8°C in its original packaging. Keep away from sources of heat and avoid excessive changes of temperature. Use until the expiry date indicated on the label. Eliminate without using if there are signs of deterioration.

SHELF LIFE

3 years.

QUALITY CONTROL

Appearance: Freeze-dried white in colour.

Microbiological Performance:

Control strains

Salmonella Enteritidis WDCM 00030 (ATCC® 13076, NCTC 12694)

Salmonella Typhimurium WDCM 00031 (ATCC® 14028, NCTC 12023)

Escherichia coli WDCM 00013 (ATCC® 25922, NCTC 12241)

Enterococcus faecalis WDCM 00009 (ATCC® 19433, NCTC 775)

Please refer to the actual batch related Certificate of Analysis (CoA).

WARNING AND PRECAUTIONS

For professional use only. Operators must be trained and have certain experience. Please read the instructions carefully before using this product. Reliability of assay results cannot be guaranteed if there are any deviations from the instructions in this document.

Consult the Safety Data Sheet (SDS) for information regarding hazards and safe handling practices.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to national and local regulations in force.

BIBLIOGRAPHY

See the references at the end of this document.

TABLE OF SYMBOLS

See the table of symbols at the end of this document.

Product	Packaging	Ref.
Novobiocin Supplement	10 vials	81021

This IFU document and the SDS are available from the online Support Center:

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