



Medical & Drug • IVD & POCT • Food Safety • Smart Wear

China Headquarter- Shenzhen

Address : R777, Hangcheng Wisdom Science Park,
Hangcheng street, Bao'an District Shenzhen 518128, China
Phone : +86 0755-23006699
Email: leo@reagen.us
Web : www.reagen.cn

USA Headquarter- California

Address : 7098 Miratech Dr Ste 110
San Diego CA 92121 USA
Phone : +1 619-3898888
Email: max@reagen.us
Web : www.reagen.us

Europe Headquarter-Norway

Address : Vollveien 7, 1433 Ås,
Norway
Phone : +47 46393984
Email: michael@reagen.us
Web : www.reagen.eu

REAGEN®

Ochratoxin A ELISA Test Kit

RNM98008

Declaration

This kit can detect Ochratoxin A in fish, Please use this extraction step to process the sample:

- 1) Weigh 2g of crushed sample into a 50mL centrifuge tube, add 10ml of 70% methanol, shake for 5 minutes, room temperature 4000 rpm centrifugal for 10 minutes;
- 2) Take 1mL of supernatant, add 1mL of 0.1M NaHCO_3 solution, and shake.
- 3) Use 50 μL of the diluted supernatant per well in the test.

Note: Dilution factor=10

Hereby declared



REAGEN