

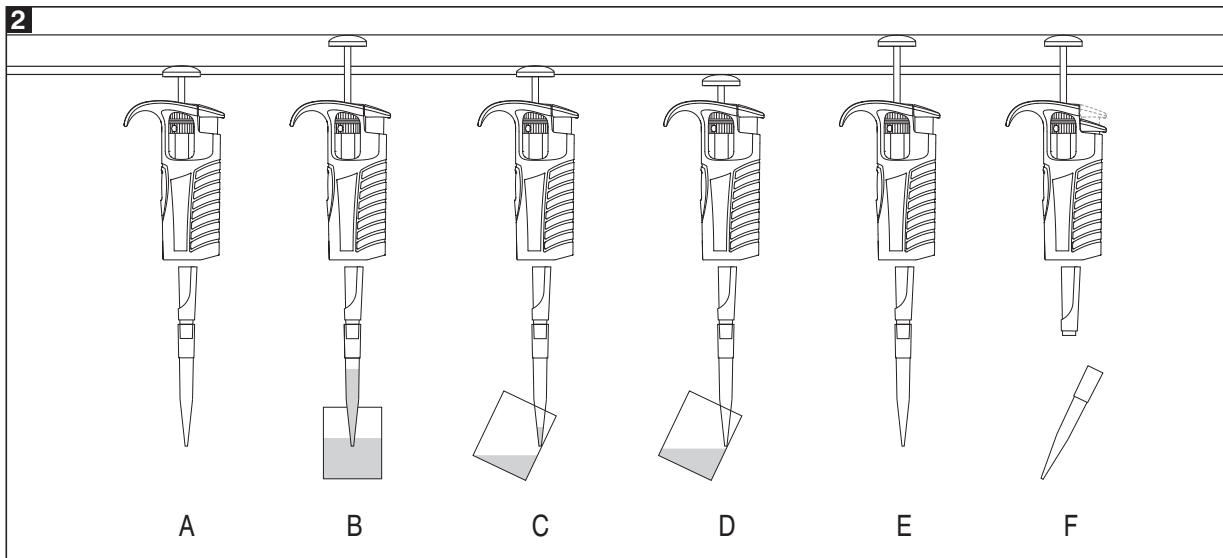
Optipette

ENGLISH	1 – 14
DEUTSCH	15 – 30
FRANÇAIS	31 – 46
ESPAÑOL	47 – 60
PORTUGUÊS	61 – 76
POLSKI	77 – 92
РУССКИЙ	93 – 108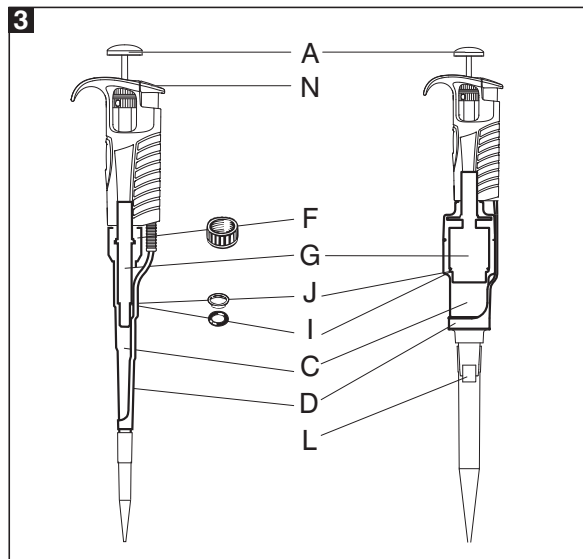

PZ HTL S.A. is in possession of a registered and certified quality management system which includes the development, production, and sales of high quality Liquid Handling products.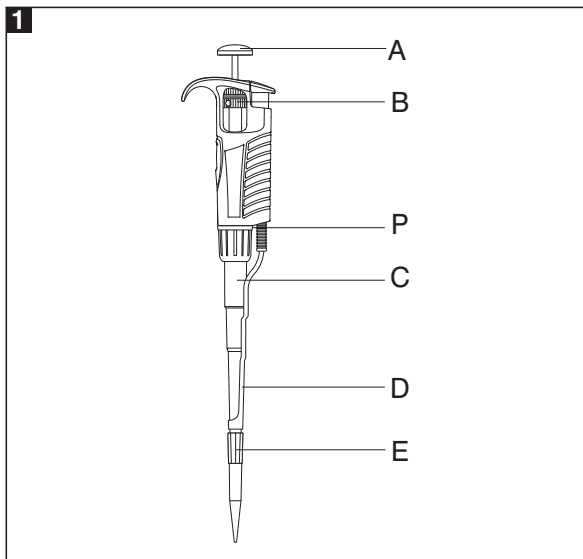

CONTENTS

1 - INTRODUCTION
2 - SETTING THE VOLUME
3 - ASPIRATING AND DISPENSING LIQUIDS
4 - PRE-RINSING
5 - DENSE AND VISCOUS LIQUIDS
6 - HTL PIPETTE TIPS
7 - RECOMMENDATIONS
8 - RECALIBRATION
9 - TROUBLESHOOTING
10 - PIPETTE KIT
11 - SPARE PARTS AND ACCESSORIES

1 - INTRODUCTION

The **OPTIPETTE** is a volumetric instrument designed to measure and transfer liquids precisely and safely. It can measure and transfer volumes from 0.1 μl to 10000 μl depending on the model.

The **OPTIPETTE** pipettes are equipped with a digital counter which shows the pipetting volume. The volume is adjusted by turning the black knurled adjustment ring, (Fig.1B), and is continuously adjustable within the volume range for the pipette. The maximum volume of the pipette is shown on the side surface of the handle.

The volume range from 0.1 μl to 10000 μl is covered by 10 pipette models.

Model	Volume range [μl]
OP2	0.1 - 2
OP10	0.5 - 10
OP20	2 - 20
OP50	5 - 50
OP100	10 - 100
OP200	20 - 200
OP250	50 - 250
OP1000	100 - 1000
OP5000	500 - 5000
OP10000	1000 - 10000
OP2, OP10	Measurement and transfer of micro-volumes, DNA sequencing and enzyme-assay applications.
OP20, OP50, OP100, OP200, OP250, OP1000	Measurement and transfer of general aqueous solution, acids and bases
OP5000, OP10000	Measurement and transfer of large volumes.

The liquid is aspirated into disposable tips attached to the pipette shaft. Disposable tips ensure maximum safety and eliminate possibility of cross contamination of the liquid samples.

To protect the user from contaminated tips the **OPTIPETTE** is equipped with a built-in tip ejector, (Fig.1D).

The construction of the ejector enables the user to set up the length. The adjustable tip ejector accommodates every variety of tips available on the market. When using narrow tubes, it may be necessary to remove the tip ejector.

Adjusting of the Tip Ejector Length

- in 2-1000 μl pipettes (Fig. 6A)

The regulated "H" spacers, allow for regulating the length of tip ejector by +1 mm or +2 mm. An "H0" spacer is inserted on a standard basis. The outside shape of the spacer identifies the adjustment.

- in 5000 and 10000 μl pipettes (Fig. 6B).

The length of tip ejector is regulated by screwing in or screwing out the tip ejector stem with a screwdriver. Turn the screwdriver counterclockwise to increase the length of tip ejector, and clockwise to reduce the length of tip ejector. The ejector can be adjusted in the range of 5 mm.

If above described method of ejector adjustment is not sufficient or the diameter of the ejector opening is too large to dispose the tip it is necessary to put the ejector cap "M" onto the ejector, (Fig. 6C).

- in 2 and 10 μl pipettes (Fig. 6D)

The 2 μl - 10 μl pipettes may require the use of an ejector cap "M" to efficiently eject certain brands of pipette tips. Simply place the cap on the bottom of the pipette shaft and slide the cap upwards until it surrounds the bottom of the tip ejector.

The **OPTIPETTE** is a high quality instrument which offers excellent accuracy and precision.

The accuracy and precision (repeatability) of liquid volume depend on the quality of tips used. The values for accuracy and precision given in the table below were obtained using **HTL** tips. Only those tips guarantee correct operation of the pipettes and ensure accuracy and precision of liquid sampling.

Model	Cat. no.	Volume [μl]	Accuracy [%]	Precision [%]	Tip μl
OP2	5601	0.2 1.0 Max 2.0	± 12.0 ± 2.7 ± 1.5	± 6.0 ± 1.3 ± 0.7	10 μl
OP10	5602	Min 0.5 5.0 Max 10.0	± 4.0 ± 1.0 ± 0.5	± 2.8 ± 0.6 ± 0.4	
OP20	5603	Min 2 10 Max 20	± 3.0 ± 1.0 ± 0.8	± 1.5 ± 0.5 ± 0.3	
OP50	5607	Min 5 25 Max 50	± 2.5 ± 1.0 ± 0.8	± 2.0 ± 0.6 ± 0.4	200 μl
OP100	5604	Min 10 50 Max 100	± 1.6 ± 0.8 ± 0.8	± 0.80 ± 0.24 ± 0.20	
OP200	5605	Min 20 100 Max 200	± 1.2 ± 0.8 ± 0.6	± 0.60 ± 0.25 ± 0.20	
OP250	5600	Min 50 125 Max 250	± 1.0 ± 0.8 ± 0.6	± 0.4 ± 0.3 ± 0.3	300 μl
OP1000	5606	Min 100 500 Max 1000	± 1.6 ± 0.7 ± 0.6	± 0.40 ± 0.20 ± 0.15	
OP5000	5608	Min 500 2500 Max 5000	± 1.2 ± 0.6 ± 0.5	± 0.50 ± 0.20 ± 0.15	
OP10000	5609	Min 1000 5000 Max 10000	± 2.5 ± 0.8 ± 0.5	± 0.6 ± 0.3 ± 0.2	10000 μl

The accuracy and precision are obtained with **HTL** tips, using a gravimetric method, performing at least 10 mea-

surements of distilled water at a temperature of $20 \pm 1^\circ\text{C}$, according to EN ISO 8655 standard.

The pipette design enables the user to perform the recalibration process according to the rules presented in section 8.


2 - SETTING THE VOLUME

The counter displays three figures to be read from top to bottom. Additional to the figures on the lower wheel are printed graduations to enable a volume setting in the range increment of each pipette model.

Examples of the meaning of the black and red figures:


Pipettes **OP2**

Red figures at the bottom = $1/100 \mu\text{l}$
Increment = $0.002 \mu\text{l}$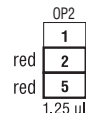

Pipettes **OP10, OP20,**

Red figures at the bottom = $1/10 \mu\text{l}$
Increment = $0.02 \mu\text{l}$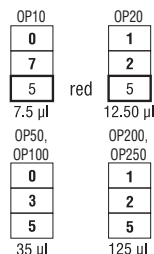


Pipettes **OP50, OP100, OP200, OP250**

Black figures only = $1 \mu\text{l}$
Increment = $0.2 \mu\text{l}$

Pipettes **OP1000, OP5000,**

Red figures at the top = ml
Increment =

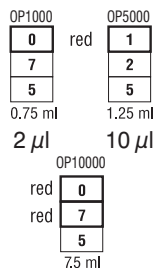Pipettes **OP10000**

Red figures at the top = ml
Increment = $20 \mu\text{l}$

The volume of the pipette is set by the black adjustment knob (Fig. 1B). To attain the maximum accuracy, set volume must be approached from a higher value by diminishing counter readings.

- If the desired volume is lower than set volume shown by the counter, the operator should turn the or the black adjustment knob (Fig. 1B) to the direction diminishing counter readings to the required volume. Before achieving the required volume slowly rotate the knob and observe carefully diminishing reading to avoid accidentally passing the setting value.
- If the desired volume is higher than set volume shown by the counter, the operator should turn the black adjustment knob (Fig. 1B) increasing the value until the lower figure wheel comes 1/3 of a turn beyond the required setting and then slowly backward until the setting reaches the desired volume. Make sure not to pass the setting value.

If the knob is accidentally turned too far, the process must be repeated. The desired volume must always be set from the higher value in the order of decreasing value.

3 - ASPIRATING AND DISPENSING LIQUIDS

Place a tip on the shaft of the pipette. See section 6 for the appropriate tip. Press the tip on firmly using a slight twisting motion to ensure a positive, airtight seal.

Important: Never aspirate liquids into the OPTIPETTE without a tip attached.

Aspiration

Press the pushbutton to the first positive stop, (Fig. 2A). Holding the pipette vertically, immerse the tip into the sample liquid. The depth to which the tip is immersed in the sample liquid depends on the model.

Model	Immersion depth (mm)
OP2	≤ 1
OP10	≤ 1
OP20, OP50, OP100	2 ÷ 3
OP200, OP250, OP1000	2 ÷ 3
OP5000	3 ÷ 6
OP10000	5 ÷ 7

Release the pushbutton slowly and smoothly to aspirate the sample, (Fig. 2B). Wait one second and then withdraw the tip from the liquid. When the pipette tip is immersed not as deeply as the recommended depth or when the pipetting pushbutton is rapidly released air may enter the disposable tip.

Avoid touching the orifice of the tip.

Dispensing

- Place the end of the tip against the inside wall of the vessel at an angle of 10 to 40 degrees.
- Press the pushbutton smoothly to the first stop, (Fig. 2C). Wait one second.
- Press the pushbutton to the second stop to expel any remaining liquid, (Fig. 2D).
- Keeping the pushbutton depressed to the very end, remove the pipette by drawing the tip against the inside surface of the receiving vessel.
- Release the pushbutton to its starting position, (Fig. 2E).
- Eject the tip by pressing the tip ejector button, (Fig. 2F). Remember to change the tip whenever a different kind of liquid is to be sampled.

Filters

Replaceable filter installed in a seat in the bottom part of the shaft is used in 5000 µl and 10000 µl pipettes (Fig. 3L). The filter prevents the aspirated liquid from entering into the shaft and thus from polluting the inside of the shaft and the piston. Using the filter is especially important when aspirating and dispensing large volumes of liquid.

If the filter becomes wet during liquid aspiration it should be replaced with a new one.

4 - PRE-RINSING

When pipetting liquids of higher viscosity or lower surface tension than water (e.g. sera or organic solvents), a film of liquid is formed on the inside wall of the pipette tip. This film can create an error. Since the film remains relatively constant in successive pipetting operations with the same tip, this error can be avoided by forming the film before transferring the first sample. This is done by aspirating a sample and dispensing it back into the same vessel. Since the film is already formed, all of the following samples will have better accuracy and repeatability.

This pre-rinsing operation should be repeated when the volume to be aspirated is changed or when a new tip is used.

5 - DENSE AND VISCOUS LIQUIDS

The **OPTIPETTE** specifications of accuracy and precision are based on pipetting distilled water. The handling of liquids with physical qualities of density, viscosity and surface tension differing extremely from water may need a gravimetrically checked compensation of the volume setting. Normally the degree of error resulting from heavy or viscous liquids is negligible if the pipetting is done slowly and carefully. It is most important to give the liquids some time to follow the change of pressure by holding the pipette tip in its position for at least 2 sec. after the aspiration and the blow out stroke.

If in extreme cases this method of operation does not result in accurate values, a compensation could be achieved as follows:

Weigh the liquid pipetted when the **OPTIPETTE** is set to the nominal value. Then calculate the set-off from the nominal value:

$$\text{Corr: val.} = 2 \times \text{nom. val.} - \frac{m}{\gamma}$$

m - weight of the sample

γ - density of liquid

Check this operation once again and correct if necessary. Note the corrected value for further pipetting the same kind of liquid.

6 - HTL PIPETTE TIPS

HTL tips are made from high performance polypropylene and their quality guarantees the precision and accuracy associated with the **OPTIPETTE**. Strict control is maintained throughout the manufacturing process to ensure the highest quality.

The accuracy and precision figures for the **OPTIPETTE** are only guaranteed when HTL tips are used. The use of inferior quality tips will seriously degrade the performance of the **OPTIPETTE**.

Tips 10

These tips are used for volumes between 0.1 μl and 10 μl . They are used with the OP2 and OP10 models.

Tips 200

These tips are used for volumes between 2 μl and 200 μl . They are used with the OP20, OP50, OP100 and OP200 models.

Tips 300

These tips are used for volumes between 50 μl and 250 μl . They are used with the OP250 models.

Tips 1000

These tips are used for volumes between 100 μl and 1000 μl . They are used with the OP1000 models.

Tips 5000

These tips are used for volumes between 500 μl and 5000 μl . They are used with the OP5000 models.

Tips 10000

These tips are used for volumes between 1000 and 10000 μl . They are used with the OP10000 models, which are equipped with a white pushbutton.

7 - RECOMMENDATIONS

Observing the following recommendations will ensure maximal possible accuracy and precision of liquid sampling.

- Make sure to operate the **OPTIPETTE** slowly and smoothly.
- The depth of immersion in the sample liquid should be the minimum necessary and should remain constant during aspiration.
- The **OPTIPETTE** should be held in a vertical position.
- Change the tip when volume setting is changed or when a different liquid is to be aspirated.
- Change the tip if a droplet remains on the end of the tip from the previous pipetting operation.
- Each new tip should be pre-rinsed with the liquid to be pipetted.
- Liquid should never enter the **OPTIPETTE** shaft. To prevent this:
 - Press and release the pushbutton slowly and smoothly.
 - Never turn the pipette upside down.
 - Never lay the pipette on its side when there is liquid in the tip.
- Never force the volume setting beyond its recommended limits.
- When pipetting liquids with temperatures different from the ambient temperature, it is recommended to pre-rinse the tip several times before use.
- Do not pipette liquids with temperatures above 70°C.

- When pipetting acids or corrosive solutions which emit vapours, it is recommended to disassemble the shaft and to rinse the piston and seal with distilled water after finishing the pipetting operation.

The pipettes are UV resistant which was confirmed by our tests. The recommended distance from the radiation source to exposed element should be not less than 50 cm. Prolonged and very intense UV exposure can cause de-coloration of pipette parts, without affecting its performance.

8 - RECALIBRATION

HTL pipettes are calibrated by gravimetric method, using HTL tips and distilled water, at the temperature $20 \pm 1^\circ\text{C}$, according to EN ISO 8655 standard.

If during pipette operation you find that the accuracy error (the difference between the real aspirated volume and the preset volume) exceeds the permissible value given in the table in section 1, the pipette recalibration procedure should be carried out.

Before starting the recalibration it is necessary to check whether the following requirements have been fulfilled during error determination:

- the ambient temperature, and the temperature of the pipette, tips and water was identical
- the density of the liquid used is close to that of distilled water
- the balance with appropriate sensitivity has been used.

Volume checked [μL]	Balance sensitivity [mg]
0.1 - 10	≤ 0.001
10 - 100	≤ 0.01
> 100	≤ 0.1

- mg/ μL conversion factor has been taken into account
- the requirements given in sections 3 and 7 have been fulfilled

If the above conditions are satisfied and the accuracy error for selected volume given in section 1 exceeds the permissible value, the pipette recalibration procedure should be carried out.

The recalibration can be performed within one full turn of the key to the right or to the left only.

Recalibration conditions:

- Ambient temperature and the temperature of the pipette, tips and liquid should be within the range $20 - 25^\circ\text{C}$ and stabilised during weighing within $\pm 0.5^\circ\text{C}$
- Measurements should be conducted using distilled water
- Balance sensitivity should be suitable for the volume to be controlled

Recalibration procedure:

- Set the dose volume depending on the pipette volume according to the following table:

Model	Range of the pipette volumes [μL]	Preset volume [μL]	Permissible volumes [μL]	Volume change ΔV for full turn of the calibration key [μL] (24 increments)
OP2	0.1 - 2	0.2	0.176 - 0.224	0.06
OP10	0.5 - 10	0.5	0.48 - 0.52	0.33
OP20	2 - 20	2	1.94 - 2.06	0.63
OP50	5 - 50	5	4.875 - 5.125	2.50
OP100	10 - 100	10	9.84 - 10.16	2.50
OP200	20 - 200	20	19.76 - 20.24	6.30
OP250	50 - 250	50	49.5 - 50.5	6.30
OP1000	100 - 1000	100	98.4 - 101.6	25.00
OP5000	500 - 5000	500	494 - 506	125.00
OP10000	1000 - 10000	1000	975 - 1025	250.00

- Perform 5 aspirations, weigh each one and calculate the average value of the aspirations.
- Calculate average aspirated volume in [μL] multiplying the average aspiration amount [mg] by the distilled water density coefficient [$\mu\text{L}/\text{mg}$], which depends on temperature and pressure according to the following table:

Temperature [$^\circ\text{C}$]	Pressure [kPa]		
	95.0	101.3	105.0
20	1.0028	1.0029	1.0029
21	1.0030	1.0031	1.0031
22	1.0032	1.0033	1.0033
23	1.0034	1.0035	1.0036
24	1.0037	1.0038	1.0038
25	1.0039	1.0040	1.0040

If the average aspirated volume exceeds the permissible value, the following should be done:

- Remove the pipetting pushbutton, (Fig. 4A),
- Holding the volume setting knob to protect it against rotation, insert the calibration key into the cuts of the calibration screw, (Fig. 4B),
- Turn the key clockwise to reduce the aspirated volume, or counterclockwise to increase the volume, (Fig. 4C). One full turn of the calibration key changes the pipette aspiration volume by the amount given in the table.
- Take out the key and fix the pipetting pushbutton (Fig. 4D).

Determine the average aspirated volume. The average volume should be within the permissible range given in the table. If the volume exceeds the values stated, the recalibration procedure should be repeated.

In case of pipetting the liquids with physical properties considerably different from those of water, follow the rules given in section 5.

More information on calibration procedure can be found on www.htl.com.pl

9 - TROUBLESHOOTING

If you notice an improper pipette operation identify the cause and eliminate the fault. Doing this, follow the instruction in the sequence provided. Replacement of elements into new ones may be required only exceptionally, and should not occur under normal pipette use.

Droplets of liquid remain in the pipette tip.

- The tip is emptied too fast.
Decrease the speed of pressing the pipette pushbutton.
- The tip wettability has increased due to extensive use.
Replace the tip with a new one.

Droplets of air appear in the liquid aspirated into the tip.

- The pipette tip immersion is too shallow.
Immerse the tip deeper according to the instructions.
- The pipette tip is incorrectly pressed onto the pipette shaft.
Press the pipette firmly.

- The tip is damaged or worn out due to extensive use.
Replace the tip with a new one.

The pipette incorrectly aspirates the liquid or liquid drops out from the tip.

- The pipette tip is incorrectly pressed onto the pipette shaft.

Press the pipette tip firmly.

- The shaft nut is loose (Fig. 3F).

Tighten the shaft nut.

- The sealing surface of the shaft is cracked or scored.
Remove the tip ejector. Unscrew the shaft nut, inspect the shaft and the piston assembly. Replace the damaged parts (see section 11). When reassembling the pipette, the nut should be hand tightened. In the models OP2, OP10 and OP20, the damage of the shaft may also cause a damage of the piston assembly. Replace the damaged parts (see section 11). When reassembling the pipette, the nut should be hand tightened.

To remove the tip ejector in models OP5000 and OP10000, remove the ejector pushbutton (Fig. 3N) and using a screwdriver unscrew the tip ejector by turning the screwdriver counterclockwise.

- Damage of the piston or seal due to prolonged use with the aggressive liquids.

Disassemble the pipette as described above. Replace the piston, seal and O-ring (see section 11). Rinse the inside of the shaft in distilled water and dry. Lubricate the seal and O-ring with the lubricant. The replacement of the piston requires conducting of calibration procedure.

Note: The parts of OP2 and OP10, pipette should be lubricated evenly with the minimum amount of lubricant.

- The pipette is reassembled improperly.
Disassemble the pipette and reassemble it, observing the proper sequence of steps (Fig. 3).
- No lubricant on the sealing elements.

Remove the tip ejector. Unscrew the shaft nut, remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts in distilled water and dry thoroughly. Lightly lubricate the inside surfaces of the seal and the O-ring with the lubricant. Reassemble the pipette in the reverse order.

- Contamination of the inside of the pipette caused by extensive aspiration of chemically aggressive liquids or because liquid got inside the pipette.

Remove the tip ejector. Unscrew the nut, remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts with distilled water and dry thoroughly. Lightly lubricate the inside surfaces of the seal and the O-ring with the lubricant. Reassemble the pipette in the reverse order.

Note: The shaft, shaft nut, ejector, piston assembly, seal and O-ring may be autoclaved at a temperature of 121°C for 20 minutes at pressure 1 bar.

The shaft of the 5000 and 10000 models should be autoclaved without the filter.

The handle and the outside of the pipette may be cleaned using a cloth damped in isopropanol.

Warning: Before using cleaning agents other than those recommended by the manufacturer, check the compatibility charts and consider chemical resistance of the following plastics: PP, PC, POM, PA, PPS, PVDF used to produce the pipette parts.

If the pipette malfunction persists after carrying out the above steps, send the pipette to your HTL service representative.

Before returning the pipette, make sure that it is free from any chemical, radioactive or microbiological contamination that might pose any danger during transportation and repairing. Clean the pipette as thoroughly as possible.

10 - PIPETTE KIT

The pipettes are delivered in the kits including:

- Pipette
- Short instruction manual
- Calibration key
- Ejector cap (for pipette models OP10000)
- Ejector spacers (for pipette models OP2-OP1000)
- CD

Note:

The following parts are not part of the standard equipment of the pipette and are available upon request:

- Ejector cap (for pipette models OP2, OP10, OP5000)
- Pipette stand
- Lubricant
- Filters (for pipette models OP5000, OP10000)
- Color identification rings (for pipette models OP2-OP1000)

11 - SPARE PARTS AND ACCESSORIES

All the spare parts and accessories indicated in Fig. 3, 4, 5 and 6:

A: Pushbutton

B: Adjustment knob

C: Shaft

D: Ejector

F: Shaft nut

G: Piston assembly

I: O-ring

J: Seal

K: Calibration key

L: Filter

M: Ejector cap

N: Ejector pushbutton

P: Identification ring

S: Pipette stand

The stand assembly diagram is shown in Fig. 5.

Spare parts and accessories can be ordered from your HTL representative (type of pipette and name of the part for this pipette should be specified).

Warning: The replacement of the piston requires conducting of calibration procedure according to section 8.

All rights reserved. Product described in this manual is subject to availability and technical modification. Errors excepted.

PZ HTL S.A. reserves the right to improve, enhance or otherwise modify its products without prior notification.

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© 2015 PZ HTL S.A.