

Single Channel Digital Pipette

Handrop® Plus



Single-channel variable volume pipette

- Ergonomic design with perfect fit in the hand, thanks to its design and non-slip bands
- Easy-to-read digit display
- Easy user calibration
- Precise volume selection mechanism activated by the top button. The selected volume is displayed digitally in an easy-to-read side window
- Automatic tip ejector
- Full volume range: 0.1 μ l to 5 ml
- Materials with excellent chemical resistance
- Autoclavable
- 1 year warranty

RECOMMENDATIONS FOR PIPETTING

- Hold the pipette upright while aspirating the liquid and submerging the tip a few millimeters into the liquid.
- Pre-rinse the tip, inhaling and expelling the liquid 5 times. It is especially important when dispensing viscous liquids and liquids of different density than water.
- Always control the movements of the pipetting button with your thumb to ensure balance.
- When pipetting liquids at different temperatures, rinse the tip several times before use.

MAINTENANCE

To obtain the best results, the cleanliness of the pipette should be checked. Special attention must be observed with the cone(s) of coupling.

The design of the pipettes allows maintenance to be carried out in the laboratory. However, we can provide you with a comprehensive repair and calibration service that includes a service report and calibration certificate.

Note: Periodically check the functioning of the pipette, e.g. every 3 months and always after any manipulation.

Pipette cleaning

To clean the pipette, use ethanol and a soft, lint-free tissue. It is recommended to clean the coupling cone regularly.