TaqMan Assays for genetic variation research

Superior performance—reliable, robust solutions

Genetic variation: decoding the blueprint for biodiversity

Research on genetic variation in animals and plants has expanded our understanding of evolution and human diseases, accelerated the pace of drug development, and helped identify and breed agricultural traits to improve the world's food and fuel supply. Researchers are looking to uncover the association between genetic makeup and phenotypes in studies focusing on single nucleotide polymorphisms (SNPs), copy number variants (CNVs), insertion/deletions (indels), and somatic mutations. A genomics revolution, fueled by advances in biotechnology tools, has significantly increased the rate at which we are able to obtain and analyze data to better understand biodiversity.

We're at the forefront of this revolution, and our reagents, Applied Biosystems™ TaqMan™ Assays, and Applied Biosystems™ platforms for genetic variation analysis, are the preeminent real-time PCR tools for variation research.

Coupled with Applied Biosystems™ capillary electrophoresis, and Ion Torrent™ DNA sequencing systems, we offer a complete solution for genetic analysis research, from discovery to confirmation.

TaqMan Assays for analyzing genetic variation

TaqMan Assays comprise preoptimized PCR primer pairs and one or two probes (depending on product family) for allelic discrimination or quantitative real-time PCR (qPCR). Each assay contains:

- An unlabeled PCR primer pair
- An Applied Biosystems™ TaqMan™ probe with a FAM™ or VIC™ dye label on the 5′ end, and a minor groove binder (MGB) and nonfluorescent quencher (NFQ) on the 3′ end

TaqMan Assays are used to amplify and detect specific variants in target genomic DNA (gDNA). Figure 1 depicts the Applied Biosystems™ TaqMan™ SNP Genotyping Assay process. Real-time PCR using TaqMan Assays is based on the 5′ nuclease activity of Taq DNA polymerase.

Here's how it works:

1. TaqMan probes hybridize to the target DNA between the two unlabeled PCR primers. Signal from the fluorescent dye on the 5′ end of a TaqMan probe is quenched by the NFQ on its 3′ end through fluorescence resonance energy transfer (FRET).
2. During PCR, Taq polymerase extends the unlabeled primers using the template strand as a guide.
3. When the polymerase reaches the TaqMan probe, it cleaves the molecule, separating the dye from the quencher. The qPCR instrument detects fluorescence from the unquenched FAM or VIC dye.

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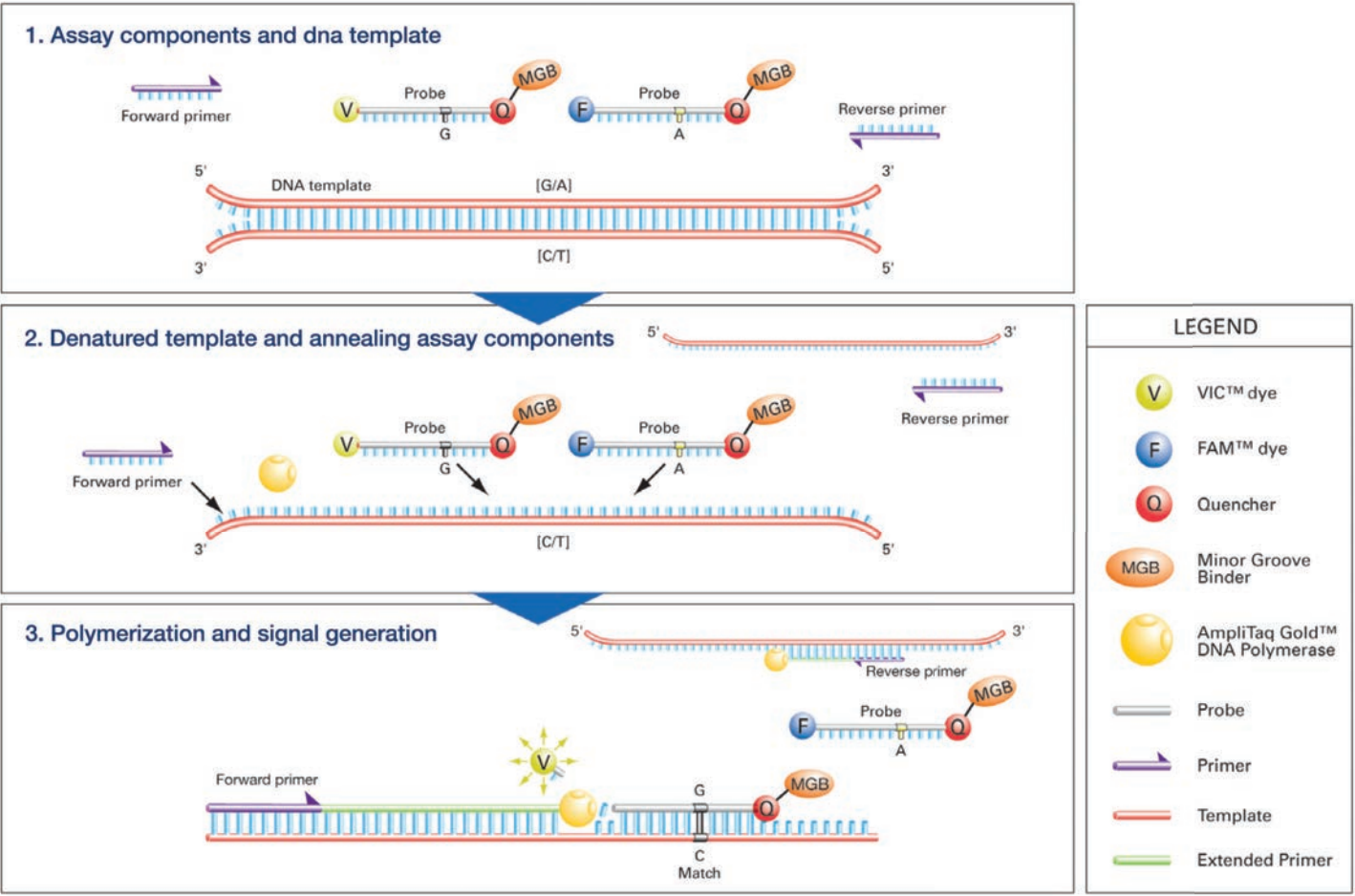
Figure 1. The TaqMan SNP Genotyping Assay. (1) The four TaqMan SNP Genotyping Assay components and the target DNA template with the SNP alleles (in brackets). (2) The denatured DNA target and annealing of the assay components. (3) Signal generation leading to specific allele detection.

TaqMan SNP Genotyping Assays

- **Better allelic discrimination**—TaqMan probes incorporate 3' MGB technology to stabilize the probe-template complex
- **Minimize failures**—TaqMan SNP Genotyping Assays are subject to a robust design pipeline, and functional QC testing for human assays on 20 gDNA samples
- **Full-coverage assay pool**—over 7 million human SNP assays (including 160,000 validated assays tested on four ethnic populations of 45 gDNAs each) and over 10,000 mouse SNP assays
- **Simplicity**—all probes and primers are contained in a single tube: no need to optimize probe, primer, salt concentrations, or temperature; all assays use universal PCR conditions
- **Integrated run and analysis solutions**—Applied Biosystems™ instruments and associated software help you move easily from run to results

SNPs are heritable single-base pair variations that occur throughout an organism's genome. SNPs comprise the most common form of genetic variation, with some estimates of SNPs in a given human genome numbering more than 10 million. SNP genotyping plays a central role in characterizing individuals and populations, studying disease traits in humans and other organisms, and identifying genes responsible for advantageous crop traits.

TaqMan SNP Genotyping Assays provide a highly flexible technology for detection of polymorphisms within any genome. TaqMan Assays have a simple workflow and provide a quick way to generate genotyping data (Figure 2). Based on powerful TaqMan chemistry and robust probe and primer designs, and coupled to dependable Applied Biosystems instruments and software, these made-to-order assays produce high-confidence results. TaqMan Assays are ideal for genotyping applications including association studies, candidate region or gene analysis, and fine-mapping studies.

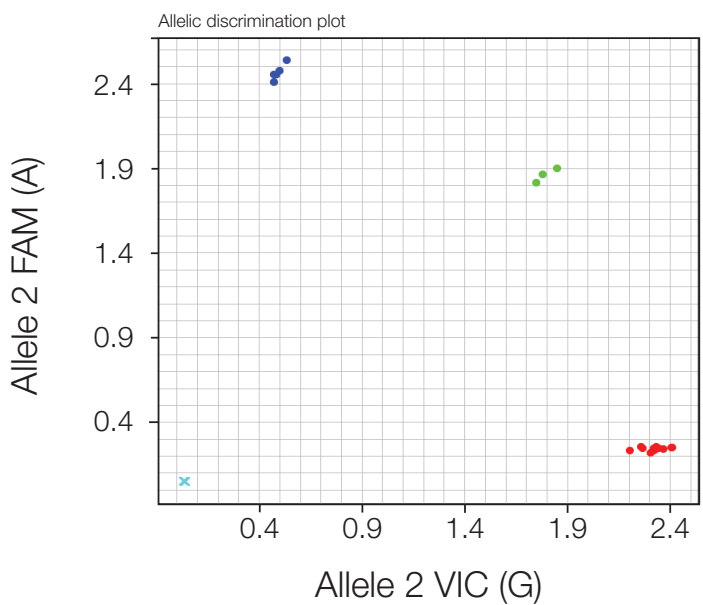
Figure 2. A three-cluster allelic discrimination plot generated with TaqMan SNP Genotyping Assay, C___1202883_20 (rs1801133) for *MTHFR* gene.

Easy online ordering

Predesigned TaqMan SNP Genotyping Assays

Find predefined assays using our new TaqMan Assay search tool at thermofisher.com/ordertaqman

- Easy-to-use interface with a powerful, logical search engine
- Search by keyword (gene, SNP ID) or genomic location
- Filter by SNP type (e.g., missense mutation, intronic, UTR)
- View results on a genome alignment map for easy selection

Custom TaqMan SNP Genotyping Assays

Can't find your assay in our predefined assay collection? Try designing a custom assay using our Applied Biosystems™ Custom TaqMan™ Assay Design Tool at thermofisher.com/snpcadt

- Manually enter your own custom target sequences or upload a file for batch design
- Enter custom primers and probes you have already designed to have us manufacture a ready-to-use assay for you

Simple workflow for quick results

TaqMan SNP Genotyping Assays constitute the simplest SNP genotyping technology available. We deliver your ready-to-use SNP genotyping assay in your choice of format: single-tube, 96- or 384-well plate (custom plating service), or Applied Biosystems™ TaqMan™ OpenArray™ plate (Figure 3). The rest is easy. Just combine the assay with Applied Biosystems™ TaqMan™ Genotyping Master Mix or TaqMan™ Universal PCR Master Mix and your purified DNA sample. There is no need to optimize probe, primer, salt concentrations, or temperature, because all assays use universal reagent concentrations and thermal cycling conditions. After generating an endpoint read using a thermal cycler or real-time PCR instrument, no transfers, washes, or additional reagents are required, and the plate remains sealed; just read the plate and analyze the genotypes. This helps reduce the chance of contamination, sample mix-ups, and sample loss. The simplicity of the chemistry allows you to easily automate the reaction for massively parallel genotyping studies, readily increasing the number of assays, number of samples, or both. Additionally, the analysis software allows you to auto-call genotypes, minimizing manual effort.

Simple data analysis

Applied Biosystems™ TaqMan™ Genotyper Software is a great resource for fast and accurate genotype calling. It is a free SNP genotyping data analysis tool for use with TaqMan SNP Genotyping Assays performed in 48-, 96-, or 384-well microtiter plates or OpenArray plates.

TaqMan Genotyper Software can be downloaded at thermofisher.com/taqmangenotyper

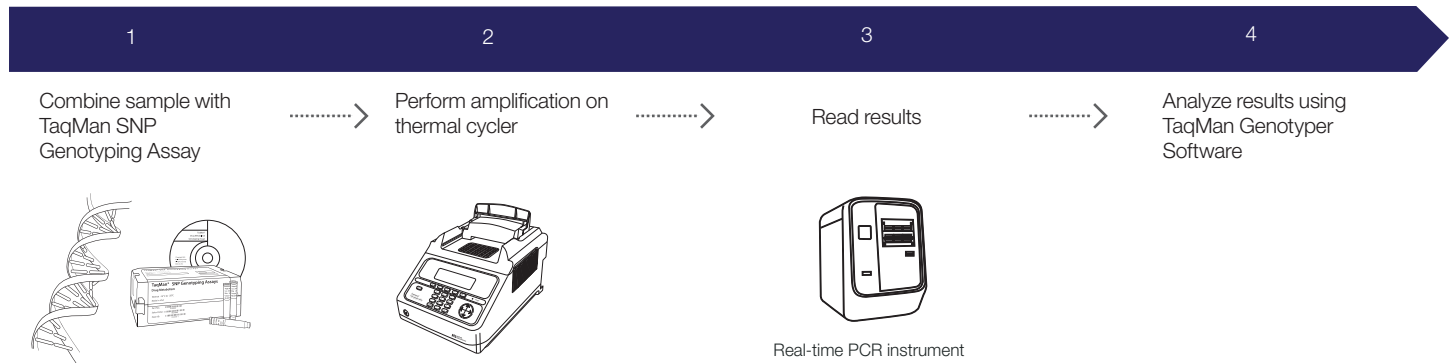
Figure 3. Workflow for TaqMan SNP Genotyping Assays.

Predesigned TaqMan SNP Genotyping Assays

Compatible Applied Biosystems™ TaqMan™ Master Mix and sample prep reagents have been developed to work in conjunction with TaqMan SNP Genotyping Assays to ensure high-quality results.

- TaqMan Genotyping Master Mix
- Applied Biosystems™ TaqMan™ Sample-to-SNP™ Kit
- Applied Biosystems™ TaqMan™ GTXpress™ Master Mix
- Applied Biosystems™ TaqMan™ Universal Master Mix II

The choice of which master mix to use depends on your sample type (tissue, blood, plant, etc.), sample preparation method (purified DNA or crude lysate), and use of fast or standard PCR

cycling. For more information, go to thermofisher.com/taqmansnp

Ordering information

	Number of SNPs	Number of 5 µL rxns (384-well plate)	Number of 25 µL rxns (96-well plate)	Assay mix formulation	Assay type	Human assays (Cat. No.)	Nonhuman assays (Cat. No.)
Predesigned TaqMan SNP Genotyping Assays for Human and Mouse							
Small-scale	>7 million	1,500	300	40X	Made-to-order	4351379	4351384*
Medium-scale	>7 million	5,000	1,000	40X	Made-to-order	4351376	4351382*
Large-scale	>7 million	12,000	2,400	80X	Made-to-order	4351374	4351380*
Custom TaqMan SNP Genotyping Assays							
Small-scale	∞	1,500	300	40X	Made-to-order	4331349	4332077
Medium-scale	∞	5,000	1,000	40X	Made-to-order	4332072	4332075
Large-scale	∞	12,000	2,400	80X		4332073	4332076
TaqMan Drug Metabolism Genotyping Assays							
Small-scale	2,700	750	150	20X	Inventoried	4362691	N/A

*Over 10,000 mouse assays available.

All assays are quality-control tested using a mass spectrometer to verify sequence and yield. In addition, all human (predesigned and custom) TaqMan SNP Genotyping Assays receive a genomic functional test on first synthesis. The subsequent syntheses of already-tested human assays and all nonhuman assays receive a fill volume check and mass spectrometry. All assays have a VIC dye–labeled probe, a FAM dye–labeled probe, and two target-specific primers.

Go to thermofisher.com/taqmansnp to order.

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- **Excellent ADME panel coverage**—target polymorphisms in 221 genes encoding drug metabolism enzymes and associated transport proteins
- **Simple protocol**—all assays in the collection are run under the same PCR conditions, and specific allele detection is achieved with the Applied Biosystems™ TaqMan™ 5´ nuclease chemistry
- **Detects multiple polymorphisms**—detect SNPs, insertion/deletions (indels), and multinucleotide polymorphisms (MNPs)
- **Rapid receipt of order**—performance-tested assays are already in inventory, ready to ship to you.
- **Assays match databases**—assays are aligned with allele nomenclature from public allele nomenclature sites

Pharmacogenetics is the study of how a person's genetic makeup affects how he or she responds to drugs. This research offers the promise of providing information that will not only allow current drugs to be dosed and delivered more effectively but also allow the development of drugs that are specifically tailored to treat an individual.

We offer 2,700 unique Applied Biosystems™ TaqMan™ Drug Metabolism Genotyping Assays for detecting polymorphisms in 221 genes that code for various drug metabolism enzymes (DMEs) and associated transport proteins. Polymorphisms

associated with these genes may influence the rate of drug metabolism within individuals, potentially affecting drug efficacy and the occurrence of side effects (Figure 4). The complex nature of these genes have had limited research conducted because few technologies and products could effectively characterize these polymorphisms. All of the assays in this collection target potentially causative polymorphisms, including those within regulatory elements, coding regions, and associated splice junctions.

TaqMan SNP Genotyping Assay technology delivers superior specificity
Each TaqMan Drug Metabolism Genotyping Assay contains two allele-specific probes and a primer pair to detect the specific SNP target. Both the probes and primers uniquely align within the genome, enabling the TaqMan genotyping technology to provide superior specificity. It is this specificity that allows these assays to detect targets residing in highly homologous gene families that may include pseudogenes.

TaqMan Drug Metabolism Genotyping Assays were developed using a high level of bioinformatics and wet-lab stringency. The assays were designed with information from several public SNP databases, including recognized public allele nomenclature sites. All assays have passed performance tests involving 180 unique DNA samples from four different populations.

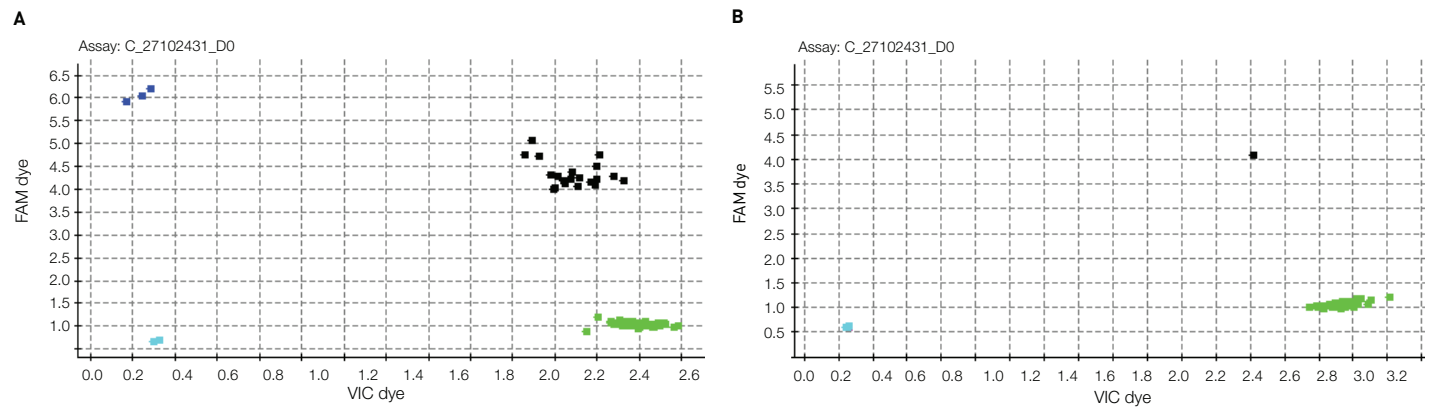
Figure 4. Allelic discrimination plots for the C__27102431_D0 assay run on (A) 45 each African-American and Caucasian, and (B) 45 each Chinese and Japanese gDNA samples. C__27102431_D0 targets the CYP2D6*4,g.1846G>A polymorphism, which encodes an mRNA splicing defect that results in a nonfunctional CYP2D6 protein. If an individual carries two nonfunctional CYP2D6 alleles, they will have the poor metabolizer (PM) phenotype and the metabolism of numerous drugs will be impacted.

Markers relevant for drug metabolism

The Applied Biosystems™ TaqMan™ DME Assay PharmaADME Core Marker Set contains a predefined group of TaqMan Drug Metabolism Genotyping and Applied Biosystems™ TaqMan™ Copy Number Assays, providing over 95% coverage of core markers in 33 ADME genes identified by the PharmaADME consortium.

This assay set greatly simplifies the study of these key putative functional genetic ADME variants and consists of:

- 164 DME assays for SNP and indel polymorphisms
- 14 copy number assays for copy number and hybrid gene variants

Assay sets are delivered in individual tubes, providing the flexibility to select a subset of assays or the entire PharmaADME Core Marker Set.

DME Assay Index

A DME Assay Index is also available with all drug metabolism assays. This file lists each assay along with context sequence, location on the NCBI assembly, the refSNP number (from dbSNP), and the common allele nomenclature from a public allele nomenclature site, when available.

Quick delivery, convenient format

For fast delivery, all assays in this collection have been manufactured and placed into inventory and are ready to ship at ambient temperature. Like other TaqMan SNP Genotyping Assays, these single-tube products consist of two allele-specific TaqMan MGB probes (labeled with either VIC or FAM dye) and two locus-specific primers. TaqMan Drug Metabolism Genotyping Assays are supplied as single tubes and in 96- and 384-well plates (custom plating service). Additionally, all products are formulated for the small-scale reaction size: a 20X single-tube assay, supporting 750 reactions at a 5 µL reaction size.

Optimized supporting reagents

Compatible TaqMan Master Mix and sample preparation reagents have been developed to work in conjunction with TaqMan Drug Metabolism Genotyping Assays to ensure high-quality results:

- TaqMan Genotyping Master Mix
- TaqMan Universal Master Mix II

Additional information about TaqMan Drug Metabolism Genotyping Assays, including links to the PharmaADME Core Marker Set and the DME Assay Index, can be found at thermofisher.com/taqmandme

TaqMan Copy Number Assays

- **Gold standard technology**—extraordinary accuracy and reliability; performance guaranteed for all predesigned assays**
- **Results in hours**—simplest method available to study CNV
- **Scalable solution**—automated workflow offers optimum platform for high-throughput validation of copy number changes
- **Comprehensive assay collection**—predesigned assays for human, mouse, and common vector marker/reporter genes
- **Option for custom assays**—Custom Plus and Custom TaqMan Assays for user-defined targets of interest

CNV, initially defined as variation in copy number of segments of DNA ≥1 kb in size, between individuals, is found in all humans as well as other animals and plants.

CNV affects a significant portion of the genome (approximately 12% of the human genome) and includes deletions, duplications, and other complex genotyping patterns. These CNVs can influence gene expression and be associated with specific phenotypes and diseases, as observed in microdeletion and microduplication syndromes.

Superior chemistry and streamlined methods offer reliable results

TaqMan Copy Number Assays combine Applied Biosystems™ TaqMan™ Assay chemistry with Applied Biosystems™ real-time PCR instruments to form a method for obtaining specific, reproducible, and easy-to-interpret copy number results (Figure 5). TaqMan Copy Number Assays are an ideal validation tool for microarray or next-generation sequencing follow-up studies and can be used to find specific targets. The workflow can be automated so that several hundred to thousands of samples can be processed in a single day.

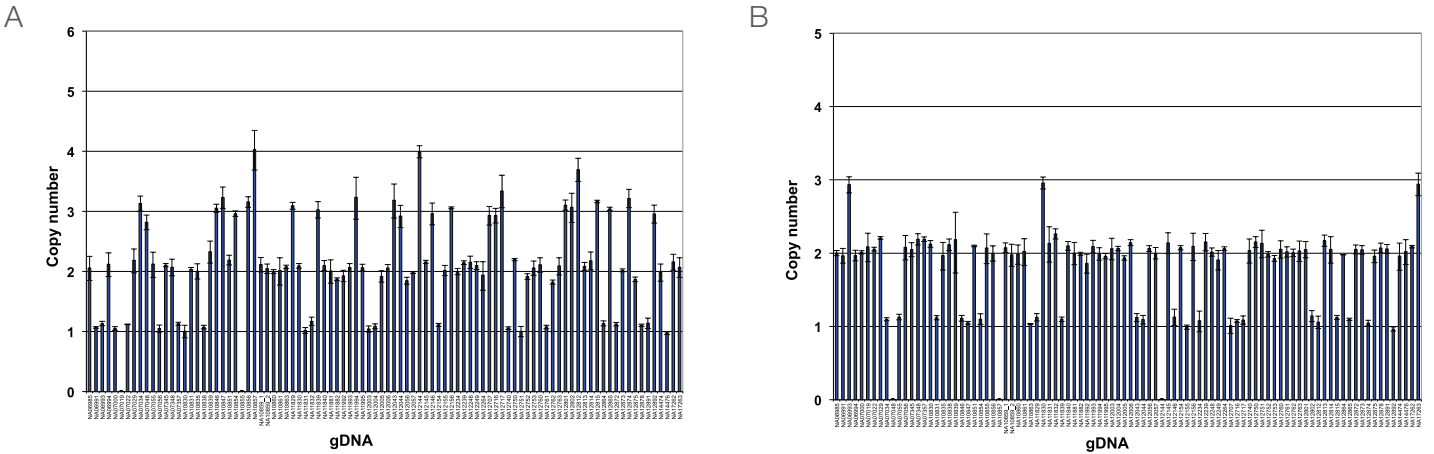
Figure 5. High specificity of TaqMan Copy Number Assays. C4A and C4B represent two isoforms of the C4 gene family. The sequences for these two genes differ in only 5 bases, but the encoded C4A and C4B proteins are functionally different. Differential detection of (A) C4A and (B) C4B is very challenging. Shown are TaqMan Copy Number Assays for C4A and C4B with the HAPMAP CEU sample set. Distinct copy number changes are observed. (JPT/CHB and YRI data not shown.)

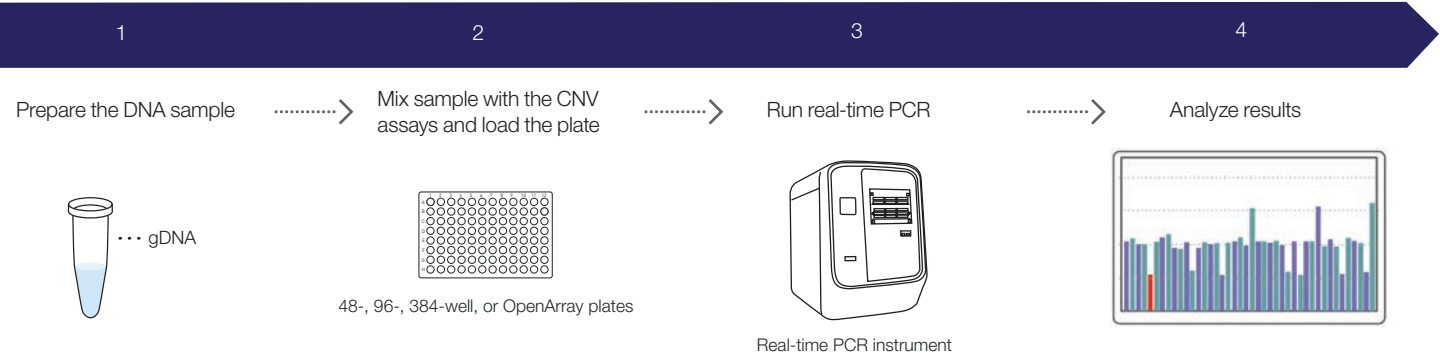
Figure 6. Workflow for TaqMan Copy Number Assays.

TaqMan Copy Number Assays

TaqMan Copy Number Assays include predesigned collections for both human and mouse genomes. The human collection includes more than 1.6 million assays for genome-wide coverage. The mouse predesigned collection includes more than 180,000 assays targeting gene exons. Predesigned assays to common vector marker and reporter genes are also available for transgenic studies.

Find predesigned assays using our online TaqMan Assay search tool at thermofisher.com/cnv

Applied Biosystems™ Custom Plus TaqMan™ Copy Number Assays are an optimal solution for studying variation in human and mouse genomic regions of interest for which a predesigned assay is not available. Custom Plus assays use the same bioinformatics pipeline used to manufacture predesigned TaqMan Copy Number Assays (which includes premasking of

SNPs and repetitive sequences and assay genome uniqueness checks) and can be generated for high-quality genomic targets of interest using the online Applied Biosystems™ GeneAssist™ Copy Number Assay Tool. Standard Custom TaqMan Copy Number Assays are an option for additional targets of interest. Unlike Custom Plus assays, standard Custom assay designs do not go through premasking or genome quality checks, but can be compared with the human or mouse reference assays for compatibility in duplex reactions.

Two Applied Biosystems™ TaqMan™ Copy Number Reference Assays are available for copy number analysis in both human and mouse species. Note that the reference assays are species-specific.

Feature	Predesigned TaqMan Copy Number Assay	Custom Plus TaqMan Copy Number Assay	Custom TaqMan Copy Number Assay
Designed using copy number–specific algorithm optimized for performance	✓	✓	✓
Availability limited to human and mouse assays	✓	✓	
Contains TaqMan FAM dye–labeled MGB probes and two unlabeled PCR primers	✓	✓	✓
Targets undergo SNP and repetitive sequence masking	✓	✓	
Genome specificity check	✓	✓	
Reference assay compatibility check	✓	✓ (optional)	✓
Assay sequences provided			✓
Assay context sequences and genome location provided	✓	✓	

A simple CNV analysis workflow

TaqMan Copy Number Assays have one of the simplest workflows of all currently available CNV analysis methods (Figure 6). The test assay (FAM dye–labeled), the reference assay (VIC dye–labeled), your sample DNA, and TaqMan Master Mix (TaqMan Genotyping Master Mix is recommended, with TaqMan Universal Master Mix II and Applied Biosystems™ TaqMan™ Gene Expression Master Mix also being compatible) are combined and then run on an Applied Biosystems real-time PCR system using standard TaqMan Assay PCR conditions. On average, setup to primary analysis takes only 3–4 hours (including a ~2 hour PCR run).

Analysis tools and methods

TaqMan Copy Number Assays are supplied in single tubes, or the assays can be custom-plated in 96- and 384-well plates. The assay reactions are run on a real-time PCR instrument, and the data are analyzed using Applied Biosystems™ CopyCaller™ Software.

Additional information on TaqMan Copy Number Assays, as well as links to CopyCaller Software and the GeneAssist Copy Number Assay Tool, can be found at thermofisher.com/cnv

	Number of 10 µL rxns (384-well plate)	Number of 20 µL rxns (96-well plate)	Assay mix formulation	Assay type	Cat. No.
Predesigned TaqMan Copy Number Assays					
Small-scale	720	360	20X	Made-to-order	4400291
Medium-scale	1,500	750	20X	Made-to-order	4400292
Large-scale	5,800	2,900	60X	Made-to-order	4400293
Custom Plus TaqMan Copy Number Assays					
Small-scale	720	360	20X	Made-to-order	4442487
Medium-scale	1,500	750	20X	Made-to-order	4442520
Large-scale	5,800	2,900	60X	Made-to-order	4442488
Custom TaqMan Copy Number Assays					
Small-scale	720	360	20X	Made-to-order	4400294
Medium-scale	1,500	750	20X	Made-to-order	4400295
Large-scale	5,800	2,900	60X	Made-to-order	4400296
TaqMan Copy Number Reference Assays (Human)					
RNase P	1,500	750	20X (1 tube)	Inventoried	4403326
RNase P	6,000	3,000	20X (4 tubes)	Inventoried	4403328
TERT	1,500	750	20X (1 tube)	Inventoried	4403316
TERT	6,000	3,000	20X (4 tubes)	Inventoried	4403315
TaqMan Copy Number Reference Assays (Mouse)					
Tfrc	1,500	750	20X (1 tube)	Inventoried	4458366
Tfrc	6,000	3,000	20X (4 tubes)	Inventoried	4458367
Tert	1,500	750	20X (1 tube)	Inventoried	4458368
Tert	6,000	3,000	20X (4 tubes)	Inventoried	4458369

Looking for a different formulation, scale, or label? The TaqMan Custom Assay and Oligo Service can accommodate special requests. To learn more, email specialoligos@thermofisher.com or contact your local sales representative.

Go to thermofisher.com/cnv to order.

TaqMan Mutation Detection Assays for somatic mutation detection

- **High specificity**—mutant allele detection is based on an allele-specific primer, while wild type background is suppressed by the proprietary MGB blocker oligonucleotide
- **High sensitivity**—assays can detect down to 0.1% mutant molecules in a background of wild type DNA, as demonstrated in spiking experiments (Figure 8)
- **Detect multiple types of mutations**—detect single- and multiple-nucleotide mutations and insertion/deletions (indels)
- **Wide dynamic range and excellent PCR efficiency**—assays demonstrate at least 4 logs of dynamic range and an average efficiency of 100% ± 10%
- **Fast, simple workflow**—like other TaqMan Assays, typically require 3 hours from sample to results, with minimum hands-on time

Somatic mutations can be present at low levels against a high background of wild type sequences, and methods used to detect and characterize these mutations in tumor specimens need to be highly sensitive and accurate. Methods that are commonly used include gene sequencing (including

pyrosequencing and traditional Sanger sequencing) and real-time PCR. Applied Biosystems™ TaqMan™ Mutation Detection Assays were designed based on a novel competitive allele-specific Applied Biosystems™ TaqMan™ (castPCR™) technology (Figure 7), which combines allele-specific TaqMan qPCR with an allele-specific MGB blocker oligonucleotide to effectively suppress nonspecific amplification of the off-target allele. These assays target mutations in 45 genes implicated in a number of cancer models:

ABL1, AKT1, ALK, APC, ATM, BRAF, CDH1, CDKN2A, CSF1R, CTNNB1, EGFR, ERBB2, ERBB4, FBXW7, FGFR1, FGFR2, FGFR3, FLT3, GNAS, HNF1A, HRAS, IDH1, JAK2, JAK3, KDR, KIT, KRAS, MET, MLH1, MPL, NOTCH1, NPM1, NRAS, PDGFRA, PIK3CA, PTEN, PTPN11, RB1, RET, SMAD4, SMARCB1, SMO, STK11, TP53, VHL

TaqMan Mutation Detection Assays
TaqMan Mutation Detection Assays contain mutant allele assays, which specifically detect one or more mutant alleles, and corresponding gene reference assays, which detect mutation-free regions of the genes in which the target mutations reside.

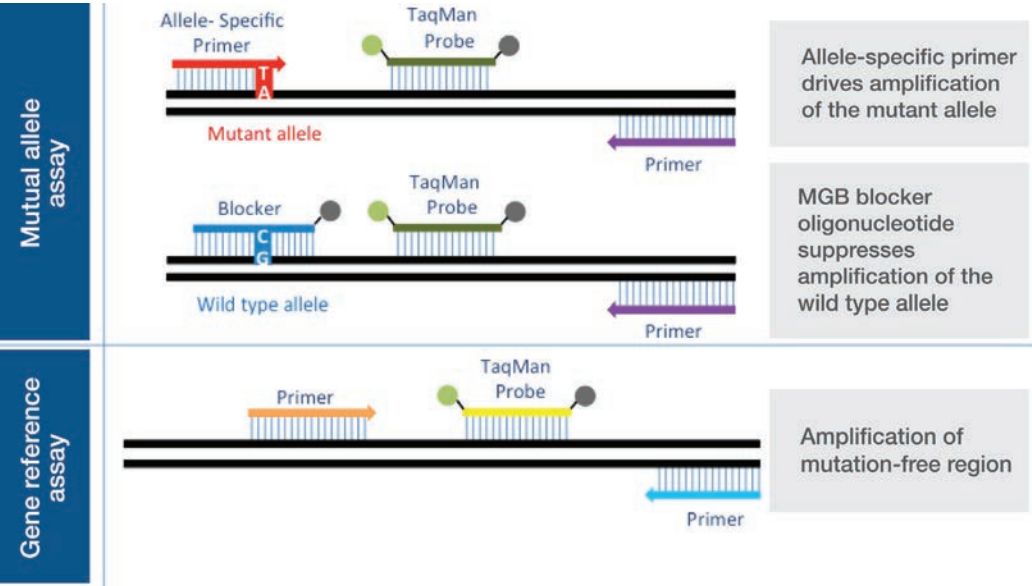
Figure 7. TaqMan Mutation Detection Assay types.

Two experiment types

Mutation detection analysis requires two types of experiments:

Detection ΔC_t cutoff determination

A mutant allele assay and corresponding gene reference assay are run on three or more wild type gDNA samples that are from the same sample type as the test samples (e.g., gDNA from FFPE tissue samples). The ΔC_t value is calculated for the mutant allele assay/gene reference assay pair, for each sample. The average ΔC_t for all samples is then calculated and is used to derive the detection ΔC_t cutoff value for the mutant allele assay.

Mutation detection

A test sample is run with one or more mutant allele assays and a corresponding gene reference assay. The ΔC_t for the mutant allele assay/gene reference assay pair is calculated, and this value is compared to the previously determined detection ΔC_t cutoff value to determine the sample's mutation status.

Simple workflow

Purified gDNA, extracted from a sample of unknown mutation status, is run with one or more mutant allele assays and corresponding gene reference assays. For each real-time PCR, the gDNA is combined with:

- A TaqMan Mutation Detection Assay
- TaqMan Genotyping Master Mix
- (Optional) Applied Biosystems™ TaqMan Mutation Detection IPC Reagent Kit—to distinguish true target negatives from PCR failure or inhibition

Reactions are run on a real-time PCR system using a universal thermal cycling protocol for mutation detection. After the run, the real-time PCR system analysis software determines the C_t for each TaqMan Mutation Detection Assay and (optional) IPC reagent reactions. Real-time results can be exported as files that can be opened in free Applied Biosystems™ Mutation Detector™ Software.

Ordering information

Product	Size	Assay type	Cat. No.
TaqMan Mutation Detection Assays	150 μ L, 10X	Inventoried	4465804
TaqMan Mutation Detection Reference Assays	150 μ L, 10X	Inventoried	4465807
TaqMan EGFR Exon 19 Deletions Assay	150 μ L, 10X	Inventoried	4465805
TaqMan Mutation Detection IPC Reagent Kit	1 kit	Inventoried	4467538

New assays for other cancer gene mutation targets will continually be released.

Go to thermofisher.com/castpqr for the most current list.

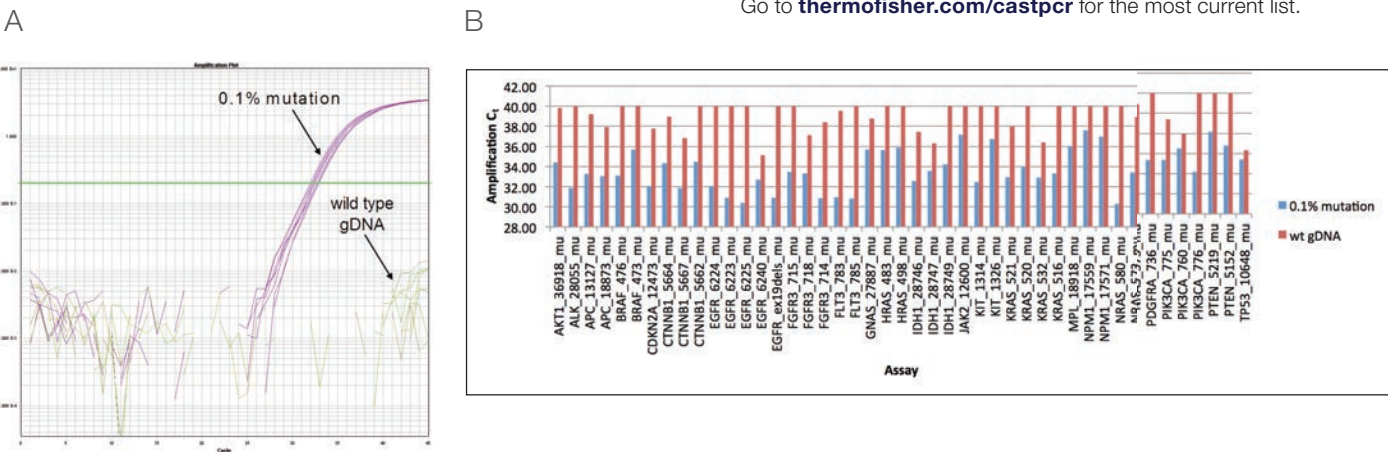
Figure 8. C_t differences between 0.1% mutation samples and wild type gDNA in TaqMan Mutation Detection Assays. For each mutant allele assay, 0.1% mutant allele samples were obtained by spiking cell line wild type gDNA (30 ng, ~10,000 copies) with 10 copies of mutant allele synthetic templates. (A) Example of amplification plot for KRAS_522_mu assay run on a 0.1% mutation sample and a wild type gDNA sample (30 ng gDNA). (B) For a majority of the assays, there is a significant difference in amplification C_t values between the 0.1% mutant allele sample and wild type gDNA (P value < 0.05).

TaqMan genotyping reagents for optimal performances

TaqMan Sample-to-SNP Kit

The TaqMan Sample-to-SNP Kit takes you from biological sample to results typically in less than an hour, without isolating DNA. The kit consists of two parts: the Applied Biosystems™ DNA Extract All Reagents and the TaqMan GTXpress Master Mix. The DNA All Lysis Reagents reduce prolonged procedures for the release of real-time PCR-ready DNA to a 5-minute protocol. They are compatible with a wide variety of samples ranging from blood to buccal swabs to animal and plant tissues. DNA extracted with DNA Extract All Reagents can be used with TaqMan SNP Genotyping Assays (not recommended for other TaqMan Assays).

TaqMan master mixes

TaqMan master mixes contain buffer, dNTPs, passive reference dye, thermostable hot-start DNA polymerase, and other components, and are provided in a convenient single-vial format. They are formulated to provide optimal results for TaqMan Assays.

- **TaqMan Genotyping Master Mix**—the TaqMan Genotyping Master Mix is optimized for end-point fluorescence detection in SNP genotyping applications in standard mode; the TaqMan Genotyping Master Mix provides excellent pre- and post-PCR stability for high-throughput setup and analysis
- **TaqMan GTXpress Master Mix**—the TaqMan GTXpress Master Mix is designed to deliver accurate genotyping results with robust performance in less than 50 minutes; the TaqMan GTXpress Master Mix is also available as part of the TaqMan Sample-to-SNP Kit

Ordering information and assay compatibility

	TaqMan Genotyping Master Mix	TaqMan GTXpress Master Mix
Cat. No. (size)	4371355 (10 mL) [†]	4401892 (10 mL)
TaqMan SNP Genotyping Assays	††	††
TaqMan Drug Metabolism Genotyping Assays	††	+
TaqMan Copy Number Assays	††	–
TaqMan Mutation Detection Assays for somatic mutation detection	††	–

[†]Other pack sizes are available.

^{††}Thermo Fisher Scientific validated: We have performed extensive testing and optimization.

+Thermo Fisher Scientific demonstrated: Limited testing has been performed. We cannot guarantee optimal performance for all TaqMan Assays.

–Not recommended.

Quality service and support at every step of your workflow

From manufacturing to follow-up—consistent reliability

TaqMan Assays are designed, manufactured, packaged, tested, and shipped using the highest-quality materials and methods. Furthermore, they are backed by our worldwide technical support teams.

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TaqMan Assays are manufactured in-house at our ISO 13485–certified manufacturing facilities and are never outsourced.

Comprehensive worldwide support

Whether you need help finding a TaqMan Assay for your target, deciding which format best suits your needs, placing your order through our online ordering system, or setting up your reactions, our sales and technical support staff are here to help.

Sales support

Your sales representative can help you find Web and print resources to help you choose the right TaqMan Assay products for your genetic variation research. For more demanding projects, she or he can also involve our technical sales specialists, who have more in-depth knowledge of TaqMan Assay technology and our relevant supporting reagents and instruments.

Technical support

If you have questions about how to use TaqMan Assays or how to analyze results, go to **thermofisher.com/support** to contact our technical support specialists. These agents are skilled in experimental planning and design, are expert troubleshooters, and are familiar with a wide variety of applications that use TaqMan Assays.

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TaqMan Assay type	Estimated delivery time (business days/weeks)
Inventoried (in stock)	1–4 days
Made-to-order/Custom TaqMan Assays (manufactured when order is placed)	5–12 days
TaqMan Custom Plating Service (configure 96- or 384-well plates with any TaqMan assays)	2–5 weeks

**The TaqMan Assays QPCR Guarantee

We stand behind every predesigned TaqMan Assay you buy. We’re committed to helping you achieve your research goals and believe our predesigned TaqMan Assays establish the benchmark for high-quality and easy-to-use real-time PCR products. If you are not satisfied with the performance of a predesigned TaqMan Assay, we’ll replace it at no cost or credit your account. For more information, and full terms and conditions of the guarantee, go to **thermofisher.com/taqmanguarantee**

appliedbiosystems

Find out more at thermofisher.com/taqman

ThermoFisher
S C I E N T I F I C

For Research Use Only. Not for use in diagnostic procedures. © 2015 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. TaqMan is a trademark of Roche Molecular Systems, Inc., used under permission and license. **CO37435 0915**

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RNeasy® Mini, Midi and Maxi Kits

For purification of total RNA from cells, tissues, and yeast, and for RNA cleanup*

RNeasy Mini, Midi and Maxi Kits provide a fast and convenient procedure for RNA purification, delivering high-quality RNA for use in many downstream applications. RNA is isolated on the silica membrane in trusted RNeasy spin columns, which have binding capacities of 100 µg (mini), 1 mg (midi) and 6 mg (maxi) RNA.

Benefits of RNeasy Mini, Midi and Maxi Kits:

- Fast procedure delivering high-quality total RNA in minutes
- Ready-to-use RNA for use in many downstream applications
- Consistent RNA yields from small to large samples
- No phenol/chloroform extraction or LiCl or ethanol precipitation

Fast, reproducible RNA purification

RNA purification with the RNeasy Mini Kit takes just 20 minutes, and can be performed using a microcentrifuge or QIAvac vacuum manifold or fully automated on the QIAcube®. With RNeasy Midi and Maxi Kits, RNA is purified in less than 1 hour using a centrifuge. All kits deliver reproducible yields of intact RNA with high RIN values and A_{260}/A_{280} ratios of 1.9–2.1 (Figures 1 and 3). Safe and convenient analysis of RNA integrity can be achieved using the QIAxcel® system (Figure 3A), which analyzes up to 96 samples on ready-to-use gel cartridges.

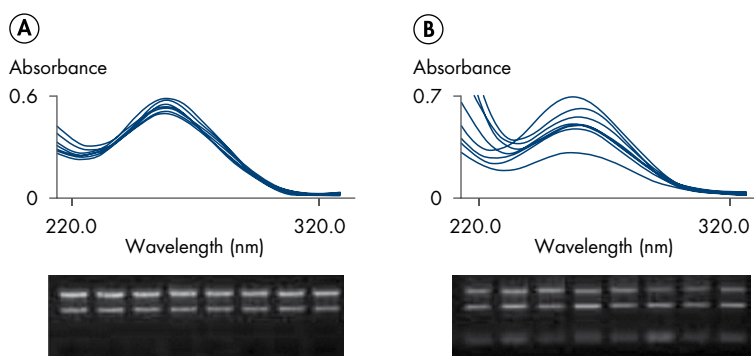
Figure 1. Reproducible RNA yields with RNeasy Kits. RNA was purified from eight HeLa cell samples (10^6 cells each) using either the RNeasy Mini Kit (**RNeasy**) or acid-phenol-extraction (**Acid Phenol**). **A** UV spectrophotometry of samples diluted in 10 mM Tris-Cl, pH 7.5. **B** Formaldehyde agarose gel analysis of 10 µg of each sample, based on A_{260} readings in **A**.

* For difficult-to-lyse tissues, specialized RNeasy Kits are available. For details, visit www.qiagen.com/RNA.

RNeasy Procedure

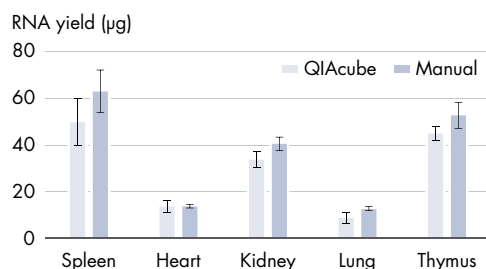
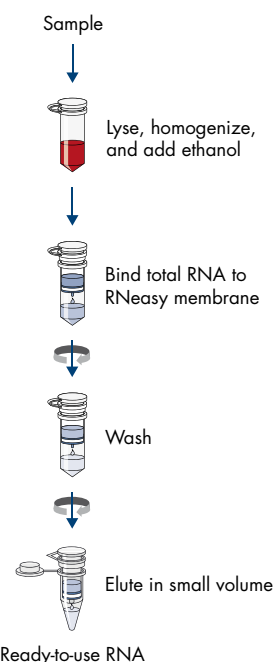
Figure 2. High RNA yields from various tissues. Rat tissues (15 mg) were stabilized in RNAlater® RNA Stabilization Reagent and disrupted using the TissueLyser II. RNA was purified using the RNeasy Mini Kit, either manually (**Manual**) or on the QIAcube (**QIAcube**).

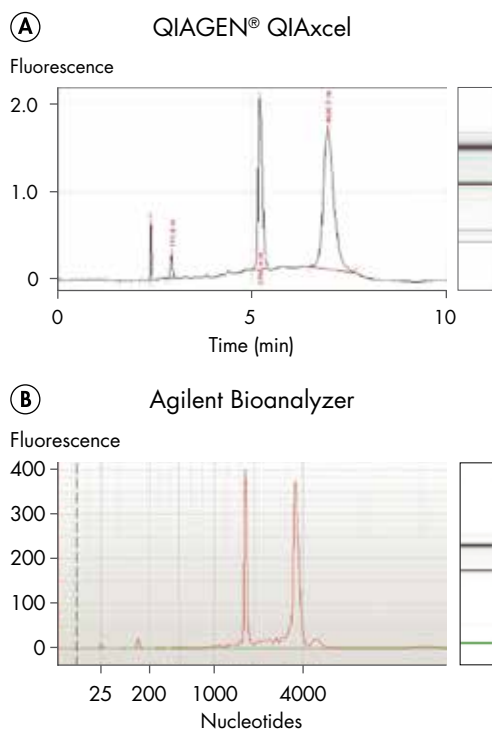
Figure 3. Highly intact RNA. RNA was purified from Jurkat cells using the RNeasy Mini Kit. The purified RNA was analyzed on the **A** QIAxcel system (ratio of 28S to 18S rRNA: 1.55) and **B** Agilent® 2100 Bioanalyzer (ratio of 28S to 18S rRNA: 1.7). A high RNA Integrity Number (RIN) of 9.6 was obtained, indicating highly intact RNA.

High performance in downstream applications

Ready-to-use RNA can be purified from a variety of sample types using RNeasy Kits, and yields are comparable when purification is either carried out manually or automated on the QIAcube (Figure 2). The innovative QIAcube automates the binding, washing, and elution of RNA, and can process up to 12 samples per run. RNA purified using RNeasy Kits is longer than 200 nucleotides and performs well in a wide range of applications, including RT-PCR and real-time RT-PCR (Figure 4).

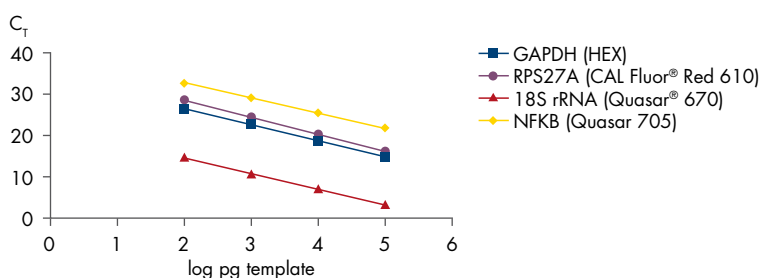
Figure 4. Reliable multiplex real-time RT-PCR. RNA was purified from HeLa cells using the RNeasy Mini Kit. 4-plex, real-time one-step RT-PCR was then performed on the Rotor-Gene® Q using the Rotor-Gene Multiplex RT-PCR Kit and self-designed TaqMan® assays for the indicated targets. The plots of C_t value versus log template amount were parallel, indicating all 4 targets were amplified with the same high efficiency.

Ordering Information

Product	Contents	Cat. no.
RNeasy Mini Kit (50)*	50 Spin Columns, Collection Tubes, Reagents, Buffers	74104
RNeasy Midi Kit (10)*	10 Spin Columns, Collection Tubes, Reagents, Buffers	75142
RNeasy Maxi Kit (12)*	12 Spin Columns, Collection Tubes, Reagents, Buffers	75162

* Mini and midi kits are available in larger kit sizes; for details, visit www.qiagen.com. Midi and maxi kits require use of a centrifuge capable of attaining 3000–5000 x g and equipped with a swing-out rotor for 15 ml (midi) or 50 ml (maxi) centrifuge tubes.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual. QIAGEN kit handbooks and user manuals are available at www.qiagen.com or can be requested from QIAGEN Technical Services or your local distributor. “RNA/ater®” is a trademark of AMBION, Inc., Austin, Texas and is covered by various U.S. and foreign patents.

Trademarks: QIAGEN®, Sample to Insight®, QIAcube®, QIAxcel®, RNeasy®, Rotor-Gene® (QIAGEN Group); Agilent® (Agilent Technologies, Inc.); CAL Fluor®, Quasar® (Biosearch Technologies, Inc.); TaqMan® (Roche Group).
© 2015 QIAGEN, all rights reserved. PROM-2155-002

Ordering www.qiagen.com/contact | Technical Support support.qiagen.com | Website www.qiagen.com

FastAP™ Thermosensitive Alkaline Phosphatase

Catalog Number EF0651, EF0652, EF0654

Pub. No. MAN0012876 Rev. C.00

WARNING! Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Sheets (SDSs) are available from thermofisher.com/support.

Contents and storage

Cat. No.	Contents	Amount	Storage
EF0651	FastAP™ Thermosensitive Alkaline Phosphatase	1000 U for 1000 reactions, 1 U/μL	-25 °C to -15 °C
	10X FastAP™ Buffer	2 x 1.5 mL	
EF0652	FastAP™ Thermosensitive Alkaline Phosphatase	5 x 1000 U for 5000 reactions, 1 U/μL	
	10X FastAP™ Buffer	10 x 1.5 mL	
EF0654	FastAP™ Thermosensitive Alkaline Phosphatase	300 U for 300 reactions, 1 U/μL	
	10X FastAP™ Buffer	1.5 mL	

BSA included

Description

Thermo Scientific™ FastAP™ Thermosensitive Alkaline Phosphatase catalyzes the release of 5'- and 3'- phosphate groups from DNA, RNA and nucleotides. This enzyme also removes phosphate groups from proteins. FastAP™ is a novel alkaline phosphatase, which is active in all Thermo Scientific restriction enzyme buffers as well as in PCR buffers. It dephosphorylates all types of DNA ends in 10 min at 37 °C. The enzyme is inactivated in 5 min at 75 °C. Therefore, removal of alkaline phosphatase is not required prior to ligation.

Applications

- Dephosphorylation of cloning vector DNA to prevent recircularization during ligation.
- Simultaneous digestion and dephosphorylation of vector DNA.
- PCR product clean-up: nucleotide degradation prior to sequencing of PCR product.
- Dephosphorylation of nucleic acid 5'-termini prior to labeling with T4 Polynucleotide Kinase.
- Other applications where dephosphorylation of DNA and RNA substrates is necessary.
- Protein dephosphorylation.

Source

E.coli cells with a cloned bacterial AP gene.

Definition of Activity Unit

One unit is the amount of the enzyme required to dephosphorylate 5'-termini of 1 μg of linearized pUC57 DNA in 10 min at 37 °C in FastAP™ buffer.

Storage Buffer

The enzyme is supplied in: 20 mM HEPES-NaOH (pH 7.4), 1 mM MgCl₂, 0.1 mM ZnCl₂, 0.1 % (v/v) Triton X-100 and 50 % (v/v) glycerol.

10X FastAP Buffer

100 mM Tris-HCl (pH 8.0 at 37 °C), 50 mM MgCl₂, 1 M KCl, 0.2 % Triton X-100 and 1 mg/mL BSA.

Inhibition and Inactivation

- Inhibitors: metal chelators.
- Inactivated by heating at 75 °C for 5 min.

Note

- Binding of FastAP™ Thermosensitive Alkaline Phosphatase to DNA may result in a band shift in agarose gels. To avoid this, incubate samples with 6X Loading Dye & SDS Solution (#R1151) at 65 °C for 10 min and chill on ice prior to electrophoresis.
- FastAP™ Thermosensitive Alkaline Phosphatase is active in all restriction enzyme buffers and may be added directly to digest DNA. Heat inactivation of the restriction enzyme before dephosphorylation reaction is not necessary.

Protocol for fast simultaneous plasmid vector linearization and dephosphorylation

1. Prepare the following reaction mixture containing:

Components	Volume
Plasmid DNA	1 µg
10X Thermo Scientific™ FastDigest™ Buffer	2 µL
FastDigest™ Restriction Enzyme	1 µL
FastAP™ Thermosensitive Alkaline Phosphatase	1 µL
Water, nuclease-free (#R0581)	to 20 µL
Total volume	20 µL

2. Mix thoroughly, spin briefly and incubate at 37 °C for 10 min.

3. Stop reactions by heating at 65 °C for 15 min or at 80 °C for 20 min (if restriction enzyme is not inactivated at 65 °C).

Note. For FastDigest™ SphI (PaeI) (#FD0604), simultaneous digestion and dephosphorylation is not recommended. Perform digestion, spin column purification and then dephosphorylation.

Protocol for nucleic acid dephosphorylation

This protocol is suitable for removal of 3' and 5' -phosphate groups from DNA and RNA.

1. Prepare the following reaction mixture:

Components	Volume
Linear DNA (~3 kb plasmid)	1 µg (~1 pmol termini)
10X reaction buffer for AP used in reaction	2 µL
FastAP™ Thermosensitive Alkaline Phosphatase	1 µL (1 U)
Water, nuclease-free (#R0581)	to 20 µL
Total volume	20 µL

2. Mix thoroughly, spin briefly and incubate 10 min at 37 °C.

3. Stop reaction by heating for 5 min at 75 °C.

Note. For efficient dephosphorylation plasmid DNA should be free of RNA and genomic DNA.

Protocol for dephosphorylation of proteins

Reaction mixture: 1X FastAP™ reaction buffer, 0.1-0.2 mg/mL of phosphoprotein, 10 U of FastAP™ Thermosensitive Alkaline Phosphatase. Incubate at 37 °C for 1 h.

For example: If you are doing a 20 µL reaction setup you need 2 µL 10X FastAP™ buffer, 2-4 µg of protein (to be in the range of 0.1-0.2 mg/mL) and 10 U of FastAP™ Thermosensitive Alkaline Phosphatase (1 U/µL).

Note

- The reaction can be stopped by addition of a final concentration of 50 mM EDTA (#R1021) or by addition of a final concentration of 10 mM sodium orthovanadate (Na₃VO₄).
- The optimal incubation time and the enzyme concentration must be determined experimentally for each substrate.

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

Thermo Fisher Scientific Baltics UAB | V.A. Graiciuno 8, LT-02241 Vilnius, Lithuania
For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2020 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

Buffered Saline Solutions

For various human clinical samples

Pub. No. MAN0018562 Rev. 3.0

 **WARNING!** Read the Safety Data Sheets (SDSs) and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Safety Data Sheets (SDSs) are available from [thermofisher.com/support](https://www.thermofisher.com/support).

Intended Use

For *in vitro* diagnostic use

Buffered saline solutions are isotonic saline solutions used to maintain pH and osmotic balance as well as provide cells with water and essential inorganic ions. Buffered saline solutions are used in a variety of laboratory applications, such as washing cells, transporting cells, diluent of human biological samples, or for reagent preparation. Thermo Fisher Scientific offers a wide range of buffered saline solutions including those with and without; calcium chloride, magnesium chloride, magnesium sulfate, and phenol red.

Buffered saline solutions are for professional use. They are used in medical laboratories by personnel who have received specialized education and training with regard to procedures utilizing In Vitro Diagnostic products. IVD products of this type are not intended as sole determinant in a diagnostic situation. Test results are interpreted by a healthcare professional as part of the clinical management of a patient.

Principle and explanation of procedure

Isotonic solutions are generally utilized to maintain cells for a short period time in a viable condition while the cells are manipulated outside of their regular growth environment. A physiological pH (6.8 to 7.4) value can be important to maintain viability when investigating human biological samples in an *in vitro* diagnostic application. Isotonic solutions can also be used for rinsing and washing steps in various diagnostic procedures (1-4).

Contents and storage

All quality control testing results are reported on lot-specific Certificate of Analysis available on our website: [thermofisher.com](https://www.thermofisher.com).

Product	Cat. No.	Storage	Shelf life ^[1]
Phosphate Buffered Saline (PBS) 7.4 (1X)	10010001 ^[2] 10010002 ^[2]	15°C to 30°C	12 months
Phosphate Buffered Saline (PBS) 7.4 (1X)	10010023 ^[3] 10010031 ^[2] 10010049 ^[3] 10010072 ^[2]	15°C to 30°C	24 months
Phosphate Buffered Saline (PBS) 7.4 (1X)	10010015 ^[4]	15°C to 30°C	36 months
Phosphate Buffered Saline (PBS) 7.2 (1X)	20012019 ^[4]	15°C to 30°C	36 months
Phosphate Buffered Saline (PBS) 7.2 (1X)	20012027 ^[2] 20012043 ^[2] 20012050 ^[2] 20012099 ^[2]	15°C to 30°C	24 months
Phosphate Buffered Saline (PBS) 7.4 (10X)	70011036 ^[4] 70011044 ^[2] 70011069 ^[2]	15°C to 30°C	24 months
Phosphate Buffered Saline (PBS) 7.2 (10X)	70013016 ^[4]	15°C to 30°C	36 months

For In Vitro Diagnostic Use.

Product	Cat. No.	Storage	Shelf life ^[1]
Phosphate Buffered Saline (PBS) 7.2 (10X)	70013032 ^[2] 70013073 ^[2]	15°C to 30°C	24 months
Dulbecco's Phosphate Buffered Saline (DPBS) (1X) [+] Calcium Chloride [+] Magnesium Chloride	14040083 ^[4] 14040091 ^[4] 14040141 ^[2] 14040133 ^[2] 14040117 ^[3] 14040097 ^[2] 14040182 ^[2] 14040216 ^[3]	2°C to 8°C	36 months
Dulbecco's Phosphate Buffered Saline (DPBS) (1X) [-] Calcium Chloride [-] Magnesium Chloride	14190086 ^[4] 14190094 ^[4] 14190144 ^[3] 14190136 ^[3] 14190230 ^[4] 14190250 ^[3] 14190235 ^[3]	15°C to 30°C	36 months
Dulbecco's Phosphate Buffered Saline (DPBS) (1X) [-] Calcium Chloride [-] Magnesium Chloride	14190342 ^[3] 14190359 ^[2] 14190367 ^[3]	15°C to 30°C	12 months
Dulbecco's Phosphate Buffered Saline (DPBS) (1X) [-] Calcium Chloride [-] Magnesium Chloride	14190240 ^[4]	15°C to 30°C	24 months
Dulbecco's Phosphate Buffered Saline (DPBS) (10X) [-] Calcium Chloride [-] Magnesium Chloride	14200059 ^[4] 14200067 ^[4] 14200075 ^[2] 14200166 ^[2]	15°C to 30°C	36 months
Dulbecco's Phosphate Buffered Saline (DPBS) (10X) [+] Calcium Chloride [+] Magnesium Chloride	14080048 ^[4] 14080055 ^[2] 14080098 ^[2]	2°C to 8°C	24 months
Hanks' Balanced Salt Solution (HBSS) (1X) [+] Calcium Chloride [+] Magnesium Chloride [-] Phenol Red	14025050 ^[4] 14025092 ^[2] 14025076 ^[3] 14025134 ^[2] 14025126 ^[2]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (1X) [+] Calcium Chloride [+] Magnesium Chloride [+] Phenol Red	24020083 ^[4] 24020091 ^[4] 24020117 ^[2] 24020198 ^[2]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (1X) [-] Calcium Chloride [-] Magnesium Chloride [-] Magnesium Sulfate [+] Phenol Red	14170120 ^[2] 14170112 ^[2] 14170097 ^[2] 14170161 ^[2]	15°C to 30°C	36 months

Product	Cat. No.	Storage	Shelf life ^[1]
Hanks' Balanced Salt Solution (HBSS) (1X) [-] Calcium Chloride [-] Magnesium Chloride [-] Magnesium Sulfate [-] Phenol Red	14175046 ^[4] 14175053 ^[4] 14175095 ^[2] 14175079 ^[2] 14175103 ^[2] 14175145 ^[2]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (1X) [-] Calcium Chloride [-] Magnesium Chloride [+] Phenol Red [+] Sodium Bicarbonate	14170070 ^[4] 14170088 ^[4]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (10X) [-] Sodium Bicarbonate	14060040 ^[4]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (10X) [-] Calcium Chloride [-] Magnesium Chloride	14180046 ^[4]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (10X) [-] Calcium Chloride [-] Magnesium Chloride [-] Sodium Bicarbonate [-] Phenol Red	14185045 ^[4]	15°C to 30°C	24 months
Hanks' Balanced Salt Solution (HBSS) (10X) [-] Sodium Bicarbonate [-] Phenol Red	14065056 ^[2] 14065098 ^[2]	15°C to 30°C	24 months
Hanks' Balanced Salt Solution (HBSS) (10X) [-] Sodium Bicarbonate [-] Phenol Red	14065049 ^[4]	15°C to 30°C	36 months
Hanks' Balanced Salt Solution (HBSS) (10X) [-] Calcium Chloride [-] Magnesium Chloride [-] Magnesium Sulfate	14185052 ^[2] 14185098 ^[2]	15°C to 30°C	24 months
Earle's Balanced Salt Solution (EBSS) (1X) [-] Calcium Chloride [-] Magnesium Chloride [-] Phenol Red	14155048 ^[4] 14155063 ^[2] 14155098 ^[2]	15°C to 30°C	36 months
BME, Basal Medium (Eagle) (1X) [+] Earle's Salts [+] Phenol Red [-] L-glutamine [-] HEPES	41010026 ^[4]	2°C to 8°C Protect from light	12 months
HEPES Buffer Solution 1 M	15630049 ^[4] 15630056 ^[4] 15630106 ^[2] 15630130 ^[2]	2°C to 8°C	24 months

^[1] Shelf life is determined from Date of Manufacture. Do not use beyond the labelled expiration date.

Precautions

Do not use the product if packaging, including bottles and vials, have been compromised and/or show evidence of microbial contamination, cloudy appearance, discoloration, drying, cracking, or other signs of deterioration.

 **CAUTION!** Human samples are potentially biohazardous. Follow standard precautions for handling, storage and disposal.

 **WARNING!** Do not use for injection or infusion! Please report any serious incidents in relation to the device to the manufacturer and the Competent Authority of the EU Member State in which the user and/or patient is established.

- Once opened, use Buffered Saline Solutions within 14 days for maximal growth performance.
- Avoid repeated warming/cooling and prolonged exposure to light.
- Do not use beyond labeled expiration date.
- All solutions that come into contact with clinical samples must be sterile. Always use proper aseptic techniques and work inside a laminar flow hood. Consult our [Gibco Cell Culture Basics](#) for aseptic handling.

Test protocol

Buffered saline solutions have many different uses in diagnostic protocols such as cell washing, diluents or as transport media. Refer to the appropriate procedures. Several procedures and applications are provided in the references.

Cellular suspensions prepared in a buffered saline solution should not be stored longer than several hours as cellular viability may decrease. Therefore, the suitability of the buffered saline solution for a specific cell type should be tested prior to use.

Quality control

Standard evaluations for cell culture media are pH, osmolality, endotoxins and sterility testing for liquid products. All quality control testing results are reported on lot specific Certificate of Analysis available on our website: thermofisher.com.

References

1. Lennette EH, Halonen P and Murphy FA . Laboratory Diagnosis of Infectious Disease - Principles and Practices (1988).Springer, New York, p.43.
2. WHO Manual for the laboratory diagnosis and virological surveillance of influenza, 2011.
3. Winn, W. C., & Koneman, E. W. (2006). Koneman's color atlas and textbook of diagnostic microbiology (6th ed.). Philadelphia: Lippincott Williams & Wilkins.
4. WHO Guidelines on the Establishment of Virology Laboratories in Developing Countries, 2008.

Labeling symbols

The symbols present on the product label are explained in the following table.

	MANUFACTURER		USE BY
	IN VITRO DIAGNOSTIC MEDICAL DEVICE		CONSULT INSTRUCTIONS FOR USE
	CATALOG NUMBER		CAUTION, CONSULT ACCOMPANYING DOCUMENTS
	BATCH CODE		UPPER AND LOWER LIMITS OF TEMPERATURE
	Sterilized using aseptic processing technique		PROTECT FROM LIGHT
	European Mark of Conformity		AUTHORISED REPRESENTATIVE IN THE EUROPEAN COMMUNITY

Limited product warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale at www.thermofisher.com/us/en/home/global/terms-and-conditions.html. If you have any questions, please contact Life Technologies at www.thermofisher.com/support.

Manufacturer:
Life Technologies Corporation |
3175 Staley Road |
Grand Island, New York 14072 USA

Life Technologies Europe B.V.
Kwartsweg 2, 2665 NN Bleiswijk
The Netherlands

Manufacturer:
Life Technologies™ Ltd. |
3 Fountain Drive, Inchinnan Business Park |
Paisley PA49RF, Scotland, United Kingdom |
Tel: +44 (0)141 81416305

Manufacturer:
Dual manufactured products

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0018562

Revision	Date	Description
3.0	30 June 2021	Added SKU 15630106 and 15630130 to HEPES Buffer Solution 1 M
2.0	28 July 2020	Changed the superscript (2) for SKU 14025076 to a superscript (3) to indicate that the product is now dual manufactured in GI and Inchinnan.
1.0	12 November 2019	New document

Important Licensing Information: These products may be covered by one or more Limited Use Label Licenses. By use of these products, you accept the terms and conditions of all applicable Limited Use Label Licenses.

©2021 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

Accurate and sensitive somatic mutation detection powered by castPCR™ technology

TaqMan® Mutation Detection Assays

- High specificity—mutant allele detection is based on an allele-specific primer, while wild type background is suppressed by the proprietary MGB blocker oligonucleotide
- High sensitivity—assays can detect down to 0.1% mutation in a background of wild type DNA, as demonstrated in spiking experiments
- Wide dynamic range and excellent PCR efficiency—assays demonstrate at least 4 logs of dynamic range and an average PCR efficiency of 100% ± 10%
- Fast, simple workflow—like other TaqMan® Assays, typically requires 3 hours from sample to results, with minimum hands-on time

Cancer research samples often contain rare somatic mutations within a high background of normal wild type DNA. Many mutation detection methods compatible with tumor specimens, including gene sequencing and real-time PCR, have been reported in the literature and are commercially available. However, commercially available kits have various limitations in terms of sensitivity, specificity, cost, workflow, and turnaround time. We have developed sensitive and easy-to-use TaqMan® Mutation Detection Assays to accurately assess mutation status. TaqMan® Mutation Detection Assays were designed based on the novel competitive allele-specific TaqMan® PCR (castPCR™) technology, which combines allele-specific TaqMan® qPCR with allele-specific MGB blocker oligonucleotides that effectively suppress nonspecific amplification from the off-target allele.

Currently, the assay portfolio covers key somatic mutations identified in various cancer genes including, but not limited to, *KRAS*, *BRAF*, *HRAS*, *NRAS*, *EGFR*, *PIK3CA*, *KIT*, *PTEN*, and *TP53* genes, which have been implicated in many types of cancer. These mutations were selected from the comprehensive Sanger COSMIC database for somatic mutations. The target selection was based on frequency of occurrence and input from leading cancer researchers. We will continually add more mutation assays to cover additional cancer gene mutations. For the most updated list of available assays, refer to the TaqMan® Mutation Detection Assay index file at lifetechnologies.com/castpcr.

About the assays

TaqMan® Mutation Detection Assays contain mutant allele assays, which specifically detect one or more mutant alleles, and corresponding gene reference assays, which detect mutation-free regions of the genes in which the target mutations reside (Figure 1). The validated assay set additionally includes corresponding wild type allele assays (not described here; refer to the TaqMan® Mutation Detection Assay protocol for further information).

Two experiment types

Two types of experiments are required for mutation detection analysis:

1. Detection ΔC_t cutoff determination

A mutant allele assay and corresponding gene reference assay are run on three or more wild type gDNA samples that are from the same sample type as the test samples

[e.g., gDNA from FFPE tissue samples, Figure 2). ΔC_t values are calculated for each sample run with a mutant allele assay/gene reference assay pair. The average ΔC_t value for all samples is then calculated and is used to derive the detection ΔC_t cutoff value for the mutant allele assay.

2. Mutation detection

A test sample is run with one or more mutant allele assays and a corresponding gene reference assay (Figure 2). The ΔC_t value for the mutant allele assay/gene reference assay pair is calculated, and this value is compared to the previously determined detection ΔC_t cutoff value to determine the sample mutation status.

Optional use of internal positive control (IPC)

You can duplex the IPC reagents with any TaqMan® Mutation Detection Assay to distinguish true target negatives from PCR failure or inhibition (Figure 3).

Figure 1. TaqMan® Mutation Detection Assay types.

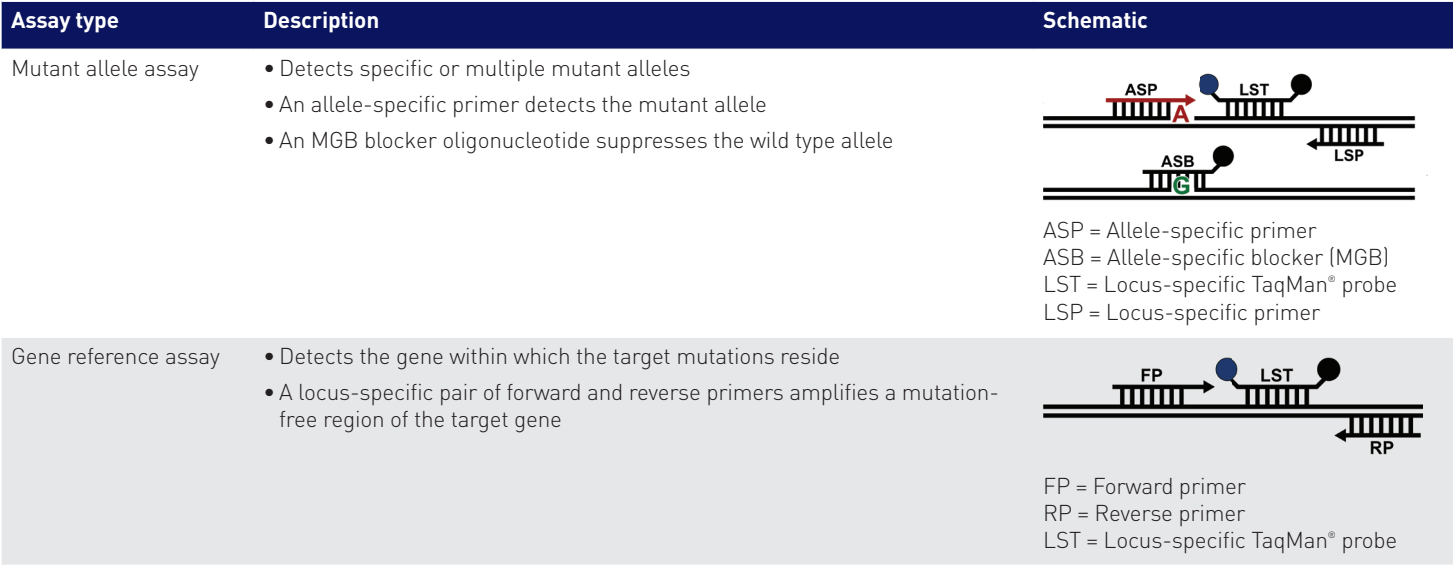
Figure 2. Gene reference and mutant allele assays are run with a genomic DNA sample to determine the mutation status of each target mutation within the cancer gene.


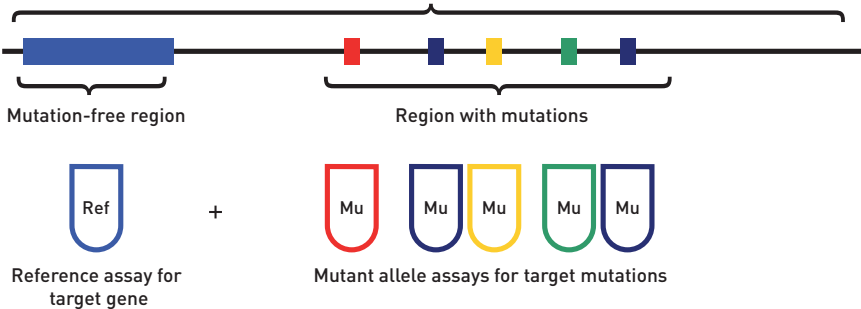
Figure 3. Internal positive controls. The TaqMan® Mutation Detection IPC Reagent Kit is a set of optional internal positive control reagents that can be duplexed with any TaqMan® Mutation Detection Assay to provide a positive PCR control result. The IPC reagents can distinguish a mutation target negative result from a PCR failure result.

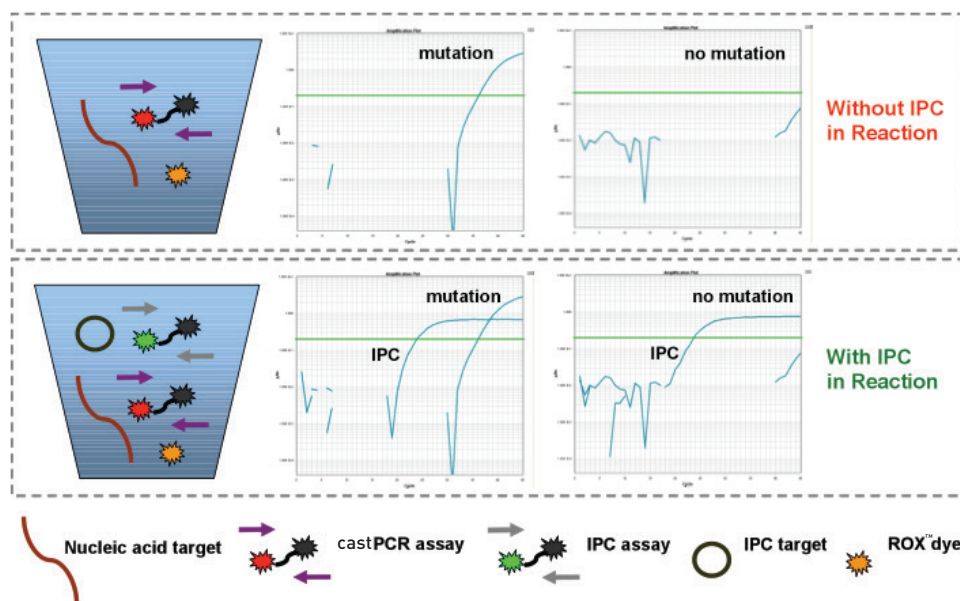Procedure

Purified gDNA, extracted from a sample with an unknown mutation status, is run with one or more mutant allele assays and the corresponding gene reference assay. For each real-time PCR reaction, the gDNA is combined with:

- **A TaqMan® Mutation Detection Assay**—contains two primers and a FAM™ dye–labeled MGB probe to detect a mutant allele or reference gene target. Mutant allele assays also contain an MGB oligonucleotide blocker.
- **TaqMan® Genotyping Master Mix**—contains AmpliTaq Gold® DNA Polymerase UP (Ultra Pure), dNTPs, and buffer
- **(Optional) TaqMan® Mutation Detection IPC Reagent Kit**—contains an internal positive control (IPC) template, two primers, and a VIC® dye–labeled TAMRA™ probe. It can be used to distinguish true target negatives from PCR failure or inhibition.

Reactions are run on a real-time PCR system, using a universal mutation detection thermal cycling protocol. After the run, the real-time PCR system's analysis software determines the C_t values for each TaqMan® Mutation Detection Assay and (optional) IPC reagent reactions. Real-time results export files can be opened in the free Mutation Detector™ Software for post-PCR data analysis. The C_t difference between each mutant allele assay and reference assay is calculated. This ΔC_t value, which represents the quantity of a specific mutant allele detected in a sample, is used to determine sample mutation status by comparison to a previously determined detection ΔC_t cutoff value. You can search for, or download a list of, currently available TaqMan® Mutation Detection Assays at lifetechnologies.com/castpcr.


Note: All TaqMan® Mutation Detection Assays have undergone extensive testing to ensure high sensitivity and specificity. The first set of released assays, covering 14 *KRAS*, 29 *EGFR*, and the *BRAF* V600E mutations, underwent additional testing, including determination of: the inherent amplification efficiency difference between mutant allele assays and corresponding reference assays, to enable quantitative analysis of percent mutation in a sample; and assay detection ΔC_t cutoff values using spiked cell line gDNA samples.

Assay performance

Specificity

Mutant allele detection is based on an allele-specific primer, while the wild type allele background is suppressed by the proprietary MGB blocker oligonucleotide. Assays can detect down to 0.1% mutant allele in the presence of a wild type allele background (Figure 4).

Figure 4. C_t difference between 0.1% mutation samples and wild type gDNA. For each assay, 0.1% mutant allele samples were obtained by spiking 10 copies of mutant allele synthetic templates into 10,000 copies of cell line wild type gDNA. **(A)** Example of amplification plot for KRAS_522_mu assay on 0.1% mutant allele sample and wild type gDNA. **(B)** There is a significant difference in amplification C_t values between the 0.1% mutant allele sample and wild type gDNA (*P* value < 0.05 for 46 out of 48 assays in the example graph).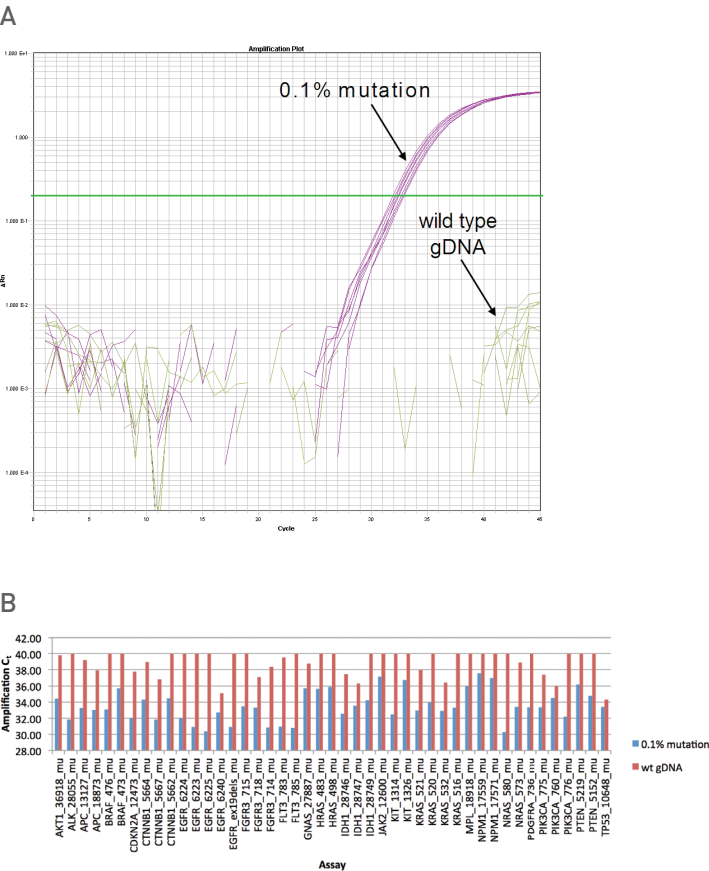

High sensitivity

TaqMan® Mutation Detection Assays can detect as few as 1–5 mutant copies in up to one million copies of wild type background. Assay sensitivity is demonstrated using synthetic template spiking experiments (Figure 5 and 6).

Wide dynamic range and excellent PCR efficiency

Assays demonstrate up to 7 logs of dynamic range and an average PCR efficiency of 100% ± 10% (Figure 6).

Figure 5. Assay sensitivity and selectivity. For every single assay, the sensitivity and selectivity were analyzed through synthetic template spiking experiments. 10 copies to 10⁵ copies of mutant allele synthetic template were spiked into a constant background of 10⁵ copies of wild type cell line genomic DNA. For a subset of the assays, 1 copy to 10⁶ copies of mutant allele synthetic template were spiked into a constant background of 10⁶ copies of wild type allele synthetic template. In the example shown, the BRAF_476_mu assay can detect 1 copy of mutant allele in a background of 10⁶ copies of wild type allele.


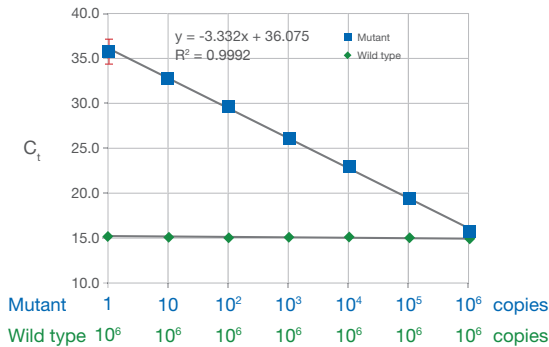
Figure 6. Assay dynamic range. Each assay was tested with 10⁵ copies to 10 copies of synthetic template within a constant background of 10⁵ copies of wild type genomic DNA. A subset of the assays was tested with 10⁷ copies to 1 copy of synthetic template within a constant background of 10⁷ copies of wild type allele synthetic template. In the example shown, the KRAS_532_mu assay has 7 logs of dynamic range, with an average PCR efficiency of 100% ± 10%.

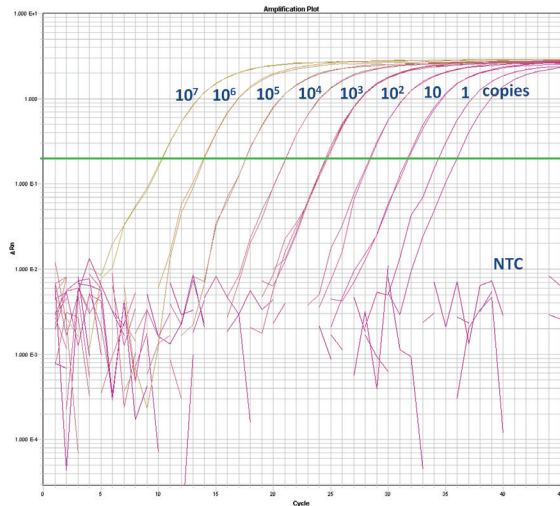Accuracy and reproducibility

Assays demonstrate excellent reproducibility and accurate quantification (Table 1).

Sample type compatibility

The assays can be used with gDNA samples extracted from FFPE tissues, fresh frozen tissues, and cell lines.

Data analysis software

For data analysis, Mutation Detector™ Software allows users to determine the mutation status and quantify the % mutation of their samples from TaqMan® Mutation Detection Assay data collected on the Applied Biosystems® ViiA™ 7, 7900HT, 7500, 7500 Fast, and StepOnePlus™ Real-Time PCR Systems (Table 2).

Table 1. Accuracy and reproducibility. Selected assays were tested in gDNA spiking experiments. In the example shown, G12C mutant cell line gDNA was spiked into wild type cell line gDNA at percentages ranging from 100% to 0.1%. The measured percent mutation was averaged from three experiment runs. The measured percent mutation is highly concordant with the expected percent mutation ($R^2 = 0.9997$). Accurate and precise quantification (CV < 20%) is obtained among the replicate runs when the target allele copy number is >30.

Copy number, target mutant allele	Expected (%)	Measured (%)	CV (%)
3,000	100.0	100.0	0.0
1,500	50.0	48.9	2.2
750	25.0	23.3	3.8
375	12.5	11.2	7.8
188	6.3	5.7	7.5
90	3.0	2.6	9.0
30	1.0	0.8	17.0
15	0.5	0.4	26.0
3	0.1	0.1	23.0

Table 2. Instrument compatibility.

Applied Biosystems® real-time PCR system	Block module	Software version
StepOnePlus™ system	Fast 96-Well Block Module	StepOne™ Software v2.X
7500 system	Standard 96-Well Block Module	SDS v1.X and v2.X
7500 Fast system	Fast 96-Well Block Module	SDS v1.X and v2.X
7900HT Fast system	Standard 96-Well Block Module, Fast 96-Well Block Module, 384-Well Block Module	SDS v2.X
ViiA™ 7 system	Standard 96-Well Block Module, Fast 96-Well Block Module, 384-Well Block Module	ViiA™ 7 Software v1.X
QuantStudio® 12K Flex system	Standard 96-Well Block Module, Fast 96-Well Block Module, 384-Well Block Module	QuantStudio® Software v1.0

Ordering information

Product	Quantity	Cat. No.
TaqMan® Mutation Detection Assays	150 µL, 10X	4465804
TaqMan® Mutation Detection Reference Assays	150 µL, 10X	4465807
TaqMan® EGFR Exon 19 Deletions Assay	150 µL, 10X	4465805
TaqMan® Mutation Detection IPC Reagent Kit	1 kit	4467538

For more information and full terms of the TaqMan® Assays QPCR Guarantee, go to lifetechnologies.com/taqmanguarantee

A Thermo Fisher Scientific Brand

appliedbiosystems

TaqMan Gene Expression Assay solutions

Proven performance for fast, reliable results

ThermoFisher
SCIENTIFIC

The leader in gene expression analysis

We are the leader in gene expression analysis, providing world-class sample preparation with Applied Biosystems™ technologies, real-time PCR using Applied Biosystems™ TaqMan™ or Applied Biosystems™ SYBR™ Green chemistry, and industry-leading real-time PCR instruments and data analysis software.

Applied Biosystems™ TaqMan™ assay technology is the gold standard in performance, quality, and content for gene expression analysis. Developed using long-standing bioinformatic expertise in primer and probe design, and stringent testing across applications and integrated platforms, TaqMan Assays provide you with the most reliable and robust real-time PCR solutions.

With over one and a half million predesigned and preoptimized assays across a growing list of model species, a wide range of formats to scale to your needs, and a robust manufacturing quality system, we have a complete suite of solutions that will enable you to get fast, reliable, and accurate gene expression results.

Contents

TaqMan Gene Expression Assays	4
The largest selection of predesigned assays	6
Proven performance	8
Flexible formats	12
Complementary reagents	16
Support at every step of your workflow	18

TaqMan Gene Expression Assays

Proven 5' nuclease-based real-time PCR chemistry

Get results you can trust

TaqMan Gene Expression Assays are referenced in tens of thousands of publications and are considered the gold standard for gene expression quantification by scientists around the world.

TaqMan Gene Expression Assays are based on 5' nuclease chemistry, and each assay contains the primer and probe set for your target of interest. Here's how an assay works (Figures 1 - 3):


1. At the start of the real-time PCR reaction, the temperature is raised to denature the double-stranded cDNA. During this step, the signal from the fluorescent dye on the 5' end of the Applied Biosystems™ TaqMan™ probe is quenched by the MGB–nonfluorescent quencher on the 3' end of the probe.

2. In the next step, the reaction temperature is lowered to allow the primers and probe to anneal to their specific target sequences.


3. Taq polymerase synthesizes a complementary DNA strand using the unlabeled primers and template. When the polymerase reaches the TaqMan probe, its endogenous 5' nuclease activity cleaves the probe, separating the dye from the quencher.

With each cycle of PCR, more dye molecules are released, resulting in an increase in fluorescence intensity proportional to the amount of amplicon synthesized.


1. Assay components and DNA template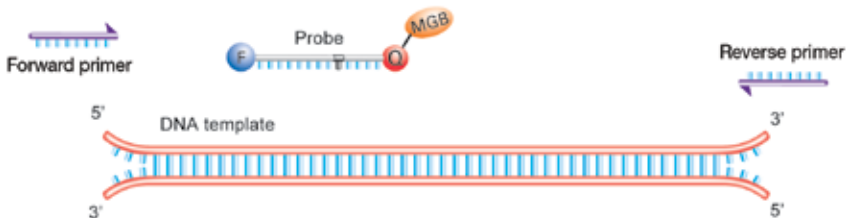

2. Denatured template and annealing assay components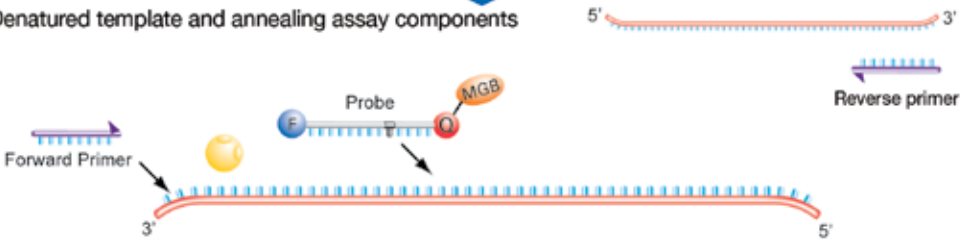

3. Polymerization and signal generation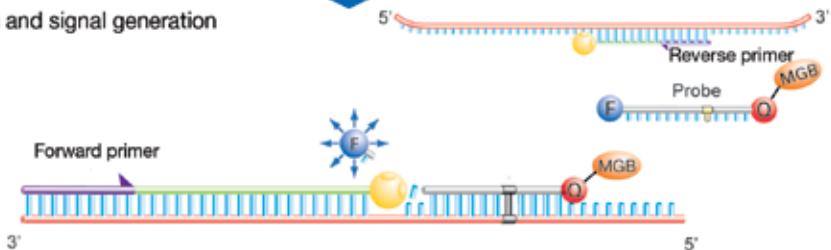

Legend


- F Applied Biosystems™ FAM™ or Vic™ dye
- Q Quencher
- MGB Minor groove binder
- Applied Biosystems™ AmpliTaq Gold™ DNA Polymerase
- Probe
- Primer
- Template

Prepare the cDNA sample

Prepare the total RNA

Perform reverse transcription


Evaluate the cDNA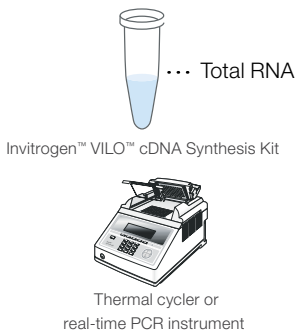

Prepare the reaction mix and load the plate

Prepare the PCR reaction mix

Load the plate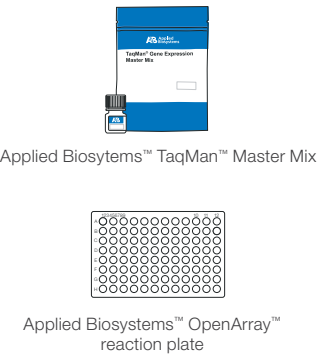

Run the real-time PCR reaction


Create the plate document/experiment

Run the plate

Analyze the results

Refer to the user guide for your real-time PCR instrument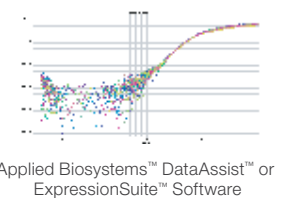

The largest selection of predesigned assays

Spend time on results, not assay design and optimization

With TaqMan predesigned assays, spend your time generating results, not designing and optimizing assays.

- Detect virtually any gene product—more than 1.5 million predesigned assays, and custom design for everything else
- Assays for nearly every human, mouse, and rat gene in the RefSeq database
- Available for 25 species, and some pathogens
- Assays for multiple locations per transcript and across nearly every exon junction in human
- Strain-neutral assays for mouse and rat

To learn more and order, go to thermofisher.com/taqmangex

- Not finding what you’re looking for in our predesigned assay collection? The Applied Biosystems™ Custom TaqMan™ Assay Design Tool lets you design and order a TaqMan Assay to detect any gene from any organism. Design and order your assays at thermofisher.com/cadt Custom TaqMan Assays are typically delivered in 5–12 business days.
- Also, try Applied Biosystems™ TaqMan™ Endogenous Controls—a collection of TaqMan Assays targeting commonly used control gene products for sample input normalization in real-time PCR.

Predesigned TaqMan Gene Expression Assays (as of November 2015)


Species	Number of assays	Gene coverage (%)*
Human (<i>H. sapiens</i>)	205,707	99.8%
Mouse (<i>M. musculus</i>)	176,510	99.5%
Chinese hamster (<i>C. griseus</i>)	154,743	88.2%
Rat (<i>R. norvegicus</i>)	146,589	89.2%
Cow (<i>B. taurus</i>)	103,562	99.6%
Rice (<i>O. sativa</i>)	99,822	95.6%
Arabidopsis (<i>A. thaliana</i>)	97,879	93.8%
Nematode (<i>C. elegans</i>)	92,687	95.1%
Rhesus monkey (<i>M. mulatta</i>)	69,310	55.8%
Zebrafish (<i>D. rerio</i>)	63,712	77.3%
Frog (<i>X. tropicalis</i>)	56,764	87.3%
Dog (<i>C. familiaris</i>)	55,558	64.3%
Chicken (<i>G. gallus</i>)	48,432	85.1%
Fruit fly (<i>D. melanogaster</i>)	41,607	94.0%
Sweet corn (<i>Z. Mays</i>)	38,493	59.5%
Cynomolgus monkey (<i>M. fascicularis</i>)	37,652	80.5%
Pig (<i>S. scrofa</i>)	16,247	90.3%
Fission yeast (<i>S. pombe</i>)	6,538	94.3%
Rabbit (<i>O. cuniculus</i>)	5,927	80.9%
Baker's yeast (<i>S. cerevisiae</i>)	5,524	93.4%
Horse (<i>E. caballus</i>)	3,891	72.8%
Soybean (<i>G. max</i>)	3,456	13.5%
Guinea pig (<i>C. porcellus</i>)	2,037	64.3%
Grape (<i>V. vinifera</i>)	965	25.3%
Wheat (<i>T. aestivum</i>)	760	43.6%
Summary	1,534,372	81.1%, 25 species

*Percent coverage refers to genes in the RefSeq database.

There are multiple assays for my gene product. How do I choose the right one?

Genomic alignment maps on our website make it easy to see exactly what gene products are detected and how they align to the genomic locus. The top of the map shows the target gene. Below it, all TaqMan Gene Expression Assays for target gene products are shown relative to the genomic locus map. The known transcripts from the locus are shown below, with their RefSeq accession numbers.

- A. Gene symbol
- B. Alignment of TaqMan amplicons to the gene. Hover over an assay to see its name and assay number as well as the transcripts it detects. Click on an assay to open an assay details pane for more information and to add the assay to your shopping cart.
- C. Assays providing the best coverage are marked with a star symbol.
- D. Narrow your results by specifying the type of assay you need.
- E. All RefSeq transcripts that map to the gene locus, showing exon usage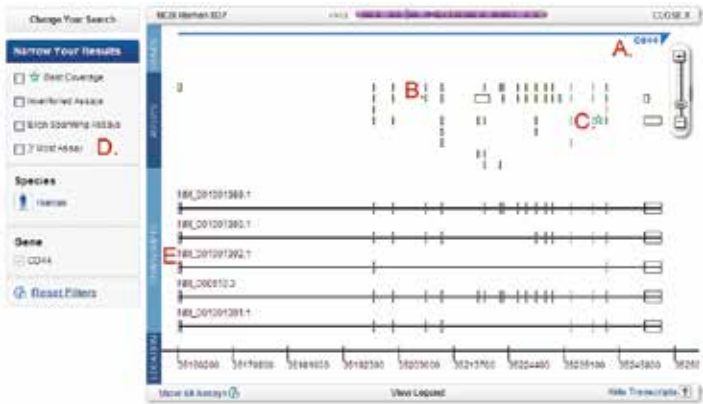

The TaqMan Assays qPCR guarantee

We stand behind every predesigned TaqMan Assay. We are committed to helping you achieve your research goals and believe our predesigned TaqMan primer and probe sets establish the benchmark for high-quality and easy-to-use real-time PCR products.

We want you to be happy with your purchase and confident in the genomic tools we provide. Therefore, we guarantee every TaqMan Assay in terms of:

- **Quality**—high-quality manufacturing for reproducible results from lot to lot
- **Performance**—superior sensitivity, specificity, and accuracy
- **Content**—the largest collection of primer and probe sets using the world’s best and most extensively validated assay design pipeline
- **Results**—enables you to obtain data you can trust

If you are not satisfied with the performance of a predesigned TaqMan Assay, we’ll replace it at no cost or credit your account. For more information, and to see the full terms and conditions of the guarantee, go to thermofisher.com/taqmanguarantee

Proven performance

Reliable reagents for confidence in your results

TaqMan MGB probes bind more tightly—shorter, more specific probes

TaqMan probes include an MGB moiety at the 3' end that increases the T_m of the probe and stabilizes probe–target hybrids. This means that TaqMan probes can be significantly shorter than traditional probes, providing better sequence discrimination and flexibility to accommodate more targets.


Nonfluorescent quencher (NFQ) maximizes sensitivity

TaqMan probes incorporate an NFQ to absorb (quench) signal from the fluorescent FAM or VIC dye label at the other end of the probe. The properties of the NFQ combined with the short length of MGB probes result in lower background signal than with non-MGB/NFQ probes. Lower background noise results in increased sensitivity and precision in your data.

TaqMan probe outperforms non-MGB probe in real-time PCR

Input	C_t		Standard deviation	
	TaqMan Assay	Non-MGB assay	TaqMan Assay	Non-MGB assay
10 ng	9.72	13.35	0.02	0.15
1 ng	13.36	16.82	0.04	0.18
0.1 ng	16.76	20.23	0.07	0.13
10^{-2} ng	20.19	23.72	0.04	0.13
10^{-3} ng	23.64	27.31	0.03	0.10
10^{-4} ng	27.01	30.66	0.04	0.12
10^{-5} ng	30.24	32.82	0.13	0.19

Figure 2. TaqMan probes provide better sensitivity and precision. Comparison of two 5' nuclease PCR assays for 18S rRNA. Ten-fold dilutions of Universal Human Reference RNA (10 – 10^{-5} ng) were prepared and analyzed in 11 replicate real-time PCR reactions using either the TaqMan Gene Expression Assay (FAM dye–labeled, with NFQ) or the non-MGB assay (FAM dye–labeled, with BHQ). Real-time PCR was run according to the respective manufacturers' recommended conditions. Across a 6-log range of input template, the TaqMan Assay displayed earlier C_t values and better reproducibility across all data points. In addition, the TaqMan Assay had higher signal and lower background, resulting in better sensitivity and higher precision.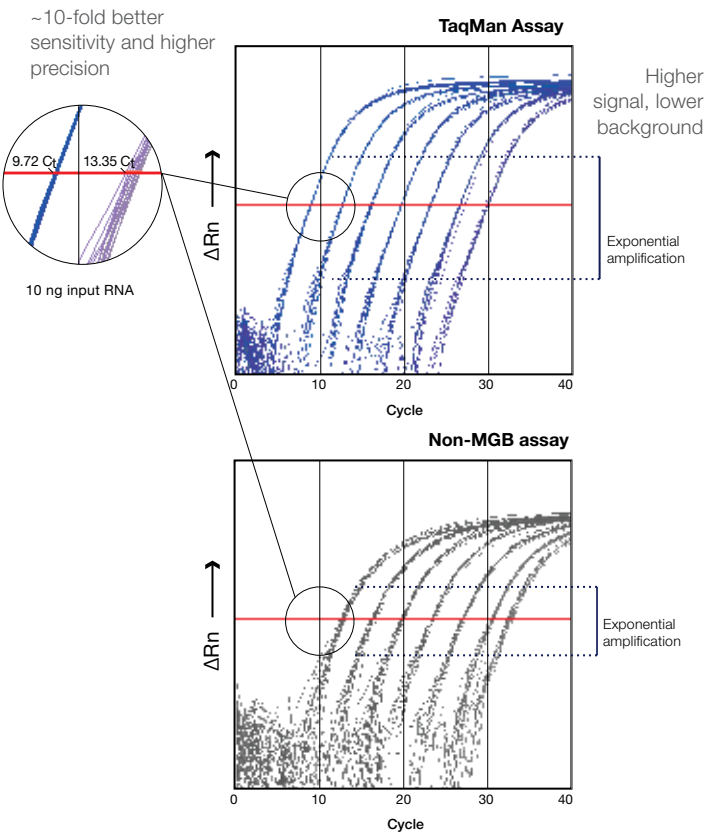

- **Specificity:** Advanced primer/probe sequence selection criteria plus MGB probe enhancement deliver the specificity and reproducibility you need for confidence in your results. Your results are generated from amplification of the intended target, not from nonspecific dye binding or amplification of closely related genes or pseudogenes.
- **Sensitivity:** The NFQ on TaqMan probes minimizes background, and intelligent PCR primer and probe design maximizes amplification efficiency. Get better sensitivity and accuracy—reliably detect targets present at 10 or fewer copies.
- **Reproducibility:** Accurately reproduce results from well to well, day to day, and lab to lab—even across manufacturing lots.
- **Wide dynamic range:** Detect from a handful to millions of target molecules with the same reaction setup. Capture the full spectrum of expression variability in virtually any experimental scenario.
- **High amplification efficiency:** All TaqMan Gene Expression Assays have a PCR efficiency of 100% ($\pm 10\%$). Use the comparative C_t ($\Delta\Delta C_t$) method of quantification confidently.
- **Ease of use:** All assays use a single, universal thermal cycling profile. Run any assay combination on a single plate. Avoid instrument-programming errors.
- **Comprehensive assay information:** Genomic mapping data are provided prior to purchase.

Detect as few as 10 target molecules with high sensitivity and large dynamic range

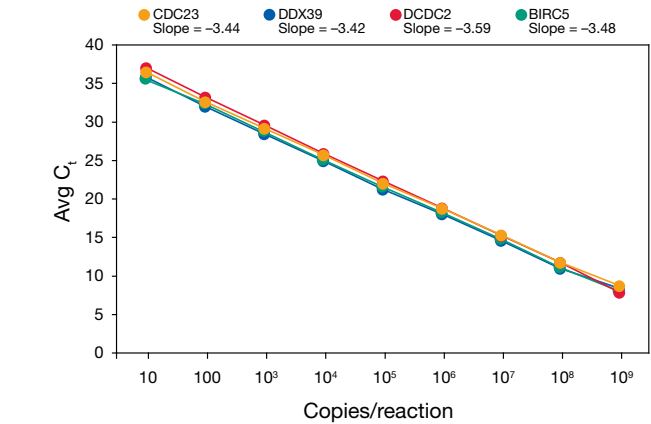
Figure 3. Sensitivity and wide dynamic range. Sequential 10-fold dilutions of synthetic sense RNA corresponding to 4 gene products—CDC23, DDX39, DCDC2, and BIRC5—were added to a background of yeast RNA to evaluate the sensitivity and dynamic range of TaqMan Gene Expression Assays. Samples containing 50 to 5×10^9 target molecules were reverse transcribed, and 20% of each RT reaction was used in quadruplicate PCR reactions using TaqMan Gene Expression Master Mix. Reactions containing as few as 10 copies were detected ($C_t \sim 35$).

Reproducible quantification with virtually 100% amplification efficiency

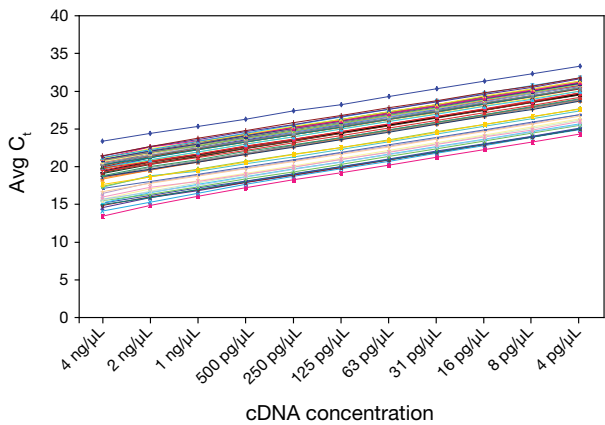
Figure 4. Reliable performance and wide dynamic range. TaqMan Gene Expression Assays were used to analyze expression of 60 targets across a 2-fold dilution series of universal reference cDNA, from 4 ng/ μ L to 4 pg/ μ L. The average slope of the lines is 1.02. TaqMan Assays exhibit virtually 100% amplification efficiency at each cycle of PCR: each target molecule is copied, doubling the fluorescence signal.

Specificity for your mRNA target

TaqMan Assay design helps ensure target mRNA specificity: readily distinguish even highly homologous sequences

Specificity is built into the TaqMan Assay design pipeline. As a result, assays detect only their intended targets. Even TaqMan Gene Expression Assays for members of highly homologous gene families typically amplify their targets with C_t values at least 10 cycles earlier than the closest homolog, or with at least 1,000-fold discrimination if equal numbers of the two targets are present.

TaqMan Gene Expression Assays are designed to detect only their intended targets, easily discriminating among highly homologous sequences.

HOX gene family members HOXA10, HOXC10, and HOXD10 share ~80% sequence homology

HOXA10	AATTGGCTGACAGCAAAGAGCGGAAGGAAGAAGAGGTGCCCTATACTAAACACCAGACGCTGGAATTGGAGAAAGAATTTCTGTTCAATATGTATTTGACGCGAGAGCGCCGCTGG		
HOXC10	T T G A A A T T A G G C G T T C A A T A T G T A T T G A C G C G A G A G C C C G C T G G		
HOXD10	T T C A A A G T T A G A T A A A C T C C C C C		
Gene	RefSeq ID	TaqMan Assay ID	Homology
HOXA10	NM_018951.3	Hs00172012_m1	—
HOXC10	NM_017409.3	Hs00213579_m1	81%
HOXD10	NM_002148.3	Hs00157974_m1	79%

Clear gene expression results for HOX gene family members

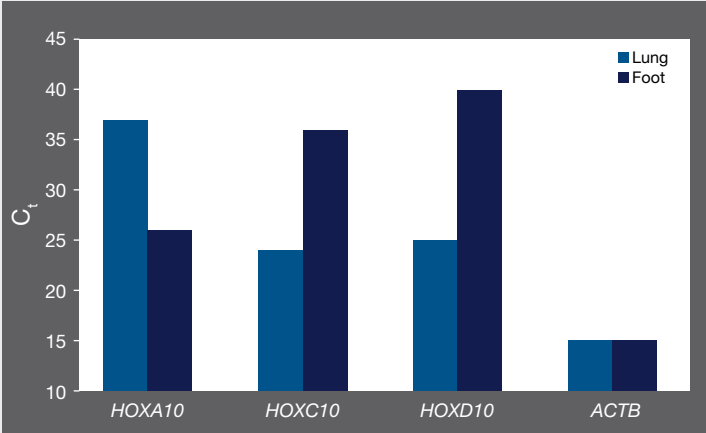
Figure 5. TaqMan Gene Expression Assays detect only their intended targets, even among the highly homologous HOX gene family members. In vertebrates, as in *Drosophila*, location-appropriate expression of members of the HOX gene family is essential for normal embryogenesis. Tissue-specific expression of 3 closely related HOX genes, comparable to published data, was easily detected using TaqMan Gene Expression Assays.

Advanced bioinformatics

TaqMan Gene Expression Assays are designed using our sophisticated design pipeline that has been stringently validated by functionally testing more than 18,000 assays (a statistically significant subset). Since then, our customers have consistently confirmed through their own validation experiments that TaqMan Gene Expression Assays enable reliable, reproducible results.

This process is used to design all TaqMan Gene Expression Assays, including inventoried assays, made-to-order assays, and Applied Biosystems™ Custom Plus assays. We offer ~73,000 inventoried assays and over 1.5 million made-to-order assays, which are manufactured when an order is placed. Applied Biosystems™ Custom Plus TaqMan™ RNA Assays are ideal for newly identified genes and specific splice variants, and offer the same performance as predesigned TaqMan Assays.

TaqMan Assay design and manufacture


- Target selection**
mRNA sequences (NCBI)

Preprocessing
–Map to genome
–Mask SNPs, repeats, and discrepancies
–Identify exon–exon junction

Assay design
Thermodynamic and chemistry parameters
–Balance T_m for universal thermal cycling
–Avoid secondary structure, optimize GC content
–Optimize amplicon size
–Eliminate primer-dimer formation

In silico QC
–Score assays for target specificity
–Score assays for genome specificity

Assay selection
High-quality TaqMan Gene Expression Assays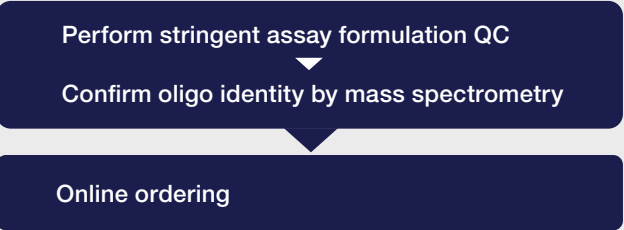

Flexible formats

A variety of formats for different research needs


Configurations to fit your research goals

Are you analyzing hundreds (or thousands) of samples, and expression from a handful of genes? Or does your research involve a few samples that need to be analyzed for a long list of mRNA targets? No matter what experiment you are performing, there is a TaqMan Gene Expression Assay format and real-time PCR instrument for your research needs.

Single tubes

- Low entry price
- Flexible
- Run on any real-time PCR instrument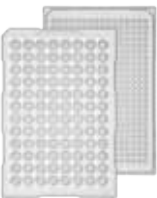

96- or 384-well plates

- Optimal for small to medium projects
- Balances flexibility with streamlined reaction setup
- Run on any 96- or 384-well real-time PCR instrument

384-well microfluidic cards

- Low cost per reaction
- Optimal for medium to large projects
- Run on Applied Biosystems™ QuantStudio™ 7 & 12K Flex, ViiA™ 7, and 7900HT Real-Time PCR Systems

OpenArray plates

- Lowest cost for large projects
- Ultimate throughput
- Run on QuantStudio 12K Flex Real-Time PCR System

TaqMan Gene Expression Assays (single tubes)

Predesigned assays come in four different sizes so that you can order only the number of assays appropriate for your research. In addition, for made-to-order assays in small, medium, and large sizes, you can choose FAM or VIC dye labeling, and non-primer-limited or primer-limited formulation. (Extra small assays are only available with FAM dye labels.)

For more information, go to thermofisher.com/allgenes

Size	No. of reactions*	Concentration	Reporter dye	Cat. No.
Extra small (inventoried)†	75	20X	FAM	4453320
Extra small (made-to-order)‡	75	20X	FAM	4448892
Small (inventoried)†	250	20X	FAM	4331182
Small (made-to-order)‡	360	20X	FAM or VIC	4351372, 4448489 (VIC) 4448484 (VIC-PL**)
Medium (made-to-order)‡	750	20X	FAM or VIC	4351370, 4448490 (VIC) 4448485 (VIC-PL**)
Large (made-to-order)‡	2,900	60X	FAM or VIC	4351368, 4448491 (VIC) 4448486 (VIC-PL**)

* Reaction number is based on 20 µL reaction size.
** Primer-limited.
† Inventoried assays are typically delivered in 1–4 business days.
‡ Made-to-order assays are typically delivered in 5–12 business days.

Applied Biosystems™ TaqMan™ Arrays: 96-well plates or 384-well microfluidic cards

- Configure a Custom TaqMan Array containing inventoried predesigned assays, or select from our gene signature assay collections
- TaqMan Gene Expression Assays are loaded into one of two TaqMan Array formats: 96-well plates (Fast or standard) or 384-well microfluidic cards

(To include made-to-order or custom assays on your plate or card, order using our Applied Biosystems™ TaqMan™ Custom Plating Service, or contact your sales representative for other options.)

Custom TaqMan Array 96-well plates

- Choose any inventoried TaqMan Gene Expression Assay
- 6-plate minimum order
- Choose standard (20 µL rxn) or Fast (10 µL rxn) format

Typically delivered in 4–14 business days

To learn more and order, go to thermofisher.com/arrayplates

Assays + controls	Assay replicates	Samples per plate	Name	Cat. No. (standard)	Cat. No. (Fast)
95 + 1*	1	1	Format 96	4391524	4413255
92 + 4**	1	1	Format 96 +	4391525	4413256
47 + 1*	2	1–2	Format 48	4391526	4413257
44 + 4**	2	1–2	Format 48 +	4391527	4413258
31 + 1*	3	1–3	Format 32	4391528	4413259
28 + 4**	3	1–3	Format 32 +	4391529	4413260
15 + 1	6	1–6	Format 16	4413264	4413261
12 + 4	6	1–6	Format 16 +	4413265	4413262
7 + 1	12	1–12	Format 8	4413266	4413263

* Available with one manufacturing control assay for 18S ribosomal RNA. These formats are required for plates with assays for rhesus, canine, or a mixture of species.
** Includes the manufacturing control assay for 18S ribosomal RNA, plus assays for 3 additional candidate endogenous control genes: *GAPDH*, *HPRT1*, and *GUSB*, appropriate for human, mouse, or rat sample analysis.

Custom TaqMan Array 384-well microfluidic cards

- Choose any inventoried TaqMan Gene Expression Assays
- 10-card minimum order
- Run on the QuantStudio 7 & 12K Flex, ViiA 7, and 7900HT Fast Real-Time PCR Systems
- No robotics required: cards have 8 sample-loading ports, each connected to 48 wells containing dried-down TaqMan Assays
- 1 µL reactions (2 µL including channel filling and overage)
- Typically delivered in 3–4 weeks

To learn more and order, go to thermofisher.com/arraycards

Assays + controls*	Assay replicates	Samples per card	Name	Cat. No.
11 + 1	4	8	Format 12	4342247
15 + 1	3	8	Format 16	4346798
23 + 1	2 (or 4)	8 (or 4)	Format 24	4342249
31 + 1	3	4	Format 32	4346799
47 + 1	1 (or 2)	8 (or 4)	Format 48	4342253
63 + 1	3	2	Format 64	4346800
95 + 1	1 (or 2)	4 (or 2)	Format 96a	4342259
95 + 1	2 (or 4)	2 (or 1)	Format 96b	4342261
191 + 1	2	1	Format 192	4346802
380 + 4	1	1	Format 384	4342265

* These arrays are available with one manufacturing control assay for 18S ribosomal RNA.

Applied Biosystems™ TaqMan™ Array Gene Signature Plates and Cards

- Predesigned, preloaded TaqMan Assays for gene products specific to pathways, biomarkers, or disease target classes to facilitate drug discovery and disease research
- Endogenous control panels are also available to identify the best housekeeping gene products for your research
- Gene signature plates are typically delivered in 5–10 business days, and gene signature cards in 1–4 business days

Here is a sampling of what’s available:

- Apoptosis
- Endogenous controls
- Cancer
- Immune system and inflammation
- Cell cycle proliferation and regulation
- Neurology
- Development and stem cells
- Signal transduction
- ECM matrix and adhesion
- Toxicology and drug metabolism


To see the complete collection of 96-well gene signature plates, go to thermofisher.com/signatureplates
To see the collection of 384-well gene signature microfluidic cards, go to thermofisher.com/signaturecards

OpenArray Real-Time PCR Plates

- TaqMan Assays loaded and dried down into the 3,072 through-holes on OpenArray Real-Time PCR Plates
- Process up to 576 samples to obtain over 43,000 data points, with a single operator in an 8-hour day, without the use of robotics
- For use with the QuantStudio 12K Flex Real-Time System with an Applied Biosystems™ OpenArray™ block configuration and supporting reagent kits only
- OpenArray plates with inventoried assays are typically delivered in 4–5 weeks, and within 5–6 weeks for custom assays

To learn more about OpenArray technology on the QuantStudio 12K Flex system,go to thermofisher.com/openarray

Assays + controls	Assay replicates	Samples per plate	Name	Cat. No.
18	3	Up to 48	Format 18	4471124
56	1	Up to 48	Format 56	4471125
112	1	Up to 24	Format 112	4471126
168	1	Up to 16	Format 168	4471127
224	1	Up to 12	Format 224	4471128

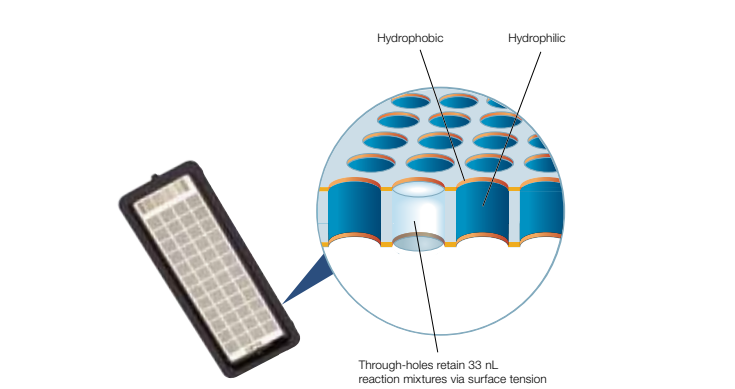
TaqMan Custom Plating Service: 96- or 384-well plates

- Configure 96- or 384-well plates with any TaqMan Gene Expression Assays, including custom assays designed to your target sequences and made-to-order assays.
- Set up custom configurations of any TaqMan Assays, including inventoried, made-to-order, custom, or Custom Plus gene expression assays or custom TaqMan probes and primers
 - Choose 96- or 384-well plate, and Fast or standard format
 - Receive in dried-down or liquid formulation
 - Typically delivered in 2–5 weeks

Complementary reagents

Everything you need for reliable results

We provide everything you need for real-time PCR analysis, starting with isolating RNA from virtually any sample type, to reverse transcription into cDNA, optional preamplification to stretch small samples for analysis of many gene products, and of course, real-time PCR data analysis.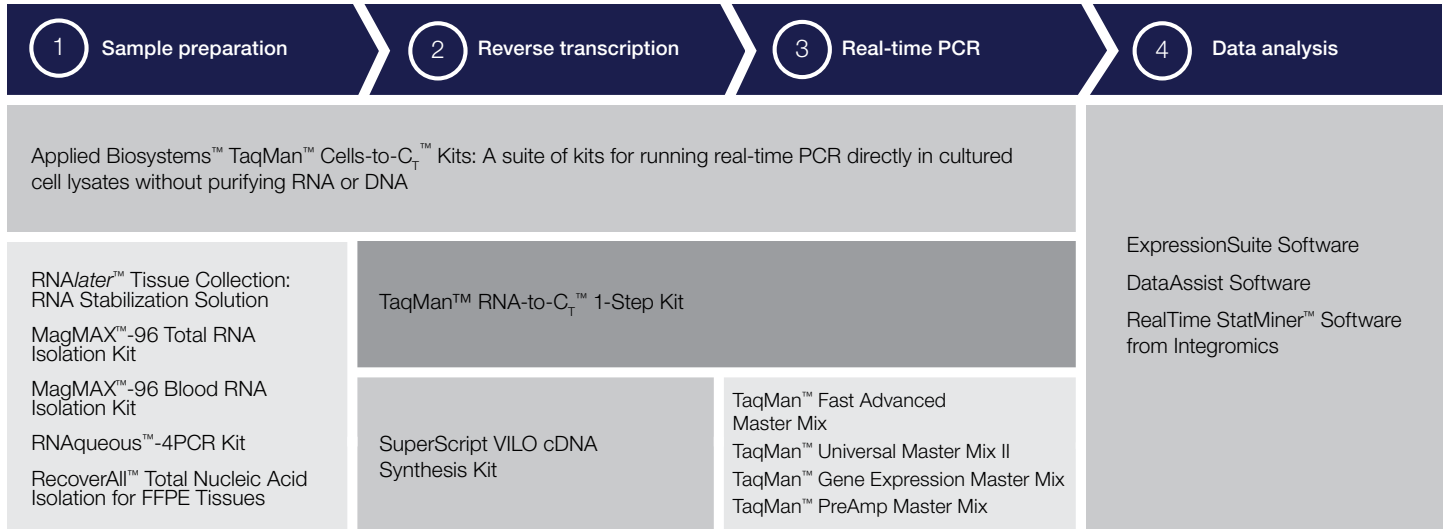

Support at every step of your workflow

Consistent reliability from manufacturing to follow-up

Quality manufacturing and stringent quality control

TaqMan Assays are manufactured in-house under rigorous quality processes at our ISO 13485–certified manufacturing facilities, and are never outsourced.

Comprehensive worldwide support

Whether you need help finding a TaqMan Assay for your target, deciding which format best suits your needs, placing your order through our online ordering system, or setting up your reactions, our global sales and technical support teams are here to help.

Technical support

If you have questions about how to use TaqMan Assays or how to analyze results, call or email our technical support specialists. These scientists are skilled in experimental planning and design, are expert troubleshooters, and are familiar with a wide variety of applications that use TaqMan Assays.

Rapid delivery

We continually strive to minimize delivery time on TaqMan Assay products. To that end, we have implemented streamlined order processing systems that interface with our new manufacturing facilities to help reduce delivery times.

Everything you need to meet the MIQE guidelines for peer-reviewed publications

The Minimum Information for Publication of Quantitative Real-Time PCR Experiments (MIQE) guidelines, published by Bustin et al. in *Clinical Chemistry* (April 2009), are meant to ensure that real-time PCR experiments are meaningful, accurate, and reproducible. We support this initiative and commend the MIQE scientists for their leadership.

We provide the following for easier adherence to these guidelines:

- **TaqMan Assay annotation**—Information requested under the real-time PCR target, oligonucleotide, and protocol sections of the guidelines is provided in your assay shipment and on our website. All biologically relevant information is available, including assay location, transcripts detected, and amplicon size. Protocols with recommended reagents and reaction conditions are also available on our website.

- **Publications**—There are >9,900 peer-reviewed publications that cite TaqMan Assays, so including the TaqMan Assay ID in lieu of sequences is sufficient and widely accepted.
- **Instrument software**—Applied Biosystems™ instrument software reports C_t values for quantification. The C_t can be used to generate standard curves, determine slope, and derive R² values. To help adhere to the MIQE guidelines, the term quantification cycle (C_q) may be used directly in place of C_t.
- **Data analysis**—We offer data analysis software, including ExpressionSuite and DataAssist Software; simple-to-use tools for calculating relative gene expression using statistical analysis and visualization; and RealTime StatMiner Software (Integromics) for additional statistical analysis workflows.

applied
biosystems

Find out more at **[thermofisher/allgenes](https://thermofisher.com/allgenes)**

ThermoFisher
S C I E N T I F I C

For Research Use Only. Not for use in diagnostic procedures. TaqMan and AmpliTaq Gold are registered trademarks of Roche Molecular Systems, Inc., used under permission and license. RealTime StatMiner is a trademark of Integromics, S.L. **CO020031 1215**

Real-time PCR master mixes and instrument compatibility

Applications	TaqMan® and TaqPath real-time PCR master mixes								One-step real-time RT-PCR master mixes					
	Gene expression/ small RNA	Genotyping/ CNV	Gene expression/ genotyping	Gene expression/ genotyping	Gene expression	Genotyping/ SNP	Genotyping/ SNP	Gene expression	Gene expression/ pathogen detection	Gene expression/ pathogen detection	Gene expression/ pathogen detection	Gene expression for 1-step and multiplex	Gene expression/ pathogen detection	Gene expression/ pathogen detection
	TaqMan Fast Advanced Master Mix	TaqPath ProAmp Master Mix	TaqMan Universal PCR Master Mix	TaqMan Universal Master Mix II	TaqMan Gene Expression Master Mix	TaqMan Genotyping Master Mix	TaqMan GTXpress Master Mix	TaqMan Fast Universal PCR Master Mix	TaqMan RNA-to-C _T 1-Step Kit	Power SYBR Green RNA-to-C _T 1-Step Kit	TaqMan Fast Virus 1-Step Master Mix	TaqPath 1-Step Multiplex Master Mix	SuperScript III Platinum One-Step qRT-PCR Kit	SuperScript III Platinum One-Step qRT-PCR Kit w/ ROX
	4444556 4444557 4444963 4444964 4444965 4444558	A32704, A30865, A30866, A30871, A30867, A30872, A32705, A30868, A30869, A30873, A30870, A30874	4304437, 4305719, 4318157, 4326708, 4364338, 4364340, 4324018, 4324020, 4326614, 4364341, 4364343	4440043, 4440040, 4440047, 4440048, 4440049, 4440041, 4440042, 4440038, 4440044, 4440045, 4440046, 4440039	4370048 4369016 4369514 4369510 4369542 4370074	4371353 4371355 4381656 4371357 4381657	4403311 4401892 4401890 4401857	4352042 4364103 4366072 4366073 4367846	4392653 4392938 4392656	4391178 4389986	4444432 4444434 4444436	A28521, A28522, A28523, A28525, A28526, A28527	11732020 11732088	11745500 11745100
Instruments														
Analytik Jena qTOWER	√	√	√	√	√	√	√	√	√	√	√	√	√	
Applied Biosystems 7500	√	√	√	√	√	√	√	√	√	√	√	√	√	
Applied Biosystems 7500 Fast	√	√	√	√	√	√	√	√	√	√	√	√	√	
Applied Biosystems 7300	√	√	√	√	√	√	√	√	√	√	√	√*	√	√
Applied Biosystems 7900HT	√	√	√	√	√	√	√	√	√	√	√	√*	√	√
Applied Biosystems QuantStudio 12K Flex	√	√	√	√	√	√	√	√	√	√	√	√	√	
Applied Biosystems QuantStudio 6	√	√	√	√	√	√	√	√	√	√	√	√	√	
Applied Biosystems QuantStudio 3/5	√	√*	√	√	√	√	√	√	√	√	√	√*	√	
Applied Biosystems QuantStudio 7	√	√	√	√	√	√	√	√	√	√	√	√	√	
Applied Biosystems StepOne	√	*	√	√	√	√	√	√	√	√	√	*	√	√
Applied Biosystems StepOnePlus	√	*	√	√	√	√	√	√	√	√	√	*	√	√
Applied Biosystems ViiA 7	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad CFX384	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad CFX96	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad IQ5	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad MiniOpticon	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad/MJ Chromo4	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad/MJ Opticon	√	√	√	√	√	√	√	√	√	√	√	√	√	
Bio-Rad/MJ Opticon 2	√	√	√	√	√	√	√	√	√	√	√	√	√	
Eppendorf Mastercycler ep realplex	√	√	√	√	√	√	√	√	√	√	√	√	√	
Qiagen/Corbett Rotor-Gene 6000	√	√	√	√	√	√	√	√	√	√	√	√	√	
Qiagen/Corbett Rotor-Gene Q	√	√	√	√	√	√	√	√	√	√	√	√	√	
Qiagen/Corbett Rotor-Gene 3000	√	√	√	√	√	√	√	√	√	√	√	√	√	
Roche LightCycler 480	√	√	√	√	√	√	√	√	√	√	√	√	√	
Roche LightCycler Nano	◊	◊	◊	◊	◊	◊	◊	◊	◊	◊	◊	◊	◊	
Agilent/Stratagene MX3000P	√	√	√	√	√	√	√	√	√	√	√	√	√	
Agilent/Stratagene MX3005P	√	√	√	√	√	√	√	√	√	√	√	√	√	
Agilent/Stratagene MX4000	√	√	√	√	√	√	√	√	√	√	√	√	√	
TaKaRa T800	◊	◊	◊	◊	◊	◊	◊	◊	◊	√	◊	◊	◊	

√ – Indicates preferred kit and tested on this instrument.

◊ – Indicates kit should work, but has not been tested on this instrument.

* – Not compatible with multiplexing. The QuantStudio 3 system can be used for limited multiplexing; it is not compatible with Mustang Purple or Cy®5 dyes. The QuantStudio 5 system is compatible with multiplexing.

Real-time PCR master mixes and instrument compatibility

SYBR Green real-time PCR master mixes						
Applications	Gene expression	Gene expression	Gene expression	Gene expression	Gene expression	Gene expression
	SYBR Select Master Mix	Fast SYBR Green Master Mix	Power SYBR Green PCR Master Mix	SYBR Green PCR Master Mix	SYBR GreenER qPCR SuperMix Universal	PowerUp SYBR Green Master Mix
Instruments	4472903, 4472908, 4472913, 4472918, 4472919, 4472920	4385610, 4385612, 4385616, 4385617, 4385618, 4385614	4367659, 4367660, 4368577, 4368702, 4368706, 4368708	4344463, 4309155, 4364344, 4364346, 4334973, 4312704	1176202K, 11762100, 11762500	A25742, A25743, A25776, A25777, A25778, A25779, A25780, A25918
Applied Biosystems 7500	✓	✓	✓	✓	✓	✓
Applied Biosystems 7500 Fast	✓	✓	✓	✓	✓	✓
Applied Biosystems 7300		✓	✓	✓	✓	✓
Applied Biosystems 7900HT		✓	✓	✓	✓	✓
Applied Biosystems QuantStudio 12K Flex	✓	✓	✓	✓	✓	✓
Applied Biosystems QuantStudio 6	✓	✓	✓	✓	✓	✓
Applied Biosystems QuantStudio 3/5	✓	✓	✓	✓	✓	✓
Applied Biosystems QuantStudio 7	✓	✓	✓	✓	✓	✓
Applied Biosystems StepOne	✓*	✓	✓	✓	✓	✓
Applied Biosystems StepOnePlus	✓*	✓	✓	✓	✓	✓
Applied Biosystems ViiA 7	✓	✓	✓	✓	✓	✓
Bio-Rad CFX384	✓	✓	✓	◇	◇	✓
Bio-Rad CFX96	✓	✓	✓	◇	◇	✓
Bio-Rad iQ5	✓	✓	✓	✓		✓
Bio-Rad MiniOpticon	✓	✓	✓	✓	✓	✓
Bio-Rad/MJ Chromo4	✓	✓	✓	✓	✓	✓
Bio-Rad/MJ Opticon	✓	✓	✓	✓	✓	✓
Bio-Rad/MJ Opticon 2	✓	✓	✓	✓	✓	✓
Eppendorf Mastercycler ep realplex		✓	✓	✓		✓
Qiagen/Corbett Rotor-Gene 6000	✓*	✓	✓	✓	✓	✓
Qiagen/Corbett Rotor-Gene Q	✓*	✓	✓	✓	◇	✓
Qiagen/Corbett Rotor-Gene 3000	✓*	✓	✓	✓	◇	✓
Roche LightCycler 480	✓	✓	✓	✓		✓
Roche LightCycler Nano		◇	◇	◇		◇
Agilent/Stratagene MX3000P	✓	✓	✓	✓	✓	✓
Agilent/Stratagene MX3005P	✓	✓	✓	✓	✓	✓
Agilent/Stratagene MX4000	✓	✓	✓	✓	✓	✓
TaKaRa T800		✓	✓	✓		✓

✓ – Indicates preferred kit and tested on this instrument.

✓* – Ensure primer concentration guide is followed (<200 nM for standard mode and 300–400 nM for fast mode). For StepOne/StepOnePlus System, install software version 2.3 or higher update.

◇ – Indicates kit should work, but has not been tested, on this instrument.

Find out more at thermofisher.com/mastermixes

ThermoFisher
SCIENTIFIC

For Research Use Only. Not for use in diagnostic procedures. © 2017 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. TaqMan is a registered trademark of Roche Molecular Systems, Inc., used under permission and license. Analytik Jena and qTOWER are trademarks of Analytik Jena AG. Bio-Rad, CFX384, CFX96, iQ, MiniOpticon, Chroma4, and Opticon are trademarks of Bio-Rad Laboratories, Inc. Eppendorf, Mastercycler, and realplex are trademarks of Eppendorf AG. Qiagen and Rotor-Gene are trademarks of Qiagen GmbH. Roche, LightCycler, and LightCycler Nano are trademarks of Roche Diagnostics, GmbH. Stratagene, MX3000P, MX3005P, and MX4000 are trademarks of Agilent Technologies, Inc. TaKaRa is a trademark of TaKaRa Holdings Japan. Cy is a trademark of GE Healthcare. **COL13556 0317**

TaqMan Fast Advanced Master Mix

Performance superior to standard master mixes in less than half the time

Features and benefits

- **Best-in-class performance**—superior sensitivity, accuracy, dynamic range, and specificity compared to standard mixes in standard mode
- **Engineered for enhanced benchtop stability**—stable at room temperature for up to 72 hours in preassembled reactions
- **Optimized for multiplexing**—validated for duplexing with exogenous and endogenous internal positive control assays
- **Reduced run times**—optimized on fast instruments and for fast cycling conditions on standard instruments
- **Seamless transition through the workflow**—validated with Applied Biosystems™ TaqMan® Assays for gene expression and microRNAs and Applied Biosystems™ TaqMan® Array Microfluidic Cards

Applied Biosystems™ TaqMan® Fast Advanced Master Mix has been designed to perform better than standard master mixes (Figure 1), requiring shorter run times (<40 minutes) and delivering superior results.

Our best-in-class gene expression master mix employs Applied Biosystems™ AmpliTaq™ Fast DNA Polymerase, which has been engineered for enhanced stability, allowing your preassembled reactions to be left at room temperature for up to 72 hours without impacting performance. The formulation has been optimized for duplex PCR with both endogenous and exogenous control assays, enabling you to run a control in every well to further increase confidence in your results.

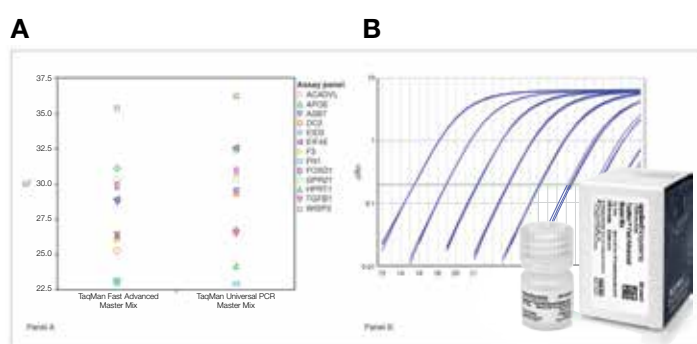
Figure 1. Performance of TaqMan Fast Advanced Master Mix vs. Applied Biosystems™ TaqMan® Universal PCR Master Mix. (A) Comparison of C_t values across a panel of 13 TaqMan Gene Expression Assays. **(B)** Representative amplification plot from real-time PCR of a dilution series of human cDNA amplified in 4 replicate reactions using the Applied Biosystems™ 7500 Fast Real-Time PCR System and the FN1 TaqMan Gene Expression Assay.

For maximum flexibility, TaqMan Fast Advanced Master Mix has also been optimized for use on both fast instruments and fast PCR cycling conditions on standard instruments. TaqMan Fast Advanced Master Mix has been rigorously tested and optimized to help ensure success with all TaqMan gene expression and microRNA assays, enabling a seamless transition through your workflow.

Table 1. Dynamic range comparison between TaqMan Fast Advanced Master Mix and products from other leading suppliers. Comparison of detection range (in number of logarithmic units) across a panel of various TaqMan Gene Expression Assays. The range of detection must have PCR efficiency between 85% and 115% and R² values ≥0.98. Each master mix was tested using cDNA template and run according to the manufacturers' respective recommended protocols. Reactions (6 replicates) were run on the Applied Biosystems™ 7900HT Fast Real-Time PCR System.

Assay	Assay type	TaqMan Fast Advanced Master Mix	Roche FastStart Reagent	Qiagen QuantiTect Reagent	Qiagen QuantiFast Reagent	Bio-Rad iTaq Supermix	Bio-Rad iTaq Fast Supermix	Orders of magnitude	Final (ng/μL)
<i>APOA1</i>	Good Fast	7	5	5	5	5	5	7	0.00001
<i>APOA1 (FAM)/ GAPDH (VIC)</i>	Good Fast	7	4	4	5	5	5	6	0.0001
<i>APOA1 (FAM)/ GAPDH (VIC)</i>	Housekeeping	7	7	7	7	7	7	5	0.001
<i>UBC</i>	Housekeeping	6	4	4	5	5	5	4	0.01
<i>HIST1H3F</i>	LenAmpLong	5	3	3	3	3	3	3	0.1
<i>TXNDC</i>	GCampLow, PrimerLong	5	2	2	3	3	3	2	1
<i>FOXD1</i>	GCampHigh	4	2	2	2	2	2	1	10
<i>GPR34</i>	GCProbeLow, Low dRn	3	1	2	2	2	2		
<i>WISP</i>	HighProbeTm	2	0	0	1	1	1		

Best-in-class performance

TaqMan Fast Advanced Master Mix was designed to perform better than current standard master mixes. Our master mix was benchmarked against the leading suppliers' standard and fast master mixes to demonstrate our superior sensitivity, accuracy, dynamic range, and specificity.

The unparalleled dynamic range of TaqMan Fast Advanced Master Mix is shown in Table 1. These results demonstrate the ability of the master mix to provide dependable target quantitation over a wider dynamic range compared to leading suppliers' standard and fast master mixes. For a variety of assays, TaqMan Fast Advanced Master Mix was capable of detection across 2 additional orders of magnitude when run under identical conditions.

Benchtop stability for high-throughput handling and convenience

TaqMan Fast Advanced Master Mix has been engineered to retain its high level of performance in preassembled reactions for up to 72 hours. If you use high-throughput liquid handling systems, the stability of this mix helps to ensure that the results on the first plate will mimic those of the last plate. For less extreme throughput needs, the enhanced stability of this master mix provides overall added convenience to your workflow, as you are no longer constrained to immediately running your plates upon assembly.

Figure 2 shows an assay that was run upon assembly (time 0) and after 72 hours of incubation at 30°C, simulating the most extreme room temperature scenario. The results

after 72 hours show excellent PCR efficiency and R² values, almost identical to those at time 0, as well as a ΔC_t between time 0 and 72 hours of less than 1.


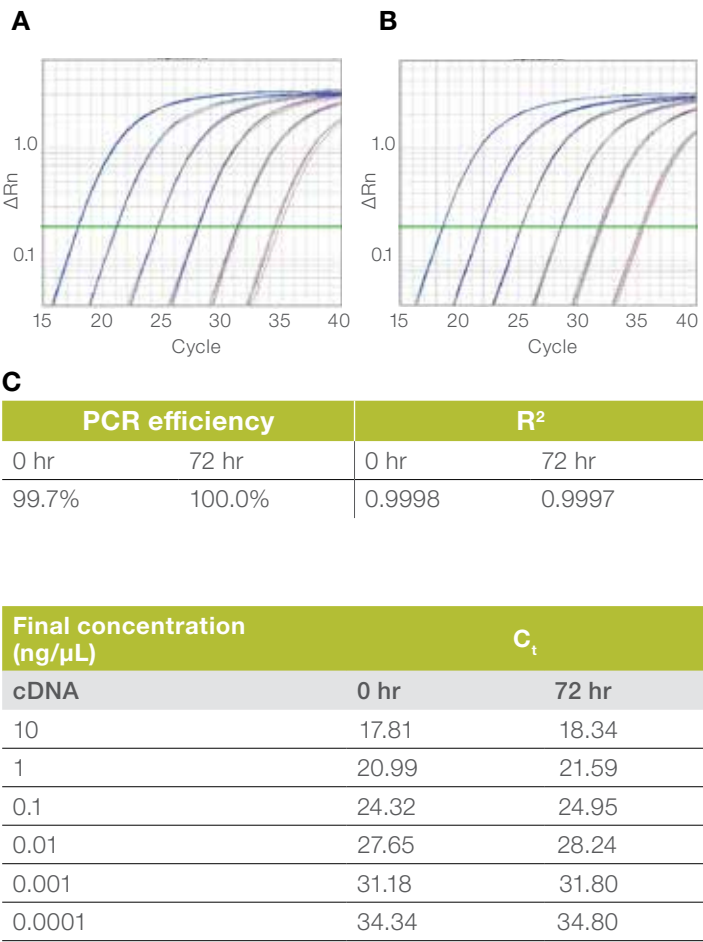
Figure 2. Benchtop stability of TaqMan Fast Advanced Master Mix. This B2M TaqMan Gene Expression Assay was run **(A)** upon assembly (time 0) and **(B)** after 72 hr of incubation at 30°C. **(C)** The results after 72 hr show excellent PCR efficiency, R² values, and C_t values when compared to time 0.

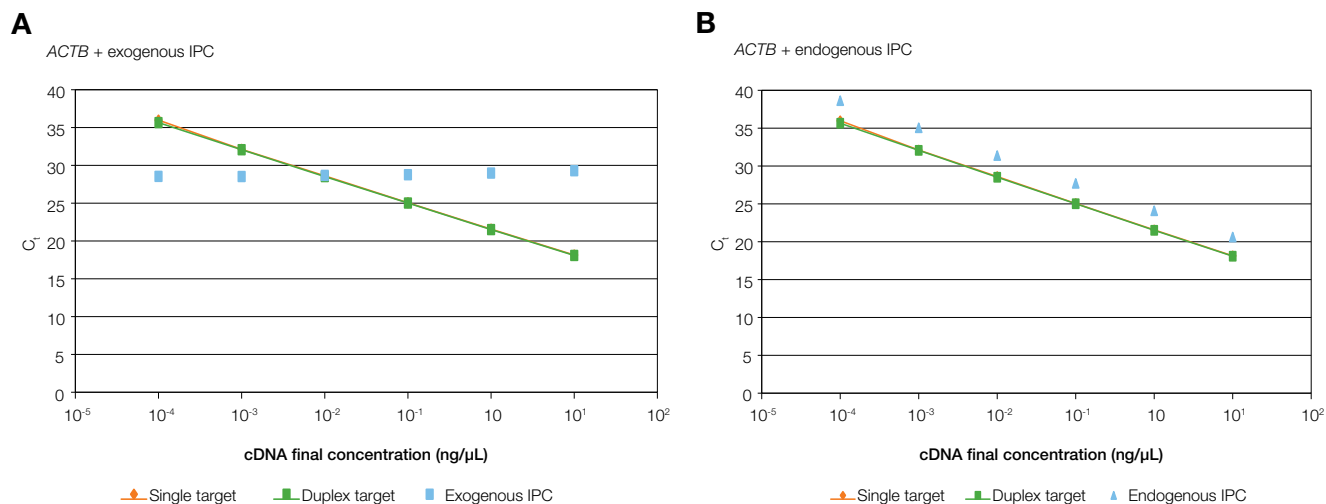Exogenous duplex

	PCR efficiency	R ²
Single target	92.9%	0.999
Duplex target	95.1%	1.000

Endogenous duplex

	PCR efficiency	R ²
Single target	92.9%	0.999
Duplex target	92.6%	1.000

Figure 3. TaqMan Fast Advanced Master Mix is optimized for multiplexing with exogenous or endogenous control assays. Results are shown for *ACTB* (β -actin gene), which was serially diluted and amplified in single-target reactions and duplex reactions. The duplex reactions included the single target *ACTB* and either **(A)** a constant quantity of exogenous target or **(B)** a relative quantity of endogenous target.

Optimized for multiplexing

We realize that confidence is paramount when it comes to your results. For added confidence in your results, TaqMan Fast Advanced Master Mix was designed to help deliver accurate results for duplex reactions using an internal positive control (IPC). Figure 3 shows results for the experimental target gene *ACTB* (β -actin), which was serially diluted and amplified in both single-target reactions and duplex reactions. The duplex reactions included the single target *ACTB* and either a constant quantity of exogenous target (Figure 3A) or a relative quantity of endogenous target (Figure 3B). TaqMan Fast Advanced Master Mix succeeded in providing nearly identical PCR efficiency, R², and C_t values for *ACTB* in both simplex and duplex environments.

Validated for microRNA assays

TaqMan Fast Advanced Master Mix has been validated for multiple real-time PCR applications, including microRNA assays. The formulation provides high specificity and a large dynamic range, the two most critical performance attributes that define successful results when working with microRNAs. The data in Figure 4 demonstrate excellent PCR linearity over a range of inputs, covering 6 orders of magnitude.

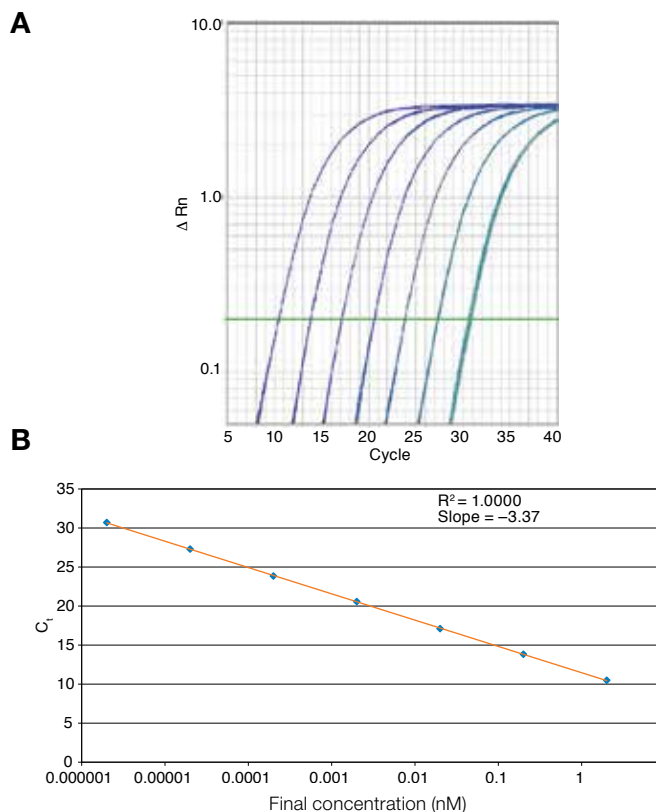
Figure 4. miRNA linear target amplification over a dynamic range of 6 orders of magnitude of input. **(A)** Amplification plot and **(B)** standard curve from real-time PCR of a dilution series of a synthetic target amplified in 4 replicate reactions using the 7900HT Fast Real-Time PCR System and the Let-7c TaqMan MicroRNA Assay.

Reduced run times on standard instrumentation

TaqMan Fast Advanced Master Mix has been optimized for use with both fast and standard instrumentation, enabling researchers who currently own standard instruments to realize the performance benefits and time savings this mix provides. Figure 5 showcases the impressive results achieved when using TaqMan Fast Advanced Master Mix under fast thermal cycling conditions on the Applied Biosystems™ QuantStudio™ 12K Flex Real-Time PCR System. The mix has been tested with all Applied Biosystems™ standard real-time PCR instrumentation, including the QuantStudio, 7900HT, 7500, and 7300 systems, to enable success whether or not you own a fast-enabled instrument.

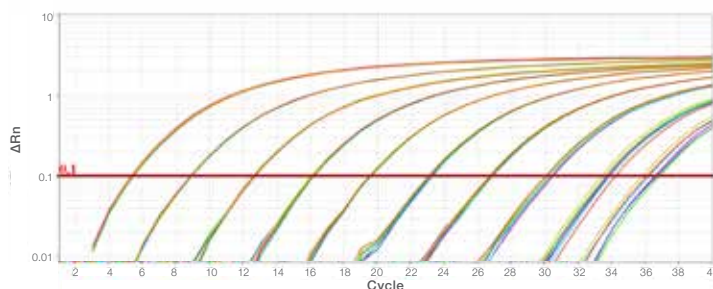
Figure 5. Results on the QuantStudio 12K Flex Real-Time PCR System using TaqMan Fast Advanced Master Mix. Amplification plot from real-time PCR of a dilution series of human cDNA amplified in 8 replicate reactions using the Eukaryotic 18S rRNA TaqMan Gene Expression Assay and the QuantStudio 12K Flex Real-Time PCR System.

Ordering information

Size	Quantity	No. of 20 μ L rxns	Cat. No.
TaqMan Fast Advanced Master Mix			
Mini Pack	1 x 1 mL	100	4444556
1 Pack	1 x 5 mL	500	4444557
2 Pack	2 x 5 mL	1,000	4444963
5 Pack	5 x 5 mL	2,500	4444964
10 Pack	10 x 5 mL	5,000	4444965
Bulk Pack	1 x 50 mL	5,000	4444558

Find out more at thermofisher.com/mastermixes

ThermoFisher
SCIENTIFIC

TaqMan Genotyping Master Mix


Real-time PCR master mix tailored for SNP genotyping studies

Tailored for unrivaled cluster resolution for unambiguous single-nucleotide polymorphism (SNP) allelic discrimination, the Applied Biosystems™ TaqMan™ Genotyping Master Mix is optimized for genotyping applications, including:

- Candidate gene studies
- Drug target validation
- Disease association studies
- Population genetics
- Linkage mapping
- Agricultural applications
- Copy number variation analysis

Introduction

TaqMan Genotyping Master Mix is designed to deliver reliable, cost-effective SNP detection for accurate and reproducible allelic discrimination. The master mix optimizes the preferential binding of the allele-specific probe, providing exceptional separation and clustering of alleles and consistently strong fluorescent signals. Powered with the highly purified Applied Biosystems™ AmpliTaq Gold™ DNA Polymerase, UP (Ultra Pure), TaqMan Genotyping Master Mix can replace Applied Biosystems™ TaqMan™ Universal PCR Master Mix in existing SNP genotyping protocols using the same reaction setup and thermal cycling conditions.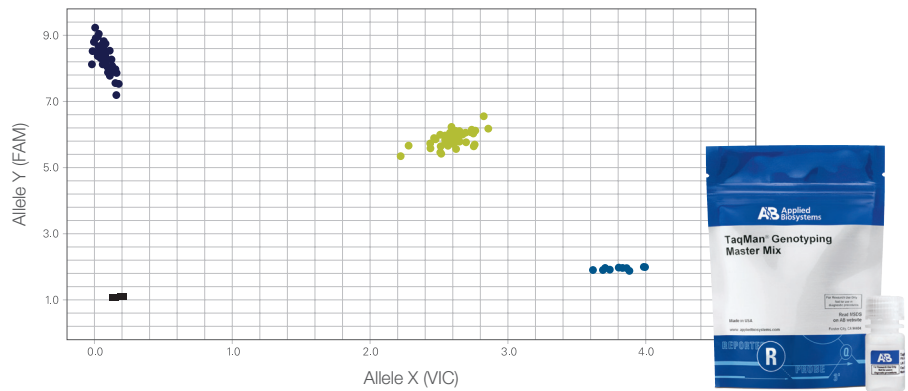

Benefits

- Specifically formulated for endpoint fluorescent detection of SNPs and insertions/deletions
- Discrete clusters and high call rates for accurate and reproducible allelic discrimination
- Reliable discrimination of SNPs in difficult targets
- Excellent room-temperature stability for flexible pre- and post-PCR setup and analysis
- Universal thermal cycling conditions for consistent results
- Validated for use with Applied Biosystems™ TaqMan™ SNP Genotyping Assays, TaqMan™ Copy Number Assays, and TaqMan™ Mutation Detection Assays

Optimized formulation for exceptional performance

TaqMan Genotyping Master Mix is a convenient 2X mix for TaqMan probe-based genotyping reactions. It includes the following components:

- AmpliTaq Gold DNA Polymerase, UP (Ultra Pure), a highly purified DNA polymerase. This hot-start enzyme is inactive at room temperature, so reactions can be set up on the benchtop. The enzyme is activated during thermal cycling.
- Optimized components including buffer and dNTPs for consistent, reliable genotypes
- Passive internal reference based on proprietary ROX™ dye for precise data analysis

Setting a new standard for allelic discrimination

For clear genotyping results, each allele-specific TaqMan™ probe must yield bright and consistent fluorescent signals to provide discrete clusters that are widely separated, indicating excellent specificity. The performance of TaqMan Genotyping Master Mix was tested using 3 ng samples of human genomic DNA (gDNA) and a validated SNP assay to genotype dbSNP rs2293052 in the gene *NOS1*. The resulting cluster plot (Figure 1) shows strong fluorescent signals for each allele and clear separation between the three clusters—easily discriminating the two homozygous and one heterozygous genotypes. In a comparison against five commercially available mixes, TaqMan Genotyping Master Mix shows the highest average call rate (Figure 2). Tight, well-separated clusters for each genotype provide exceptional call rates and, most importantly, accurate and efficient SNP analysis.

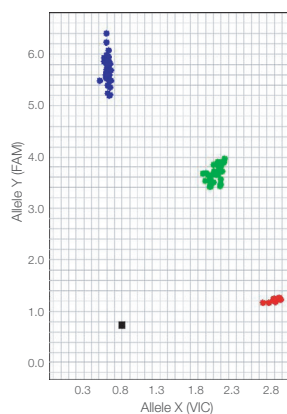
Figure 1. TaqMan Genotyping Master Mix provides bright fluorescence signals for discrete, well-separated allelic clusters. Cluster plot of 94 gDNA samples and two no-template controls genotyped using Applied Biosystems™ TaqMan™ SNP Genotyping Assay C__15969983_10, with PCR performed on the Applied Biosystems™ GeneAmp™ PCR System 9700 and allelic discrimination on the 7900HT Fast Real-Time PCR System.

Consistent performance—even with difficult templates

TaqMan Genotyping Master Mix offers unambiguous allelic discrimination even for the most challenging assays. For example, GC-rich targets can present amplification challenges that reduce SNP detection because of persistent secondary structure. Human gDNA samples were genotyped for a SNP in a GC-rich region using a TaqMan SNP Assay to genotype dbSNP rs12214 in the cathepsin D gene. As shown in Figure 3, TaqMan Genotyping Master Mix yields brighter fluorescent signals, tighter clusters, and more accurate allele calling compared to a mix from supplier “S”. These data demonstrate that TaqMan Genotyping Master Mix provides higher call rates for reliable SNP genotyping in difficult targets, eliminating the need to retest uncalled samples.

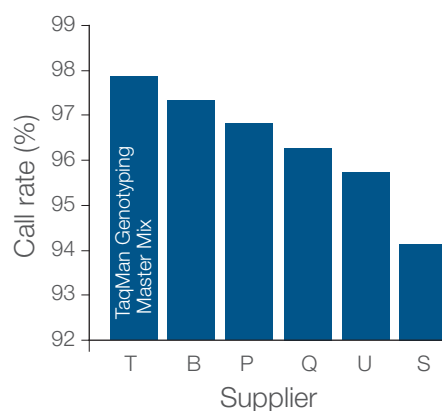
Figure 2. TaqMan Genotyping Master Mix provides the highest SNP call rates, outperforming other master mixes. Average call rates for 94 gDNA samples and two no-template controls genotyped using Applied Biosystems™ TaqMan™ SNP Genotyping Assay C__27102425_10, with PCR performed on the GeneAmp PCR System 9700 and allelic discrimination on the 7900HT Fast Real-Time PCR System.

Copy number variation applications

Copy number variation is an important polymorphism in the human genome that can be associated with certain genomic disorders as well as some simple genetic and complex diseases. TaqMan Genotyping Master Mix, used with TaqMan Copy Number Assays, provides relative quantitation of an experimental gene compared to a reference gene in a duplex PCR. Between 1 and 3 copies of *CYP2D6*, the gene for a drug-metabolizing enzyme, were detected for 92 human gDNA samples when the samples were amplified using TaqMan Genotyping Master Mix (Figure 4).

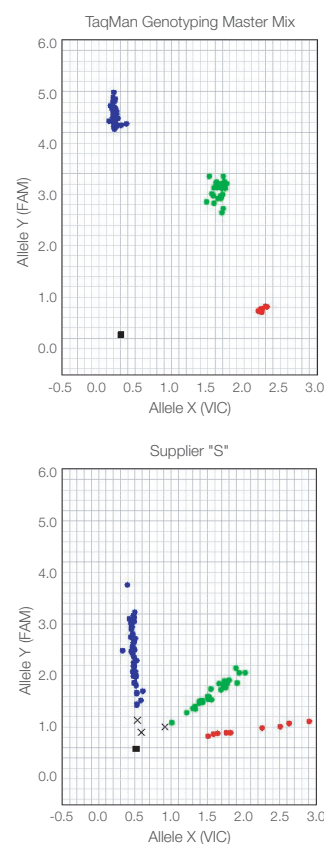
Figure 3. Consistent, reliable SNP detection in a GC-rich region using TaqMan Genotyping Master Mix. Genotyping assays were compared using TaqMan Genotyping Master Mix and a PCR master mix from supplier “S” on a set of 94 human gDNA samples (3 ng) and two no-template controls, using Applied Biosystems™ TaqMan™ SNP Assay C__12050942_10. PCR was performed on the GeneAmp PCR System 9700 and allelic discrimination on the 7900HT Fast Real-Time PCR System.

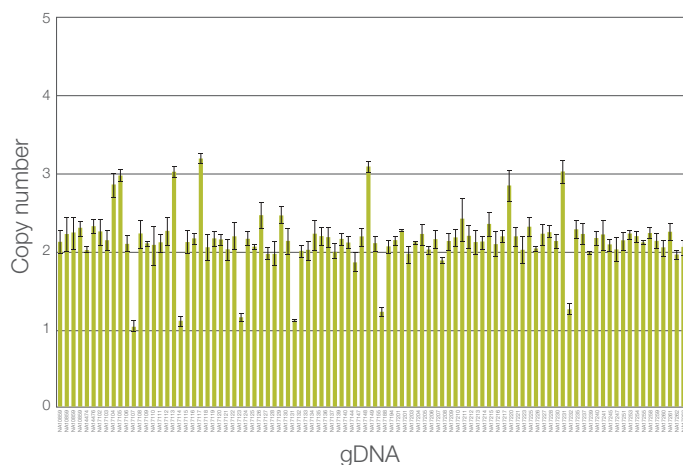
Figure 4. TaqMan Genotyping Master Mix is used for amplification in TaqMan Copy Number Assays. TaqMan Genotyping Master Mix is used with a TaqMan Copy Number Assay designed to target the *CYP2D6* gene, determining the copy number for this target in 92 gDNA samples. The RNase P reference gene is present in two copies per diploid genome or one copy per haploid genome.

Somatic mutation detection applications

TaqMan Mutation Detection Assays can detect somatic mutations in genes that are associated with cancer from different sample types, such as cell lines, formalin-fixed, paraffin-embedded tissue samples, and fresh frozen tissue samples. TaqMan Genotyping Master Mix, combined with TaqMan Mutation Detection Assays, which use competitive allele-specific Applied Biosystems™ TaqMan™ PCR (castPCR™) technology, can help detect rare amounts of mutated DNA in a sample that contains large amounts of normal, wild type DNA.

Pre- and post-PCR stability

Benchtop stability of real-time PCR mixes provides the flexibility to perform experiments over multiple days. To demonstrate the stability of TaqMan Genotyping Master Mix, both pre- and post-PCR storage conditions were tested to determine the effects on genotyping data. To evaluate pre-PCR stability, reactions were set up at room temperature (24°C), stored in the dark for up to three days, thermal-cycled for PCR, and read for endpoint fluorescence to assign alleles. To assess post-PCR stability, PCR was conducted immediately after reaction setup, but reactions were left on the bench for up to three days before measuring endpoint fluorescence for allelic discrimination. Even after three days at room temperature, either before or after PCR, TaqMan Genotyping Master Mix yielded tight clusters and reproducible results (Figure 5). The excellent benchtop stability of TaqMan Genotyping Master Mix gives ample flexibility for experimental setup and sample processing.

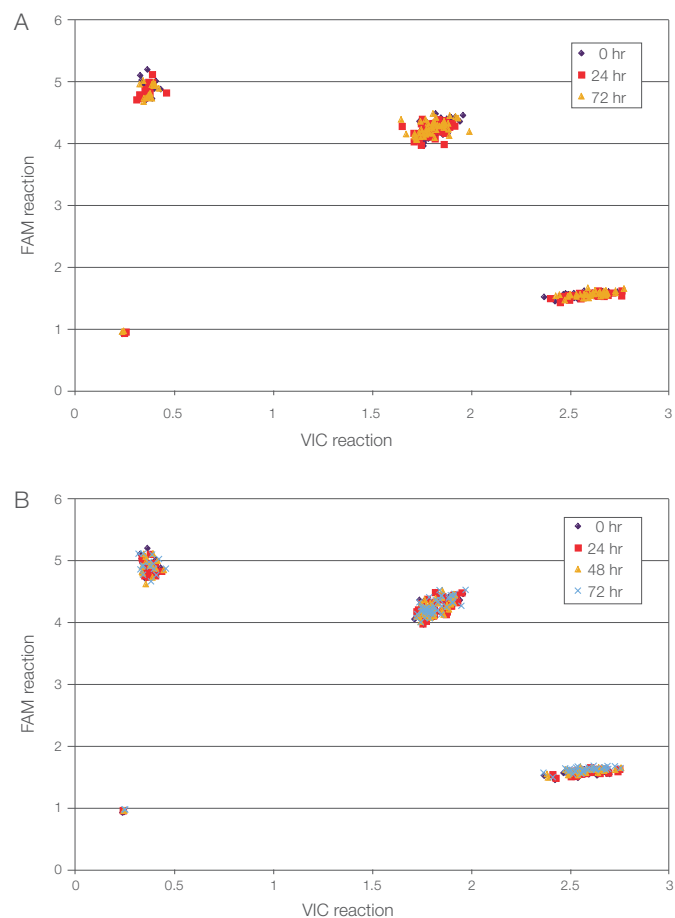
Figure 5. TaqMan Genotyping Master Mix provides pre- and post-PCR stability for up to 3 days. Genotyping reactions were set up using 94 gDNA samples and two no-template controls with TaqMan Genotyping Master Mix and TaqMan SNP Assay C____2188620_10. Reactions were left on the bench either (A) before or (B) after thermal cycling for the indicated amounts of time. PCR was conducted on the GeneAmp PCR System 9700 and allelic discrimination on the 7900HT Fast Real-Time PCR System.

Conclusion

TaqMan Genotyping Master Mix:

- Demonstrates extremely reliable allelic discrimination for SNP genotyping, with discrete clusters for high call rates even with challenging targets
- Provides reliable quantitation of DNA copy number when used with TaqMan Copy Number Assays
- Offers robust benchtop stability at room temperature, pre- and post-PCR, and consistent results across multiple instruments over multiple days to meet all throughput needs
- Complements TaqMan Mutation Detection Assays to provide high specificity and sensitivity for mutant allele detection

Instruments and assays compatible with TaqMan Genotyping Master Mix (standard thermal cycling mode)

Instruments and assays
QuantStudio 3/5/6/7/12 Real-Time PCR Systems
Applied Biosystems™ StepOne™ and StepOnePlus™ Real-Time PCR Systems
Applied Biosystems™ 7000, 7300, 7500, 7500 Fast, and 7900HT Fast Real-Time PCR Systems
Applied Biosystems™ Veriti™ Thermal Cyclers
GeneAmp PCR System 9700
Applied Biosystems™ 9800 Fast Thermal Cycler
TaqMan SNP Genotyping Assays
TaqMan Drug Metabolism Genotyping Assays
TaqMan Copy Number Assays
TaqMan Mutation Detection Assays
Applied Biosystems™ Custom TaqMan™ SNP Genotyping Assays
21 CFR Part 11 compliance module

Ordering information

Product	Unit size	Reactions*	Cat. No.
TaqMan Genotyping Master Mix			
Mini pack	1 mL tube	40	4371353
1-pack	10 mL bottle	400	4371355
2-pack	2 x 10 mL bottles	800	4381656
Single bulk pack	50 mL bottle	2,000	4371357
Multi-bulk pack	2 x 50 mL bottles	4,000	4381657
Quick Reference Card	1 card	—	4371130
Protocol	1 protocol	—	4371131

* Assumes 50 µL reaction volume; consult protocol for other recommended reaction volumes.

Find out more at thermofisher.com/taqmanmm

ThermoFisher
SCIENTIFIC

TaqMan SNP Genotyping Assays

Applied Biosystems™ TaqMan™ SNP Genotyping Assays from Thermo Fisher Scientific provide a highly flexible technology for detection of polymorphisms within any genome. Applied Biosystems™ TaqMan™ Assays have the simplest workflow available and are the quickest way to generate genotyping data. Based on powerful Applied Biosystems™ TaqMan™ probe and primer chemistry and designs, and coupled to dependable Applied Biosystems™ instruments and software, these made-to-order assays produce high-confidence results. TaqMan Assays are ideal for genotyping applications, including screening, association, candidate region, candidate gene, and fine-mapping studies.

Content-rich marker selection tools simplify study design and help you select from a library of human and mouse assays. This library includes over 7 million genome-wide human assays (of which 3.5 million are HapMap SNP-based assays, 160,000 are validated assays, and over 950,000 are coding region assays) and 10,000 mouse assays. We also offer 2,700 inventoried drug metabolism genotyping assays. Additionally, with Applied Biosystems™ Custom TaqMan™ SNP Genotyping Assays you can confidentially submit target SNP sequences for any genome to create your own assays. Let TaqMan SNP Genotyping Assays accelerate the pace of your discovery by eliminating time-consuming experimental design and optimization.

Powerful, proven chemistry

Whether your genotyping studies require targeted detection of essential SNPs, or the flexibility to choose SNPs for mapping, TaqMan SNP Genotyping Assays are the technology of choice. Proven TaqMan

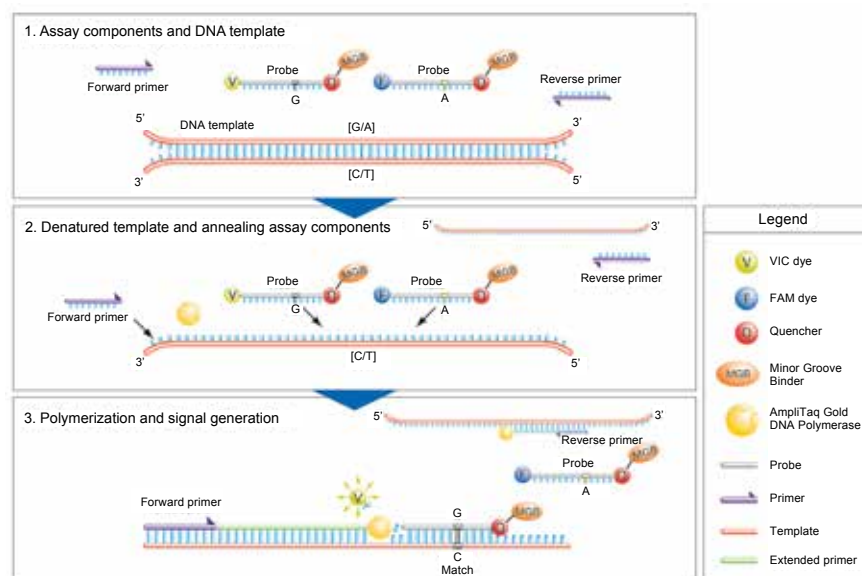
Figure 1. Allelic discrimination is achieved by the selective annealing of TaqMan MGB probes.

probes, which incorporate minor groove binder (MGB) technology at the 3' end, deliver superior allelic discrimination. The MGB molecule binds to the minor groove of the DNA helix, improving hybridization-based assays by stabilizing the MGB probe-template complex. This increased binding stability permits the use of probes as short as 13 bases for improved mismatch discrimination and greater flexibility when designing assays for difficult or variable sequences. In addition to SNP detection, TaqMan probes can be designed to detect multiple nucleotide polymorphisms (MNPs) and insertion/deletions (indels).

Detection is achieved with proven 5' nuclease chemistry by means of exonuclease cleavage of an allele-specific 5' dye label, which generates the permanent assay signal (Figure 1). All MGB probes include a nonfluorescent

quencher (NFQ) that virtually eliminates the background fluorescence associated with traditional quenchers, and provides a greater signal-to-noise ratio for superior assay sensitivity.

TaqMan SNP Genotyping Assays collection

TaqMan SNP Genotyping Assays are the world's largest collection of single-tube, ready-to-use SNP assays available. The TaqMan SNP Genotyping Assays library consists of two collections of human assays and one of mouse assays, and can be supplemented with assays designed using our Custom TaqMan SNP Genotyping Assays Service.

Over 7 million human SNP genotyping assays

This assay group contains over 7 million genome-wide SNPs, providing unprecedented marker coverage. Included in this collection are 160,000 validated assays that have approximately 10 kb spacing across gene regions. These assays were subjected to an extensive minor allele frequency test in 2–4 ethnic populations (45 individual samples per ethnic group) and as a result, offer the highest success rate. Also included are over 600,000 assays for the detection of nonsynonymous SNPs in coding regions, including many putative functional SNPs. Visit thermofisher.com/taqmansnp for more information.

Over 10,000 mouse SNP genotyping assays

The Applied Biosystems™ Mouse TaqMan™ Predesigned SNP Genotyping Assays collection consists of over 10,000 assays, and can be supplemented with assays designed using our Custom TaqMan SNP Genotyping Assays Service.

TaqMan Drug Metabolism Genotyping Assays

The collection of Applied Biosystems™ TaqMan™ Drug Metabolism Genotyping Assays includes 2,700 assays that target high-value polymorphisms in 221 drug metabolism genes. These assays have proven performance in four different ethnic population samples, consisting of 45 individuals each. To enable easy identification, these assays have been mapped to the common public allele nomenclature websites where possible. Visit thermofisher.com/taqmandme for more information.

All TaqMan SNP Genotyping Assays are generated using next-generation algorithms from the Thermo Fisher Scientific bioinformatics pipeline. For all predesigned assays, bioinformatics evaluation of target SNP sequences includes the masking of adjacent SNPs and ambiguous bases so that assay design and subsequent performance is not affected by the poor quality of the underlying sequence. Lastly,

the assay designs are aligned to the human genome using BLAST to ensure that each assay binds uniquely to the intended polymorphism. As the Custom TaqMan SNP Genotyping Assay Service is confidential and secure, you simply perform your own bioinformatics analysis prior to submitting your sequence for assay design.

Custom assay service for any possible SNP

Custom TaqMan SNP Genotyping Assays can be developed for any SNP in any organism. This service can generate assays for the detection of SNPs, MNPs, indels of up to 6 bases, or QSY™-labeled probes for multiplexing SNP assays.

Custom TaqMan SNP Genotyping Assays provide you with a complete service that includes secure and confidential ordering, assay design and manufacturing, and quality-control testing for synthesis accuracy and formulation completeness. Additionally, custom human assays are subjected to a functional test on 20 unique DNA samples.

Use the free Applied Biosystems™ Custom TaqMan™ Assay Design Tool to input and submit your sequence for assay design. This easy-to-use online resource lets you quickly submit your sequence information and start the ordering process securely and confidentially. Access the Custom TaqMan Assay Design Tool at thermofisher.com/snpcadt

Quality design and manufacturing

Probes and primers used in TaqMan SNP Genotyping Assays are designed using our rigorous bioinformatics pipeline. This proprietary group of algorithms has generated millions of TaqMan Assay designs by utilizing heuristic design rules deduced from both manufacturing and assay performance data. All assays are designed to perform under universal reaction conditions, as calculated probe and primer melting temperatures are consistent and include contributions from associated probe conjugates (i.e., dyes and MGB).

After manufacturing, assay components undergo extensive laboratory testing at our state-of-the-art manufacturing facility. Quality-control testing includes mass spectrometry for sequence verification and formulation assessments of probe and primer concentrations. Additionally, all human SNP genotyping assays are functionally tested to ensure allelic discrimination.

Simple workflow for quick results

TaqMan SNP Genotyping Assays constitute the simplest SNP genotyping technology available. We deliver your ready-to-use SNP genotyping assay at ambient temperature in a convenient, single-tube format. The rest is easy. Just combine the assay with Applied Biosystems™ TaqMan™ Genotyping Master Mix or TaqMan™ Universal PCR Master Mix and your purified DNA sample (Figure 2). There is no need to optimize probe, primer, salt concentrations, or temperature because all assays use universal reagent concentrations and thermal cycling conditions.

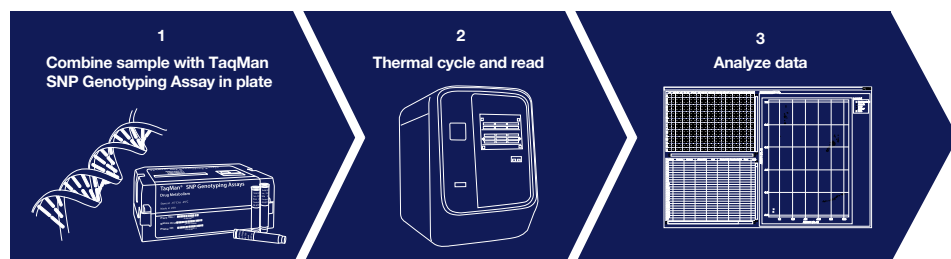
Figure 2. A simple workflow and reliable instruments combine to generate fast, high-confidence results.

After generating an endpoint read using a thermal cycler or real-time PCR instrument, no transfers, washes, or additional reagents are required, and the plate remains sealed; just read the plate and analyze the genotypes. This reduces the chance of contamination, sample mix-up, and sample loss. The simplicity of the chemistry allows you to easily automate the reaction for massively parallel genotyping studies, readily increasing the number of assays, number of samples, or both. Additionally, the analysis software allows you to auto-call genotypes, minimizing manual intervention.

Reliable real-time PCR platforms

A suite of superior Applied Biosystems instrument platforms is available for processing and analyzing TaqMan SNP Genotyping Assays (Table 1). These instruments, which meet all throughput needs and budgets, include the 7500, 7500 Fast, 7900HT Fast, ViiA™ 7, StepOne™, and StepOnePlus™ Real-Time PCR Systems, and the QuantStudio™ 3, 5, 6, 7, and 12K Flex Real-Time PCR Systems (Figure 3). Following PCR amplification, an endpoint read can be performed on any Applied Biosystems real-time PCR system. All of these dependable instruments offer the advanced multicolor detection capabilities required for highly accurate and reproducible allelic discrimination assays.

Data analysis software

The sophisticated SDS software package provided with all Applied Biosystems real-time PCR systems facilitates experimental setup, data collection, and assay performance analysis. The SDS software uses an advanced multicomponent algorithm to calculate the distinct signal contribution of each allele of a marker from the fluorescence measurements of each sample well during the assay plate read. The multicomponent data collected from the plate read are stored as SDS files, ready for genotype determination by the SDS software or optional Applied Biosystems™ TaqMan™ Genotyper Software (Figure 4).

Table 1. Applied Biosystems instrument capacities.

Instrument	Capacity
7500/7500 Fast Real-Time PCR System	96-well block (standard or Fast)
7900HT Fast Real-Time PCR System	96- and 384-well blocks (standard or Fast)
ViiA 7 Real-Time PCR System	96-well (standard or Fast), 384-well, and TaqMan Array Card blocks
StepOne Real-Time PCR System	48-well block (standard or Fast)
StepOnePlus Real-Time PCR System	96-well block (standard or Fast)
QuantStudio 3 Real-Time PCR System	96-well block (standard or Fast)
QuantStudio 5 Real-Time PCR System	96-well (standard or Fast) and 384-well blocks
QuantStudio 6 Real-Time PCR System	96-well (standard or Fast) and 384-well blocks
QuantStudio 7 Real-Time PCR System	96-well (standard or Fast), 384-well, and TaqMan Array Card blocks
QuantStudio 12K Flex Real-Time PCR System	96-well (standard or Fast), 384-well, TaqMan Array Card, and OpenArray plate blocks

Figure 3. The QuantStudio 5 Real-Time PCR System (left) and the QuantStudio 12K Flex Real-Time PCR System (right), which offers the highest throughput of all Applied Biosystems real-time PCR instruments.

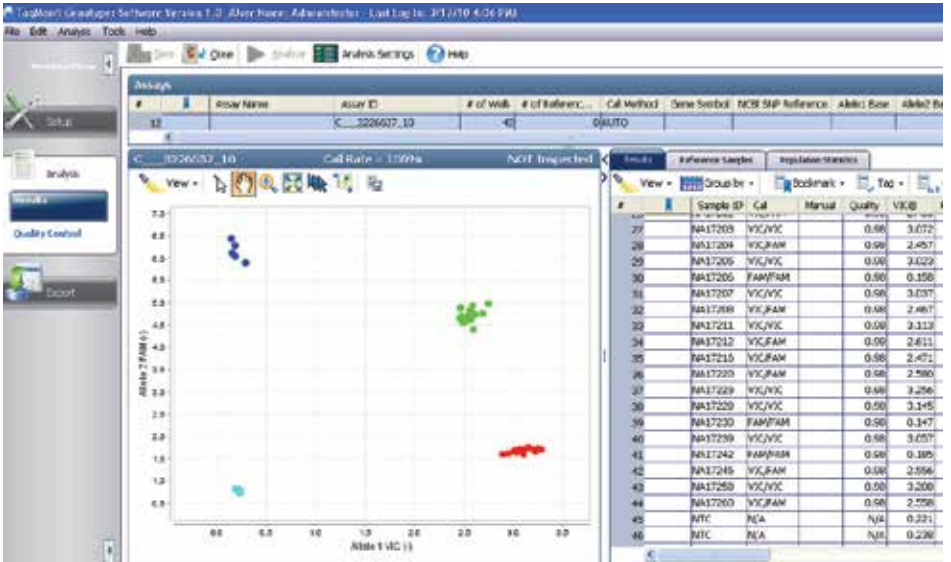
Figure 4. TaqMan Genotyper Software automatically determines sample genotypes and displays data.

TaqMan Genotyper Software is a great resource for fast and accurate genotype calling. It is a free SNP genotyping data analysis tool for use with TaqMan SNP Genotyping Assays performed in 48-, 96-, or 384-well microtiter plates or on Applied Biosystems™ TaqMan™ OpenArray™ Plates. It has a state-of-the-art genotype-calling algorithm, an intuitive user interface, and enhanced study-based analysis features. The software enables multi-plate data analysis for high-throughput workflows and improved accuracy in genotype calling; versatile export features and comprehensive quality-control features facilitate streamlining of the entire workflow. TaqMan Genotyper Software can be downloaded at thermofisher.com/taqmangenotyper

Simple ordering

Selecting and ordering TaqMan SNP Genotyping Assays is as simple as “point and click.” Use SNPbrowser Software to select the most informative SNPs for your genotyping studies. As you identify SNPs of interest, simply upload your selected TaqMan SNP Genotyping Assays to our online ordering tool.

The TaqMan Assay online ordering tool (Figure 5) enables you to search, select, and order from our catalog of over 7 million made-to-order predesigned TaqMan SNP Genotyping Assays. You can search for SNPs using any of several criteria: National Center for Biotechnology Information (NCBI) gene ID, NCBI SNP reference ID (rs#), or gene symbol. You can further refine your search by using SNP type (i.e., intragenic, 5’ or 3’ UTR, chromosome, etc.).

Our Custom TaqMan SNP Genotyping Assays supply you with SNPs that are not available from our predesigned assay collection, including those from any nonhuman organism. This service designs assays for all possible SNP, MNP, and indel targets but without the up-front bioinformatic preparation used for the predesigned made-to-order assays. Our complementary Custom TaqMan Assay Design Tool conveniently formats your target sequence for submission to our manufacturing facilities. To order custom assays, simply prepare your target sequence according to the Design and Ordering Guide, and upload your submission file at thermofisher.com/snpcaadt

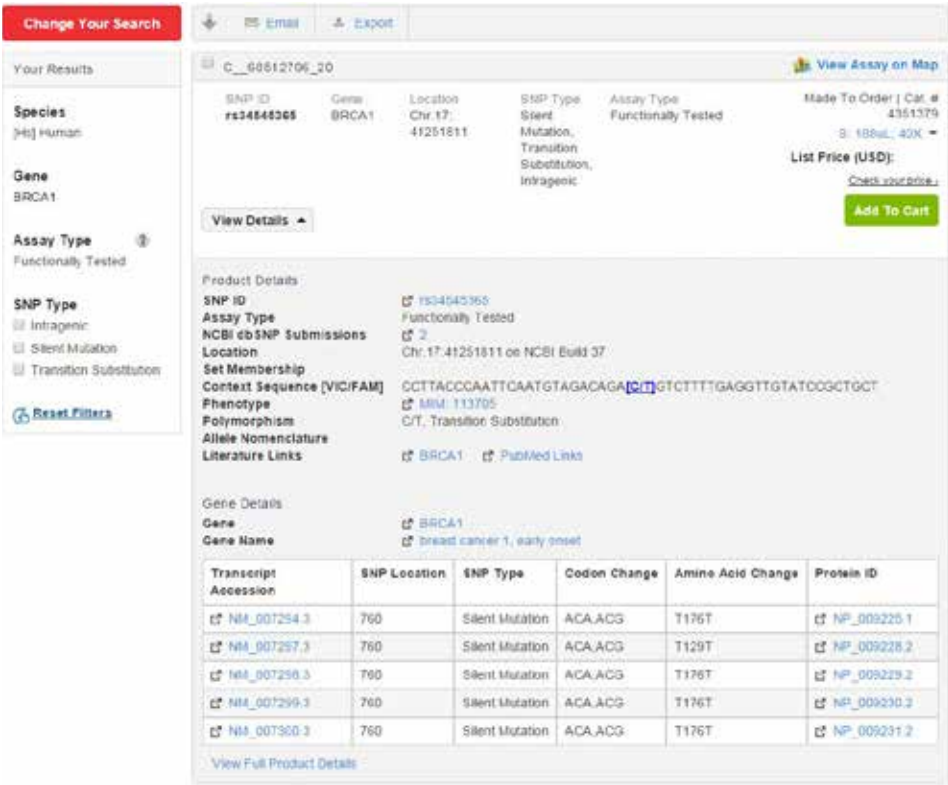
Figure 5. Our TaqMan Assay search and order tool makes online ordering easy. For convenient online ordering and multiple search options for all our genotyping assays, including keyword, batch, and location searches, visit thermofisher.com/taqmansnp

Ordering information

Size	Human assays (Cat. No.)	Nonhuman assays (Cat. No.)	Number of SNPs	No. of 5 µL reactions (384-well plate)	No. of 25 µL reactions (96-well plate)	Assay mix formulation	Assay type
Predesigned TaqMan SNP Genotyping Assays for Human and Mouse							
Small	4351379	4351384*	>4.5 million	1,500	300	40X	made-to-order
Medium	4351376	4351382*	>4.5 million	5,000	1,000	40X	made-to-order
Large	4351374	4351380*	>4.5 million	12,000	2,400	80X	made-to-order
Custom TaqMan SNP Genotyping Assays							
Small	4331349	4332077	∞	1,500	300	40X	made-to-order
Medium	4332072	4332075	∞	5,000	1,000	40X	made-to-order
Large	4332073	4332076	∞	12,000	2,400	80X	made-to-order
TaqMan Drug Metabolism Genotyping Assays							
Small	4362691	NA	2,700	750	150	20X	inventoried

All assays are quality-control tested using a mass spectrometer to verify sequence and yield. All assays have a VIC™ dye–labeled probe, a FAM™ dye–labeled probe, and two target-specific primers. All assays, excluding Custom TaqMan SNP Genotyping Assays, undergo bioinformatics evaluation of target SNP sequences.
Functional testing against 20 unique genomic DNA samples is performed on all custom and predesigned made-to-order human TaqMan SNP Genotyping Assays. Validation testing against four populations (45 samples/population) was performed on all 160,000 validated TaqMan SNP Genotyping Assays, and all TaqMan Drug Metabolism Genotyping Assays.
* Over 10,000 mouse assays available.

applied
biosystems

Find out more at thermofisher.com/taqmansnp

ThermoFisher
S C I E N T I F I C

For Research Use Only. Not for use in diagnostic procedures. © 2016 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified. TaqMan is a trademark of Roche Molecular Systems, Inc., used under permission and license. **CO020271 0116**

Understanding the Differences in PCR Plastics Format

Publication Number MAN0025596 Revision A.0

■ Description	1
■ PCR plate formats	1
■ PCR tube formats	6
■ PCR sealing closures	8
■ How to use the PCR Plastics Selection Tool	9
■ Quality	13


Description

The items listed throughout this user bulletin are representative of available plastics and other consumables. Choosing the correct plastic for your instrument is important for ensuring success. Use the qPCR and PCR Plastics Selection Tool to find the right plastic for you:

<http://www.thermofisher.com/findplastics>.

PCR plate formats

Number of wells	
96 wells	384 wells
	
Most commonly used.	For use with high-throughput automation platforms.

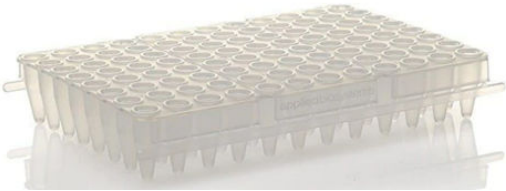Number of wells, segmented ^[1,2]	
8 wells	24 wells
	
A segmented 96-well plate that can be broken into 8-tube strips.	A plate provided with 24 wells.
32 wells	48 wells
	
A segmented 96-well plate that can be broken into three 32-well plates.	A plate provided with 48 wells.

[1] 8 and 32-well PCR plates often come as segmented 96-well plates that can easily be broken apart.

[2] Convenient format allows for a lower number of samples.

Material	
Polypropylene frame and wells	Polycarbonate frame with polypropylene wells
	
• Can withstand rapid changes in temperature. • Minimizes absorption of reaction components.	Same polypropylene features but also contains a rigid frame compatible with high-throughput robotic applications.

Frame color

Available in various colors to provide visual organization and identification of samples.

Well color ^[1]		
Clear or colored	White	Frosted
		
Recommended for end-point PCR applications.	- Ensures the highest level of sensitivity for qPCR reactions. - Recommended for low copy number samples.	Alternative to white plastic for qPCR if the fluorescence detector is oversaturated.


^[1] Refer to qPCR instrument manufacturer for recommendations on well color for proper plate calibration.

Skirt type		
Skirted	Semi-skirted	Non-skirted
		
- Full height panel surrounding the edge of the plate. - Fits securely over a thermal cycler with a raised block. - Enhanced mechanical strength for use with robotic platforms.	- Short panel around the edge of the plate. - Adequate support during pipetting. - Mechanical strength for robotic handling.	- No panel surrounding the edge of the plate. - Fits the block of most thermal cyclers and real-time PCR instruments. - Not suitable for robotic applications.

Automation compatible
• Skirted and semi-skirted plates provide a necessary side surface for grippers to grasp. • Polycarbonate-framed plates are able to withstand the forces exerted by grippers.

Thermal block fit		
96-well 0.2 mL (standard profile)	96-well 0.1 mL (low profile/fast)	384-well
		
• Most commonly used. • 0.3 mL maximum well volume.	• Reduced height reduces evaporation and enhances thermal conductivity. • 0.2 mL maximum well volume.	• Ideal for high-throughput applications. • 40 µL maximum well volume.

Alphanumeric labeling	
Printed	Engraved/molded
	
High contrast, easy to read.	May improve sealing of the outer edges.

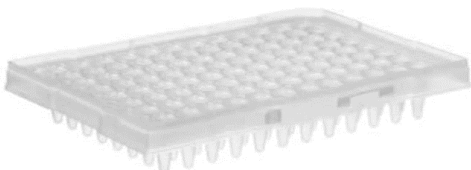Deck type	
Flat	Raised
	
• Universal fit with most thermal cyclers. • Facilitates sealing and handling.	• Required for some thermal cycler lids and instruments. • Balances lid pressure without the need for adapters.

Frame notch

A modified corner of a PCR plate that some instruments require for optimal fit.

Barcoding

Allows for easy recording and tracking of samples in high-throughput experiments.

PCR tube formats

To find the correct PCR plate or tube for your instrument use the qPCR and PCR Plastics Selection Tool: <http://www.thermofisher.com/findplastics>.

Tube format	
Single	Strip tubes
	
Individual PCR tubes with separate or attached caps provide the flexibility to set up the exact number of reactions to run.	Offered in 8 or 12 tube strips with separate or attached caps.

Well profile	
0.1 mL (low profile)	0.2 mL (standard profile)
	
- Fits 0.1 mL thermal blocks. - Reduced height minimizes air space above the reaction to reduce evaporation and enhance thermal conductivity.	- Fits 0.2 mL thermal blocks. - Most commonly used.

Tube color

Available in various colors to provide visual organization and identification of samples.

Cap type	
Domed	Flat
	
• For end-point PCR applications. • May be required for instruments with concave lids.	• For qPCR applications, use optical or ultra-clear caps. • May be required for instruments with flat lids.

Cap format ^[1]	
Separate	Attached
	
Strip caps can be used with standard or low profile PCR strip tubes and plates.	Can open and close tubes independently to prevent sample contamination.

^[1] To ensure compatibility select strip tubes and caps from the same brand, Applied Biosystems™ or Thermo Scientific™

PCR sealing closures

To find the correct PCR plate or tube for your instrument use the qPCR and PCR Plastics Selection Tool: <http://www.thermofisher.com/findplastics>.


Closure type		
Sealing films	Sealing foils	Strip caps
		
Creates a tight seal over wells to prevent evaporation and cross-contamination. • Clear adhesive films—For PCR, economical and sticky. • Optical sealing films—For qPCR applications, non-tacky pressure sensitive adhesive.	Creates a tight seal over wells to prevent evaporation and cross-contamination. • For use with light-sensitive samples. • Pierceable with pipette tip for sample transfer.	Seals securely with minimal pressure. Can be used with strip tubes or PCR plates. • Domed caps—For PCR. • Flat caps—For qPCR applications, maximizes the passage of fluorescence signals and minimize distortion.

How to use the PCR Plastics Selection Tool

1. Go to <http://www.thermofisher.com/findplastics>.

qPCR and PCR Plastics Selection Tool

Find the right fit, **fast**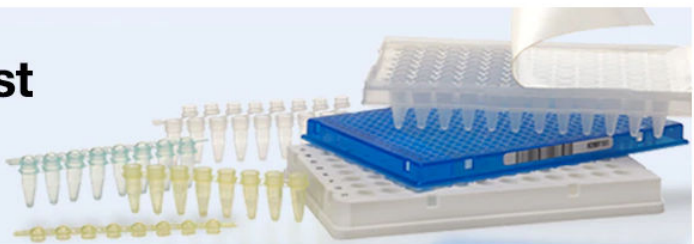

Select your real-time PCR (qPCR) or thermal cycler (PCR) instrument from the drop-down menu, and easily find the plates, tube strips, single tubes, caps, seals, adhesive films, and accessories compatible with your instrument.

1. Manufacturer

2. Model

3. Block Format

Please choose a manufacturer, model, and block format first.

2. Select the instrument manufacturer.

1. Manufacturer

2. Model

3. Block Format

Thermo Fisher Scientific / Applied Biosystems
Bio-Rad
Agilent
Analytik Jena / Biometra
Axygen
Bioer / Bulldog
Bioneer

3. Select the instrument model.


1. Manufacturer

2. Model

3. Block Format


Applied Biosystems MiniAmp
Applied Biosystems MiniAmp Plus
Applied Biosystems ProFlex
Applied Biosystems QuantStudio 12k
Applied Biosystems Quant Studio 3

4. Select the block format.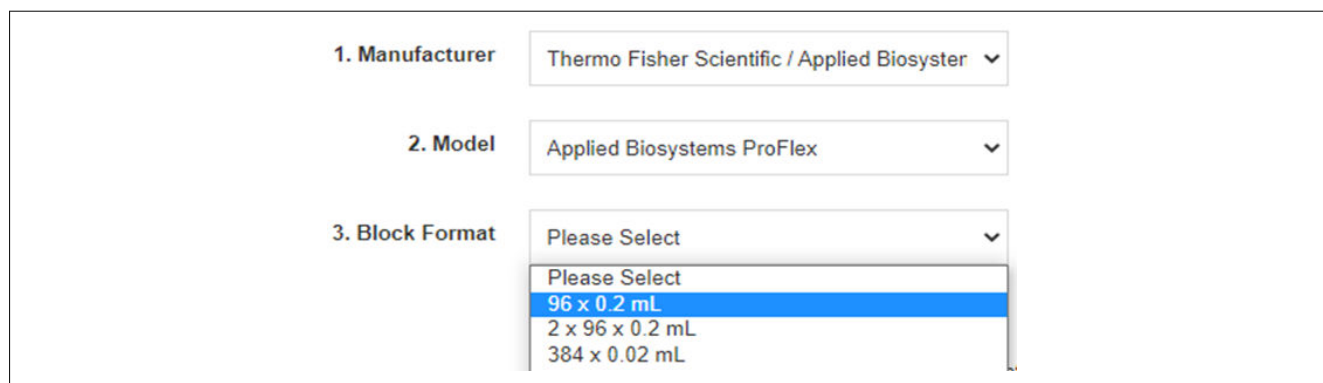

The screenshot shows a web interface with three dropdown menus. The first menu, labeled '1. Manufacturer', is set to 'Thermo Fisher Scientific / Applied Biosystem'. The second menu, labeled '2. Model', is set to 'Applied Biosystems ProFlex'. The third menu, labeled '3. Block Format', is open, showing a list of options: 'Please Select', '96 x 0.2 mL', '2 x 96 x 0.2 mL', and '384 x 0.02 mL'. The '96 x 0.2 mL' option is highlighted in blue.

5. Click **Display Results**.

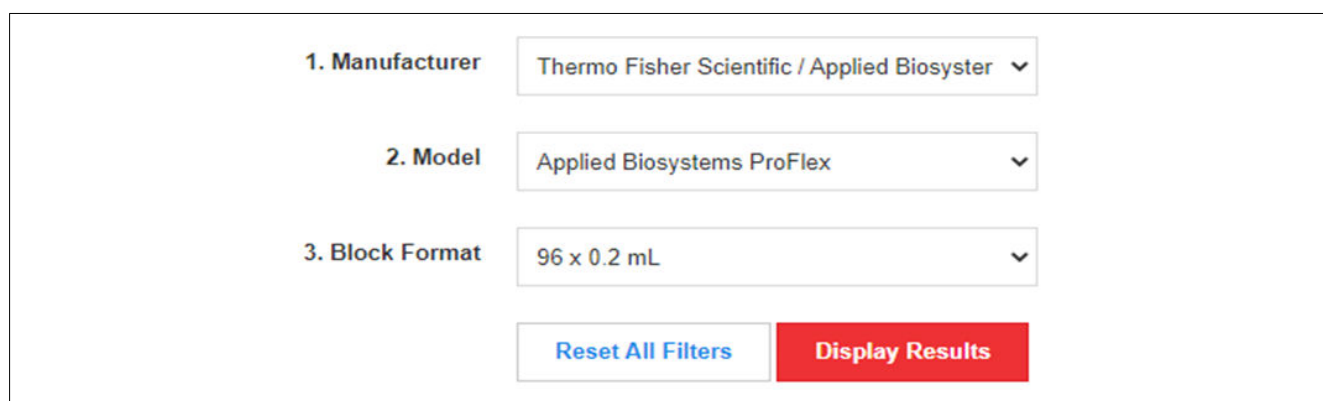The screenshot shows the same web interface as before, but the 'Block Format' dropdown menu is now closed and set to '96 x 0.2 mL'. Below the dropdowns are two buttons: a blue 'Reset All Filters' button and a red 'Display Results' button.

6. Filter results using features in left-hand column.

Consumable Type Please Select ▼ Number of Wells Please Select ▼ Plate Skirt Type Please Select ▼ Well Profile Please Select ▼ Color Please Select ▼ Well Color Please Select ▼ Barcoded Please Select ▼ Automation Compatible Please Select ▼ Reset filters	132 results found <table border="1"> <thead> <tr> <th>Catalog #</th> <th>Product Name</th> <th>Product Size</th> </tr> </thead> <tbody> <tr> <td>403012</td> <td>MicroAmp™ Optical 96-Well Reaction Plate with Barcode & Optical Caps</td> <td>20 plates</td> </tr> <tr> <td>4306311</td> <td>MicroAmp™ Clear Adhesive Film</td> <td>100 seals</td> </tr> <tr> <td>4306737</td> <td>MicroAmp™ Optical 96-Well Reaction Plate with Barcode</td> <td>20 plates</td> </tr> <tr> <td>4311971</td> <td>MicroAmp™ Optical Adhesive Film</td> <td>100 seals</td> </tr> <tr> <td>4312063</td> <td>MicroAmp™ Splash-Free 96-Well Base</td> <td>10 bases</td> </tr> <tr> <td>4313950</td> <td>MicroAmp™ Multi Removal Tool</td> <td>1 tool</td> </tr> <tr> <td>4314320</td> <td>MicroAmp™ Optical 96-Well Reaction Plate with Barcode & Optical Adhesive Films</td> <td>100 plates</td> </tr> <tr> <td>4316567</td> <td>MicroAmp™ Optical 8-Tube Strip, 0.2 mL</td> <td>125 strips</td> </tr> <tr> <td>4316813</td> <td>MicroAmp™ Optical 96-Well Reaction Plate</td> <td>500 plates</td> </tr> <tr> <td>4323032</td> <td>MicroAmp™ Optical 8-Cap Strips</td> <td>300 strips</td> </tr> <tr> <td>4326659</td> <td>MicroAmp™ Optical 96-Well Reaction Plate with Barcode</td> <td>500 plates</td> </tr> <tr> <td>4333183</td> <td>MicroAmp™ Adhesive Film Applicator</td> <td>5 applicators</td> </tr> <tr> <td>4360954</td> <td>MicroAmp™ Optical Adhesive Film</td> <td>25 seals</td> </tr> </tbody> </table>	Catalog #	Product Name	Product Size	403012	MicroAmp™ Optical 96-Well Reaction Plate with Barcode & Optical Caps	20 plates	4306311	MicroAmp™ Clear Adhesive Film	100 seals	4306737	MicroAmp™ Optical 96-Well Reaction Plate with Barcode	20 plates	4311971	MicroAmp™ Optical Adhesive Film	100 seals	4312063	MicroAmp™ Splash-Free 96-Well Base	10 bases	4313950	MicroAmp™ Multi Removal Tool	1 tool	4314320	MicroAmp™ Optical 96-Well Reaction Plate with Barcode & Optical Adhesive Films	100 plates	4316567	MicroAmp™ Optical 8-Tube Strip, 0.2 mL	125 strips	4316813	MicroAmp™ Optical 96-Well Reaction Plate	500 plates	4323032	MicroAmp™ Optical 8-Cap Strips	300 strips	4326659	MicroAmp™ Optical 96-Well Reaction Plate with Barcode	500 plates	4333183	MicroAmp™ Adhesive Film Applicator	5 applicators	4360954	MicroAmp™ Optical Adhesive Film	25 seals
Catalog #	Product Name	Product Size																																									
403012	MicroAmp™ Optical 96-Well Reaction Plate with Barcode & Optical Caps	20 plates																																									
4306311	MicroAmp™ Clear Adhesive Film	100 seals																																									
4306737	MicroAmp™ Optical 96-Well Reaction Plate with Barcode	20 plates																																									
4311971	MicroAmp™ Optical Adhesive Film	100 seals																																									
4312063	MicroAmp™ Splash-Free 96-Well Base	10 bases																																									
4313950	MicroAmp™ Multi Removal Tool	1 tool																																									
4314320	MicroAmp™ Optical 96-Well Reaction Plate with Barcode & Optical Adhesive Films	100 plates																																									
4316567	MicroAmp™ Optical 8-Tube Strip, 0.2 mL	125 strips																																									
4316813	MicroAmp™ Optical 96-Well Reaction Plate	500 plates																																									
4323032	MicroAmp™ Optical 8-Cap Strips	300 strips																																									
4326659	MicroAmp™ Optical 96-Well Reaction Plate with Barcode	500 plates																																									
4333183	MicroAmp™ Adhesive Film Applicator	5 applicators																																									
4360954	MicroAmp™ Optical Adhesive Film	25 seals																																									

Note: To return to the full list of result, click **Reset Filters**.

Understanding the filter options

Consumable Type

- Choose between plate, tube strip, individual tube, cap strip, seal, and accessory.

Number of Wells

- Choose 8 or 12 for strip tubes and caps.
- Choose 8, 24, 32, 48, 96, or 384 for plates.

Plate Skirt Type

- Choose skirted, semi-skirted, or non-skirted.

Well Profile

- Choose between standard or low profile/fast.

Color

- Choose from the available plate colors. Common options include clear, white, black, and single or assorted colors.

Well Color

- Choose from the available colors. Common options include clear, white, black, single and assorted colors.

Barcoded

- Narrow options based on whether or not they have a barcode.

Automation Compatible

- Narrow options based on whether or not they are designed for use with automated robotic applications.

Quality

Production

- The production process is carried out in a Class 100,000 cleanroom under ISO 13485 or ISO 9001 guidelines.

Materials

- All PCR plastics are manufactured using medical-grade virgin polypropylene.

Design

- Made with polished tools designed to produce chemical-free and ultra-smooth well surfaces that help prevent inhibition and binding.

QC testing

- Each well is visually inspected and electrostatically tested for defects; samples of each lot are run through a PCR cycling test to check for sealing and contaminants.

For descriptions of symbols on product labels or product documents, go to thermofisher.com/symbols-definition.

The information in this guide is subject to change without notice.

DISCLAIMER: TO THE EXTENT ALLOWED BY LAW, THERMO FISHER SCIENTIFIC INC. AND/OR ITS AFFILIATE(S) WILL NOT BE LIABLE FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE, OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING YOUR USE OF IT.

Revision history: Pub. No. MAN0025596

Revision	Date	Description
A.0	7 September 2021	Baseline for this revision history.

Important Licensing Information: This product may be covered by one or more Limited Use Label Licenses. By use of this product, you accept the terms and conditions of all applicable Limited Use Label Licenses.

TRADEMARKS: All trademarks are the property of Thermo Fisher Scientific and its subsidiaries unless otherwise specified.

©2021 Thermo Fisher Scientific Inc. All rights reserved.

MicroAmp™ plastic consumables compatibility chart for Applied Biosystems™ endpoint PCR systems and genetic analyzers

		3 x 32-well		96-well 0.2 mL			96-well 0.1 mL		384-well		Genetic analyzers	
		ProFlex™	ProFlex, SimpliAmp™, VeritiPro™, Veriti™, MiniAmp™ Plus, MiniAmp		2720	9700	Veriti	ProFlex, Veriti , VeritiPro™		9700	310	3130, 3130xl, 3500, 3500xL, 3730, 3730xl
Product	Cat. No.											
96-well 0.2 mL reaction plates												
Optical 96-Well Plate	N8010560, 4316813		•	•	•							•
Optical 96-Well Plate with Barcode	4306737, 4326659		•	•	•							•
96-Well Plate with Barcode & Optical Caps	403012		•	•	•							
Optical 96-Well Plate with Barcode & Optical Adhesive Films	4314320		•	•	•							
EnduraPlate Optical 96-Well Clear Plate	A36924		•	•	•							•
EnduraPlate Optical 96-Well Clear Plate with Barcode*	4483354, 4483352		•	•	•							•
TriFlex 3 x 32-Well Reaction Plate	A32810, A32811	•	•	•	•							
96-well 0.1 mL reaction plates												
Fast Optical 96-Well Plate, 0.1 mL	4346907						•				•	•
Fast Optical 96-Well Plate with Barcode, 0.1 mL	4346906, 4366932						•				•	•
EnduraPlate Optical 96-Well Fast Plate	A36930						•				•	•
EnduraPlate Optical 96-Well Fast Clear Plate with Barcode*	4483485, 4483494						•				•	•
384-well reaction plates												
Optical 384-Well Plate	4343370							•	•			•
Optical 384-Well Plate with Barcode	4309849, 4326270, 4343814							•	•			•
EnduraPlate Optical 384-Well Plate	A36931							•	•			•
EnduraPlate Optical 384-Well Clear Plate with Barcode*	4483285, 4483273							•	•			•
Strip tubes and caps												
Fast 8-Tube Strip, 0.1 mL	4358293		•	•	•	•	•					
Optical 8-Tube Strip with Attached Optical Caps, 0.2 mL	A30588	•	•	•	•	•						
8-Tube Strip with Attached Domed Caps, 0.2 mL	A30589	•	•	•	•	•						
8-Tube Strip, 0.2 mL*	N8010580	•	•	•	•	•					•	
Optical 8-Tube Strip, 0.2 mL	4316567	•	•	•	•	•						
8-Cap Strip*	N8010535, N8011535	•	•	•	•	•	•					
Optical 8-Cap Strip	4323032	•	•	•	•	•	•					
12-Cap Strip*	N8010534, N8011534		•	•	•	•	•					
Single tubes												
Fast Reaction Tube with Cap, 0.1 mL	4358297, 4358293						•					
Reaction Tube with Cap, 0.2 mL*	N8010540, N8010612, N8011540	•	•	•	•	•						
Reaction Tube without Cap, 0.2 mL*	N8010533, N8011533	•	•	•	•	•						
Optical Tube without Cap, 0.2 mL	N8010933	•	•	•	•	•						
Seals and covers												
Clear Adhesive Film	4306311		•	•	•	•	•	•	•			
Optical Adhesive Film	4360954, 4311971		•	•	•	•	•	•	•			
96-Well Full Plate Cover	N8010550				•	•						
32-Well Clear Adhesive Film	A32812	•	•	•	•	•						
Accessories												
Splash-Free 96-Well Base	4312063		•	•	•	•	•					
96-Well Support Base	4379590		•	•	•	•	•					•
96-Well Base	N8010531		•	•	•	•						
96-Well Reaction Tube/Tray/Retainer Set, 0.2 mL	403083, 403086			•	•	•						

* Multiple colors are available.

Note: Experiments using one or two 8-tube strips with attached caps require blank tube strips to balance lid pressure on the block or the use of the Applied Biosystems™ MicroAmp™ 96-Well Tray/Retainer Set (Cat. No. 4381850)—bottom part of the tray *only*. For use with the 96-well block of the ProFlex, SimpliAmp, Veriti, VeritiPro, MiniAmp Plus, and MiniAmp thermal cyclers.

Find out more at thermofisher.com/pcrplastics

MicroAmp™ plastic consumables compatibility chart for Applied Biosystems™ real-time PCR systems

Product	Cat. No.	48-well	96-well 0.2 mL			96-well 0.1 mL			384-well
		StepOne™	7000	7300, 7500	QuantStudio™ 3, 5, 6, 6 Pro, 7, 7 Pro, 12K; ViiA™ 7; 7900HT	StepOnePlus™	7500	QuantStudio 3, 5, 6, 6 Pro, 7, 7 Pro, 12K; ViiA 7; 7900HT**	QuantStudio 5, 6, 6 Pro, 7, 7 Pro, 12K; ViiA 7; 7900HT
96-well 0.2 mL reaction plates									
Optical 96-Well Plate	N8010560, 4316813		•	•	•				
Optical 96-Well Plate with Barcode	4306737, 4326659		•	•	•				
Optical 96-Well Plate with Barcode & Optical Caps	403012		•	•	•				
Optical 96-Well Plate with Barcode & Optical Adhesive Films	4314320		•	•	•				
EnduraPlate Optical 96-Well Clear Plate	A36924			•	•				
EnduraPlate Optical 96-Well Clear Plate with Barcode*	4483354, 4483352			•	•				
96-well 0.1 mL reaction plates									
Fast Optical 96-Well Plate, 0.1 mL	4346907					•	•	•	
Fast Optical 96-Well Plate with Barcode, 0.1 mL	4346906, 4366932					•	•	•	
EnduraPlate Optical 96-Well Fast Plate	A36930					•	•	•	
EnduraPlate Optical 96-Well Fast Clear Plate with Barcode*	4483485, 4483494					•	•	•	
384-well reaction plates									
Optical 384-Well Plate	4343370								•
Optical 384-Well Plate with Barcode	4309849, 4326270, 4343814								•
EnduraPlate Optical 384-Well Plate	A36931								•
EnduraPlate Optical 384-Well Clear Plate with Barcode*	4483285, 4483273								•
48-well reaction plates									
Fast Optical 48-Well Plate	4375816	•							
Strip tubes and caps									
Fast 8-Tube Strip, 0.1 mL	4358293	•				•	•	•	
Optical 8-Tube Strip with Attached Optical Caps, 0.2 mL	A30588		•	•	•				
Optical 8-Tube Strip, 0.2 mL	4316567		•	•	•				
Optical 8-Cap Strip	4323032	•	•	•	•	•	•	•	
Single tubes and caps									
Fast Reaction Tube with Cap, 0.1 mL	4358297	•				•		•	
Optical Tube without Cap, 0.2 mL	N8010933		•	•					
Seals and covers									
Optical Adhesive Film	4360954, 4311971		•	•	•	•	•	•	•
48-Well Optical Adhesive Film	4375323	•							
Reaction trays									
96-Well Tray/Retainer Set	403081		•						
Fast 48-Well Tray	4375282	•							
96-Well Tray for VeriFlex Blocks	4379983					•			
Accessories									
Splash-Free 96-Well Base	4312063		•	•	•	•	•	•	
96-Well Support Base	4379590		•	•	•	•	•	•	
96-Well Base	N8010531		•	•	•	•	•	•	

* Multiple colors are available.

** 7900HT (0.2 mL and 0.1 mL) instruments are not recommended for use with individual tubes or tube strips.

Note: Experiments using one or two 8-tube strips with attached caps require blank tube strips to balance lid pressure on the block or the use of the MicroAmp 96-Well Tray/Retainer Set (Cat. No. 4381850)—bottom part of the tray *only*. For use with the 96-well block of the 7000, 7300, 7500, and ViiA 7 systems, and the QuantStudio 3, 5, 6, 6 Pro, 7, 7 Pro, and 12K instruments.

Find out more at thermofisher.com/pcrplastics

MSC-3000, Centrifuge/Vortex Multispin

DESCRIPTION

Centrifuge/vortex Multi-Spin **MSC-3000** is product of many years evolution of Spin-Mix-Spin technology that is intended for collecting micro volumes of reagents on the microtube's bottom (first centrifugation spin), following mixing (mix) and collecting the reagents again from the walls and cap of the microtube (second spin). We named this repetitive algorithm of operation that is aimed at reducing the mistakes during sample preparation for PCR analysis a "sms-algorithm".

Multi-Spin is a fully automatic device for reproducing sms-algorithm for 12 tubes at one time, thus saving time considerably. A must-have instrument for PCR and DNA analyses laboratory.

Multi Spin is four devices combined in one:

1. Centrifuge — Maximum RCF: up to $800 \times g$
2. Vortex (3 mixing modes — soft, medium, hard; regulated time; Vortexing regulation timer 1–20 sec)
3. Centrifuge/Vortex;
4. SMS-cycler for realization of the "sms-algorithm".

Table to the right compares FVL-2400N, **MSC-3000** and MSC-6000

Multi-Spin allows considerable time saving compared to Combi-Spin by automatically performing cycling program of sample mixing and spinning according to the set spin-mix-spin cycle for 12 microtubes simultaneously.

SPECIFICATIONS

Speed control range	1000–3500 rpm (increment 100 rpm)
RCF max.	$800 \times g$
Spin timer	1 s – 99 min
Vortexing intensity	Soft, medium, hard
Vortexing time	0–20 s (increment 1 s)
SMS-cycle regulation	1 - 999 cycles
Timer sound signal	+
Display	LCD, 2 x 16 signs
Safety	Autostop at open lid
Overall dimensions (W×D×H)	190x235x125 mm
Weight	2.1 kg
Input current/power consumption	12 V, 11 W (0.9 A)
External power supply	Input AC 100–240 V; 50/60 Hz; Output DC 12 V

CAT. NUMBER

Including rotors R-1.5, R-0.5/0.2	Including rotors R-1.5, R-0.5/0.2
BS-010205-AAN	230VAC 50/60Hz Euro plug
BS-010205-AAK	230VAC 50/60Hz UK plug, 230VAC 50/60Hz AU plug, 100VAC 50/60Hz US plug, 120VAC 60Hz US plug
BS-010205-EK	IQ OQ document
BS-010205-GK	PQ document

R-0.5/0.2
BS-010205-BK
rotor

Rotor for 12 x 0.5 ml and 12 x 0.2 ml microtest tubes

R-1.5
BS-010205-AK
rotor

Rotor for 12 x 1.5/2 ml microtest tubes


R-2/0.5
BS-010205-CK
rotor

Rotor for 8 x 2/1.5 ml and 8 x 0.5 ml microtest tubes

R-2/0.5/0.2
BS-010205-DK
rotor

Rotor for 6x2/1.5 ml + 6x0.5 ml + 6x0.2 ml microtest tubes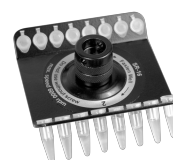

SR-16
BS-010202-AK
rotor

Rotor for 2 x 8-section 0.2 ml microtube strips

SR-32
BS-010205-FK
rotor

Rotor for 4 x 8-section 0.2 ml microtube strips

* not compatible with Combi-Spins produced before 2015

Bio RS-24, Mini-rotator

DESCRIPTION

Mini-rotator **Bio RS-24** provides vertical rotation of the platform. The rotator is an ideal instrument for preventing blood coagulation in tubes and for fulfilment of procedures of biological components extraction.

The device is simple to operate; it is designed as a low cost solution.

Rotator can be used in cold rooms or incubators, operating at ambient temperature range +4°C to +40°C.

PRS - series platforms are equipped with universal rubber clamps for different size tube fixation;
PRSC - series platforms have metal clamps able to hold heavier solutions (e.g. soil, sand)

SPECIFICATIONS

Speed control range	5-30 RPM
Digital time setting	1 min–24 hrs / non–stop
Vertical rotation movement	overhead, 360°
Maximum continuous operation time	8 hrs
Overall dimensions (W×D×H)	325×190×155 mm
Weight	1.4 kg
Input current/power consumption	12 V, 110 mA/1.3 W
External power supply	Input AC 100–240 V 50/60 Hz; Output DC 12 V

CAT. NUMBER

Including PRS-22 platform	Including PRS-22 platform
BS-010133-AAG	230VAC 50/60Hz Euro plug
BS-010133-AAK	230VAC 50/60Hz UK plug, 230VAC 50/60Hz AU plug, 100VAC 50/60Hz US plug, 120VAC 60Hz US plug
BS-010133-BK	IQ OQ document
BS-010133-CK	PQ document

ACCESSORIES

PRS-22
BS-010117-FK
platform

Can accommodate 22 tubes, 10-16 mm diameter (1.5 ml - 15 ml tubes).

PRSC-18
BS-010117-EK
platform

Can accommodate 18 tubes, 16 mm diameter (15 ml).

PRS-4/12
BS-010117-AK
platform

Can accommodate 4 tubes, 20-30 mm diameter (up to 50ml tubes) and 12 tubes, 10-16 mm diameter (1.5 ml - 15 ml tubes).