CERTIFIED COMPANY UNI EN ISO 9001:2008 & UNI CEI EN ISO 13485:2012

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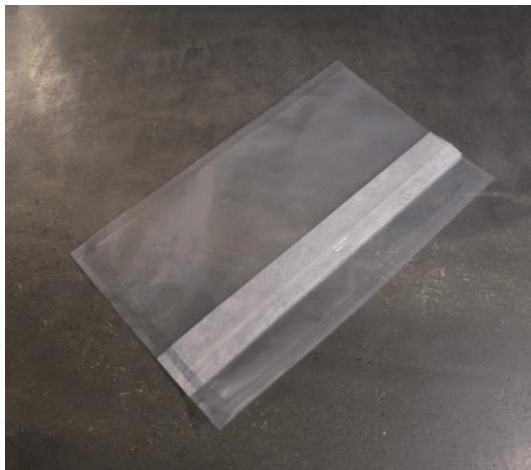
#### **Sterile bags with filter for Stomacher® blenders**

Rigid and transparent polyethylene (PE) bags, sterile R. Bags are tested by various parameters, in detail by opening capacity, sealing at the base and dimensions.

Bags available in two different type:

- > with full surface filter (ideal for pasty sample)
- > with lateral filter (ideal for fibrous sample)

Suitable for foodstuff in accordance with applicable law. Latex free device.

In compliance with ISO 4833-2 new standards

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*Intended purpose: "GENERAL LABORATORY USE". Ideal for sampling, mincing and homogenisation of foodstuff or similar. PRODUCT NOT SUBJECT TO CE MARKING.*

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*With full-surface calibrated microperforated filter for sample homogenization and automatic filtration, particularly suitable for pasty/viscous samples (chocolate, cakes, cheese...).*

*Filter porosity 280 micron. Sterile.*

*Flora is extracted from the sample and goes through the filter during blending. You can then pipet behind the filter with no residue or particles of the sample.*

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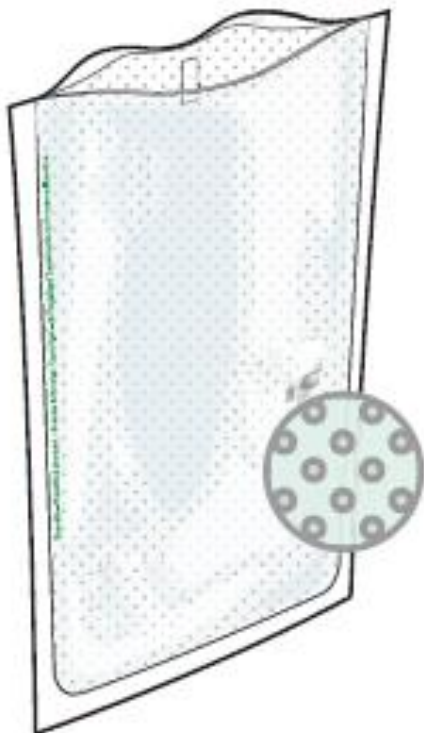
**With internal side non-woven filter** for sample homogenization and automatic filtration, particularly suitable for fibrous samples and for pipetting. Filter porosity inferior to 250 micron. Sterile

Flora is extracted from the sample and goes through the filter during blending. You can then pipet behind the filter with no residue or particles of the sample.

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### Particularly suitable for pipetting

#### INSTANT FILTRATION

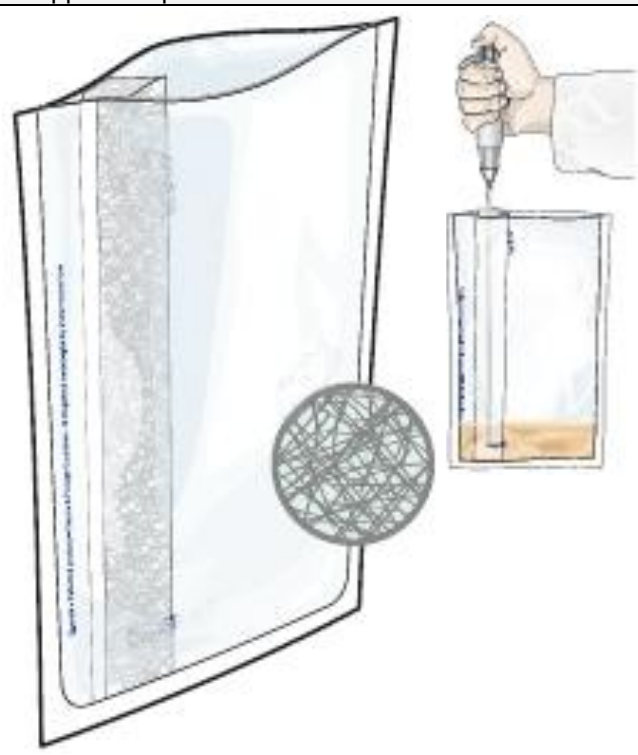
- Filtration during homogenization
- No need to wait for debris sedimentation
- Particle-free solution, no debris on the Petri dish: quick colony counting
- Precise results and excellent reproducibility

#### NO CROSS-CONTAMINATION

- Sterile bags
- Same bag is used to homogenize, filter and pipet
- No contact between the sample and the blender during homogenization

#### USER-FRIENDLY

- Adapted to all samples
- Compatible with any lab blender
- Approved for food contact

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Here is a comparison of a sample\* homogenized in a bag without a filter and in a bag with filter. The bacterial concentration remains the same after filtration

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\* *Lactobacillus analysis in a cheese sample, plated on MRS agar incubated at 30°C for 48h.*

## SACCHETTI IDONEI PER / BAGS ADAPTED FOR

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**Food and agricultural microbiological analyses** (Ex: meat, cooked meat, fish, prepared dishes, fruits, vegetables)

**Pharmacological microbiological analyses** (Ex: pills, tablets, blisters, powders)

**Environmental microbiological analyses** (Ex: plants, soil, filtration membranes, tissues)

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*Keep out of flame or heat sources which might damage the product*

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*Do not use after expiry date or if packing is opened*

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*Do not re-use: Disposable device*

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*Do not vary the intended purpose of the product*

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*Store in dry place, Temperature range: min -10°C max +50°C*

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*Disposal: use appropriate personal protective equipment and act according to applicable regulations*

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*Before use with particular substances, check the resistance / compatibility chart on our catalogue*

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