**Labkem cleaner M66 Alkaline liquid detergent AUX**  
**Detergente Labkem cleaner M66 liquido alcalino AUX**  
**Détergent Labkem cleaner M66 liquide alcalin AUX**

Taric code: 34022090

H315-H319 P260-P280-P302+P352-P332+P313-P305+P351+P338-P337+P313

## SPECIFICATIONS:

*Concentrated alkaline liquid cleaning agent for manual use, soaking and ultrasonic cleaning of glassware and laboratory equipment made of stainless steel, porcelain, ceramic and plastic.*

*LABKEM CLEANER M 66 is biodegradable and presents high penetrating wetting power. It efficiently replaces corrosive acidic mixtures and hazardous solvents.*

*LABKEM CLEANER M 66 is used to remove grease, oils, blood, cell tissue, chemicals, distillation residues and silicon oils.*

### DIRECTIONS FOR USE:

Prepare a 2% solution in cold, warm or hot water (optimal temperature 50°C).

Immerse the items to be cleaned and ensure a mechanical action with a brush.

The use of an ultrasonic cleaner improves the cleaning and reduces contact time.

For difficult residues, increase the temperature and/or the detergent concentration.

Use cold water for blood and proteins residues and gradually raise the temperature.

Rinse promptly two time with deionized water and ensure adequate drying.

LABKEM CLEANER M 66 rinses away easily to give reliable results and no interfering residues.

Avoid contact with soft metals (aluminum, zinc and alloys)

Alkaline detergents are not recommended for cleaning plastics made of polycarbonates or polystyrene.

### APPLICATIONS:

Manual, soaking, Ultrasonic cleaner

### INGREDIENTS:

Alkalis, surfactants, chlorinated agents

### PHYSICO-CHEMICAL PROPERTIES:

Appearance .....	Light yellow
Odor .....	Characteristic smell
pH (pure) .....	12.75 - 13.75
pH (2% v/v) .....	11.20 - 12.20
Density at 20°C .....	1.112 - 1.118 g/cm <sup>3</sup>
Dry residue .....	16.5 - 17.5 %

#### Note:

Material Safety Data Sheets and Certificates of Analysis are downloadable free of charge at [www.labbox.com](http://www.labbox.com). Disclaimer: labkem chemical products are exclusively for laboratory use and by instructed and experienced users. These chemical products may be dangerous, so please follow instructions mentioned on the Material Safety Data Sheets. Due to shipping restrictions or local/national regulations, all products may not be available in all countries or for some particular customers.

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H315-H319 P280-P302+P352-P332+P313-P305+P351+P338-P337+P313

## SPECIFICATIONS:

*Phosphate free concentrated liquid alkaline cleaning agent.*

*Recommended for manual use, soaking and ultrasonic cleaning of glassware and laboratory equipment made of stainless steel, porcelain, ceramic and plastic.*

*LABKEM CLEANER M 67 is used in water analysis and control laboratories.*

*It efficiently removes grease, oils, blood, cell tissue, chemicals and biological residues.*

### DIRECTIONS FOR USE:

Prepare a 2% solution in cold, warm or hot water (optimal temperature 50°C).

Immerse the items to be cleaned and ensure a mechanical action with a brush.

The use of an ultrasonic cleaner improves the cleaning and reduces contact time.

For hard to remove residues, increase the temperature and/or the detergent concentration.

Use cold water for blood and proteins residues and gradually raise the temperature.

Rinse promptly two time with deionized water and ensure adequate drying.

LABKEM CLEANER M 67 rinses away easily to give reliable results and no interfering residues.

Avoid contact with soft metals (aluminum, zinc and alloys)

Alkaline detergents are not recommended for cleaning plastics made of polycarbonates or polystyrene.

### APPLICATIONS:

Manual, soaking, Ultrasonic cleaner

### INGREDIENTS:

Alkalis, surfactants, carbonates

### PHYSICO-CHEMICAL PROPERTIES:

Appearance .....	Light yellow to yellow
Odor .....	Characteristic smell
pH (pure) .....	12.70 - 13.20
pH (2% v/v) .....	11.60 - 12.20
Density at 20°C .....	1.072 - 1.082 g/cm <sup>3</sup>
Dry residue .....	12.0 - 14.0 %

#### Note:

Material Safety Data Sheets and Certificates of Analysis are downloadable free of charge at [www.labbox.com](http://www.labbox.com). Disclaimer: labkem chemical products are exclusively for laboratory use and by instructed and experienced users. These chemical products may be dangerous, so please follow instructions mentioned on the Material Safety Data Sheets. Due to shipping restrictions or local/national regulations, all products may not be available in all countries or for some particular customers.

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H319 P280-P305+P351+P338-P337-P313

## SPECIFICATIONS:

*Neutral Concentrated detergent for gentle cleaning of sensitive glassware and materials.*

*Suitable for manual, soaking and ultrasonic cleaning.*

*LABKEM CLEANER M 68 ensures complete removal of organic soils, oils, blood and biological residues on sensitive glassware and laboratory equipment made of soft metal like aluminium, zinc and alloys.*

### DIRECTIONS FOR USE:

Prepare a 5% solution in warm or hot water (50-60°C).

Immerse the items to be cleaned and ensure a mechanical action with a brush.

The use of an ultrasonic cleaner improves the cleaning and reduces contact time.

Rinse promptly two time with deionized water and ensure adequate drying.

LABKEM CLEANER M 68 rinses away easily to give reliable results and no interfering residues.

Compatible with all materials

### APPLICATIONS:

Manual, soaking, Ultrasonic cleaner

### INGREDIENTS:

Surfactants, complexing agents

### PHYSICO-CHEMICAL PROPERTIES:

Appearance .....	Yellowish translucent
Odor .....	Characteristic smell
pH (pure) .....	6.4
pH (2% v/v) .....	7.7
Density at 20°C .....	1.07 g/cm <sup>3</sup>
Dry residue .....	13.1 %

#### Note:

Material Safety Data Sheets and Certificates of Analysis are downloadable free of charge at [www.labbox.com](http://www.labbox.com). Disclaimer: labkem chemical products are exclusively for laboratory use and by instructed and experienced users. These chemical products may be dangerous, so please follow instructions mentioned on the Material Safety Data Sheets. Due to shipping restrictions or local/national regulations, all products may not be available in all countries or for some particular customers.

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