



PROGARD[®] PACK

Cat. # PROGOTOS2 - PROGOTNP2

Replacement

- Where used
- Milli-Q[®] Integral systems
 - Elix[®] and Elix Advantage systems
 - All Essential range
-

Shipping box content

1 Progard Pack

When to replace the Progard Pack?

For Integral range:

The Progard pack should be replaced when one of the following Alert messages is displayed:

- REPLACE PROGARD AND TANK VENT FILTER IN XX DAYS, or
- REPLACE PROGARD AND TANK VENT FILTER.

For Essential range:

- When the system displays the yellow alert:



- How to proceed
- For the Integral range, the procedure is described below.
 - For the Essential range, the User Manual describes this procedure.
-



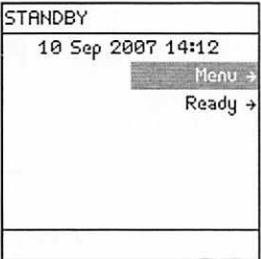
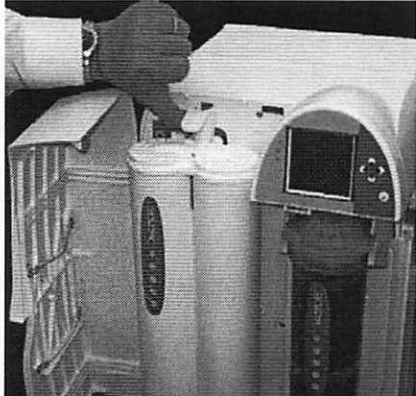
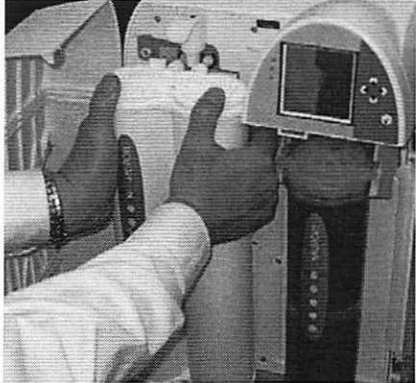
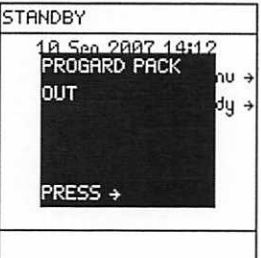
www.ServAPure.com

Order Below MSRP at
Web: www.ServAPure.com
E-mail: servapure@servapure.com
Phone: 800.338.4905 or 989.892.7745

Procedure

Removal

Follow the steps below to remove the Progard pack.

Step	Action	Diagram
1	Place the Milli-Q System into STANDBY Mode.	
2	Depressurize the system: <ul style="list-style-type: none"> • Push the Q-POD plunger all the way down and release it. • Wait until water stops flowing. • Push the Q-POD plunger all the way down and release it. 	
3	<ul style="list-style-type: none"> • Open the Milli-Q System left door. • Lift up the pack locking handle. 	
4	Remove the used Progard pack.	
5	The system indicates that the Progard pack is removed.	

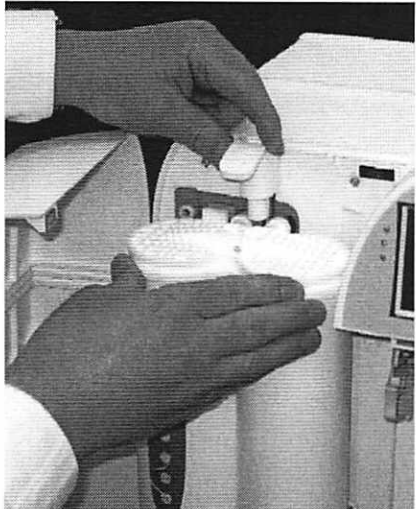
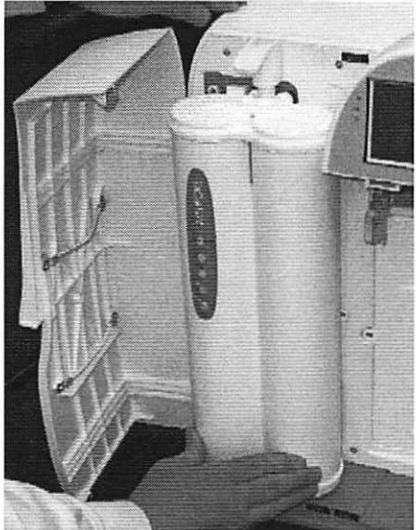
Continued on next page

Procedure, Continued

Installation

Follow the steps below to install the Progard pack.

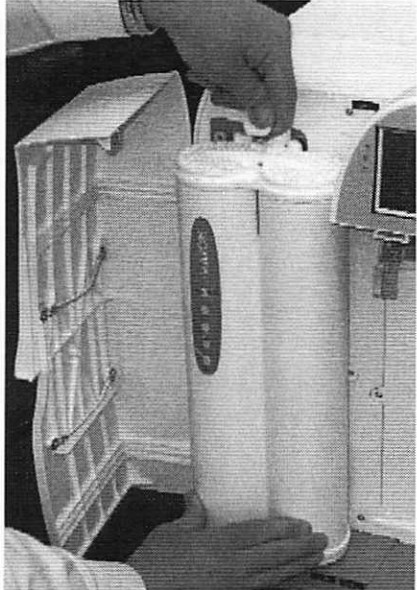

- Remove the covers on the 2 ports of the Progard pack.
- Make sure the rubber O-rings are firmly in place inside these ports.
- Wet the O-rings with water.

Step	Action	Diagram
1	<ul style="list-style-type: none">• Lift the pack locking handle up.• Push the top of the Progard pack into the ports on the Milli-Q system.	
2	Push on the bottom of the Progard pack.	

Continued on next page

Procedure, Continued

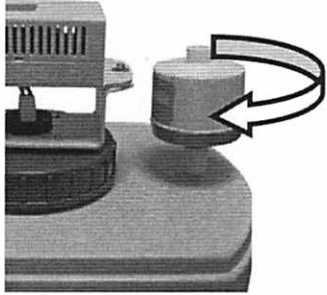
Installation
(continued)

Step	Action	Diagram								
3	<ul style="list-style-type: none"> • Push the pack locking handle down. • Close the left door. 									
4	The system indicates that a new Progard is installed.	<table border="1" style="width: 100%;"> <tr> <td style="text-align: center;">INSTALL PROGARD</td> </tr> <tr> <td>A new Progard has been installed. Catalogue N°: PR060T002 Lot N°: F6DN27324. Press → to start Progard flushing.</td> </tr> <tr> <td> </td> </tr> </table>	INSTALL PROGARD	A new Progard has been installed. Catalogue N°: PR060T002 Lot N°: F6DN27324. Press → to start Progard flushing.						
INSTALL PROGARD										
A new Progard has been installed. Catalogue N°: PR060T002 Lot N°: F6DN27324. Press → to start Progard flushing.										
5	Press  to start the Progard flush.	<table border="1" style="width: 100%;"> <tr> <td style="text-align: center;">INSTALL PROGARD</td> </tr> <tr> <td>Progard Flush procedure in progress. Remaining Time : 15 min. Press → to cancel.</td> </tr> <tr> <td> </td> </tr> </table>	INSTALL PROGARD	Progard Flush procedure in progress. Remaining Time : 15 min. Press → to cancel.						
INSTALL PROGARD										
Progard Flush procedure in progress. Remaining Time : 15 min. Press → to cancel.										
6	When the Progard Pack flush has finished, the Milli-Q System goes to READY Mode.	<table border="1" style="width: 100%;"> <tr> <td style="text-align: center;">READY</td> </tr> <tr> <td style="text-align: center;">10 Sep 2007 14:22</td> </tr> <tr> <td style="text-align: center;">Menu →</td> </tr> <tr> <td style="text-align: center;">Standby →</td> </tr> <tr> <td>Elix R : 12.5 M₂CON TC</td> </tr> <tr> <td>Elix T : 25.1°C</td> </tr> <tr> <td>Tank : 80.0 %</td> </tr> <tr> <td> </td> </tr> </table>	READY	10 Sep 2007 14:22	Menu →	Standby →	Elix R : 12.5 M ₂ CON TC	Elix T : 25.1°C	Tank : 80.0 %	
READY										
10 Sep 2007 14:22										
Menu →										
Standby →										
Elix R : 12.5 M ₂ CON TC										
Elix T : 25.1°C										
Tank : 80.0 %										

Continued on next page

Procedure, Continued

Installation
(continued)

Step	Action	Diagram
7	Unscrew and replace the vent filter.	



Order Below MSRP at
Web: www.ServAPure.com
E-mail: servapure@servapure.com
Phone: 800.338.4905 or 989.892.7745