



## McFarland Standard set

**R092**

McFarland standards are used to perform spectrophotometric comparisons of bacterial densities in water, saline or liquid growth medium. It provides laboratory guidance for the standardization of numbers of bacteria for susceptibility testing or other procedure requiring a standardization of the inoculum like growth promotion test (GPT).

### Set Contains:

R092A (Standard 0.5)- 1 tube

R092B (Standard 1)-1 tube

R092C (Standard 2)- 1 tube

R092D (Standard 3)- 1 tube

R092E (Standard 4)- 1 tube

### Directions

Prepare the inoculum of culture required for testing by using sterile saline. Match the density of the resultant suspension with the density of the desired standard. The standards must be thoroughly mixed on a vortex mixture at the time of use to obtain a uniform suspension. Adjust the density of cell suspension by adding saline if it is more turbid as compared to the desired standard or by adding culture if it is dilute. Check the density of the turbidity by determining the absorbance of 0.5 McFarland standard using a spectrophotometer with a 1 cm light path. The absorbance at 625 nm should be 0.08 to 0.10. The standards should be checked regularly to ensure the density accuracy.

### Interpretation

McFarland standards are a set of tubes with increasing concentration of Barium Sulphate suspension. The turbidity of Barium Sulphate's white precipitation is used as a point of comparison of bacterial suspensions to known bacterial turbidity.

McFarland Standard	0.5	1	2	3	4
Approximate Corresponding suspension x $10^8$ CFU/ml	1.5	3	6	9	12

**Limitation of procedure**

1. Coloured media may interfere with result interpretation and give incorrect results.
2. Bacterial suspensions of older cultures may not be comparable with expected bacterial counts.

**Storage**

Store the standards at 2-8°C, away from light after each use.

**Reference**

1. McFarland, J. 1907. Nephelometer: JAMA 14:1176-1178
2. Murry, PR; Baron, EJ; Jorgensen, JH; Landry, ML; Pfaller, MA; Manual of Clinical Microbiology 9th edition ASM press, Washington DC.

**Disclaimer :**

User must ensure of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ Publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal diagnostic or therapeutic use but for laboratory, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.