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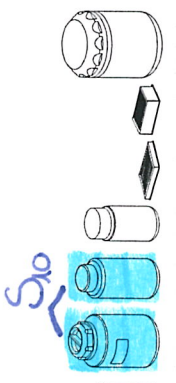
MPW-380 and MPW-380R centrifuges belong to the family of one of the largest, universal laboratory centrifuges. Thanks to the possibility of using a large number of compatible high-speed rotors [RCF above 10,000 x g for tubes up to 250 ml], low-speed horizontal rotors [RCF below 4,700 x g] and their equipment items, they are appreciated by Users of large laboratories, among others, medical, pharmaceutical, biochemical, industrial and guarantee comfort of work under time pressure. Diversity in quantity (up to 100 pieces at a time in the horizontal rotor), as well as capacity [from 0.2 ml to 500 ml in the horizontal rotor] of centrifuged tubes and centrifugation vessels during one centrifugation cycle make centrifuges being used and irreplaceable wherever the highest efficiency, professionalism, high parameters, centrifugation precision and repeatability of centrifugation effects count. One-off centrifugation of large sample volumes (up to 750 ml) definitely facilitates and speeds up the work in parasitological laboratories, in environmental protection when dewatering sewage sludge, and specialized equipment allows to perform works on testing industrial oils (ASTM system) and food oils.

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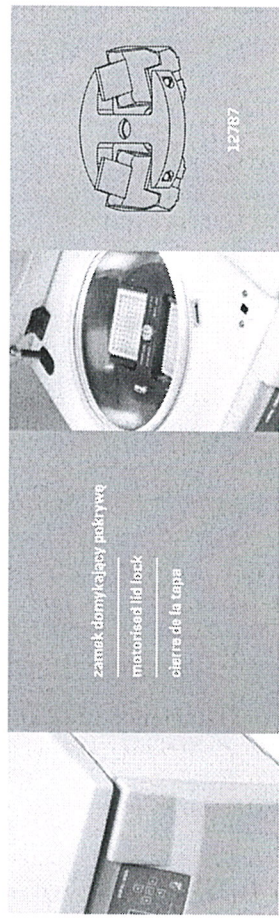
Extras Inopwra pag.3

3. Technical specification

manufacturer	"MPW MED. INSTRUMENTS" SPÓŁDZIELNIA PRACY, Boremlowska 46 Street, 04-347 Warsaw									
type	MPW - 380					MPW – 380R				
cat. number (REF)	10380/2-56	10380/1-56/100	10380/1-56/110	10380/1-56	10380/1-56/127	10380R/2-56	10380R/1-6/100	10380R/1-6/110	10380R/1-6	10380R/1-6/127
mains voltage (L1+N+PE)	230V	100V	110V	120V	127V	230V	100V	110V	120V	127V
	±10%	±5%				±10%	±5%			
mains frequency,	50/60Hz					50/60Hz	60Hz			
power consumption (max.)	800W					1400W				
current protection [A]	10A	12A				18A	20A			
cooling medium	-					R452A				
capacity (max.)	3000 ml									
speed – RPM	90 ÷ 18000 RPM (step 1 RPM)									
force – RCF	31876 x g (step 1 x g)									
kinetic energy (max.)	45000 J									
running time	00:00:01 ÷ 99:59:59 – [hours, min., sec] (step 1s) – S ₄									
time counting	since start button is pressed / since preselected speed is reached									
short-time operation mode – SHORT	yes									
continuous operation mode – HOLD	yes									
menu languages	Polish, English, German, Spanish, Italian, Portuguese, Russian, Swedish, French, Czech									
user programs	100									
adjustable temperature	-					-20 ÷ 40°C* (step 1°C)				
cooling / heating	no / no					yes / no				
guaranteed temperature with max. rotor speed	-					≤4°C				
cooling	no					yes				
cooling without centrifuging	no					yes				
acceleration (ACEL)	10 linear curves									
deceleration (DECEL)	10 linear curves									
programmable non-linear curves:										
acceleration	10									
deceleration	10									
USB communication	no									
Electromagnetic compatibility	according to EN 61326-2-6:2006									
degree of protection (according to PN-EN 60034-5:2021-01)	IP21					IP21				
height (H)	455 mm					455 mm				
width (W)	515 mm					715 mm				
depth (D)	650 mm					650 mm				
height with open cover (H _{oc})	960 mm					960 mm				
noise level	<67 dB – S ₈									
weight 230V	approx. 76 kg					approx. 114 kg				
weight 120V	approx. 81,5 kg					approx. 120 kg				

*time and possibility of obtaining a set temperature is dependent on multiple factors , including rotor type, established RPM, ambient temperature; accuracy: - ±1°C appropriate for place of temperature sensor

3.1 Environmental conditions

- The device may only be used indoors.
- The permissible ambient temperature is 2°C to 40°C.
- Maximum allowed relative humidity 80% at temperature up to 31°C decreasing linearly to 50% relative humidity at 40°C.
- The mains voltage fluctuations must not exceed ± 10% of the nominal voltage.
- Maximum altitude 2,000 m above sea level.
- Overvoltage category II.
- Pollution degree 2.

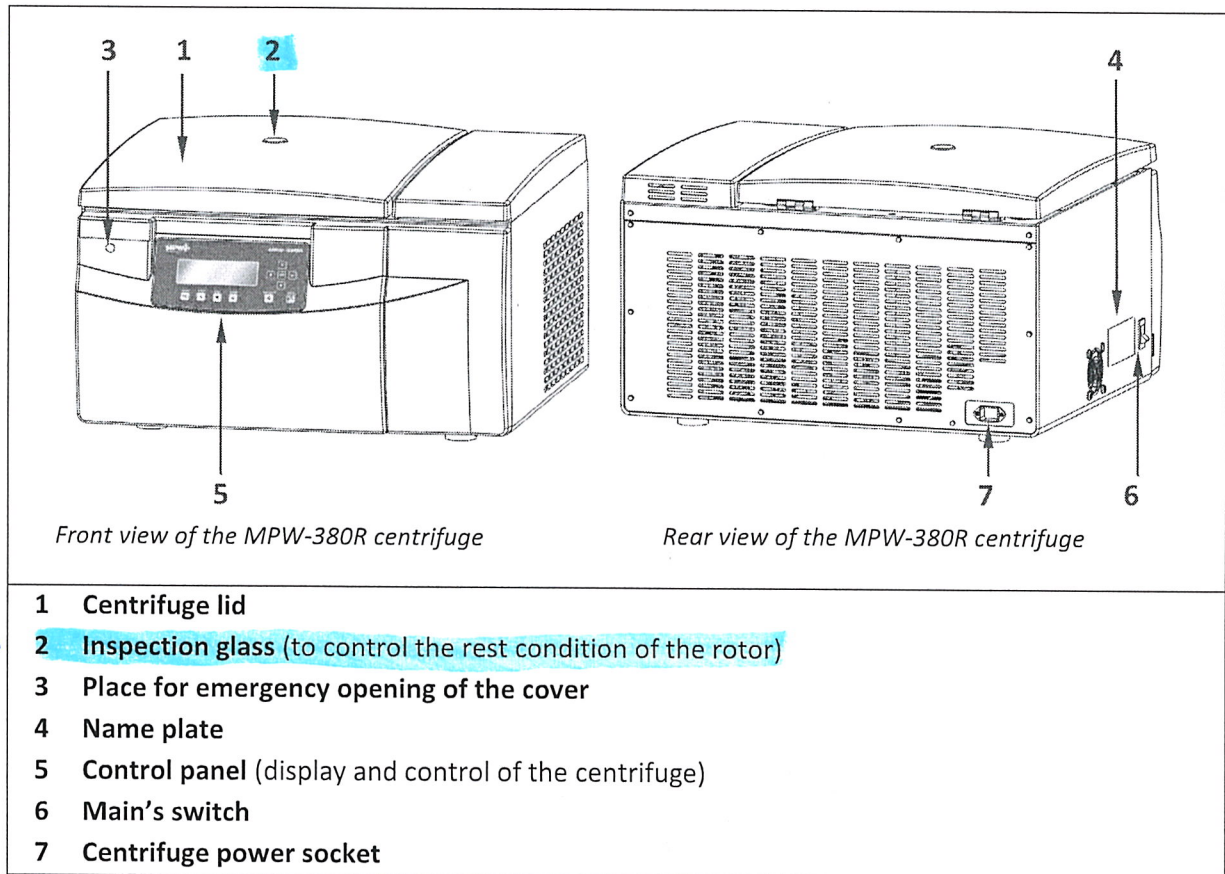
Extras manual de utilizare pag.7

6. Product description

6.1 Product Design and Appearance

A new generation of MPW MED. INSTRUMENTS laboratory centrifuges is provided with state-of-the-art microprocessor control systems, durable and quiet asynchronous brushless motors, and accessories consistent with requirements of the present-day user.

The centrifuge has rigid self-supporting structure. Housing was made of sheet aluminum, back made of steel sheet. Front and cover were made of ABS type plastic. Cover is fixed on steel hinges and from the front it is locked with two electromagnetic locks blocking possibility of opening during centrifugation. Rotation chamber casing was made of thick steel sheet. The spin chamber is made of stainless-steel sheet.

7. Centrifuging

Power switching ON/OFF is carried out with master switch situated on the left side wall of the centrifuge. All settings on the centrifuge are done by means of the control panel.

7.1 Control panel

The control panel placed on the front casing serves the purpose of controlling centrifuge operation.

	SHORT ¹	short-time centrifuging
	START	start centrifugation run
	STOP ²	end centrifugation run
	COVER	cover opening
	FAST COOL	start fast cooling mode (MPW-380R only)
	BACK/ OPTIONS	exit the current menu / enter to submenu of options (keep held down within 1 s.)
	UP	navigation in menu / increasing values
	DOWN	navigation in menu / decreasing values
	LEFT	navigation in menu
	RIGHT	navigation in menu
SET	SET	changing parameters / confirming changes

¹ the centrifuge is working as long as the key is pressed

² first-time pressing press – will make stopping centrifuging with acceleration characteristics set in the current program, second-time pressing – will make the centrifuging as fast as possible (quickest characteristic). During setting of the parameters, it serves for exiting zones on the primary screen without introducing changes.

7.2 Display

The display is located in the centre of the control panel. The main screen variants are presented below. In the user manual exemplary screens from MPW-380R are shown. For MPW-380 (without refrigeration) temperature is not shown. Blinking of field on display mean it is ready to set. Blinking of field is visualised as highlighted in the user manual.

	After switching on centrifuge, welcome screen appears. After disappearing the welcome screen, it is possible to setting up parameters.
--	---

Extras manual de utilizare pag. 20

10.1 Screen saver

Setting time of screen saver	MENU/ CONFIGURATION / SCREEN
SCREEN ↔ 1/5 ☒ SCREENSAVER: 15 min ☐ VISUAL ALARM ☒ NORMAL DISPLAY ☐ SIMPLIFIED DISPLAY	- With ▲ ▼ ◀ ▶ keys select SCREENSAVER. - Press SET and then ▼ and SET. - With ▲ ▼ keys select demanded value from 1 to 60 minutes. - Mark selection by pressing SET. - Leave the menu by pressing BACK.

10.2 Visual alarm - S6

Visual alarm	MENU/CONFIGURATION/ SCREEN
SCREEN ↔ 1/5 ☒ SCREENSAVER: 15 min ☐ VISUAL ALARM ☒ NORMAL DISPLAY ☐ SIMPLIFIED DISPLAY	- Via ▲ ▼ keys choose VISUAL ALARM - Mark it by pressing SET. VISUAL ALARM cause blinking screen after ending of centrifuging or after message occurring.

10.3 Types of main screen

To ensure optimal adaptation to the user's preferences, work is possible in two basic screen modes.

NORMAL DISPLAY - contains an expanded number of parameters visible on the display.

SIMPLIFIED DISPLAY - contains only the most important parameters visible on the display.

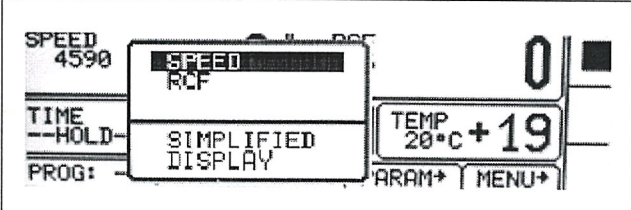
For each of the above modes, you can choose priority RPM display or RCF.

By default, the **SIMPLIFIED DISPLAY** is set

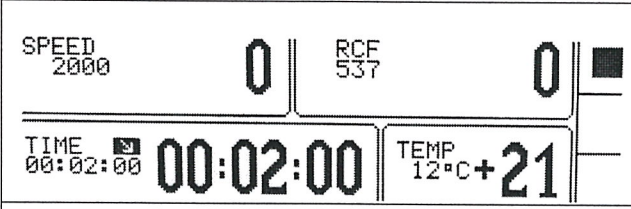
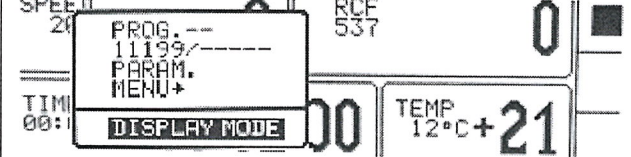
Types of main screen	
NORMAL DISPLAY	SIMPLIFIED DISPLAY
Switch between the SPEED (RPM) and RCF display priority modes	
- In the NORMAL DISPLAY mode, selecting the SPEED or RCF display mode is obtained by pressing and holding BACK for 1 sec. - then use the ▲ ▼ buttons to select the desired mode (SPEED or RCF) and press SET.	- In the SIMPLIFIED DISPLAY mode, the selection of the SPEED or RCF display mode is obtained by pressing and holding the BACK key for 1 second. - then use ▲ ▼ keys to select DISPLAY MODE, press SET, and then use ▲ ▼ keys to select the desired mode (SPEED or RCF) and press SET.

Extras manual de utilizare pag. 37

10.3.1 Switching the normal display to simplified display

	Press the BACK button for 1 sec. then:
	Via ▲▼ keys select SIMPLIFIED DISPLAY. Press SET.

10.3.2 Switching the simplified screen to normal display

	Press the BACK button for 1 sec. then:
	- Via ▲▼ keys select DISPLAY MODE (blinking). - Press SET. - Then choose NORMAL DISPLAY and press SET.

10.4 Rotating runtime

Way of time counting	MENU/CONFIGURATION/ ROTATING RUNTIME
ROTATING RUNTIME ◀▶ 2/5 ☒ COUNTING FROM PRESSING START ☐ COUNTING FROM REACHING SPEED ☐ DESCENDING ☐ ASCENDING	- Via ▲▼ choose demanded option. - Mark it by pressing SET.
Counting from: Pressing start → COUNTING SINCE ROTOR IS IDENTIFIED Reaching speed → COUNTING FROM ASSIGNED SPEED	
Presenting mode: Descending → COUNTING DOWN Ascending → COUNTING UP	

10.5 Buzzer – S6

Switching ON/OFF short audible signals accompanying every pressing of any key. Switching ON/OFF signals after centrifuging.	MENU/ CONFIGURATION /BUZZER
BUZZER ◀▶ 3/5 ☐ KEY TONE ☐ CONTINUOUS ALARM	- With ▲▼ keys select demanded option. - Mark selection by pressing SET.