AE-B Range **Automatic Autoclaves with fractionated vacuum**

Suitable to sterilize: unwrapped and wrapped solids, porous objects, and hollow ware and liquids.

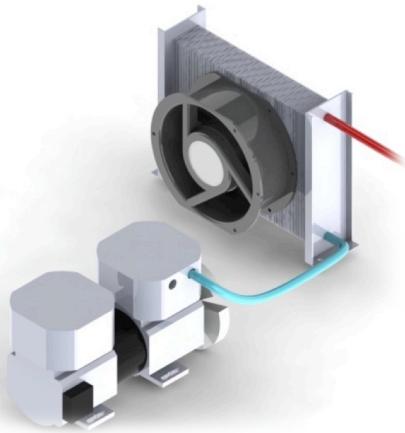
AE-B Range

Operating Principal

The purpose of the sterilization with pre-vacuum is the efficient steam penetration throughout the entire load. This performance is achieved with a vacuum pump, which removes air by fractionated pre-vacuum.

A built-in cooling system by radiator coil enables a better performance of the vacuum pump, removing air pockets from all load types.

Immediate production of steam through the built-in steam generator allows optimization of cycle times.

AE-B Range

Touch Screen Display – Brand new Touch Screen color display for an easy programming and parameters selection for sterilization process.

Main screen


Sterilization screen

Programs – Easy to operate with 12 pre-defined programs

Nº Program	Program name	Nº Pre-vacuum	Temperature (°C)	Sterilization time (min.)	Drying time (min.)
P1	Helix-Bowie&Dick Test	3	134	4'	4'
P2	Vacuum Test	1	-	-	-
P3	Porous-134	3	134	4'	10'
P4	Prion-134	3	134	18'	10'
P5	Porous-121	3	121	20'	10'
P6	Hollowware-134°C	3	134	4'	5'
P7	Hollowware -121°C	3	121	20'	5'
P8	Wrapped-134°C	1	134	4'	5'
P9	Wrapped -121°C	1	121	20'	5'
P10	Solids-134°C	1	134	4'	5'
P11	Solids -121°C	1	121	20'	5'
P12	Flash-134°C	1	134	3'	1'

Up to 50 programs user-defined with the following parameters setting:

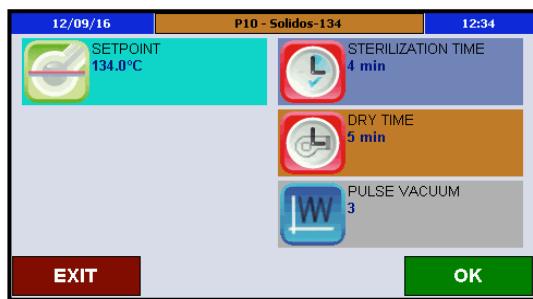- Sterilization temperature
- Sterilization time
- Drying time
- Number of pre-vacuum

Parameters setting screen

AE-B Range

Water Filling – Pure water for sterilization can be added manually into the water pan. Automatic filling is also available by connecting the device to pure mains water.

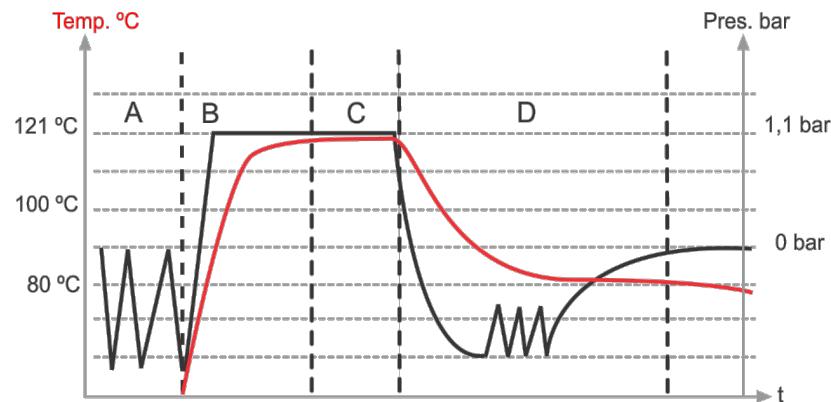
Sterilization Process

Heating Stage – Fast saturated steam production by means of a powerful built-in steam generator.

Pre-vacuum Stage – Vacuum pump with built-in cooling system by radiator coil provides a better vacuum performance.

Sterilization Control – A PT-100 temperature probe controls the chamber temperature, with the possibility of installing another flexible PT-100 temperature probe for liquid programs.

Drying Stage – Powerful heating jacket together with fractionated pre-vacuum provides a fast and intense drying.

AE-B Range

Security – All safety standards are covered with safety valve, safety thermostat and safety pressure switch, door opening prevention system in case of overpressure, open door sensor and water level detector, overheat shutoff. EU Directive and Standards:

2014/35/UE Low Voltage

2014/30/UE Electromagnetic Compatibility

2014/68/UE Pressure Equipment

EN-61010-1

EN-61010-2-040

EN-61326

AD 2000 Merkblatt

Communications and Documentation

USB Port – For electronic storage of process data to a USB flash drive.

Ethernet Port – For live monitoring with direct data storage. It is needed the customized software SOFT-SW8000.

Integrated Printer – For quality printouts after each cycle. Intervals can be adjustable. Printer has to be requested before ordering.

AE-B Range**Technical Data**

Referencia Reference Référence	Volumen total / útil Total volume / usable Volume total / utile	Dimensiones útiles cámara Usable chamber dimensions Dimensions utile chambre	EXTERIOR L x D x H mm Ø x H mm	Temp. °C	Presión máxima Max. pressure Pression max. bar	Potencia Power Puissance	Peso bruto Gross weight Poids brut Kg	Tensión Voltage Tensione	Frecuencia Frequency Fréquence Hz	Depósito Container Réservoir H ₂ O l
AE-50-B	55/50	300 x 710	505 x 580 x 1290	105-134	2,1	3600	110	230	50/60	9
AE-75-B	79/75	400 x 600	610 x 700 x 1185	105-134	2,1	3600	140	230	50/60	12
AE-110-B	115/110	400 x 850	610 x 700 x 1435	105-134	2,1	6000	180	400	50/60	12
AE-150-B	175/153	500 x 780	750 x 820 x 1400	105-134	2,1	6000	265	400	50/60	20

Consultar por otras tensiones / Ask for other voltages / Demander d'autres tensions

Electrical connection for equipment of 50 and 75 L: 230V (1P + N), 50/60 Hz.

Electrical connection for 110 and 150 L equipment: 400V (3P + N), 50/60 Hz.

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Intended use

- To lift heavy materials to put into the autoclave.
- Lift with hoist powerful system with adjustable arm.

Features

- Easy to use.
- Suitable for vertical autoclaves: 110L and 150L, consult for other models.
- Rotation up to 200 degrees.

Safety

- Emergency Stop Button.
- Motor with auto braking system.

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SET AUTOMATIC FILLING

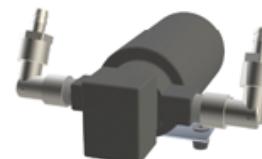
- To be installed in our facilities.

References:
KLL: for B and DRY from 28 up to 150 L
KLL-21: for AH-21B and AH-21L
KLL-MP: for MP

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FLEXIBLE "HEART" PROBE PT-100

- Recommended for sterilizing liquids in flasks of 1 litre or bigger.
- The control is done from the microprocessor.
- This accessory should be installed in factory.

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THERMAL PAPER PRINTER

- It prints program number, cycle number, temperature, date and hour of the run.
- Error messages are also printed.
- To be installed in our facilities.

References:
IT: LCD Screen
IT/TS: Touch Screen

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* Serie AE-B, AES-RFG, AE-MP

PC CONNECTION SW8000

- Communication software between the equipment and the PC.
- Connection for Ethernet.
- USB Adapter.
- Compatible with touch screen circuits.

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* Todos los modelos / All models / Tous les modèles

STERILIZATION CONTROL TAPE

- Chemical indicator for steam. The change of color indicates that the materials have been processed, without being a guarantee of proper sterilization. Class 1.
- To verify that sterilization is effective, additional methods are needed as biological indicators EN ISO 11138.
- Tape roll of 50 m x 19 mm

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* Serie AHS, AH-B, AE-B

BOWIE & DICK TEST PACK

- Class B indicator.
- Check the correct steam penetration in the porous bodies.
- Indicator printed with non-toxic inks and laminated.
- Box of 30 tests.

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* Serie AH-B

HELIX PCD TEST

- Class B indicator.
- Check the correct steam penetration in the close lumen bodies.

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