

BENCHTOP CENTRIFUGE

— ROTINA 380 | 380 R

HIGH SPEED WITH SPACE SAVING DESIGN

High performance unit supporting a wide array of accessories for multiple applications in just one rotor. Its size and performance make it the ideal bench top centrifuge for most clinical, research and industrial laboratories.

This unit is available with refrigeration and a temperature range from -20 °C to +40 °C (ROTINA 380 R).



— HIGHLIGHTS

- RPM: 50 - 15,000 min⁻¹
Adjustable in increments of 10
- RCF: 50 - 24,400
Adjustable in increments of 1
- max. capacity: 4 x 290 ml
- choice of 9 rotors
- IVDR-conform according to regulation (EU) 2017/746
- max. noise level of 54 dB(A) with rotor 1760
- easy operation with keypad and control knob
- 98 program memories for more individuality
- 9 individual acceleration and 10 deceleration stages
- model 380 R coolable from -20 to +40 °C with pre-cooling function

— FEATURES

- metal housing and lid
- viewing port in the lid
- powered one-hand lid lock
- lid dropping protection
- emergency lid lock release
- stainless steel chamber
- automatic rotor recognition
- brushless drive
- error display
- display in °C and °F possible
- imbalance switch-off
- backlit panel with actual values of all parameters
- auditory message after completion of the centrifugation run

TECHNICAL DATA

	ROTINA 380 non-refrigerated	ROTINA 380 R refrigerated
voltage *)	200 – 240 V 1 ~	200 – 240 V 1 ~
frequency	50 – 60 Hz	50 – 60 Hz
consumption	650 VA	1,300 VA
emission, immunity	EN/IEC 61326-1, class B	EN/IEC 61326-1, class B
max. capacity	4 x 290 ml	4 x 290 ml
max. RPM	15,000 min ⁻¹	15,000 min ⁻¹
max. RCF	24,400	24,400
running time	1 – 99 h: 59 min: 59 sec, ∞ continuous run, short cycle mode	1 – 99 h: 59 min: 59 sec, ∞ continuous run, short cycle mode
dimensions (WxDxH)	457x600x418 mm	457x750x418 mm
weight	approx. 51 kg	approx. 81 kg
noise level	54 dB (A) with rotor 1760	54 dB (A) with rotor 1760
temperature control, infinitely variable	-	from -20 to +40 °C
Cat. No.	1701	1706
100 – 127 V 1 ~ / 50 – 60 Hz *)	1701-01	1706-01
consumption	700 VA	1,400 VA
emission, immunity	FCC class B	FCC class B
weight	approx. 58.5 kg	approx. 88.5 kg

*) Other voltages on request.

AVAILABLE ROTORS



SWING-OUT ROTORS		angle	max. RPM	max. capacity	Cat. No.	page
	swing-out rotor, 4-place	90°	5,000 min ⁻¹	4 x 290 ml	1754	3
	swing-out rotor, 4-place	90°	4,000 min ⁻¹	4 x 290 ml	1798	5
	swing-out rotor, 6-place	90°	4,000 min ⁻¹	6 x 50 ml	1726	8
	swing-out rotor, 2-place	90°	4,000 min ⁻¹	10 plates	1760	9
	swing-out rotor, 2-place	90°	5,100 min ⁻¹	12 plates	1770	9
ANGLE ROTORS						
	angle rotor, 6-place	45°	10,000 l 11,000 min ⁻¹	6 x 94 ml	1720	10
	angle rotor, 6-place	45°	10,000 l 11,000 min ⁻¹	6 x 94 ml	1792	11
	angle rotor, 30-place	45°	15,000 min ⁻¹	30 x 2 ml	1789-A	12

SWING-OUT ROTOR, 4-PLACE | 1754

Rotor	
max. RPM max. RCF	5,000 min ⁻¹ 4,863
max. capacity	4 x 290 ml
run-up run-down, braked in sec	42 27
angle max. noise level	90° 60 dB (A)
temperature in °C ¹⁾	0
Cat. No.	1754




Bucket	
lid bioseal ²⁾	1751
Cat. No.	1752

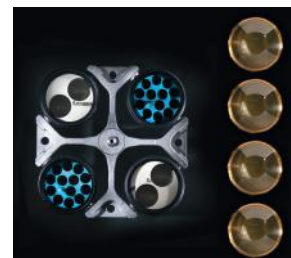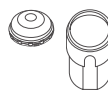

Vessels	
capacity in ml	1.5 2.0 3 4 5 6 7 9 15 20 25 45 50 94 100
Ø x L in mm	11 x 38 11 x 38 10 x 60 12 x 60 12 x 75 12 x 82 12 x 100 14 x 100 17 x 100 21 x 100 24 x 100 31 x 100 34 x 100 38 x 102 40 x 115
max. RCF ²⁾	top / bottom 3,494 / 4,779 3,494 / 4,779 4,779 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,807 4,640
radius in mm	top / bottom 125 / 171 125 / 171 171 167 167 167 167 167 167 167 167 167 167 172 166
Cat. No.	microliter tubes tubes²⁾




Adapter	
boring Ø x L in mm	11.2x39.5 11.2x39.5 11.2x39.5 12.5x56 12.5x56 12.5x56 12.5x56 17.5x62 17.5x62 25.5x81 25.5x81 35.5x81 35.5x81 38.5x80 41x97
vessels per rotor	144 144 72 96 96 96 96 52 52 24 24 12 12 8 4
Cat. No.	1761 1761 1761 1762-A 1762-A 1762-A 1762-A 1763-A 1763-A 1764 1764 1765 1765 1777 1767

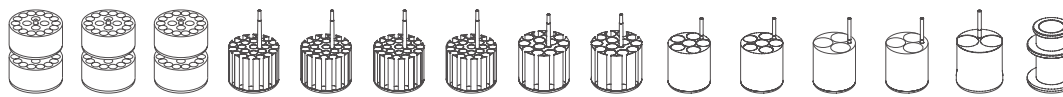

Vessels	
capacity in ml	100 250 1.1–1.4 2.6–3.4 2.7–3 4–5.5 4.5–5 4.5–5 4.9 4.9 7.5–8.5 9–10 10 1.6–5 4–7
Ø x L in mm	44 x 100 65 x 115 8 x 66 13 x 65 11 x 66 15 x 75 11 x 92 11 x 92 13 x 90 13 x 90 15 x 92 16 x 92 15 x 102 13 x 75 13 x 100
max. RCF ²⁾	4,640 4,640 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668 4,668
radius in mm	166 166 167 167 167 167 167 167 167 167 167 167 167 167 167
Cat. No.	tubes²⁾ blood collection tubes / urine tubes




Adapter	
boring Ø x L in mm	45 x 87 66 x 104.5 9 x 49 13.5 x 54 12.5 x 56 17.5 x 62 12.5 x 56 13.2 x 72 13.5 x 54 13.2 x 72 17.5 x 62 17.5 x 62 17.5 x 62 13.5 x 54 13.5 x 54
vessels per rotor	4 4 96 64 96 52 96 76 64 76 52 52 52 64 64
Cat. No.	1766 1768 1781 1783-A 1762-A 1763-A 1762-A 1787 1783-A 1787 1763-A 1763-A 1763-A 1783-A 1783-A

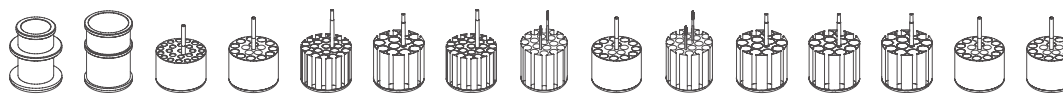



1) For cooled versions: Lowest temperature achievable with precooling and max. speed.
2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.
The max. RCF for glass tubes annotated with footnote 2) is 4,000.
5) Tested by TÜV in conformity with DIN EN 61010, section 2 - 020.


SWING-OUT ROTOR, 4-PLACE | 1754

	
Rotor	
max. RPM max. RCF	5,000 min ⁻¹ 4,863
max. capacity	4 x 290 ml
run-up run-down, braked in sec.	42 27
angle max. noise level	90° 60 dB (A)
temperature in °C ¹⁾	0
Cat. No.	1754

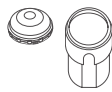	
Bucket	
lid bioseal ⁵⁾	1751
Cat. No.	1752

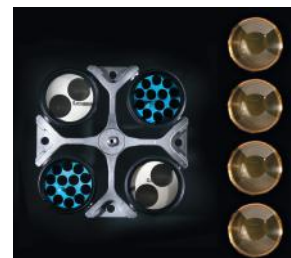

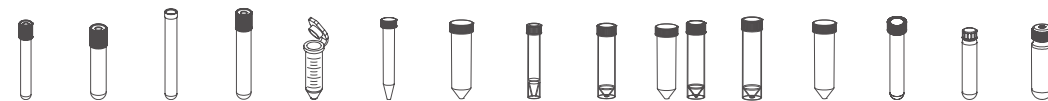	
Vessels	
capacity in ml	4 – 7 4 – 7 8 8.5 – 10 5 15 50 12 25 30 50 50 14 10 30
Ø x L in mm	13x100 16x75 16x125 16x100 17x59 17x120 29 x 115 17x100 25x90 25x110 29x115 29x115 16,5x106 16x80 26x95
max. RCF ²⁾	4,668 4,668 4,668 4,668 4,863 4,863 4,863 4,696 4,528 4,528 4,752 4,752 4,668 4,668 4,807
radius in mm	167 167 167 167 174 174 174 168 162 162 170 170 167 167 172
Cat. No.	blood collection tubes / urine tubes - tubes with screw cap

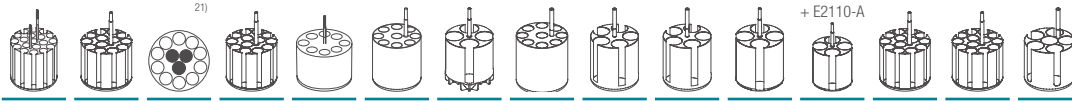	
Adapter	
boring Ø x L in mm	13.2x72 17.5x62 17.5x62 17.5x62 17x52 17x84 30x84 17x78 26.5x72 26.5x72 30x80 30x80 17.5x62 17.5x62 26.5x70
vessels per rotor	76 52 12 52 32 36 16 36 20 20 16 16 52 52 24
Cat. No.	1787 1763-A 1763-A 1763-A 1738 1771-A 1772-A 1773 1779 1779 1774-A 1774-A 1763-A 1763-A 1775

	
Vessels	
capacity in ml	50 85 94 175 175 200 225 250 290
Ø x L in mm	29x107 38x106 38x102 61 x 118 62 x 144 60 x 130 61 x 137 61 x 122 62 x 137
max. RCF ²⁾	4,752 4,807 4,807 4,863 4,863 4,863 4,863 4,863 4,863
radius in mm	170 172 172 174 174 174 174 174 174
Cat. No.	tubes with screw cap Falcon Nalgene Nunc® Falcon 5127 ²⁴⁾ - ²⁴⁾

	
Adapter	
boring Ø x L in mm	30 x 80 38.5 x 80 38.5 x 80 62 x 111 62 x 111 62 x 111 62 x 111 62 x 100 62 x 100
vessels per rotor	16 8 8 4 4 4 4 4 4
Cat. No.	1774-A 1777 1777 1782 1778 1778 1782 1769 1769

1) For cooled versions: Lowest temperature achievable with precooling and max. speed.

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

3.3) When using these tubes, bucket 1752 or 5051 cannot be closed with lid 1751 or 5053.


5) Tested by the TÜV in conformity with DIN EN 61010, section 2 - 020.

21) Adapter must be loaded as illustrated.



24) At temperatures of over +40 °C and / or when not filled to capacity, bottles may warp during centrifugation.

SWING-OUT ROTOR, 4-PLACE | 1798


Rotor	
max. RPM max. RCF	4,000 min ⁻¹ 2,898
max. capacity	4 x 100 ml
run-up run-down, braked in sec	24 17
angle max. noise level	90° 55 dB (A)
temperature in °C ¹⁾	-8
Cat. No.	1798

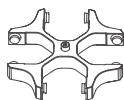


Bucket	
lid	5053
Cat. No.	5051

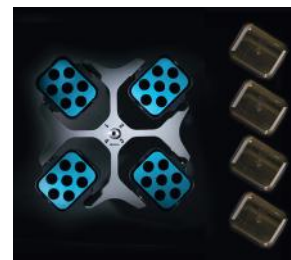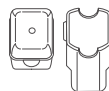

Vessels	
capacity in ml	1.5 2 1.5 2 3 5 6 7 9 15 25 50 100 100
Ø x L in mm	11x38 11x38 11x38 11x38 10x60 12x75 12x82 12x100 14x100 17x100 24x100 34x100 40x115 44x100
max. RCF ²⁾ top / bottom	1,950 / 2,826 1,950 / 2,826 2,826 2,826 2,737 2,773 2,773 2,755 2,773 2,755 2,755 2,755 2,755 2,755
radius in mm	109 / 158 109 / 158 158 158 153 155 155 154 155 154 154 154 154 154
Cat. No.	microliter tubes tubes³⁾

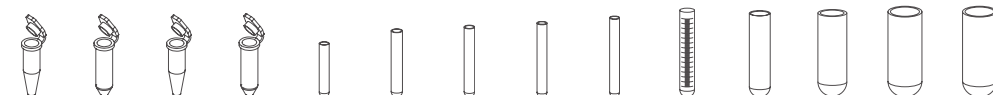


Adapter	
boring Ø x L in mm	12.5x42 12.5x42 11.5x50 11.5x50 11 x 44 12.5x42 12.5x42 12.5x86 16x50 17.5x86 26x86 36x86 42x86 45.5x86
vessels per rotor	160 160 64 64 80 80 80 80 48 48 20 8 4 4
Cat. No.	5257 5257 5281 5281 5267 5227 5227 5247¹⁵⁾ 5264 5248¹⁵⁾ 5242 5243 5249 5262

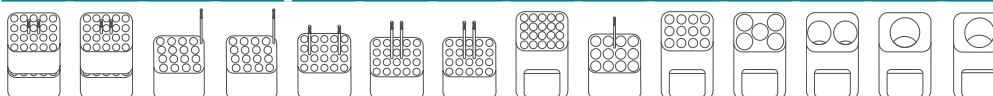


Vessels	
capacity in ml	1.1–1.4 2.6–3.4 2.7–3 4–5.5 4.5–5 4.9 7.5–8.5 9–10 10 1.6–5 4–7 4–7 8.5–10
Ø x L in mm	8x66 13x65 11 x 66 15 x 75 11 x 92 13x90 15x92 16x92 15x102 13x75 13x100 16x75 16x100
max. RCF ²⁾	2,737 2,808 2,773 2,773 2,773 2,808 2,773 2,755 2,755 2,808 2,808 2,773 2,755
radius in mm	153 157 155 155 155 157 155 154 154 157 157 155 154
Cat. No.	blood collection tubes / urine tubes



Adapter	
boring Ø x L in mm	11 x 44 13.5x52 12.5x42 16x50 12.5x42 13.5x52 16x50 17.6x86 17.6x86 13.5x52 13.5x52 16x50 17.5x86
vessels per rotor	80 48 80 48 80 48 48 44 44 48 48 48 48
Cat. No.	5267 5268 5227 5264 5227 5268 5264 5258 5258 5268 5268 5264 5248

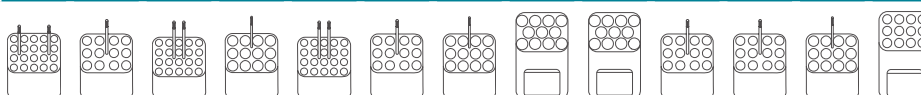

Vessels	
capacity in ml	30 50 15 12 30 50 1–8
Ø x L in mm	25x110 29x115 17x120 17x100 25x110 29x115 simple / multiple
max. RCF ²⁾ top / bottom	2,755 2,826 2,898 2,898 2,755 2,755 1,735 / 2,737
radius in mm	154 158 162 162 154 154 97 / 153
Cat. No.	tubes with screw cap cyto chambers

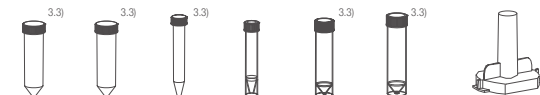


Adapter	
boring Ø x L in mm	26x86 30x86 17x90 17x90 26 x 86 36x86 -
vessels per rotor	20 8 28 28 20 8 8
Cat. No.	5266 5259 6306 6306 5266 5243 5280





- 1) For cooled versions: Lowest temperature achievable with precooling and max. speed.
- 2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.
- 3.3) When using these tubes, bucket 1752 or 5051 cannot be closed with lid 1751 or 5053.
- 15) Also available with decanting aid. (Cat. No. 5247-91 or 5248-91).


SWING-OUT ROTOR, 4-PLACE | 1798

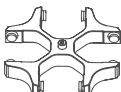	
Rotor	
max. RPM max. RCF	4,000 min ⁻¹ 3,095
max. capacity	4 x 290 ml
run-up run-down, braked in sec	24 17
angle max. noise level	90° 56 dB (A)
temperature in °C ¹⁾	-8
Cat. No.	1798


	
Bucket	
lid bioseal ⁵⁾	5093
Cat. No.	5092

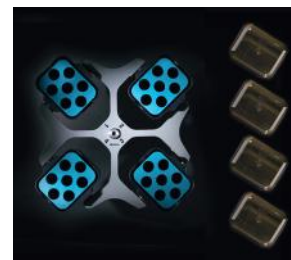


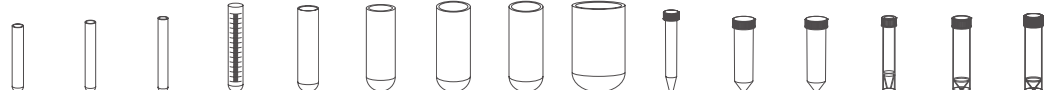
	
Vessels	
capacity in ml	5 6 7 15 25 50 100 100 250 15 30 50 12 25 30
Ø x L in mm	12x75 12x82 12x100 17x100 24x100 34x100 44x100 40x115 65x115 17x120 25x110 29x115 17x100 25x90 25x110
max. RCF ²⁾	3,005 3,005 3,005 2,952 2,898 2,952 2,952 2,952 3,095 3,095 2,898 3,095 3,005 2,826 2,898
radius in mm	168 168 168 165 162 165 165 165 173 173 162 173 168 158 162
Cat. No.	tubes ²⁾ tubes with screw cap

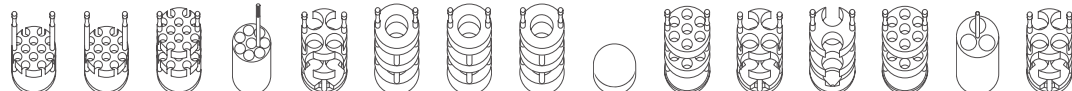	
Adapter	
boring Ø x L in mm	12.8x45 12.8x45 13.2x82 17.5x60.7 25.5x82 35.5x85 45.5x85 42x85 66x103 17x85 25.5x82 30x85 17.5x82 26x73 25.5x82
vessels per rotor	48 48 48 32 16 4 4 4 4 28 16 8 28 12 16
Cat. No.	5128 5128 5120 5136 5122 5124 5125 5126 1791 5129 5122 5123 5121 5134 5122

	
Vessels	
capacity in ml	50 10 250 290 1.1–1.4 2.6–3.4 2.7–3 4–5.5 4.5–5 4.9 7.5–8.5 9–10 10 1.6–5 4–7
Ø x L in mm	29x115 16x80 61x122 62x137 8x66 13x65 11x66 15x75 11x92 13x90 15x92 16x92 15x102 13x75 16x75
max. RCF ²⁾	3,023 2,952 3,095 3,095 2,540 2,540 2,540 2,952 3,005 2,952 2,952 2,952 2,952 2,540 2,952
radius in mm	169 165 173 173 142 142 142 165 168 165 165 165 165 142 165
Cat. No.	tubes with screw cap 5127 ²⁴⁾ - ²⁴⁾ blood collection tubes / urine tubes

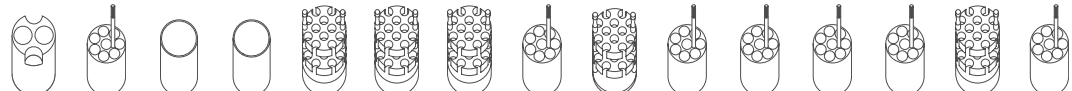	
Adapter	
boring Ø x L in mm	30x99 17.5x60.7 66x103 66x103 13.2x54.5 13.2x54.5 13.2x54.5 17.5x60.7 13.2x82 13.5x60.7 17.5x60.7 17.5x60.7 17.5x60.7 13.2x54.5 17.5x60.7
vessels per rotor	8 32 4 4 48 48 48 32 48 32 32 32 32 48 32
Cat. No.	5135 5136 6319 6319 5138 5138 5138 5136 5120 5137 5136 5136 5136 5138 5136

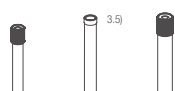	
Vessels	
capacity in ml	4–7 8 8.5–10
Ø x L in mm	13x100 16x125 16x100
max. RCF ²⁾	3,005 3,059 2,952
radius in mm	168 171 165
Cat. No.	blood collection / urine tubes

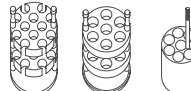	
Adapter	
boring Ø x L in mm	13.2x82 17.5x82 17.5x60.7
vessels per rotor	48 28 32
Cat. No.	5120 5121 ⁴⁾ 5136

- For cooled versions: Lowest temperature achievable with precooling and max. speed.
- Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.
The max. RCF for glass tubes annotated with footnote 2) is 4,000.
- When using these tubes, bucket 5092 cannot be closed with lid 5053.
- Please remove the spacer.
- Tested by the TÜV in conformity with DIN EN 61010, section 2 - 020.
- At temperatures of over +40 °C and / or when not filled to capacity, bottles may warp during centrifugation.

SWING-OUT ROTOR, 6-PLACE | 1726



Rotor

max. RPM max. RCF	4,000 min ⁻¹ 2,808
max. capacity	6 x 50 ml
run-up run-down, braked in sec	19 18
angle temperature in °C ¹⁾	90° -6
Cat. No.	1726



Vessels

capacity in ml	5	5	6	7	9	15	20	25	45	50	8.5 x 10	30	1-8
Ø x L in mm	12x75	12x75	12x82	12x100	14x100	17x100	21x100	24x100	31x100	34x100	16x100	26x95	simple / multiple
max. RCF ²⁾	2,576	2,522	2,665	2,665	2,665	2,665	2,719	2,808	2,719	2,808	2,665	2,808	2,003
radius in mm	144	141	149	149	149	149	152	157	152	157	149	157	112
Cat. No.	tubes²⁾											-	cyto chambers



Carrier

boring Ø x L in mm	12.5x64.4	13.5x65	12.5x71.5	12.5x71.5	14.6x74	17.6x71.5	21.5x74	26x78	32x74	35x78	17.6x71.5	26x78	-
vessels per rotor	24	102	24	24	30	24	12	12	6	6	24	12	6
Cat. No.	1369-91	1372	1369-92	1369-92	1370	1369	1346	1745	1345	1746	1369	1745	1660

CYTO

Cyto system available for this model.
More information on [page 178](#)

SWING-OUT ROTOR, 6-PLACE | 1726



Rotor

max. RPM max. RCF	4,000 min ⁻¹ 2,808
max. capacity	42 x 15 ml
run-up run-down, braked in sec	19 18
angle max. noise level	90° 54 dB (A)
temperature in °C ¹⁾	-6
Cat. No.	1726



Carrier

Cat. No.	1742
-----------------	-------------



Vessels

capacity in ml	15	2.6-3.4	4-5.5	7.5-8.5	9-10	10	1.6-5	4-7	8.5-10	15
Ø x L in mm	17x100	13x65	15x75	15x92	16x92	15x102	13x75	16x75	16x100	17x120
max. RCF ²⁾	2,808	2,683	2,683	2,808	2,808	2,808	2,683	2,683	2,808	2,808
radius in mm	157	150	150	157	157	157	150	150	157	157
Cat. No.	-²⁾	blood collection tubes / urine tubes								-



Spacer

boring Ø x L in mm	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78	17.6 x 78
vessels per rotor	42	42	42	42	42	18	42	42	42	18
Cat. No.	-	0716	0716	-	-	-	0716	0716	-	-

1) For cooled versions: Lowest temperature achievable with precooling and max. speed.


2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.
The max. RCF for glass tubes annotated with footnote 2) is 4,000.

21) Adapter must be loaded as illustrated.

SWING-OUT ROTOR, 6-PLACE | 1726

			
Rotor		Bucket	
max. RPM max. RCF	4,000 min ⁻¹ 2,325	Cat. No.	1366
max. capacity	72 x 4 ml		
run-up run-down, braked in sec.	19 18		
angle max. noise level	90° 54 dB (A)		
temperature in °C ¹⁾	-6		
Cat. No.	1726		

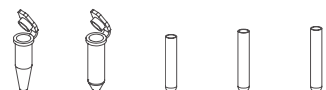



	
Vessels	
capacity in ml	1.5 2.0 1 3 4
Ø x L in mm	11 x 38 11 x 38 6 x 45 10 x 60 12 x 60
max. RCF ²⁾	2,325 2,325 2,308 2,290 2,290
radius in mm	130 130 129 128 128
Cat. No.	microliter tubes Rhesus tubes ²⁾


	
Adapter	
boring Ø x L in mm	11.5 x 38 11.5 x 38 6.5 x 23 10.5 x 23 12.5 x 44
vessels per rotor	54 54 180 72 72
Cat. No.	5277 5277 1357 1327 1326

SWING-OUT ROTOR, 6-PLACE | 1726

			
Rotor		Carrier	
max. RPM max. RCF	4,000 min ⁻¹ 2,808	Cat. No.	1741
max. capacity	60 x 9 ml		
run-up run-down, braked in sec.	19 18		
angle max. noise level	90° 54 dB (A)		
temperature in °C ¹⁾	-6		
Cat. No.	1726		

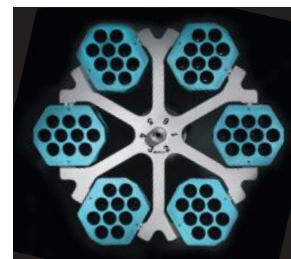

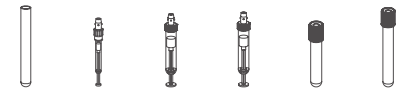	
Vessels	
capacity in ml	9 1.1 – 1.4 4.5 – 5 4.9 1.6 – 5 4 – 7
Ø x L in mm	14 x 100 8 x 66 11 x 92 13 x 90 13 x 75 13 x 100
max. RCF ²⁾	2,773 2,773 2,808 2,808 2,486 2,808
radius in mm	155 155 157 157 139 157
Cat. No.	- ²⁾ blood collection tubes / urine tubes

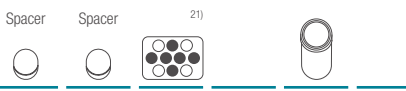
	
Adapter	
boring Ø x L in mm	14.6 x 78 14.6 x 78 14.6 x 78 14.6 x 78 14.6 x 78 14.6 x 78
vessels per rotor	60 60 36 60 60 60
Cat. No.	0701 0701 - - 6x0768 ¹⁶⁾ -

- For cooled versions: Lowest temperature achievable with precooling and max. speed.
- Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.
The max. RCF for glass tubes annotated with footnote 2) is 4,000.
- Packaging unit 10 pieces.
- Adapter must be loaded as illustrated.

SWING-OUT ROTOR, 2-PLACE | 1760

Rotor

max. RPM max. RCF	4,000 min ⁻¹ 2,397
max. capacity	10 plates
run-up run-down, braked in sec	30 23
angle max. noise level	90° 54 dB (A)
temperature in °C ¹⁾	-8
Cat. No.	1760

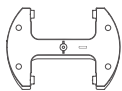

Plates

W x D x H in mm	128x86x15	128x86x17.5	128x86x22	128x86x44.5	128x86x46	128x86x83	84x59x11	124x82x20	-
capacity in ml	-	-	-	-	-	-	-	-	0.2
max. RCF ²⁾	2,397	2,397	2,397	2,397	2,397	2,397	2,397	2,397	2,397
radius in mm	134	134	134	134	134	134	134	134	134
Cat. No.	MTP	MTP	CP	DWP	MS	QP	Microtest plates	PCR plate	PCR strips




Bucket

boring Ø x L in mm	-	-	-	-	-	-	-	-	-
plates / strips per rotor	10	8	6	2	2	2	4	2	24 x 8
Cat. No.	1753-A	1753-A	1753-A	1753-A	1753-A	1753-A	1753-A	1753-A+1485	1753-A+1485

SWING-OUT ROTOR, 2-PLACE | 1770

Rotor

max. RPM max. RCF	5,100 min ⁻¹ 3,926
capacity in ml	12 plates
run-up run-down, braked in sec	65 30
angle temperature in °C ¹⁾	90° -3
Cat. No.	1770

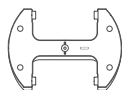

Bucket

lid bioseal ²⁾	4627
Cat. No.	4745



Plates

W x D x H in mm	128x86x15	128x86x17.5	128x86x22	128x86x44.5	128x86x46	128x86x83	84x59x11	124x82x20	-
Ø x L in mm	-	-	-	-	-	-	-	-	0,2
max. RCF ²⁾	3,926	3,926	3,926	3,926	3,926	3,926	3,926	3,926	3,926
radius in mm	135	135	135	135	135	135	135	135	135
Cat. No.	MTP	MTP	CP	DWP	MS	QP	Microtest plates	PCR plate	PCR strips



Removal frame

boring Ø x L in mm	-	-	-	-	-	-	-	-	-
plates / strips per rotor	12	10	8	2	2	2	4	2	24 x 8
Cat. No.	4626	4626	4626	4626	4626	4626	4626	4626 + 1485	4626 + 1485

1) For cooled versions: Lowest temperature achievable with precooling and max. speed.

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

ANGLE ROTOR, 6-PLACE | 1720



Rotor

max. RPM	ROTINA 380 380 R	10,000 min ⁻¹ 11,000 min ⁻¹
max. RCF		13,528 16,369
max. capacity		6 x 94 ml
run-up / run-down, braked in sec		39 / 36 45 / 44
angle max. noise level		45° 59 dB (A)
temperature in °C ¹⁾		+1
Cat. No.		1720



Vessels

capacity in ml		1.5	2.0	15	25	50	94	7.5 – 8.5	9 – 10	8.5 – 10	5	15	50	10	30	50
Ø x L in mm		11 x 38	11 x 38	17 x 100	24 x 100	34 x 100	38 x 102	15 x 92	16 x 92	16 x 100	17 x 59	17 x 120	29 x 115	16 x 80	26 x 95	29 x 107
max. RCF ²⁾	ROTINA 380	12,969	12,969	12,745	12,410	13,304	13,528	12,745	12,745	12,745	12,857	13,081	12,745	12,410	12,410	12,969
max. RCF ²⁾	ROTINA 380 R	15,692	15,692	15,422	15,016	16,098	16,369	15,422	15,422	15,422	15,557	15,828	15,422	15,016	15,016	15,692
radius in mm		116	116	114	111	119	121	114	114	114	115	117	114	111	111	116
Cat. No.		microliter tubes		tubes ²⁾				blood collection tubes /			-	tubes with screw cap				



Adapter

boring Ø x L in mm	11.4 x 39	11.4 x 39	17.5 x 92	26 x 85	35 x 89	38.4 x 89	17.5 x 92	17.5 x 92	17.5 x 92	17 x 51	17 x 106	29.8 x 97	16.5 x 74	26 x 85	29 x 92
vessels per rotor	24	24	6	6	6	6	6	6	6	6	6	6	12	6	6
Cat. No.	1449	1449	1451	1447	1463	-	1451	1451	1451	1476	1466	1454	1448	1447	1446

Vessels

capacity in ml	85	94
Ø x L in mm	38 x 106	38 x 102
max. RCF ²⁾	ROTINA 380	13,528
max. RCF ²⁾	ROTINA 380 R	16,369
radius in mm	121	121
Cat. No.	tubes with screw cap	



Adapter

boring Ø x L in mm	38.4 x 89	38.4 x 89
vessels per rotor	6	6
Cat. No.	-	-

1) For cooled versions: Lowest temperature achievable with precooling and max. speed.

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

ANGLE ROTOR, 6-PLACE | 1792

			+		
Rotor					
max. RPM	ROTINA 380 380 R	10,000 min ⁻¹ 11,000 min ⁻¹		Cat. No.	INCLUSIVE
max. RCF		13,640 16,504			
max. capacity		6 x 94 ml			
run-up run-down, braked in sec		40 / 37 48 / 44			
angle max. noise level		45° 59 dB (A)			
temperature in °C ¹⁾		+4			
Cat. No.		1792			



																
Vessels																
capacity in ml		1.5	2.0	15	25	50	94	7.5 – 8.5	9 – 10	10	8.5 – 10	5	15	50	10	30
Ø x L in mm		11 x 38	11 x 38	17 x 100	24 x 100	34 x 100	38 x 102	15 x 92	16 x 92	15 x 102	16 x 100	17 x 59	17 x 120	29 x 115	16 x 80	26 x 95
max. RCF ²⁾		ROTINA 380	13,081	12,857	12,522	13,416	13,640	12,857	12,857	12,857	12,857	12,745	13,081	13,304	12,857	12,522
max. RCF ²⁾		ROTINA 380 R	15,828	15,557	15,151	16,233	16,504	15,557	15,557	15,557	15,557	15,422	15,828	16,098	15,557	15,151
radius in mm		117	117	115	112	120	122	115	115	115	115	114	117	119	115	112
Cat. No.		microliter tubes			tubes ²⁾			blood collection tubes / urine tubes				-		tubes with screw cap		
																
Adapter																
boring Ø x L in mm		11.4x39	11.4x39	17.5x92	26x85	35x89	38.2x89.6	17.5x92	17.5x92	17.5x92	17.5x92	17x51	17x106	29.8x97	16.5x74	26x85
vessels per rotor		24	24	6	6	6	6	6	6	6	6	6	6	6	12	6
Cat. No.		1449	1449	1451	1447	1463	-	1451	1451	1451	1451	1476	1466	1454	1448	1447

			
Vessels			
capacity in ml		50	85
Ø x L in mm		29 x 107	38 x 106
max. RCF ²⁾	ROTINA 380	13,081	13,640
max. RCF ²⁾	ROTINA 380 R	15,828	16,504
radius in mm		117	122
Cat. No.		tubes with screw cap	

			
Adapter			
boring Ø x L in mm		29 x 92	38.2 x 89.6
vessels per rotor		6	6
Cat. No.		1446	-

1) For cooled versions: Lowest temperature achievable with precooling and max. speed.

2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers. The max. RCF for glass tubes annotated with footnote 2) is 4,000.

5) Tested by TÜV in conformity with DIN EN 61010, section 2 - 020.

ANGLE ROTOR, 30-PLACE | 1789-A

		+	Lid bioseal⁵⁾, phenol-resistant 	INCLUSIVE
Rotor				
max. RPM max. RCF	15,000 min ⁻¹ 24,400 min ⁻¹		Cat. No.	
max. capacity	30 x 2 ml			
run-up run-down, braked in sec	24 24			
angle max. noise level	45° 60 dB (A)			
temperature in °C ¹⁾	+4			
Cat. No.	1789-A			



							
Vessels							
capacity in ml	0.2	0.4	0.5	0.8	1.5	2	0.5
Ø x L in mm	6x18	6x45	8x30	8x45	11 x38	11x38	10.7x36
max. RCF ²⁾	24,400	24,400	24,400	24,400	24,400	24,400	23,394
radius in mm	97	97	97	97	97	97	93
Cat. No.	microliter tubes						Pediatric

								
	Adapter							
	boring Ø x L in mm	6x40	6x40	8x40	8x40	10.2x19.3	11.2x40.9	11.2x40.9
	vessels per rotor	30	30	30	30	30	30	15
Cat. No.	2024	2024	2023	2023	2031 ⁷⁾	-	0788 ¹⁷⁾	

1) For cooled versions: Lowest temperature achievable with precooling and max. speed.
2) Please note that the RCF values indicated refer only to rotor performance. The max. permissible RCF of tubes used should be verified with the individual manufacturers.
The max. RCF for glass tubes annotated with footnote 2) is 4,000.
5) Tested by the TÜV in conformity with DIN EN 61010, section 2 - 020.
7) For centrifugation at high speeds, we recommend to use form-fitting, phenol-resistant adapters. Cat. No. 2031.
17) Packaging unit 15 pieces.

CERTIFICATIONS / REGISTRATIONS

Product certification:



Product registration:



according to regulation (EU) 2017/746

Company certifications:





DOWNLOADS

[↓ Operating manual – ROTINA 380 | 380 R](#)

[↓ General Catalog](#)

DIMENSIONS

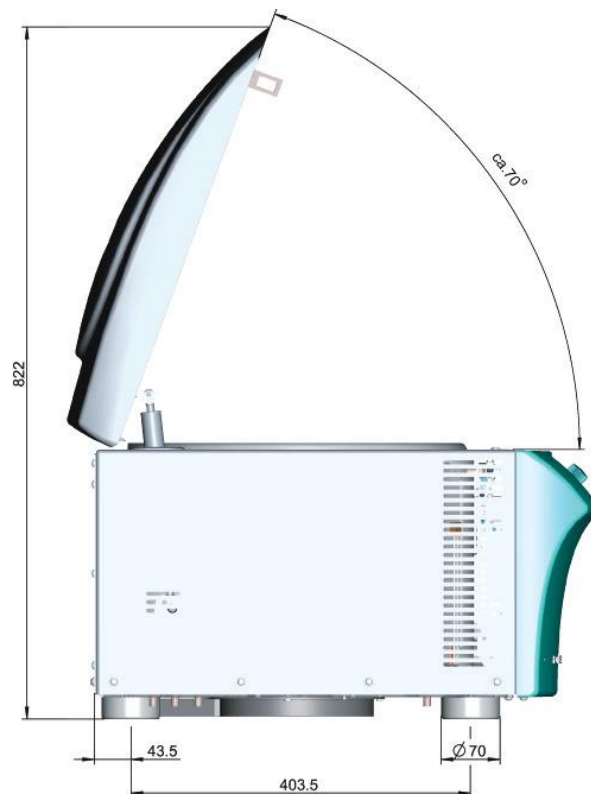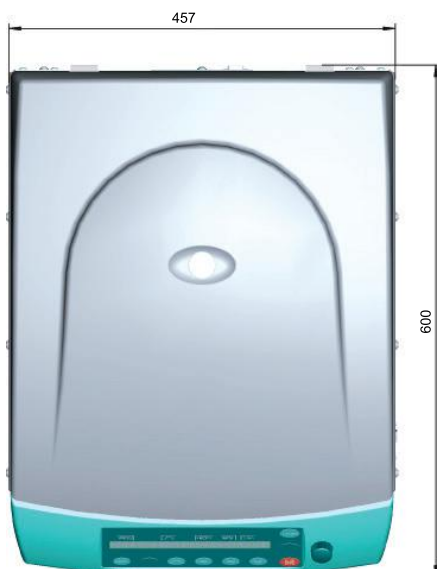

DIMENSIONS

