

OPERATING INSTRUCTIONS

LABORATORY & PHARMACEUTICAL UNIT

100/270/300/500/660/700/1400

ATEX - Approved Units

Internal storage compartment

ARCTIKO

Attention !

Atex 94/9/EC zone 2: Internal storage compartment

Note: That Zone 2 class only applies to storage compartment.- EX II - / 3 G IIA

Fan iQ-Motor

Equipment protection by type and protection “n”
(Explosive atmospheres)

EN 60079 -15:2010

All material stored in this product must be in airtight packing.

INDEX	PAGE
Before using the appliance	3
Safeguarding the environment	3
Power Failure	3
Precautions and general recommendations	3
Installing the appliance	4
Electrical connection	4
Before commissioning	5
Starting procedure	5
Starting up the appliance	5
Disconnecting the appliance	5
Control Panel.	6
Setting icons	7
Custom Settings	8
Advanced Settings	9
Service Settings	10
Advanced Service Settings	11
Status	11
Change / Reset Password	11
Download and upload Data	12
Alarms on display	13
Default settings.	14-17
Wiring Diagram	18-29

IMPORTANT

Before using the appliance

For safety reasons, the instructions should be read before the appliance is commissioned.
The operating instructions should always be available and accessible to personnel!

Low Voltage Directive 2006/95/EEC.

Electromagnetic Compatibility Directive 2004/108/EEC.
Only use the appliance for its specific purpose.

Safeguarding the environment:

Packing

The packing material is entirely recyclable, please follow the local regulation for scrapping.

Appliance

By disposal deliver the appliance to an authorised centre for recovery of refrigerant. Make the appliance unusable by cutting of the power supply cord.

Information

This appliance contains a mixture of various refrigerants. For more information please refer to the rating plate affixed to the freezer.

Power Failure:

The unit is equipped with battery back up for display and contact for remote monitoring only.
Back up time is min. 48 hours, the freezer must be connected voltage 100V, 115V or 230V, in minimum 72 hour before maximum battery is available. Please note the lifetime of the battery is estimated to be 3 years. But should be checked by authorized personnel every 6. Month.

Precautions and general recommendations:

WARNING: Keep ventilation openings, in the appliance enclosure or in the built-in structure, clear of obstruction.

WARNING: Do not use mechanical devices or other means to accelerate the defrosting process, other than those recommended by the manufacturer.

WARNING: Do not damage the refrigerant system.

WARNING: Do not use electrical appliances inside the appliance storage compartment, unless they are of the type recommended by the manufacturer.

WARNING: Observe the Mandatory sign on the refrigerator door:

NOTE ! All material stored in this cabinet must be sealed (airtight).

- Always keep the keys in separate place and out of reach of children.
- Before servicing or cleaning the appliance, unplug it from the mains or disconnect the electrical power supply.
- The power supply cable may only be replaced by an authorised person.

Installing the appliance

1. Always allow an authorised electrician to install the appliance.
2. The appliance should be located in a cool, dry place out of direct sunlight. The appliance gives off a great deal of heat when in operation. If the ambient temperature is too high there is a risk that the appliance will not operate correctly. It is recommended that the room be ventilated so that the room temperature does not exceed ambient temperature.
3. Place the appliance on a solid and flat substrate. This will eliminate any vibration and irritating noise. The appliance should be placed with at least 10cm free to the sides, at least 16cm free at the back.

Electrical Connection

The appliance meets current EU directives. Low voltage 2006/95/EEC.

Electromagnetic compatibility 2004/108/EEC.

The appliance should be given extra protection in accordance with Electricity Supply Regulations in order to protect the user against dangerous electric shocks in the event of faults.

If the switch is for three-pin plug, a three-pin plug should be used and the conductor with yellow/green insulation connected to the Earth terminal.

Data regarding voltage and absorbed power/current is given on the rating plate.

Electrical connection must be made in accordance with the local regulations.

Warning

Regulations require that this appliance be earthed:

The manufacture declines all liability for injury to persons or animals and for damage to property resulting from failure to observe the above procedures and reminders.

If the socket and plug are not of the same type either the socket or the plug must be replaced by a qualified electrician.

Under no circumstances use extension leads or adaptors.

1. The appliance should be connected by an authorised electrician or by the company that sold the appliance. Incorrect installation may damage the appliance.
2. Place the appliance on a flat floor that can bear the weight of the appliance.
3. The electrical installation should be earthed.
4. Do not use extension cables. Connect the appliance directly to a fixed installation. If the appliance is connected to a long extension cable there is a risk that the cable will become hot.
5. The appliance should be placed in a dry location. Do not allow water to come into contact with the cabinet. Water entering electrical parts may cause short-circuiting.
6. In the event of functional or electrical irregularities, please contact an authorised service engineer.

Before Commissioning

Clean the appliance both inside and out using a damp, wrung out cloth. Dry with a dry cloth.

Starting procedure

Put the plug in the socket.

The thermostat pre-set by the producer. The appliance should run for at least 6 hours before the thermostat is changed to allow for stabilisation of the cooling system.


Starting up the appliance

Set the digital thermometer at the required.

Disconnecting the appliance


Electrical disconnecting must be possible either by unplugging the appliance or by means of a double pole switch located up-line from the socket.

Control Panel (G-214 Controller)

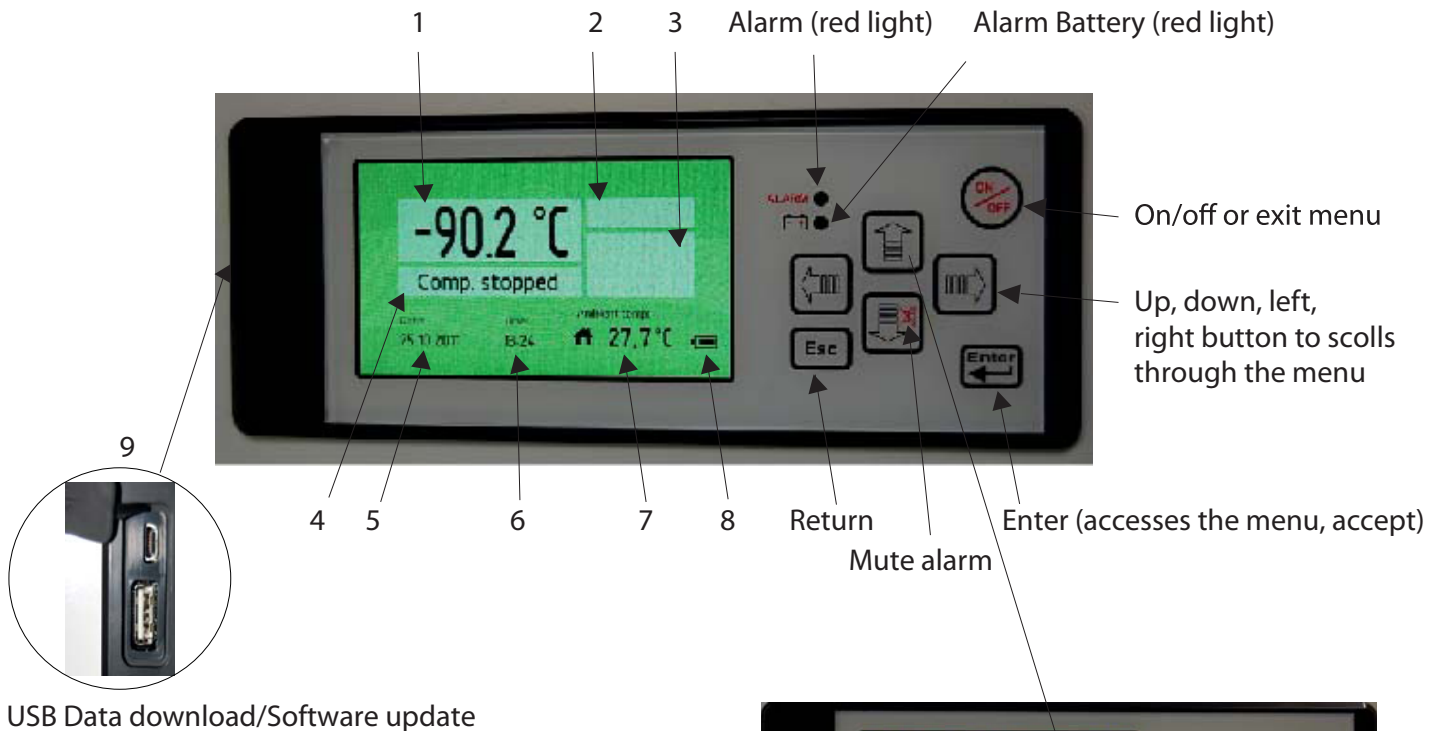

Control Panel Description

1. Temperature indicator
2. Alarm indicator
3. Alarm-icons
4. Compressor indication
5. Date indication
6. Hour indication
7. Ambient temperature indicator
8. Battery level
9. USB-Connection

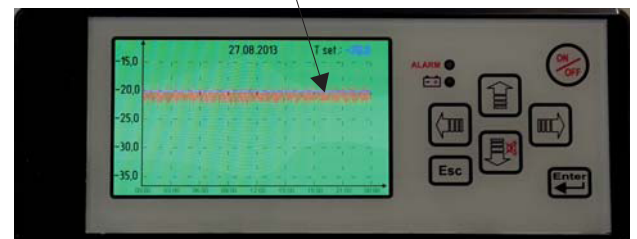

Push the **UP** button in order to generate a graph.

Push the **LEFT** button in order to return up to 10 days back.

Push the **Esc** button in order to exit the menu.

Setting icons



Custom Settings

Custom Settings



Advanced Settings

Advanced Settings



Advanced Service Settings

Advanced Service Settings



Status

Status



Change / Reset password

Change / Reset password



Custom Settings

Custom Settings



Enter password

Enter password

The menu **Custom setting** is protected with a password, which is "0000" to access the menu.



Setpoint

Under the menu **Setpoint** the temperature for the unit will be set.



Alarm Settings

Under the menu **Alarm Settings** the below menu will be available.



Select language

Under the menu **Select language** you can select the desired language.



Alarm delay

The menu **Alarm delay** is the settings of the time from an alarm will occur on the unit and until it will be shown on the display.



Door open alarm

The menu **Door open alarm** is the settings of the time from an alarm will occur on the unit and until it will be shown on the display.



High temp. alarm

The menu **High temp. alarm** is the settings of the highest temperature the device must be inside before it comes with an alarm.



Low temp. alarm

The menu **Low temp. alarm** is the settings of the lowest temperature the device must be inside before it comes with an alarm.



Probe eprom failure

The **Probe eprom failure** alarm indicates that the probe is not working.



Power failure

The **Power failure** alarm indicates that there is no power to the unit.



Logging time interval

The menu **Logging time interval** is the settings of time for the controller to log data from the unit.



Advanced Settings



Enter password

Enter password

The menu **Advanced Settings** is protected with a password which is "0000" to access the menu.



Calibration

The menu **Calibration** is the settings for offset of the temperature in the display.



Automatic defrost

The menu **Automatic defrost** is the setting of the time where a defrost automatic will be performed on the unit.



Manual defrost

Under the menu **Manual defrost** the defrost can be started immediately.



Hysteresis

The menu **Hysteresis** is the setting of the different from the setpoint where the compressor will start and stop.



Temp. range limits

The menu **Temp. range limits** is the settings of the maximum and minimum temperature the unit can be used for.



Service

Service

See Page 10



Password protection

The menu **Password protection** is the setting of the protection for the unit not to be turned off and is protected with a password.



Set date / time

The menu **Set date / time** is the setting of date and time.



Ambient temp. settings

The menu **Ambient temp. settings** is the setting of the ambient temperature (min / max.)



Service

Service Settings

Under the menu **Service**, there is information about the unit, which is important for service on the unit.



Compressor hours

The menu **Compressor hours** shows, how many hours the compressor has run.



Fan hours

The menu **Fan hours** shows, how many hours the fan has run.



Probe

Under the menu **Probe** the temperature for the sensor mounted in the unit is displayed.



Probe inside unit

Under the menu **Probe inside unit** the temperature of the sensor mounted inside the unit is displayed. There is mounted a sensor as standard in the unit, but it is possible to mount 3 sensors inside the device. (Controller version)



Probe Compressor

Under the item **Probe Compressor** the temperature of the sensor mounted on the compressor is displayed. (Not all models)



Probe evaporator

Under the point **Probe evaporator** the temperature of the sensor, that is located on the Evaporator, is displayed.



Advanced Service Settings

Advanced Service Settings

This menu is only used by the service department at the manufacturer.



Status

Status



Current alarms

The **Current alarms** shows the alarm, which have been on the unit with data and time.



Temp. last 24 hours

The **Temp. last 24 hours** shows the temperature of the unit the last 24 hours.



Change/Reset password

Change / Reset password



Change password

Under **Change password** it is possible to change the password for **Customer Settings**, **Advanced Settings** and **Advanced Service Settings**.



Reset password

Under **Reset password** it is possible to reset the password for the **Customer Settings**, **Advanced Settings** and **Advanced Service Settings**.
Contact the manufacturer to get the password for reset password.

Download Data



Place the USB pen drive in the connection on the display and press the button right for **read data** from the unit to the USB pen.



The **Read data** will be shown in the display when the data is loading to the USB pen drive.



The **Transfer complete** will be shown on the display when the loading is finished and the USB pen can be removed from the display.

Open the 2 files:
data00 and **param00**
on a computer in Excel or similar.


Upload Data (Only for new settings from the manufactory)



Turn-off the display before start.
Place the USB pen drive in the connection on the display and press the button left for upload parameters from the USB pen to the unit.



The **Data upload** will be shown in the display when the data is loading to the unit.

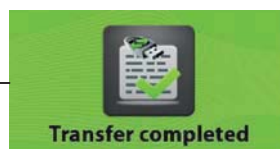

The **Transfer complete** will be shown on the display when the loading is finished and the USB pen can be removed from the display.

Alarms on display



Alarm icons



Door open alarm

The **Door open alarm** indicate, that the door is not correct closed



High temp. alarm

The **High temperature alarm** indicate, that the temperature inside the unit is higher than permitted from setting of the unit.



Low temp. alarm

The **Low temperature alarm** indicate, that the temperature inside the unit is lower than permitted from setting of the unit.



Probe eeprom failure

The **Probe eeprom failure** alarm indicate, that the probe is not working.



Power failure

The **Power failure** alarm indicate, that there is no power to the unit.



Low battery

The **Low battery** alarm indicate, that the power on the battery is to low.

Default settings LR / LF

Description	Arctiko default settings	Customer settings
Custom Settings (Password)	0000	
Set point		
Refrigerator	+4,0	
Freezer	-20,0	
Alarm Settings:		
Alarm delay		
Refrigerator	15 min	
Freezer	15 min	
Door open alarm		
Refrigerator	Enable	
Freezer	Enable	
High temp. alarm		
Refrigerator	+6	
Freezer	-15	
Low temp. alarm		
Refrigerator	+2	
Freezer	-30	
Probe/eprom failure		
Refrigerator	Enable	
Freezer	Enable	
Power failure		
Refrigerator	Enable	
Freezer	Enable	
Logging time interval		
Refrigerator	1	
Freezer	1	

Description	Arctiko default settings	Customer settings
Advanced Settings (Password)	0000	
Calibration		
Refrigerator	0,0	
Freezer	0,0	
Automatic defrost		
Refrigerator	06	
Freezer	06	
Manual defrost		
Refrigerator	OK	
Freezer	OK	
Hysteresis		
Refrigerator	2,0	
Freezer	2,0	
Temp. range limits		
Refrigerator Max.	+10,0	
Refrigerator Min.	+1,0	
Freezer Max.	-10,0	
Freezer Min.	-30,0	

Default settings LFF

Description	Arctiko default settings	Customer settings
Custom Settings (Password)	0000	
Set point		
Refrigerator	+4,0	
Freezer	-20,0	
Alarm Settings:		
Alarm delay		
Refrigerator	15 min	
Freezer	15 min	
Door open alarm		
Refrigerator	Enable	
Freezer	Enable	
High temp. alarm		
Refrigerator	+6	
Freezer	-15	
Low temp. alarm		
Refrigerator	+2	
Freezer	-30	
Probe/eprom failure		
Refrigerator	Enable	
Freezer	Enable	
Power failure		
Refrigerator	Enable	
Freezer	Enable	
Logging time interval		
Refrigerator	1	
Freezer	1	

Description	Arctiko default settings	Customer settings
Advanced Settings (Password)	0000	
Calibration		
Refrigerator	0,0	
Freezer	0,0	
Automatic defrost		
Refrigerator	06	
Freezer	06	
Manual defrost		
Refrigerator	OK	
Freezer	OK	
Hysteresis		
Refrigerator	1,0	
Freezer	1,0	
Temp. range limits		
Refrigerator Max.	+10,0	
Refrigerator Min.	+1,0	
Freezer Max.	-10,0	
Freezer Min.	-30,0	

Default settings PR / PF

Description	Arctiko default settings	Customer settings
Custom Settings (Password)	0000	
Set point		
Refrigerator	+4,0	
Freezer	-20,0	
Alarm Settings:		
Alarm delay		
Refrigerator	15 min	
Freezer	15 min	
Door open alarm		
Refrigerator	Enable	
Freezer	Enable	
High temp. alarm		
Refrigerator	+6	
Freezer	-15	
Low temp. alarm		
Refrigerator	+2	
Freezer	-30	
Probe/eprom failure		
Refrigerator	Enable	
Freezer	Enable	
Power failure		
Refrigerator	Enable	
Freezer	Enable	
Alarm log time interval		
Refrigerator	1	
Freezer	1	


Description	Arctiko default settings	Customer settings
Advanced Settings (Password)	0000	
Calibration		
Refrigerator	0,0	
Freezer	0,0	
Automatic defrost		
Refrigerator	06	
Freezer	06	
Manual defrost		
Refrigerator	OK	
Freezer	OK	
Hysteresis		
Refrigerator	2,0	
Freezer	2,0	
Temp. range limits		
Refrigerator Max.	+10,0	
Refrigerator Min.	+1,0	
Freezer Max.	-10,0	
Freezer Min.	-23,0	

Default settings BBR


Description	Arctiko default settings	Customer settings
Custom Settings (Password)	0000	
Set point		
Refrigerator	+4,0	
Alarm Settings:		
Alarm delay		
Refrigerator	5 min	
Door open alarm		
Refrigerator	Enable	
High temp. alarm		
Refrigerator	+6	
Low temp. alarm		
Refrigerator	+2	
Probe/eprom failure		
Refrigerator	Enable	
Power failure		
Refrigerator	Enable	
Logging time interval		
Refrigerator	1	

Description	Arctiko default settings	Customer settings
Advanced Settings (Password)	0000	
Calibration		
Refrigerator	0,0	
Automatic defrost		
Refrigerator	06	
Manual defrost		
Refrigerator	OK	
Hysteresis		
Refrigerator	1,0	
Temp. range limits		
Refrigerator Max.	+6,0	
Refrigerator Min.	+2,0	


Wiring diagram LR 100 / 300 500 / 700 / -ST

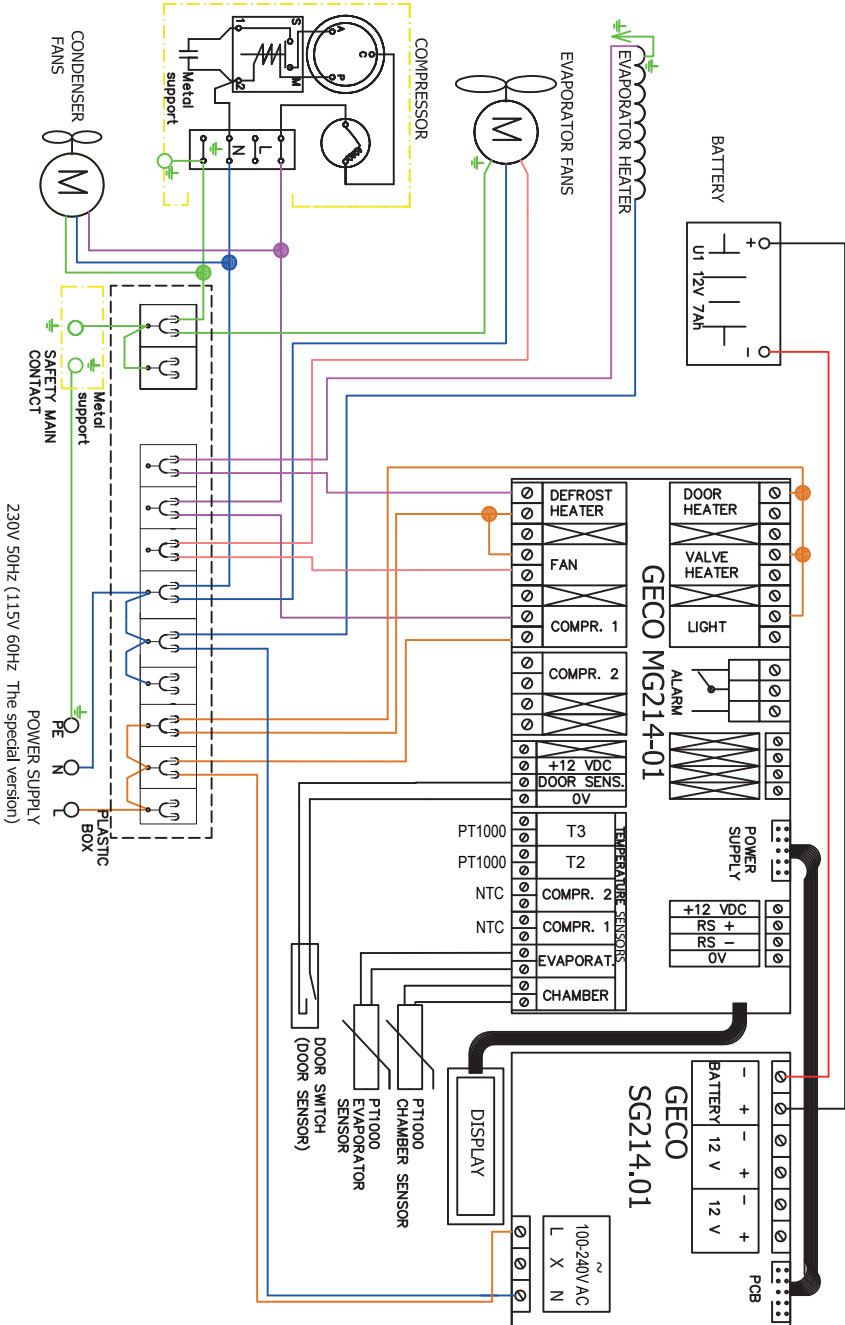


Wiring diagram LR 1400 / -ST

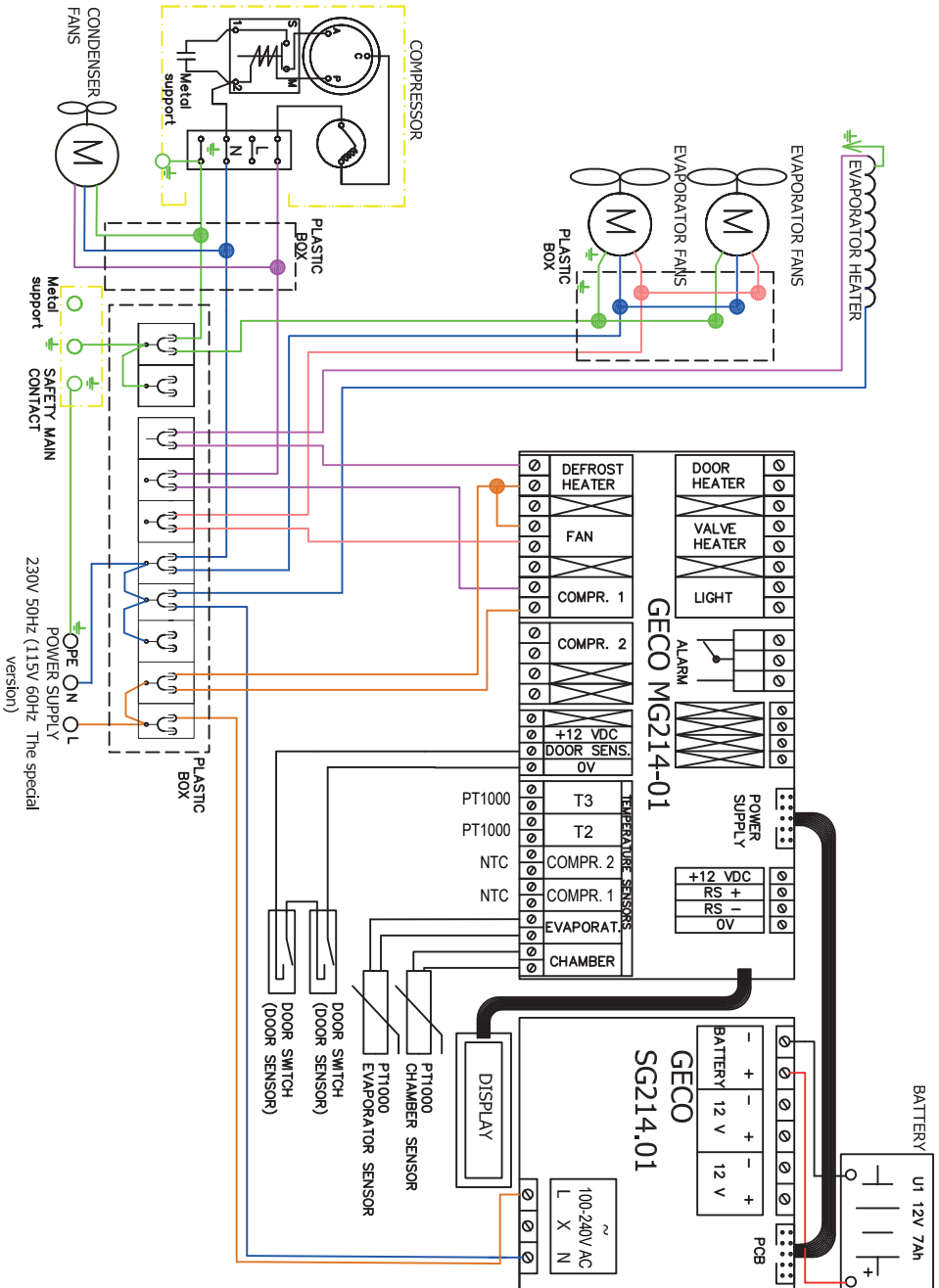


Wiring diagram LF 300 / -ST

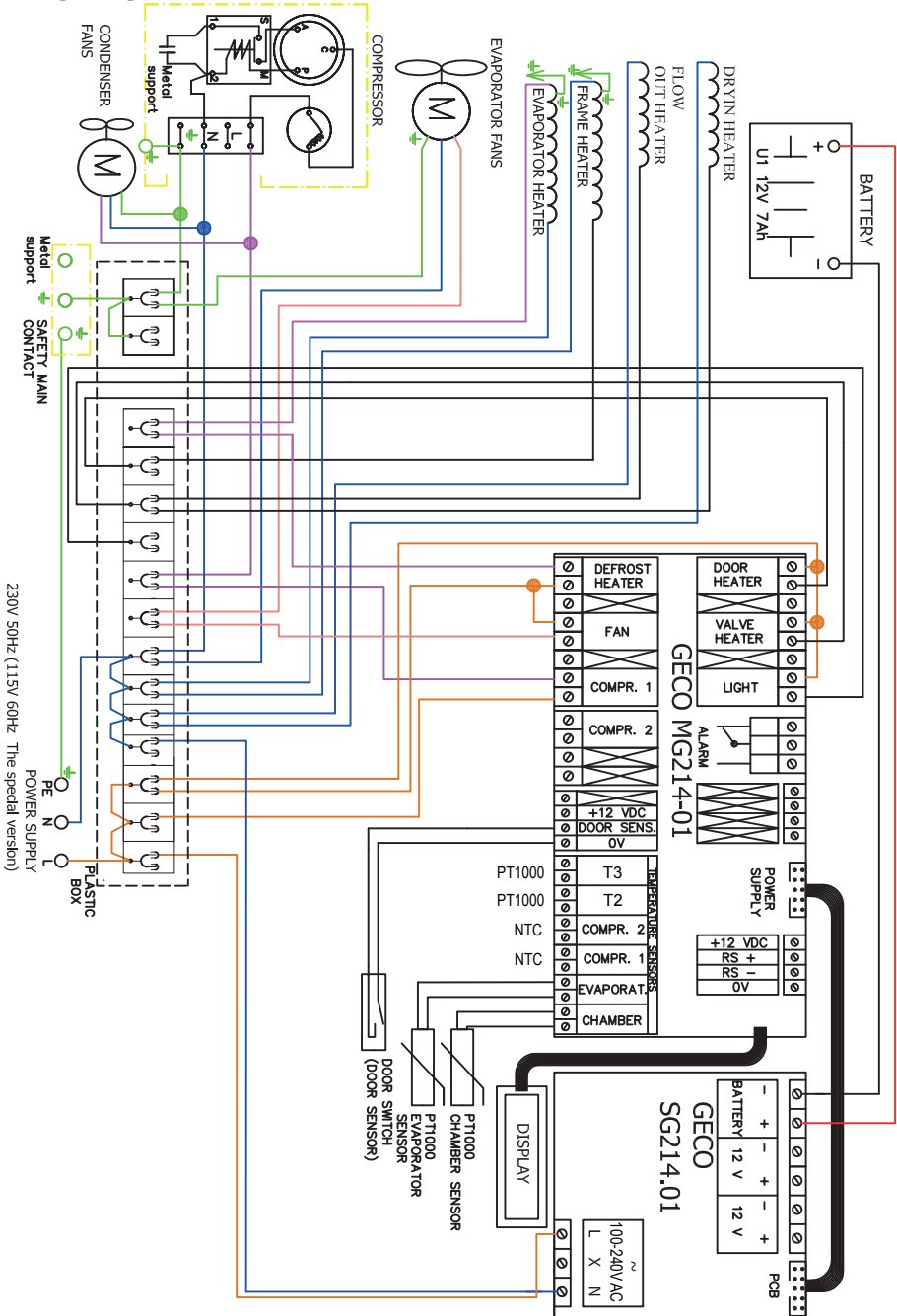


Wiring diagram LF 500 / 700 / -ST

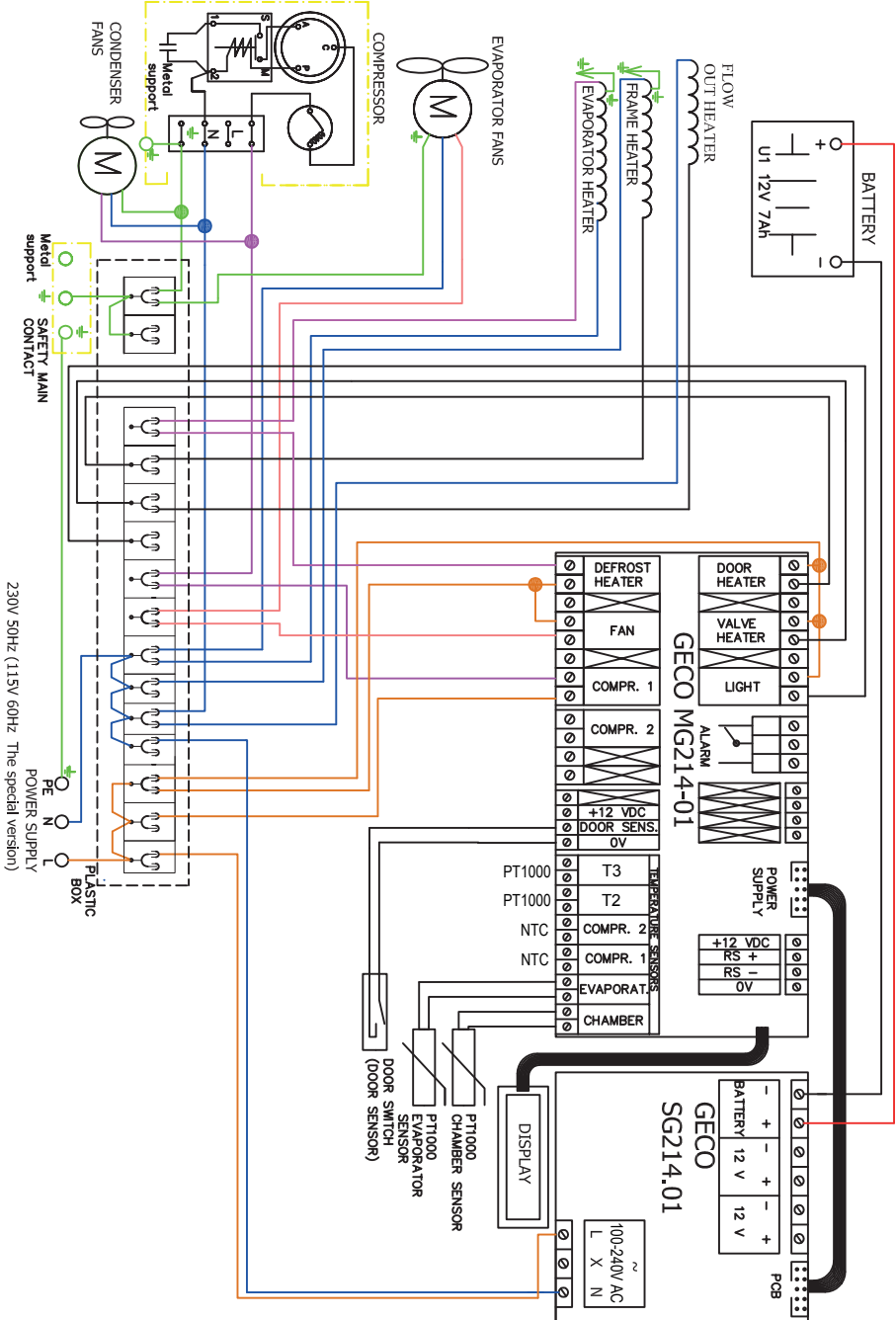


Wiring diagram LF 1400 / -ST

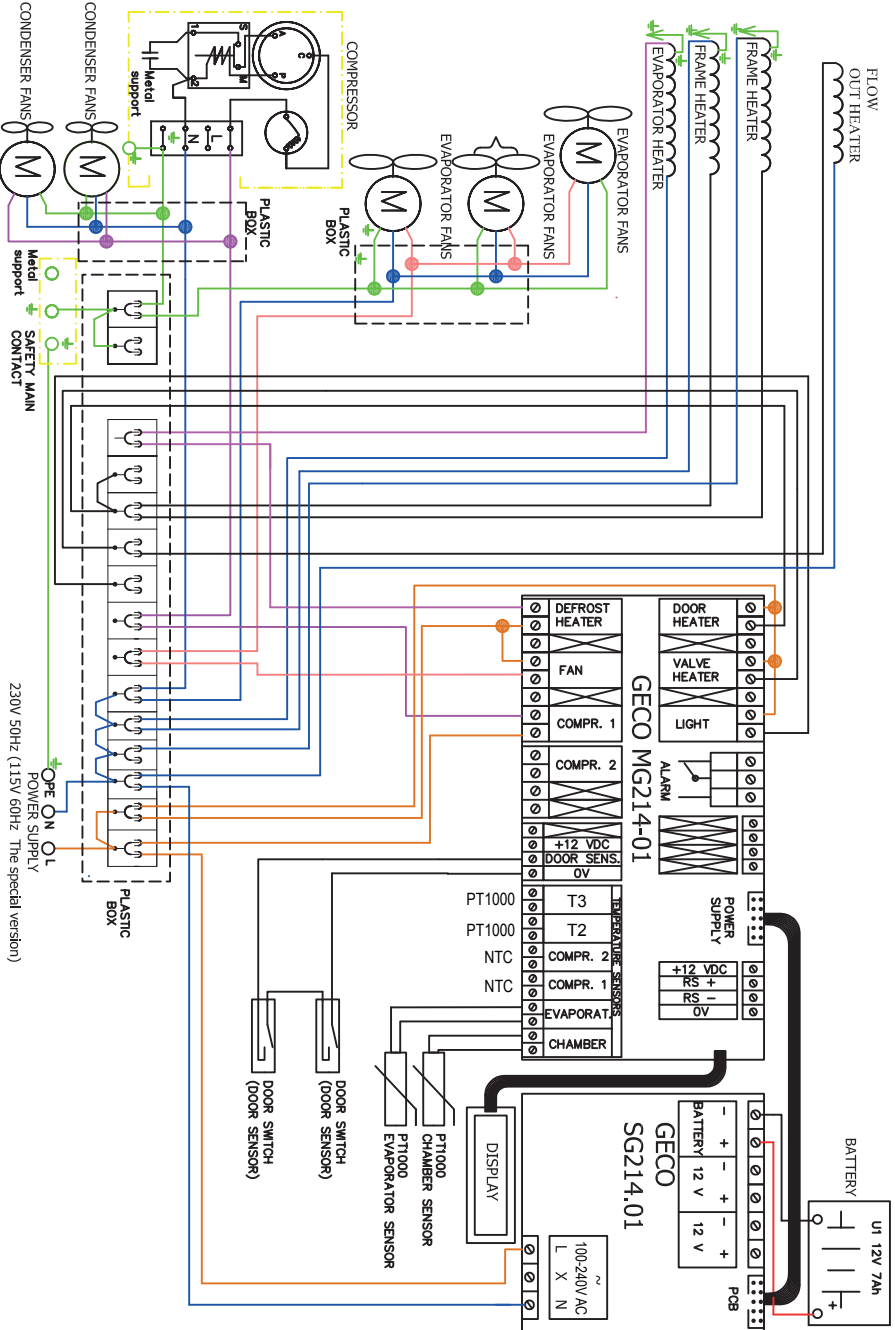


Wiring diagram LFF 270 / -ST

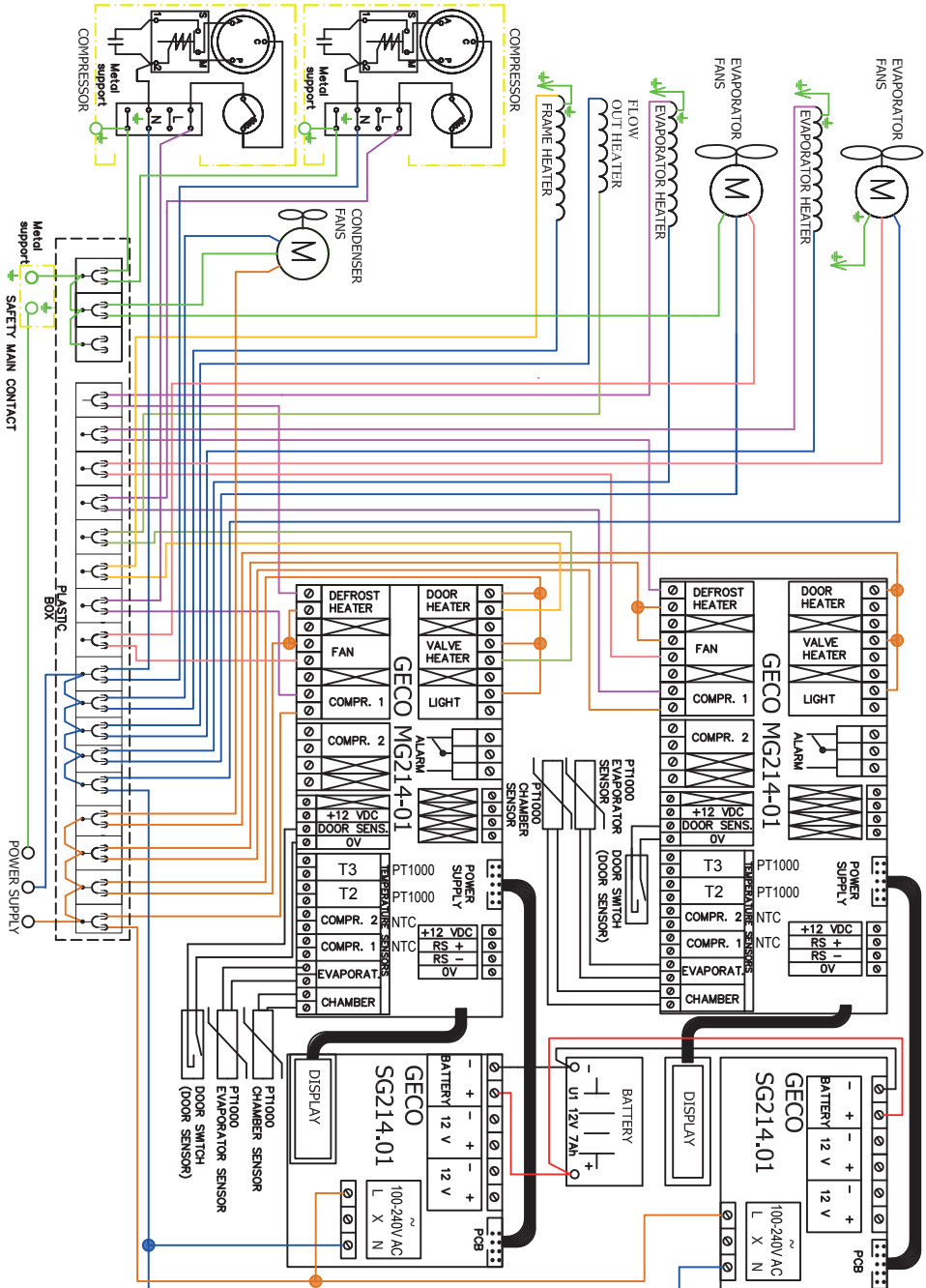


Wiring diagram LFF 660 / -ST

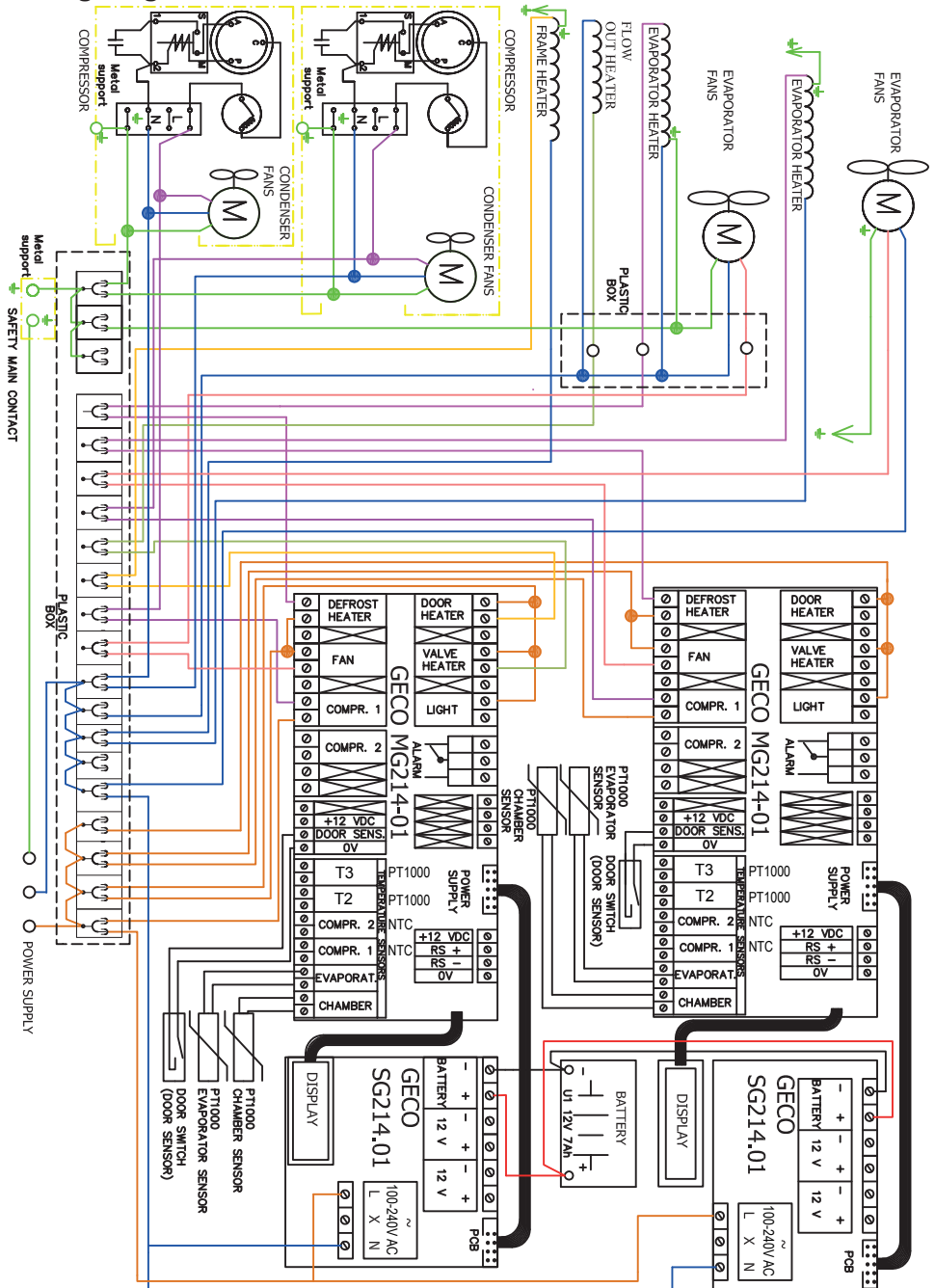

Wiring diagram BBR / PR / 300 / 500 / 700 / -ST

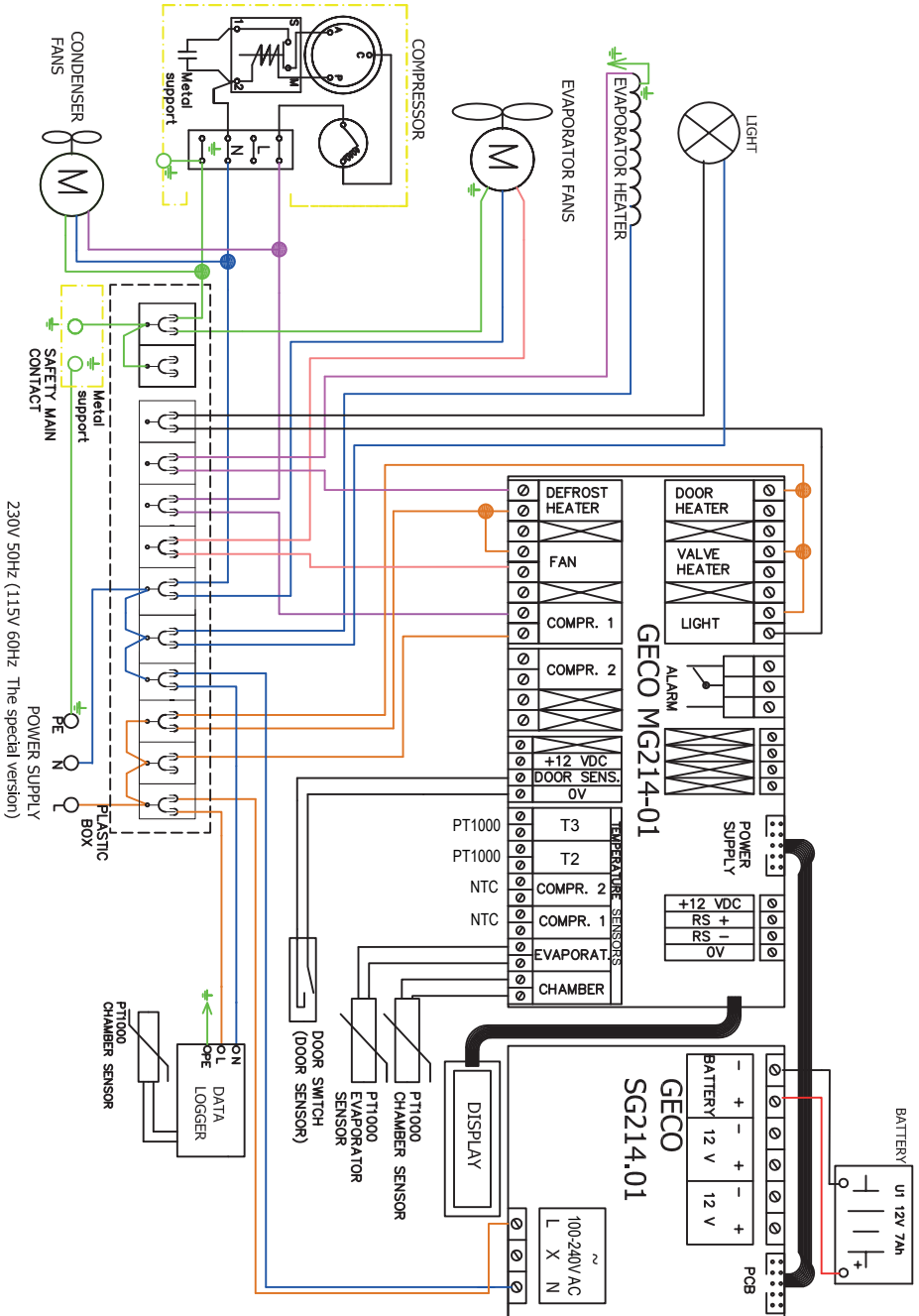


Wiring diagram BBR / PR / 1400 / -ST

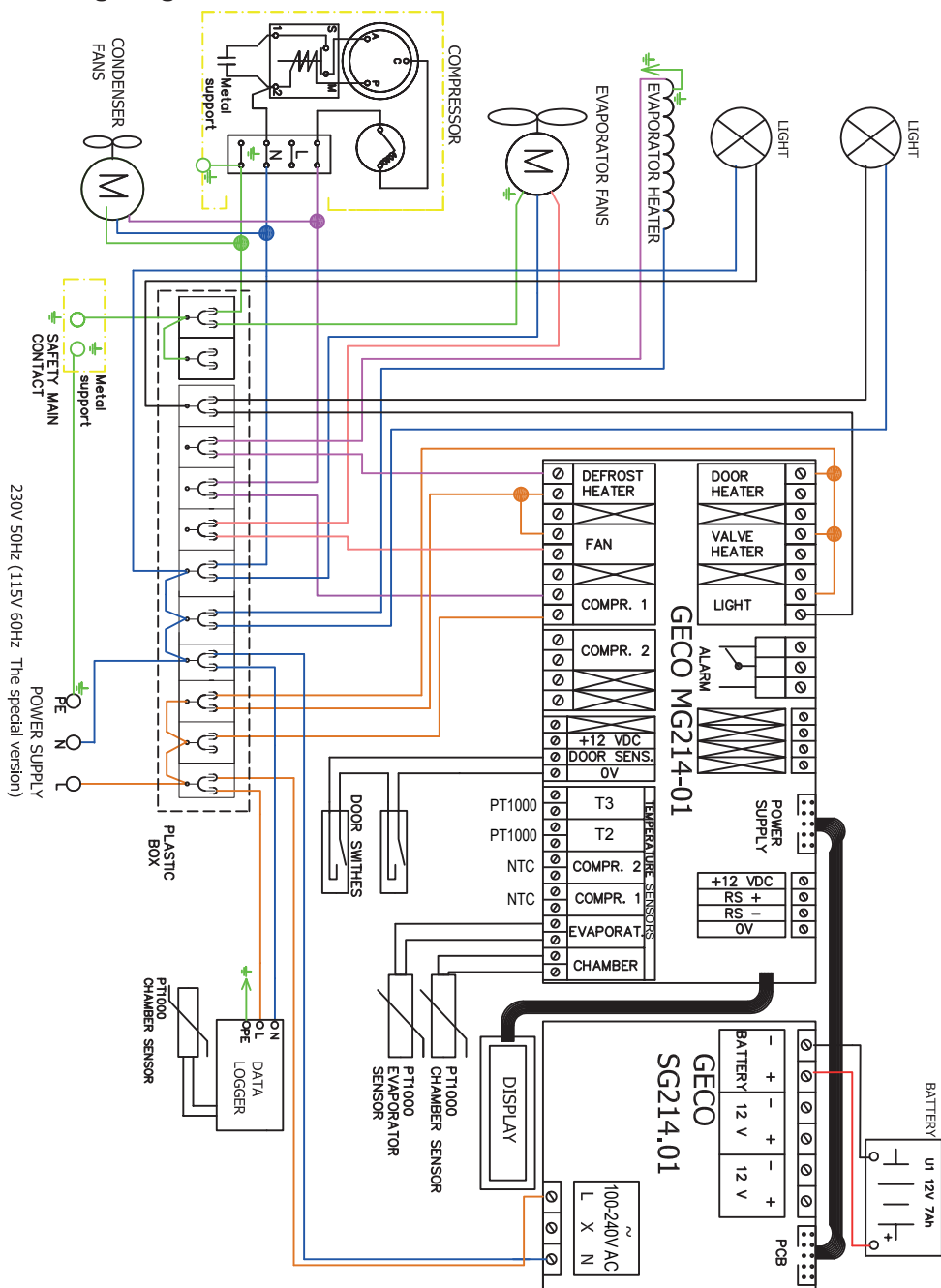




Wiring diagram PF 1400 / -ST

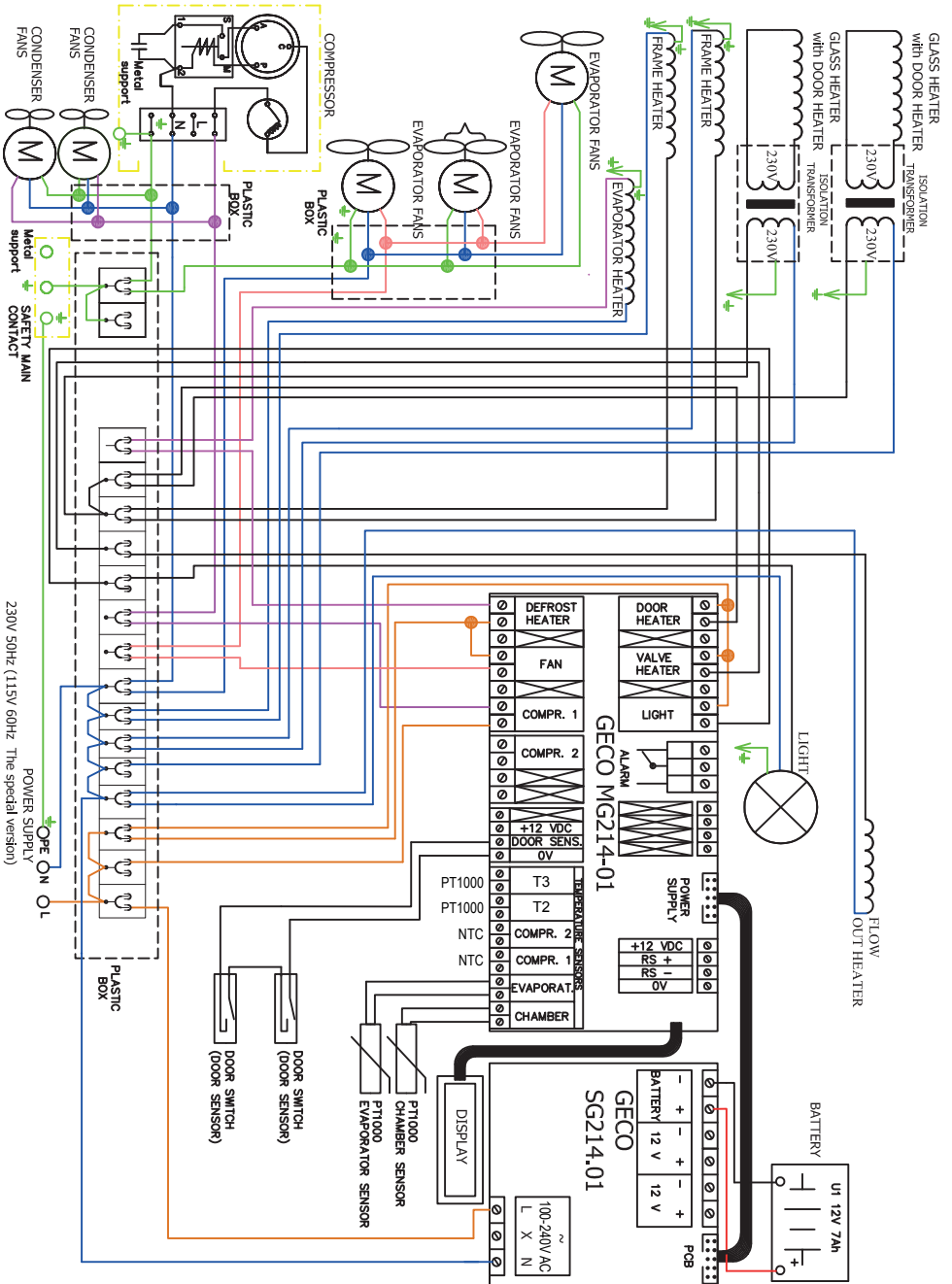

Wiring diagram BBR 100

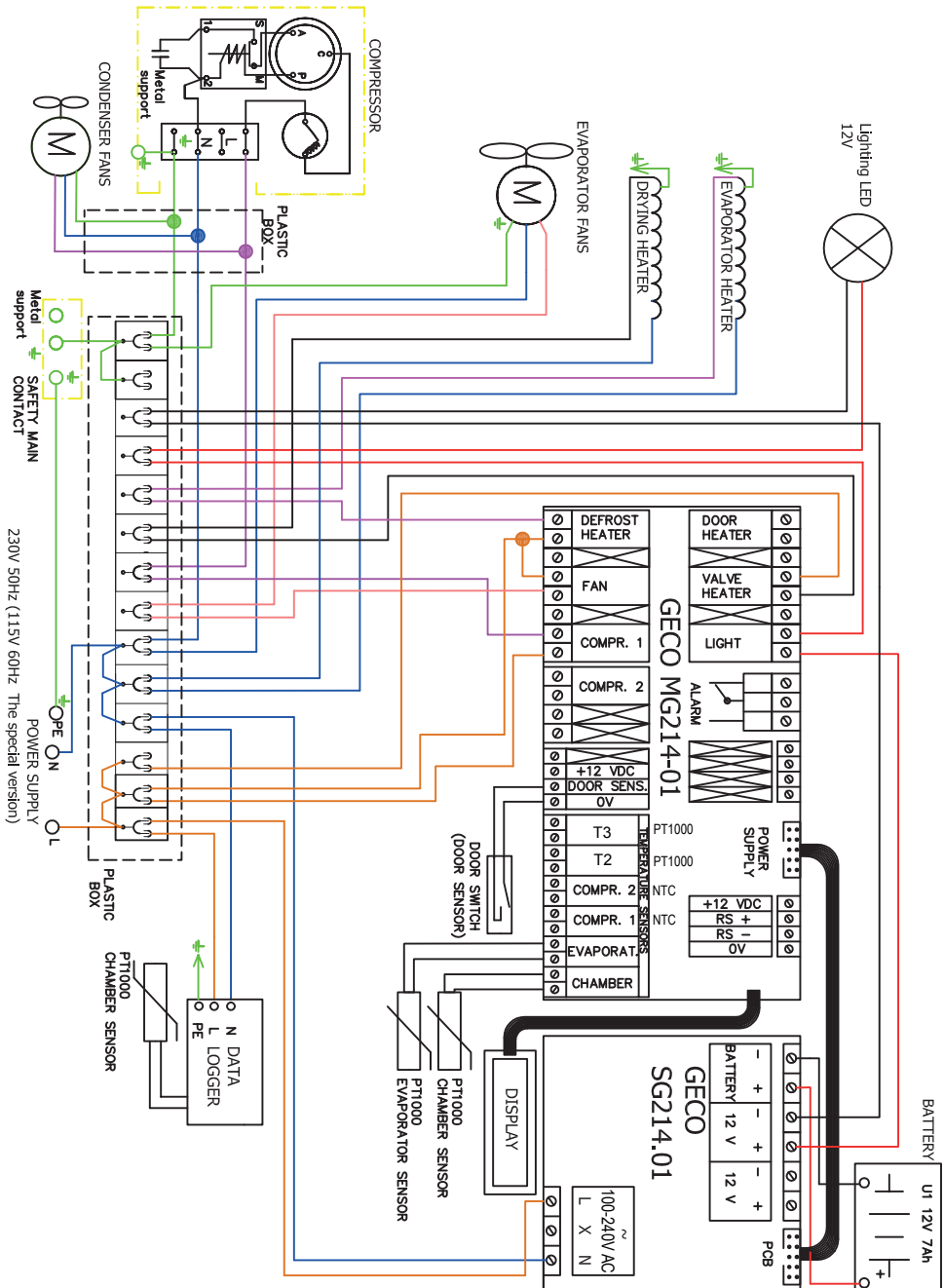

Notes:

Notes:



WWW.ARCTIKO.COM

LAMMEFJORDSVEJ 5
DK-6715 ESBJERG N
DENMARK

TEL. +45 70 20 03 28
FAX. +45 70 20 03 29

INFO@ARCTIKO.COM
WWW.ARCTIKO.COM