## Oxidase Test Stick

Rapid test for detection of cytochrome oxidase enzymatic activity.

### DESCRIPTION

Oxidase Test Stick is a diagnostic test used for differentiation and microbial identification, particularly of Gram-negative bacteria, on the basis of the presence of enzyme cytochrome oxidase.

The product matches with recommendations of EN ISO 16266 and ISO 9308-1 for detection of *Pseudomonas aeruginosa* and for confirmation of *Escherichia coli* and coliform bacteria, respectively.

### CONTENTS OF THE PACKAGES

Each package contains 50 Oxidase Test Stick.

### METHOD PRINCIPLE

Oxidase-positive bacteria produces the enzyme cytochrome oxidase (indophenol oxidase) that catalyzes the transport of electrons from donor compounds (NADH) to electron acceptors (usually oxygen).

Tetramethyl-p-phenylenediamine dihydrochloride contained in Oxidase Test Stick acts as an artificial electron donor and is oxidized by oxidase-positive bacteria forming the coloured compound indophenol blue.

### COMPOSITION

Oxidase Test Stick is made of a special paper with a zone impregnated with a solution of N,N,N',N'-tetramethyl-p-phenylenediamine dihydrochloride.

### TEST PROCEDURE

1. Allow container to come to room temperature before opening, for minimizing condensation on the stick.
2. Pick up one or more than one well isolated colony and wipe off by the zone of the stick indicated by arrows. Alternatively, transfer a drop of the suspension of the test organism to the reaction zone of the stick or dip directly the point of the stick into the microbial suspension.
3. Examine for an immediate color change (within 60 seconds) at the position of the inoculated area (NB. The usage of very dilute microbial suspensions may result in longer reactions time)

### INTERPRETING RESULTS

The development of a blue-purple color indicates a positive reaction. No color change corresponds to a negative test, i.e. the organism under investigation does not produce the enzyme cytochrome oxidase.

### LIMITATIONS

The most suitable cultures for the oxidase test are those from culture media without dyes, indicators or inhibitors. Bacterial colonies taken from media with pH values below 5.5 (e.g. after the metabolism of carbohydrates with subsequent acidification of the culture medium) can give a false negative oxidase reaction. Colonies taken from media containing nitrate may give unreliable results. Do not use steel, nichrome or iron containing loops to pick the colony. A platinum or plastic loop, or wooden applicator stick is recommended.

### STORAGE

Store at 2-8°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

### SHELF LIFE

2 years.

### QUALITY CONTROL

Control strains are indicated in the QC table.

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### PRESENTATION

	Contents	Ref.
Oxidase Test Stick	50 sticks	88029

### TABLE OF SYMBOLS

<b>LOT</b> Batch code	<b>IVD</b> <i>In vitro</i> Diagnostic Medical Device	Manufacturer	Use by	Fragile, handle with care
<b>REF</b> Catalogue number	Temperature limitation	Contains sufficient for <n> tests	Caution, consult Instruction For Use	Do not reuse

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## MUELLER HINTON AGAR

Medium for susceptibility test (Kirby-Bauer method).

### TYPICAL FORMULA (g/L)

Meat Extract	2.0
Casamino Acids, Technical	17.5
Starch	1.5
Agar	15.0
Final pH 7.3 ± 0.1	

### DESCRIPTION

MUELLER HINTON AGAR is used for antimicrobial susceptibility testing of rapidly growing aerobic microorganisms by the disk diffusion technique.

### PRINCIPLE

Casamino acids and meat extract are a source of amino acids, nitrogen, minerals, vitamins, carbon and other factors which increase the growth of microorganisms. Starch acts as a protective substance against toxic molecules which can be present in the medium. Hydrolysis of starch during sterilization supplies a little amount of glucose which represents a source of energy. Agar is the solidifying agent. Kirby-Bauer method is based on the diffusion, through the agar, of antimicrobial substances which soaks paper disks: microorganism growth shows an inhibition halo around the disk and the diameter of the halo is correlated to the Minimal Inhibiting Concentration (MIC).

### PREPARATION

Suspend 36.0 g of powder in 1 litre of distilled or deionized water. Heat to boiling and shake until completely dissolved. Sterilize in autoclave at 121°C for 15 minutes. Dispense in final containers.

### TECHNIQUE

Transfer 4-5 colonies in an appropriate broth.

Place it in a 37°C incubator until an opacity is obtained equivalent to the standard opacity of 0.5 on the MacFarland scale. Introduce a sterile swab into the inoculum and inoculate the agar passing 2 or 3 times onto the entire surface.

Press the disk containing the antimicrobial on the agar surface.

Incubate at 36±1°C for 18 hours, measure the inhibition zone with a compass and compare to the NCCLS recommended zone ranges.

### INTERPRETATION OF RESULTS

Compare obtained values of inhibition halo diameter with the values reported on NCCLS M100(M2) document.

### STORAGE

10-30°C away from light, until the expiry date on the label or until signs of deterioration or contamination are evident.

### WARNING and PRECAUTIONS

The product is not classified as hazardous by current legislation and does not contain harmful substances in concentrations of ≥1%. The product is designed for *In vitro* diagnostic use and must be used only by properly trained operators.

### DISPOSAL of WASTE

Disposal of waste must be carried out according to national and local regulations in force.

### REFERENCES

1. Bauer et al. (1966). J. Clin. Pathol. 45:493-496.
2. Mueller, J.H., and Hinton. 1941. Proc. Soc. Exp. Biol. Med. 48: 330-333.
3. NCCLS. Performance standards for susceptibility testing; Twelve Informational Supplement. NCCLS Document M100-S12, January 2002.

**Liofilchem s.r.l.**

Via Scozia-Zona industriale - 64026 Roseto degli Abruzzi Tel. +39.085.8930745 - Fax +39.085.8930330  
Web site: <http://www.liofilchem.net> E-mail: [liofilchem@liofilchem.net](mailto:liofilchem@liofilchem.net)

## PRODUCT SPECIFICATIONS

### NAME

MUELLER HINTON AGAR

### PRESENTATION

Dehydrated culture medium

### STORAGE

10-30°C

### PACKAGING

Code	Content	Packaging
610033	500 g	500 g of powder in plastic bottle
620033	100 g	100 g of powder in plastic bottle
6100335	5 kg	5 kg of powder in plastic container

### pH OF THE MEDIUM

7.3 ± 0.1

### USE

MUELLER HINTON AGAR is used for antimicrobial susceptibility testing of rapidly growing aerobic microorganisms by the disk diffusion technique.

### TECHNIQUE

Refer to technical sheet of the product.

### APPEARANCE OF THE MEDIUM

Amber medium, slightly opalescent.

### SHELF LIFE

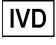
4 years

### QUALITY CONTROL

- Control of general characteristics, label and print
- Sterility control  
 7 days at 25 ± 1°C, in aerobiosis  
 7 days at 36 ± 1°C, in aerobiosis
- Microbiological control  
 Incubation conditions: 18-24 h at 36 ± 1°C

Microorganism	ATCC	Growth	Characteristics
<i>Enterococcus faecalis</i>	ATCC 29212	Good	White colonies
<i>Escherichia coli</i>	ATCC 25922	Good	Colorless colonies
<i>Proteus mirabilis</i>	ATCC 25933	Good	Colorless colonies
<i>Staphylococcus aureus</i>	ATCC 25923	Good	White colonies

### TABLE of SYMBOLS

 <b>IVD</b> In vitro Diagnostic Medical Device	 <b>LOT</b> Batch code	 Manufacturer	 Contains sufficient for <n> tests
 <b>REF</b> Catalogue number	 Temperature limitation	 Use by	 Caution, consult accompanying documents

**Liofilchem s.r.l.**

Via Scozia-Zona industriale - 64026 Roseto degli Abruzzi Tel. +39.085.8930745 - Fax +39.085.8930330  
 Web site: <http://www.liofilchem.net> E-mail: [liofilchem@liofilchem.net](mailto:liofilchem@liofilchem.net)

## Sabouraud Dextrose Agar

Medium for the cultivation and enumeration of yeasts and moulds from different materials, according to EN ISO 11133 and USP/EP/JP.

### DESCRIPTION

Sabouraud Dextrose Agar (SDA) is a non selective isolation medium used for the growth and maintenance of pathogenic and non-pathogenic fungi from clinical and nonclinical specimens. It is also used for recovery and total counting of yeasts and moulds in environmental monitoring.

This medium complies with EN ISO 11133 for microbiological examination of food, animal feed and water, where it is described as the main reference medium to carry out quantitative testing on culture media intended for fungi.

Its formula conforms to the recommendations of the harmonized method in the United States Pharmacopoeia (USP), European Pharmacopoeia (EP) and Japanese Pharmacopoeia (JP) for the microbiological examination of non sterile products. The medium is also available as gamma-irradiated triple bagged plates, particularly suitable for use in restricted areas like isolators and clean rooms.

### TYPICAL FORMULA

	(g/l)
Pancreatic Digest of Casein	5.0
Peptic Digest of Animal Tissue	5.0
Dextrose	40.0
Agar	15.0
Final pH 5.6 ± 0.2 at 25°C	

### METHOD PRINCIPLE

Pancreatic digest of casein and peptic digest of animal tissue provide amino acids, nitrogen, carbon, vitamins and minerals for organisms growth. Dextrose is an energy source. Agar is the solidifying agent. The high concentration of dextrose and the acidic pH of the medium permit selectivity of fungi.

The medium can be supplemented with chloramphenicol to increase bacterial inhibition and recovery of dermatophytes.

### PREPARATION

<u>Dehydrated medium</u>	Suspend 65 g of the powder in 1 liter of distilled or deionized water. Mix well. Heat to boil shaking frequently until completely dissolved. Sterilize in autoclave at 121°C for 15 minutes.
<u>Medium in bottles</u>	Melt the content of the bottle in a water bath at 100°C (loosing the cap partially removed) until completely dissolved. Then screw the cap and check the homogeneity of the dissolved medium, if it is the case turning the bottle upside down. Cool at 45-50°C, mix well avoiding foam formation and aseptically distribute into Petri dishes.

### TEST PROCEDURE

#### For use in medical microbiology

Streak the specimen as soon as possible after it is received in the laboratory to obtain isolated colonies. Prepared tubed slants primarily are intended for use with pure cultures for maintenance or other purposes. Incubation conditions may vary according to the type of specimen and the microorganisms being tested for.

#### For use in food, animal feed and water testing

Refer to EN ISO 11133 for specific instructions.

#### For use in industrial microbiology

##### Control of non-sterile products

Refer to the procedure described in the harmonized chapters of the Pharmacopoeia.

##### Passive Air Monitoring

Take the lid off the settle plate and leave the medium exposed to the air for a period of time no longer than 4 hours (settling plates filled with 30 ml of medium may compensate for water loss during extended incubation periods). Plates can be placed according to the 1/1/1 scheme (for 1 h, about 1 above the floor, at least 1 m from the walls or any obstacle).

##### Surfaces and Personnel Hygiene Monitoring

Take a swab sample for irregular surfaces or use the sampling template 10x10 (ref. 96762) to sample a well defined area of the test surface. Inoculate a 90 mm plate by streaking the swab over the agar surface. Furthermore, the medium is suitable for personnel hygiene monitoring to detect microbial contamination of gloves or hands e.g. in a 5-finger-print.

Incubate the plates at 20-25°C for 5-7 days or at 30-35°C for 24-48 hours.

### INTERPRETING RESULTS

Transfer of growth from slants to plated media may be required in order to obtain pure cultures of fungi. Examine for fungal colonies exhibiting typical microscopic and colonial morphology. Biochemical tests may be required for final identification.

The total combined yeasts/moulds count (TYMC) is considered to be equal to the number of CFU found per each plate. When an acceptable criterion for microbiological quality is prescribed it is interpreted as follows:

- 10<sup>1</sup> CFU: maximum acceptable count = 20;
- 10<sup>2</sup> CFU: maximum acceptable count = 200;
- 10<sup>3</sup> CFU: maximum acceptable count = 2000, and so forth.

In procedures intended for environmental and personnel hygiene monitoring, observe daily for the formation of colonies.

**APPEARANCE**

Dehydrated medium: free-flowing, homogeneous, light beige.

Prepared medium: slightly opalescent, light amber.

**STORAGE**

The powder is very hygroscopic, store the powder at 10-30°C, in a dry environment, in its original container tightly closed. Store bottles, tubes and prepared plates at 10-25°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

**SHELF LIFE**

Dehydrated medium: 4 years.

Medium in bottles: 2 years.

Medium in tubes: 1 year.

Ready-to-use plates (90 and 60 mm): 6 months.

Contact plates (55 mm): 9 months

**QUALITY CONTROL**

The medium is inoculated with the microbial strains indicated in the QC table.

Inoculum for productivity: 50-100 CFU.

Incubation conditions: 32.5 ± 2.5°C for 24-48 h (*C. albicans*) and at 22.5 ± 2.5°C for up to 5 days (all listed organisms), under aerobic atmosphere.

**QC Table.**

Microorganism		Growth
<i>Candida albicans</i>	WDCM 00054	Good
<i>Aspergillus brasiliensis</i>	WDCM 00053	Good
<i>Saccharomyces cerevisiae</i>	WDCM 00058	Good

**WARNING AND PRECAUTIONS**

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for *in vitro* diagnostic use and must be used only by properly trained operators.

**DISPOSAL OF WASTE**

Disposal of waste must be carried out according to national and local regulations in force.

**BIBLIOGRAPHY**

1. EN ISO 11133:2014+Amd1:2018. Microbiology of food, animal feed and water – Preparation, production, storage and performance testing of culture media.
2. European Pharmacopoeia 6.5 (2009) 2.6.13. Microbiological examination of non-sterile products: Test for specified microorganisms.
3. United States Pharmacopoeia 32 NF 27 (2009) <62> Microbiological examination of non-sterile products: Test for specified microorganisms.
4. Japanese Pharmacopoeia 4.05 (2008) Microbiological examination of non-sterile products: Test for specified microorganisms.
5. Sabouraud, R. (1892) Ann. Dermatol. Syphilol. 3:1061.

PRESENTATION	Category	Packaging	Ref.
Sabouraud Dextrose Agar	90 mm plates	20 plates	10035
Sabouraud Dextrose Agar	90 mm plates	100 plates	10035*
Sabouraud Dextrose Agar	90 mm plates (triple-wrapped and gamma-irradiated)	20 plates	10035S
Sabouraud Dextrose Agar	90 mm plates (triple-wrapped and gamma-irradiated, 30 ml filling volume)	20 plates	10114S f
Sabouraud Dextrose Agar	60 mm plates	20 plates	163402 f
Sabouraud Dextrose Agar	60 mm plates	450 plates	173402 f
Sabouraud Dextrose Agar	Tubes - Bottles	10 x 9 ml slant tubes	30093
Sabouraud Dextrose Agar	55 mm contact plates	20 plates	15327 f
Sabouraud Dextrose Agar	55 mm contact plates irradiated	20 plates	15327S f
Sabouraud Dextrose Agar	Tubes - Bottles	20 x 9 ml slant tubes	31093
Sabouraud Dextrose Agar	Tubes - Bottles	6 x 500 ml bottles	470040
Sabouraud Dextrose Agar	Tubes - Bottles	6 x 200 ml bottles	412280
Sabouraud Dextrose Agar	Tubes - Bottles	25 x 200 ml bottles	452280
Sabouraud Dextrose Agar	Tubes - Bottles	6 x 100 ml bottles	402280
Sabouraud Dextrose Agar	Dehydrated culture medium	500 g of powder	610103
Sabouraud Dextrose Agar	Dehydrated culture medium	100 g of powder	620103
Sabouraud Dextrose Agar	Dehydrated culture medium	5 kg of powder	610103S

f: Not CE Marked

**TABLE OF SYMBOLS**

<b>LOT</b> Batch code	<b>IVD</b> <i>In vitro</i> Diagnostic Medical Device	 Manufacturer	 Use by	 Fragile, handle with care
<b>REF</b> Catalogue number	 Temperature limitation	 Contains sufficient for <n> tests	 Caution, consult Instruction For Use	 Do not reuse

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## BILE BACTERIOLOGICAL

Bacteriological bile obtained by bile purification.

### USE

BILE BACTERIOLOGICAL is ox bile purified and dehydrated. It's a fine beige powder, easily soluble in water. It contains a mix of biliary salts and is used in media for enterobacteria, as selective agent, and for identification of enterococci.

### PHYSICO-CHEMICAL CHARACTERISTICS

	Standard	Method
Solubility in water at 5%	Complete	Eur.Ph.3rd ed.
pH of 5% solution	5.5-7.5	Eur.Ph.3rd ed.
Loss on drying	≤ 5.0%	Eur.Ph.3rd ed.
Bile acids	≥ 45.0%	Eur.Ph.10th ed.

### TECHNIQUE

Bile Bacteriological can be used as an ingredient of dehydrated culture media and need dissolution in distilled or deionized water and sterilization by autoclaving.

### QUALITY CONTROL

Dehydrated powder

Appearance: free-flowing, homogeneous.

Colour: beige.

### STORAGE

The powder is very hygroscopic: store the powder at 10-30 °C, in a dry environment, in its original container tightly closed and use it before the expiry date on the label or until signs of deterioration or contamination are evident.

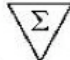
### REFERENCES

1. Cowan, S.T., Steel, K.J. (1979) Manual for the identification of medical bacteria. Edi. Ermes.

### PRESENTATION

Product	REF	$\Sigma$
BILE BACTERIOLOGICAL	611367	500 g
BILE BACTERIOLOGICAL	621367	100 g

### TABLE OF SYMBOLS

<b>LOT</b> Batch code	 Caution, consult accompanying documents	 Manufacturer	 Contains sufficient for <n> tests	<b>IVD</b> <i>In Vitro</i> Diagnostic Medical Device
<b>REF</b> Catalogue number	 Fragile, handle with care	 Use by	 Temperature limitation	 Keep away from heat source

## PEPTONE BACTERIOLOGICAL

Peptone obtained by enzymatic hydrolysis of animal tissue.

### USE

PEPTONE BACTERIOLOGICAL is an enzymatic hydrolysate of meat that supplies a limpid, colorless and very stable watery solution. It is used in the preparation of culture media as a nitrogen source readily available for bacterial growth. It is a general use very nutritive peptone, with neutral pH.

### PHYSICO-CHEMICAL CHARACTERISTICS

	Standard
Solubility in water at 2%	Complete
pH of 2% solution	7.0+/-0.5
Loss on drying	≤ 6.0%
Total nitrogen	>12.5%
α-amino nitrogen AN	3-4.5%
Ash	5.0%

### TECHNIQUE

Peptone Bacteriological can be used as an ingredient of dehydrated culture media and need dissolution in distilled or deionized water and sterilization by autoclaving.

### QUALITY CONTROL

Dehydrated powder

Appearance: free-flowing, homogeneous.

Color: cream.

### STORAGE

The powder is very hygroscopic: store the powder at 10-30 °C, in a dry environment, in its original container tightly closed and use it before the expiry date on the label or until signs of deterioration or contamination are evident.

### REFERENCES

1. Standard Methods for Examination of Water and Sewage, 15th ed., (1980).
2. J. Dairy Science, **16**: 277: (1933).

### PRESENTATION

Product	REF	$\Sigma$
PEPTONE BACTERIOLOGICAL	611701	500 g
PEPTONE BACTERIOLOGICAL	621701	100 g
PEPTONE BACTERIOLOGICAL	6117015	5 Kg

### TABLE OF SYMBOLS

<b>LOT</b> Batch code	 Caution, consult accompanying documents	 Manufacturer	 Contains sufficient for <n> tests	<b>IVD</b> <i>In Vitro</i> Diagnostic Medical Device
<b>REF</b> Catalogue number	 Fragile, handle with care	 Use by	 Temperature limitation	 Keep away from heat source

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### REFERENCIAS

- Standard Methods for Examination of Water and Sewage, 15th ed., (1980).
- J. Dairy Science, 16: 277: (1933).

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# CERTIFICATO

Nr 50 100 11497 - Rev. 02

Si attesta che / This is to certify that

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È CONFORME AI REQUISITI DELLA NORMA  
HAS BEEN FOUND TO COMPLY WITH THE REQUIREMENTS OF

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**Design and development, production and sale of in-vitro diagnostic medical devices: culture media for bacteriology, identification and susceptibility testing systems, kits for plasma protein determination (IAF 12, 29)**

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"THE VALIDITY OF THE PRESENT CERTIFICATE DEPENDS ON THE ANNUAL SURVEILLANCE EVERY 12 MONTHS AND ON THE COMPLETE REVIEW OF COMPANY'S MANAGEMENT SYSTEM AFTER THREE-YEARS"

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**Certification Mark:**

**Scope of Certificate:**

**Design and development, production and sale of in-vitro diagnostic medical devices: culture media for bacteriology, identification and susceptibility testing systems, kits for plasma protein determination. Distribution of other in-vitro diagnostic medical devices**

**Applied Standard(s):**

EN ISO 13485:2016  
Medical devices - Quality management systems - Requirements for regulatory purposes (ISO 13485:2016)  
DIN EN ISO 13485:2016

The Certification Body of TÜV SÜD Product Service GmbH certifies that the company mentioned above has established and is maintaining a quality management system, which meets the requirements of the listed standard(s). See also notes overleaf.

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---

### EC DECLARATION OF CONFORMITY

The company Liofilchem® S.r.l., registered office in Via Scozia, 64026 Roseto degli Abruzzi (TE) Italy, as a manufacturer of the *in vitro* medical-diagnostic device listed in the attached table, Revision 31.0 of 08.01.2016

hereby certifies under its own responsibility

1. that the above mentioned device complies with all the applicable provisions of Directive 98/79/EC (Annex III) and its relevant transposition into national law;
2. the above mentioned is not included in Annex II, List A and B of Directive 98/79/EC;
3. that the technical documentation referred to at Annex III of the Directive 98/79/EC is available for the national authorities in its facility and that this documentation shall be kept for 5 years after the last product has been manufactured;
4. that the manufacturing process follows suitable principles of quality assurance;
5. that, has implemented and keep up to date, a post-production surveillance system for monitoring the products;
6. that the device in question, was introduced into the market provided with CE mark.

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Rev. 31.0 del 08.01.2016

621021	HEART INFUSION BROTH
621022	MIDDLEBROOK 7H10 AGAR BASE
621210	WURTZ LACTOSE AGAR
621265	ISOSENSITEST AGAR
621367	BILE BACTERIOLOGICAL
621401	IRON SULPHITE AGAR
621402	CARY BLAIR TRANSPORT MEDIUM
621601	GLUCOSE
621618	CHROMATIC™ MH
621619	CHROMATIC™ CRE AGAR BASE
621701	PEPTONE BACTERIOLOGICAL
622202	STREPTOCOCCUS SELECTIVE AGAR
630026	LOWENSTEIN JENSEN MEDIUM w GLYCEROL 1 litre
71618	ENTEROSYSTEM 18R 20 Tests
71630	STAF SYSTEM 18 R 20 Tests
71670	COPRO SYSTEM 40 Tests
71675	COPRO SYSTEM Plus 20 Tests
71678	PATHOGENIC SYSTEM DOUBLE 40 Tests
71679	PATHOGENIC SYSTEM 20 Tests
71681	PATHOGENIC SYSTEM AST
71714	INTEGRAL SYSTEM ENTEROBATTERI 20 Tests
71718	INTEGRAL SYSTEM STAFILOCOCCI 20 Tests
71720	INTEGRAL SYSTEM STREPTOCOCCI 20 Tests
71724	INTEGRAL SYSTEM GARDNERELLA 20 TESTS
71822	INTEGRAL SYSTEM YEASTS Plus 20 Tests
72560	STREPTO SYSTEM 12 R 40 Tests
72592	MYCOPLASMA SYSTEM Plus 20 Tests
74156	A.F. GENITAL SYSTEM 20 Tests
74160	URIN SYSTEM Plus 20 Tests
74161	URIN SYSTEM Chrom 20 Tests
76010	Sensi Test gram-negative 20 Tests
76020	Sensi Test gram-positive 20 Tests
76031	SensiQuattro Gram-negative 20 Tests
76032	SensiQuattro Gram-positive 20 Tests
76033	SensiQuattro Candida EU 20 Tests
78618	ENTERO PLURI TEST 10 Tests
78619	ENTERO PLURI TEST 25 Tests
78620	OXI/FERM PLURI TEST 10 Tests
78621	OXI/FERM PLURI TEST 25 Tests
79010	Sensi Test gram-negative 4 Tests
79020	Sensi Test gram-positive 4 Tests
79031	SensiQuattro Gram-negative 4 Tests
79032	SensiQuattro Gram-positive 4 Tests
79033	SensiQuattro Candida EU 4 Tests
79156	A.F. GENITAL SYSTEM 4 Tests
79160	URIN SYSTEM Plus 4 Tests
79161	URIN SYSTEM Chrom 4 Tests
79560	STREPTO SYSTEM 12 R 8 Tests
79592	MYCOPLASMA SYSTEM Plus 4 Tests
79618	ENTEROSYSTEM 18R 4 Tests
79630	STAF SYSTEM 18 R 4 Tests
79670	COPRO SYSTEM 8 Tests
79675	COPRO SYSTEM Plus 4 Tests
79678	PATHOGENIC SYSTEM DOUBLE 8 Tests
79679	PATHOGENIC SYSTEM 4 Tests
79681	PATHOGENIC SYSTEM AST
79714	INTEGRAL SYSTEM ENTEROBATTERI 4 Tests
79718	INTEGRAL SYSTEM STAFILOCOCCI 4 Tests
79720	INTEGRAL SYSTEM STREPTOCOCCI 4 Tests
79724	INTEGRAL SYSTEM GARDNERELLA 4 Tests

79822	INTEGRAL SYSTEM YEASTS Plus 4 Tests
80009	IODINE MKTT SOLUTION 10 x 10 ml
80010	XLT 4 supplement 2 x 50 ml
80021	GLYCEROL supplement 4 x 50 ml
80022	POTASSIUM TELLURITE 1% suppl. 5 x 10 ml
80031	TWEEN 80 supplement 2 x 50 ml
80040	CHROMATIC™ SALMONELLA Supplement 2x50 ml
80047	MULLER KAUFFMANN 3X50 ml (Iodio/B.G.O.1%)
80053	VITAMIN K 1% supplement 5 x 5 ml
80056	LEGIONELLA growth supplement 10 vials
80057	H2O2 REAGENT 1 x 10 ml
80060	DECONTAM-KIT
80110	UREA 40% 6X100 ml
80219	EGG YOLK emulsion 4 x 50 ml
80252	ENTEROSYSTEM 18R REAGENT 100/200 Tests
80253	COPRO SYSTEM REAGENTS (antisera)
80257	LISTERIA SYSTEM 18R -REAG 100/200 Tests
80258	AF GENITAL SYSTEM REAGENT
80260	IDENTIF. SYSTEM-REAGENT 100/200 Tests
80271	KOVAC'S REAGENT 4x25 ml
80272	FERRIC CHLORIDE 10% 2x 25 ml
80273	NINHYDRIN 7% 10 ml
80275	MIF COLOR KIT 50 Tests
80276	ZIEHL-NEELEN 3 x 250 ml
80277	METHYLENE BLUE Solution 250 ml
80279	VASELINE OIL 4 x 50 ml
80280	V.P. TEST-Reagent 10x10ml
80281	V.P. TEST EP 10 x 10 mL
80282	Kit May-Grünwald Giemsa
80290	SAFRANIN SOLUTION 1000 ml
80291	POTASSIUM TELLURITE 3.5% suppl.5x10 ml
80292	UREA 40 % supplement 10 x 5 ml
80293	GRAM COLOR KIT 4 x 250 ml
80294	KIT COLOR ALBERT 2 x 250 ml
80295	DECOLOURIZING SOLUTION 1000 ml
80296	LUGOL PVP SOLUTION 1000 ML
80297	SAFRANIN SOLUTION 500 ml
80298	LUGOL PVP SOLUTION 250 ml
80299	CRYSTAL VIOLET SOLUTION 1000 ml
80300	TTC 1% supplement 5 x 10 ml
80350	ANTIBIOTIC TEST
80351	RAPID ANTIBIOTIC TEST 50 Tests
80380	KINYOUN COLOR KIT 2 x 250 ml
80390	FIXUR 1
80409	IODINE SOLUTION 10 x 10 ml
80410	XLT 4 SUPPLEMENT 4 x 50 ml
80422	POTASSIUM TELLURITE 1% Supplement 10 x 10 ml
80430	TTC 1% supplement 10 x 10 ml
80431	TWEEN 80 Supplement 4 x 50 ml
80453	VITAMIN K 1% SUPPLEMENT 10 x 5 ml
80491	POTASSIUM TELLURITE 3.5% Supplement 10 x 10
81001	AMPICILLIN supplement 10 vials
81002	LEGIONELLA (BMPA) supplement 10 vials
81003	BRUCELLA supplement 10 vials
81004	CAMPYLOBACTER Preston supplem 10 vials
81006	CN (Pseudomonas) supplement 10 vials
81007	CLOSTRIDIUM difficile supleme 10 vials
81008	LEGIONELLA (GVPC) supplement 10 vials
81009	IODINE solution 5 x 10 ml
81011	CLOSTRIDIUM perfringens (T.S.C.) sup.10 v.

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91200	DISC DISPENSER 8 CARTRIDGES
91203	DISC DISPENSER 6 CARTRIDGES
92000	AMOX*/SULB 2/1 AXS 0.016-256* 30 MIC Tests
920000	AMOX*/SULB 2/1 AXS 0.016-256* 100 MIC Tests
92001	RIFAMPICIN RD 0.002-32 30 MIC Tests
920010	RIFAMPICIN RD 0.002-32 100 MIC Tests
920011	RIFAMPICIN RD 0.002-32 10 MIC Tests
92002	FUSIDIC ACID FU 0.016-256 30 MIC Tests
920020	FUSIDIC ACID FU 0.016-256 100 MIC Tests
920021	FUSIDIC ACID FU 0.016-256 10 MIC Tests
92003	AMPICILLIN AMP 0.016-256 30 MIC Tests
920030	AMPICILLIN AMP 0.016-256 100 MIC Tests
920031	AMPICILLIN AMP 0.016-256 10 MIC Tests
92004	POLYMYXIN B PB 0.064-1024 30 MIC Tests
920040	POLYMYXIN B PB 0.064-1024 100 MIC Tests
920041	POLYMYXIN B PB 0.064-1024 10 MIC Tests
92005	CEFPODOXIME PX 0.016-256 30 MIC Tests
920050	CEFPODOXIME PX 0.016-256 100 MIC Tests
920051	CEFPODOXIME PX 0.016-256 10 MIC Tests
92006	CEFOTAXIME CTX 0.016-256 30 MIC Tests
920060	CEFOTAXIME CTX 0.016-256 100 MIC Tests
920061	CEFOTAXIME CTX 0.016-256 10 MIC Tests
92007	CEFOTAXIME CTX 0.002-32 30 MIC Tests
920070	CEFOTAXIME CTX 0.002-32 100 MIC Tests
920071	CEFOTAXIME CTX 0.002-32 10 MIC Tests
92008	CEFPIROME CR 0.016-256 30 MIC Tests
920080	CEFPIROME CR 0.016-256 100 MIC Tests
920081	CEFPIROME CR 0.016-256 10 MIC Tests
92009	GENTAMICIN CN 0.016-256 30 MIC Tests
920090	GENTAMICIN CN 0.016-256 100 MIC Tests
920091	GENTAMICIN CN 0.016-256 10 MIC Tests
92010	GENTAMICIN CN 0.064-1024 30 MIC Tests
920100	GENTAMICIN CN 0.064-1024 100 MIC Tests
920101	GENTAMICIN CN 0.064-1024 10 MIC Tests
92011	GATIFLOXACIN GAT 0.002-32 30 MIC Tests
920110	GATIFLOXACIN GAT 0.002-32 100 MIC Tests
920111	GATIFLOXACIN GAT 0.002-32 10 MIC Tests
92012	TEICOPLANIN TEC 0.016-256 30 MIC Tests
920120	TEICOPLANIN TEC 0.016-256 100 MIC Tests
920121	TEICOPLANIN TEC 0.016-256 10 MIC Tests
92013	ENROFLOXACIN ENR 0.002-32 30 MIC Tests
920130	ENROFLOXACIN ENR 0.002-32 100 MIC Tests
920131	ENROFLOXACIN ENR 0.002-32 10 MIC Tests
92014	SPECTINOMYCIN SPC 0.064-1024 30 MIC Tests
920140	SPECTINOMYCIN SPC 0.064-1024 100 MIC Tests
920141	SPECTINOMYCIN SPC 0.064-1024 10 MIC Tests
92015	OXACILLIN OX 0.016-256 30 MIC Tests
920150	OXACILLIN OX 0.016-256 100 MIC Tests
920151	OXACILLIN OX 0.016-256 10 MIC Tests
92016	CEFTIZOXIME CZX 0.016-256 30 MIC Tests
920160	CEFTIZOXIME CZX 0.016-256 100 MIC Tests
920161	CEFTIZOXIME CZX 0.016-256 10 MIC Tests
92017	MECILLINAM MEC 0.002-32 30 MIC Tests
920170	MECILLINAM MEC 0.002-32 100 MIC Tests
920171	MECILLINAM MEC 0.002-32 10 MIC Tests
92018	AMIKACIN AK 0.016-256 30 MIC Tests
920180	AMIKACIN AK 0.016-256 100 MIC Tests
920181	AMIKACIN AK 0.016-256 10 MIC Tests
92019	BACITRACIN BA 0.016-256 30 MIC Tests
920190	BACITRACIN BA 0.016-256 100 MIC Tests

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96021	SALMONELLA PARATYPHI C TOTALE 20 ml
96022	BRUCELLA MELITENSIS 20 ml
96023	BRUCELLA SUIS 20 ml
96031	SALMONELLA TYPHI H SLIDE 5 ml
96032	SALMONELLA TYPHI O SLIDE 5 ml
96033	SALMONELLA TYPHI TOTALE 5 ml SLIDE
96034	SALMONELLA PARATYPHI AH SLIDE 5 ml
96035	SALMONELLA PARATYPHI AO 5 ml SLIDE
96036	SALMONELLA PARATYPHI A TOTALE 5ml SLIDE
96037	SALMONELLA PARATYPHI BH 5 ml SLIDE
96038	SALMONELLA PARATYPHI BO 5 ml SLIDE
96039	SALMONELLA PARATYPHI B TOTALE 5ml SLIDE
96040	SALMONELLA PARATYPHI CH 5 ml SLIDE
96041	SALMONELLA PARATYPHI CO 5 ml SLIDE
96042	SALMONELLA PARATYPHI C TOTALE 5 ml SLIDE
96043	BRUCELLA TOTALE SLIDE 5 ml SLIDE
96044	BRUCELLA ABORTUS 5 ml SLIDE
96045	BRUCELLA MELITENSIS SLIDE 5 ml
96046	BRUCELLA BENGAL ROSE SLIDE 5 ml
96047	PROTEUS OX2 5 ml SLIDE
96048	PROTEUS OX19 5 ml SLIDE
96049	PROTEUS OXK 5 ml SLIDE
96093	CONTROLLO NEGATIVO/NEGATIVE CONTROL 0.5ml
96096	POSITIVE CONTROL FOR SALMONELLA 0.5ml
96097	POSITIVE CONTROL FOR PROTEUS 0.5ml
96098	POSITIVE CONTROL FOR BRUCELLA 0.5ml
96142	Legionella Latex Kit
96143	CAMPYLOBACTER LATEX KIT
96144	CLOSTRIDIUM DIFFICILE LATEX KIT
96148	SHIGELLA ANTISERUM
96150	E. COLI O157 LATEX KIT
96151	SALMONELLA LATEX KIT
96153	STREPTO B LATEX KIT
96154	STREPTO A LATEX KIT
96155	BENCE JONES LATEX TEST
96316	Clostridium difficile GDH Card
96317	Clostridium Difficile Toxin A+B Card
96318	Giardia Card
96319	Listeria Monocytogenes Card
96320	Salmonella Ag Card
96321	O157 E.coli Card
96401	ONE STEP AMP DRMG SCREEN 20 CARDS
96404	ONE STEP COC DRMG SCREEN
96405	ONE STEP THC DRMG SCREEN
96406	ONE STEP M-AMP DRMG SCREEN 20 CARDS ONE STEP BRUPRENORPHINE DRMG SCREEN 20 CARDS
96411	FECAL OCCULT BLOOD CARD
96415/20	FECAL OCCULT BLOOD CARD
96418	STREPTO A CARD 30 CARDS
96441	Gonorrhea Ag Card
96442	Gardnerella Vaginalis Card
96443	Trichomonas Vaginalis Card
96444	B.J. Free Kappa/Lambda Dipstick
96455	H.PYLORI CARD 20 CARD
96460	HCG URINE/SERUM CARD 50 CARD
96461	HCG URINE/SERUM CARD 100 CARD
96462	MICROALBUMIN CARD URINE 20 Cards
96465	AFP -ALFA FETO CARD 20 CARDS
96468	TUBERCOLOSI CARD 20 CARDS
96480	IgE TOTAL CARD
96485	CEA CARD 20 Cards

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Rev. 31.0 del 08.01.2016

96487	MYOGLOBIN
96488	TROPONIN 20 CARDS
96490	FERRITIN CARD
96495	SIFILIDE CARD 20 CARDS
96498	IM MONONUCLEOSIS INFECTION 20 CARDS
96590	URINE STRIP
96900	GIOTTO READER
96909	BIOMIC V3
96914	BIOMIC V3 AST
96915	BIOMIC V3 ID
96916	BIOMIC V3 CC
96919	AST Software
96931	ID Software
96932	CC Software
96933	Micropiastre 96 pozzetti Software

97800	ROTASTICK ONE STEP KIT 20 Tests
97801	RSV STICK ONE STEP 20 Tests
97802	ROTA/ADENO COMBI STICK ONE STEP 20 Tests
97803	H.PYLORI FECAL Ag ONE STEP 20 Tests
97805	STREP B STICK ONE STEP ASSAY 20 Tests
97807	ADENOSTICK ONE STEP ASSAY 20 Tests
9999	Blank Discs
99003	KPC&MBL disc kit (acc. to EUCAST)
99004	ESBL disc kit (acc. to EUCAST)
99005	ESBL disc kit (acc. to CLSI)
99006	ESBL (Chromos. Ind. AmpC) disc kit (acc. to EUCAST)
99007	KPC&MBL&OXA-48 disc kit (acc. to EUCAST)
99008	ESBL+AmpC screen disc kit
99009	AmpC disc kit

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