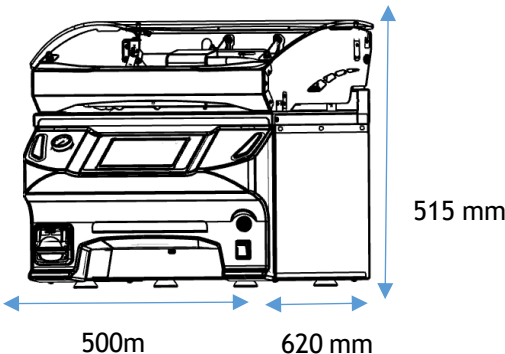


JUNE 2022

Installation Requirements for a MEDIAWEL 10 Media preparator


Thank you for purchasing the MEDIAWEL 30.

In order to optimize and perform an efficient, trouble free installation, please read carefully, fill in the here below form and return to our service team.

Access Requirements / Space Requirements		
Points to be checked:	Response:	Comments:
<p>Is the place of usage accessible through the doors?</p>  <p>500mm 620 mm 515 mm</p>		
<p>Is there enough space on the site and will it be freed at the day of the installation?</p>		
<p>In the case of an upstairs installation:</p> <ul style="list-style-type: none"> - Is there an available lift? - Should there be a handling operation to plan? 		

Mains Requirements		
Points to be checked:	Response:	Comments:
<p>Is there an electricity supply with 230 +/- 10% V mono-phase & neutral + ground, 16 amp, available at less than 2 meters from the place of usage?</p>		

Water System Requirements		
Points to be checked:	Response:	Comments:
Is the inlet water between 2 and 4 bars?		
Is the inlet water temperature between 5 and 20°C		
Is the water hardness less or more than 16°TH?		
Is the water tap located, less than 2 meters from the place of usage?		
Is the tap equipped with a barb hose to receive a flexible tubing (12.7 mm inner diameter)?		
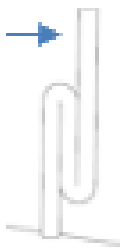
<i>Water evacuation</i>		
<i>Points to be checked:</i>	<i>Response:</i>	<i>Comments:</i>
<p><i>Is there a water evacuation point, located at less than 2 meters from the place of usage?</i> <i>4 meters are delivered with the unit</i> MEDW4012 = WATER SUPPLY HOSE 4 meters MEDW2027 = WATER SUPPLY HOSE 20 meters</p>		
<p><i>Is this evacuation, prepared to resist to temperatures up to 120°C (copper or CPVC)?</i></p>		
<p><i>Is this evacuation equipped with a U-bend / siphon?</i> <i>If yes, remove it.</i></p> 		

<i>Culture Media Dispensing</i>		
<i>Points to be checked:</i>	<i>Response:</i>	<i>Comments:</i>
<p><i>What is the concentration of salt per liters (of your saliest medium)</i> >10 Gr/L = MEDW2039 ANTI CORROSIVE SYSTEM</p>		
<p><i>Will the dispensing of the culture media take place in the same room of the media preparation?</i> MEDW2002 QUICK CONNECTOR MEDW2101 TROLLEY</p>		
<p><i>Do you want to store it under the bench</i> MEDW2002 QUICK CONNECTOR MEDW2101 TROLLEY</p>		



Media list	
Media name	Volume (maximum per batch)