(IT)

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(EN)

DESCRIPTION OF THE DEVICE:

Sterile bag for the cryopreservation of stem cells in liquid hydrogen and/or nitrogen vapours. Class II* medical device pursuant to Medical Devices Directive 93/42/EEC and subsequent amendments thereto.
CONTENT:
Each device is wrapped in a double paper blister and mated medical enclosure containing:
A primary EVA bag equipped with filling SET and sampling ports.
A secondary bag which serves as a protective wrapping for the freezing process.

VOLUMES AVAILABLE:

REF.	NOMINAL VOLUME	FILLING VOLUME*	REF.	NOMINAL VOLUME	FILLING VOLUME*
SH-6B	6 mL	3-6 mL	SH-250B	250 mL	30-70 mL
SH-25B	25 mL	20-25 mL	SH-500B	500 mL	55-100 mL
SH-50B	50 mL	10-30 mL	SH-750B	750 mL	80-180 mL

* The filling volumes specified must be adhered to, otherwise the user shall be responsible for validating different utilisation conditions.

UTILISATION:

Comply rigorously with the instructions given in the utilisation technique manual received with the first delivery.
To be used exclusively by specialist personnel.
Admissible utilisation temperatures: from -196°C to +37°C.

SPECIAL WARNINGS AND UTILISATION PRECAUTIONS:

Use only if the packaging is intact and the device is undamaged.
Use immediately after opening the package and removing the protective enclosure.
Use before the expiry data specified.
The product must be disposed of according to the applicable national regulations via the appropriate sanitary bins.

STORAGE AND TRACEABILITY:

Store in a dry and clean environment, free of biological contaminants ad/or chemical vapours; do not expose to direct light, Validity: 5 years from date of sterilisation.
Batch number and expiry date are printed on the back of the blister. The batch number also appears on the primary bag.

PACKAGING:

Cardboard boxes containing an envelope made of LDPE with 10 pieces (see "CONTENT").

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Paolo Gobbi Frattini S.r.l.